


Oxygenates in Gasoline 2310-0098

Technical Overview

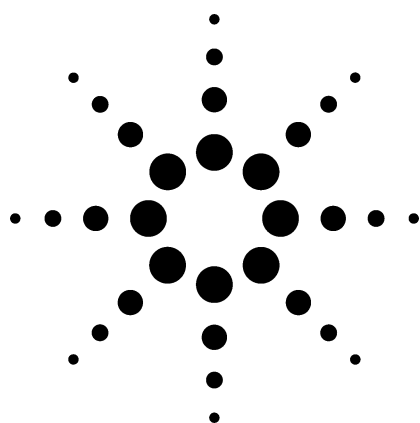

Application Highlights


- Sample is injected onto a TCEP capillary column. A 'heart' cutting technique is used to switch the oxygenates of interest onto a DB-1 capillary column where the oxygenate will separate from any co-eluting hydrocarbons. This analyzer follows the ISO E13132 method.
- Analysis time is approximately 30 minutes.

Configurations


- GC must have capillary inlet (112), two flame ionization detectors (FIDs) (211) and pneumatics control module (PCM) module (309). An autosampler and ChemStation are recommended,
- Special includes columns, Deans switching hardware, checkout samples for valve timing, factory checkout with chromatograms, method supplied on disk for ChemStation and user manual.

For More Information

For more information on our products and services, visit our Web site at www.agilent.com/chem.

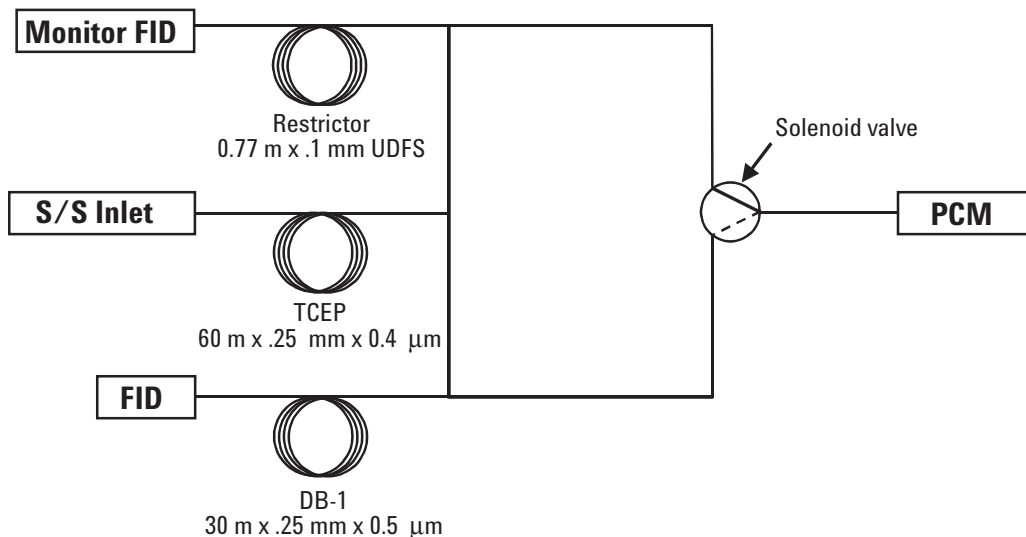

Agilent Technologies

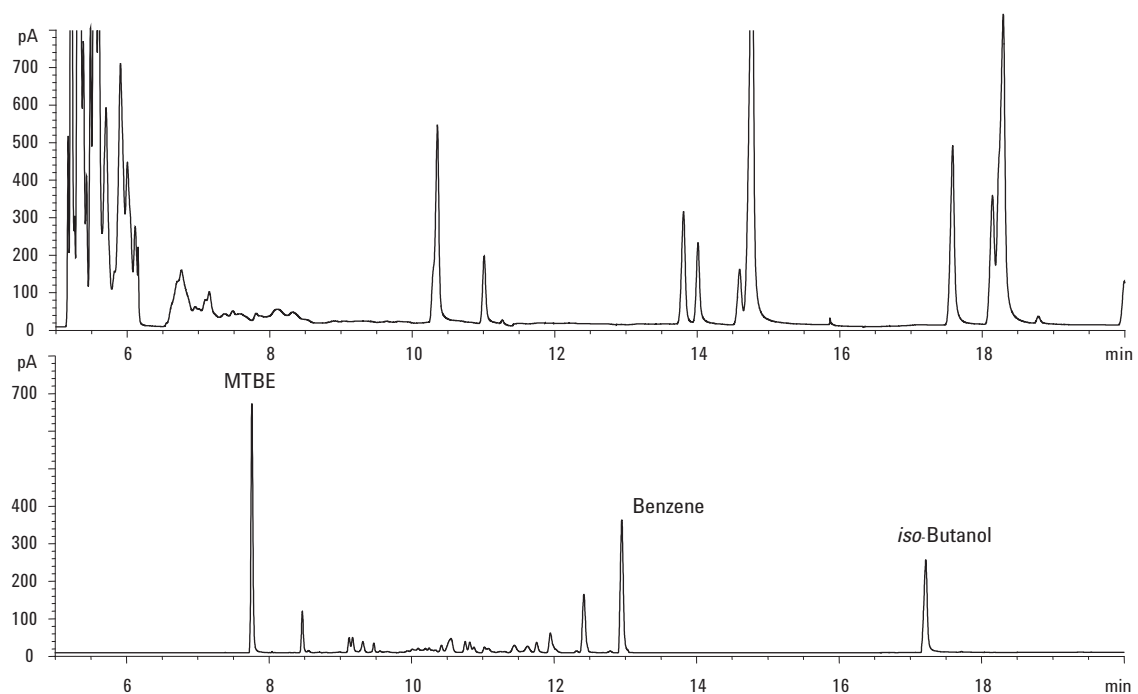

Chromatogram shows cuts for MTBE, benzene and *iso*-Butanol from gasoline.

Agilent shall not be liable for errors contained herein or for incidental or consequential damages in connection with the furnishing, performance, or use of this material.

Information, descriptions, and specifications in this publication are subject to change without notice.

© Agilent Technologies, Inc. 2003

Printed in the USA
May 9, 2003
5988-9302EN



Agilent Technologies