AAVPrime™ AAV-qPCR 滴度检测试剂盒

Cat. Nos. AA301 / AA302

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Component	Volume	Concentration	Shipping	Storage
DNase I (RNase-free)	1x20 µl 2x20 µl	2,000 units/ml	Ice pack	-20°C, Stable for at least 12 months
DNase I buffer (10x)	1x20 µl 2x20 µl		Ice pack	-20°C, Stable for at least 12 months
AAV Lysis buffer	1x200 µl 2x200 µl		Ice pack	-20°C, Stable for at least 12 months
AAV-TaqMan qPCR mix (2x)	1x500 µl 2x500 µl		Ice pack	-20°C, Stable for at least 12 months
AAV-qPCR primer mix	1x100 µl 2x100 µl	2.5 µM	Ice pack	-20°C, Stable for at least 12 months
AAV-Probe	1x100 µl 2x100 µl	2.5 µM	Ice pack	-20°C, Stable for at least 12 months
AAV-qPCR standard (DNA)	1x25 µl 2x25 µl	1x10 ⁹ copies/µl	Ice pack	-20°C, Stable for at least 12 months
ddH ₂ O	1x1 ml 2x1 ml		Ice pack	Room temperature, Stable for at least 12 months

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Cycle	Steps	Temperature	Duration	Read
1	Denaturation	95 °C	10 min	off
35	Denaturation	95°C	10 sec	off
	Annealing	60 °C	20 sec	off
	Extension	72 °C	25 sec	on

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$$\text{稀释系数} = \frac{\text{DNase reaction volume}(\mu\text{l})}{\text{Original sample volume}(\mu\text{l})} \times \frac{\text{Lysis reaction volume}(\mu\text{l})}{\text{Volume of sample use in lysis reaction}(\mu\text{l})} \times \frac{1000\mu\text{l/ml}}{\text{Volume added to PCR well}(\mu\text{l})}$$

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- DNase reaction volume: 10 μL (根据本实验流程)
- Original sample volume: The volume of AAV particles used for DNase reaction. 8 μL (根据本实验流程)
- Lysis reaction volume: 20 μL (根据本实验流程)
- Volume of sample use in lysis reaction: 10 μL (根据本实验流程)
- Volume added to PCR well: 5 μL (根据本实验流程)

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Limited Use License and Warranty

Limited use license

The following terms and conditions apply to use of the AAV-qPCR Titration Kits (the Product). If the terms and conditions are not acceptable, the Product in its entirety must be returned to GeneCopoeia within 5 calendar days. A limited End-User license is granted to the purchaser of the Product. The Product shall be used by the purchaser for internal research purposes only. The Product is expressly not designed, intended, or warranted for use in humans or for therapeutic or diagnostic use. The Product must not be resold, repackaged or modified for resale, or used to manufacture commercial products or deliver information obtained in service without prior written consent from GeneCopoeia. Use of any part of the Product constitutes acceptance of the above terms.

Limited warranty

GeneCopoeia warrants that the Product meets the specifications described in the accompanying Product Datasheet. If it is proven to the satisfaction of GeneCopoeia that the Product fails to meet these specifications, GeneCopoeia will replace the Product. In the event a replacement cannot be provided, GeneCopoeia will provide the purchaser with a refund. This limited warranty shall not extend to anyone other than the original purchaser of the Product. Notice of nonconforming products must be made to GeneCopoeia within 30 days of receipt of the Product. GeneCopoeia's liability is expressly limited to replacement of Product or a refund limited to the actual purchase price. GeneCopoeia's liability does not extend to any damages arising from use or improper use of the Product, or losses associated with the use of additional materials or reagents. This limited warranty is the sole and exclusive warranty. GeneCopoeia does not provide any other warranties of any kind, expressed or implied, including the merchantability or fitness of the Product for a particular purpose.

GeneCopoeia is committed to providing our customers with high-quality products. If you should have any questions or concerns about any GeneCopoeia products, please contact us at 020-32068595.