



# Installation Guide for JEOL MassCenter version 1.3.4 m 1/2

1. First, make sure that East Asian language support is installed.
2. Choose the right regional and language settings. Go to Start → Control Panel → Regional and Language Options → Select English (US).

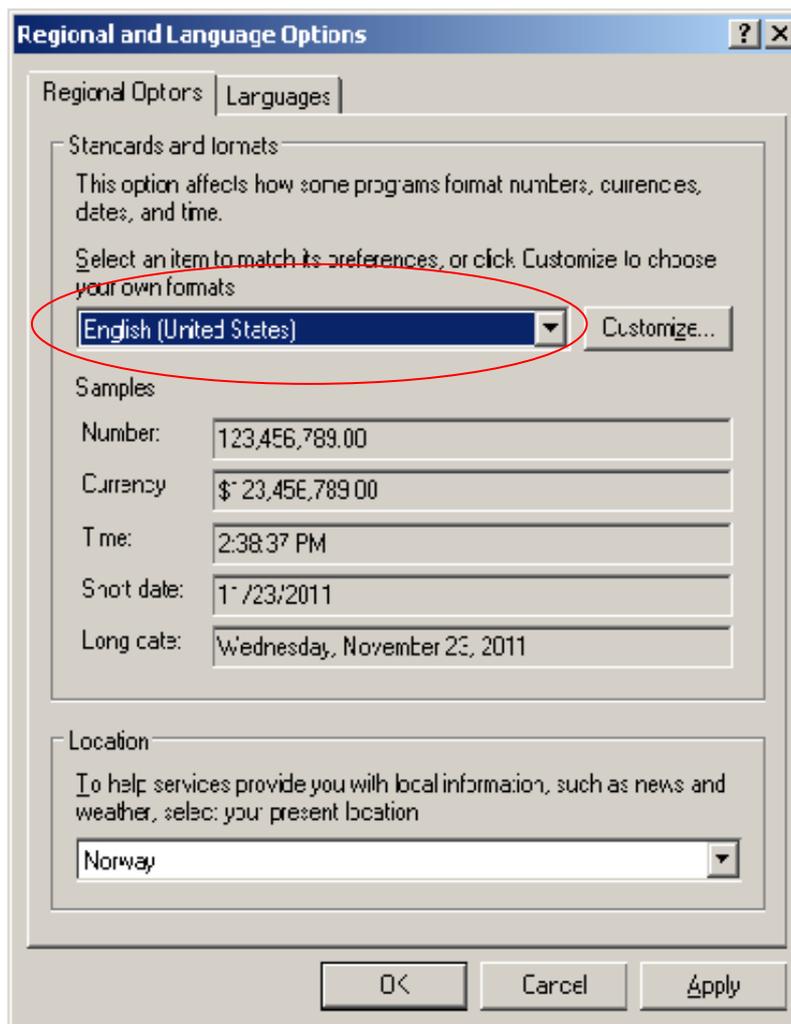

3. Then, log on to the install account on your personal computer.
4. Insert CD and start the installation of MassCenter (installation will start automatically).
5. In the step "Select MS Device Name", select *Connected device nothing*.
6. Fill in the necessary Customer Information and the CD key code;  
**MS56-1100-5013-4115**. Check mark  emulation code and click next.
7. Finish the installation and make sure that the MassCenter software icon appears on our desktop.

8. Go to C:\Program Files and mark the jeol-folder. Right mouse click and select properties. Under security give Users (KJxxxxxx\Users) Full Control.

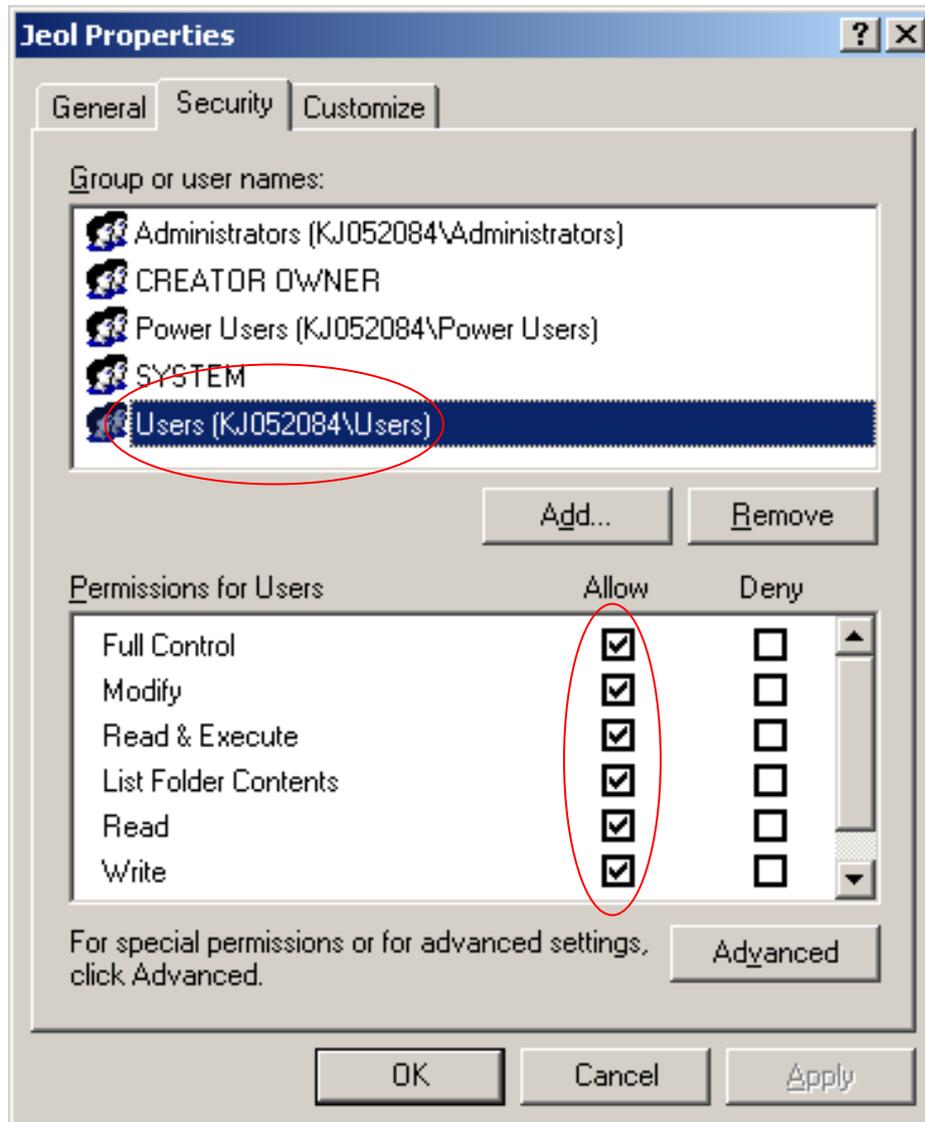

9. Go to: C:\Program Files\Jeol\MassCenter\Applications and mark the icon *prjSpectrumViewer.exe*. Right mouse click and send it as a shortcut to your desktop.
10. Do the same procedure with the icon *prjChromatogramViewer.exe*.
11. When using MassCenter you then run the program through your install account. (right mouse click the chromatogramviewer icon and run through install).



As you construct mass spectra from the total ion chromatogram spectrum viewer will start automatically)

12. To open raw data files in chromatogram viewer you need to save these files on your C:\ or D:\ disk in the correct hierarchy containing .7pj\ .7fl\ .7rw

(Which could be as the following example);

D:\MS.7pj\DART2011.7fl\Bjarte030112.7rw