The word "TECHNE" is written in a bold, sans-serif font and is enclosed within a horizontal oval border.

TECHNE

TC-512

OPERATOR'S MANUAL

CONTENTS

	PAGE
INTRODUCTION	2
SAFETY AND INSTALLATION INFORMATION	3
Operator Safety	3
Warning	3
Unpacking	3
Installation	3
Replacement Cable	4
Working Conditions	4
Guarantee	5
SÉCURITÉ ET CONSIGNES D'INSTALLATION	6
Sécurité de l'opérateur	6
Avertissement	6
Déballage	6
Installation	6
Câble de rechange	7
Conditions de travail	7
Garantie	8
SICHERHEITS - UND INSTALLATIONSINFORMATIONEN	9
Sicherheit des Bedienpersonals	9
Achtung	9
Auspacken	9
Inbetriebnahme	9
Ersatzkabel	10
Einsatzbedingungen	10
Garantie	11
INFORMACIÓN DE SEGURIDAD E INSTALACIÓN	12
Seguridad del operario	12
Advertencia	12
Desembalaje	12
Instalación	12
Cable de repuesto	13
Condiciones de trabajo	13
Garantía	14
INFORMAZIONI SULLA SICUREZZA E L'INSTALLAZIONE	15
Sicurezza dell'operatore	15
Avvertenza	15
Disimballaggio	15
Installazione	15
Cavo di ricambio	16
Condizioni di esercizio	16
Garanzia	17

INSTRUMENT CONNECTIONS	18
CONTACT INFORMATION	19
TC-512 SPECIFICATION	20
TC-512 OPERATION AND MAINTENANCE	22
Gradient	22
Adjustable Heated Lid	23
Memory Card	24
Adjustable Screen	25
Consumables	25
PC Connections	25
Cleaning the TC-512	26
Removing the Block	26
PROGRAMMING THE TC-512	28
Start up Menu	28
The System User Menu	29
Writing and Editing Programs	30
Option 1: Copying an Existing Template	31
Option 2: Editing and Saving an Existing Template	33
Option 3: Creating a New Program	34
Inserting a Segment	35
Inserting Cycle Numbers	36
Inserting Incremental or Decremental Time or Temperature	36
Inserting a Pause	38
Saving Changes to a Program	38
RUNNING A PROGRAM	38
Program Defaults	41
Data Logs	43
Printing a Data Log	43
Gradient Calculator	45
SYSTEM ADMINISTRATION	46
Program Passwords	47
System Administration Password	47
System Test	48
Controller Reset	48
Card Utilities	48
System Configuration	49
FREQUENTLY ASKED QUESTIONS	51
ADDITIONAL INFORMATION	52
Fault Finding	52
Fuses	52
Accessories	52
PC Accessories	53
Replacement Parts	53


INTRODUCTION

The TC-512 is a gradient thermal cycler, designed to accurately control the temperatures of samples between 4°C and 99°C. It has many scientific applications including DNA amplification, PCR¹, reverse transcription and sequencing.

The TC-512 is programmed via a touch screen LCD graphics display. The user is able to input various programs and parameters, including:

- Specified temperatures in °C
- Times for which the sample is to be held at the specified temperature
- Heating and cooling rates between the specified temperatures in °C/sec
- Gradient of temperatures across the thermal cycling block
- Incremental time or temperature increases or decreases when a cycle is repeated

Before using the TC-512 make sure you have read this manual carefully. If there is any doubt relating to the proper use of this equipment, the staff at Bibby Scientific Ltd or your supplier will be happy to assist you.

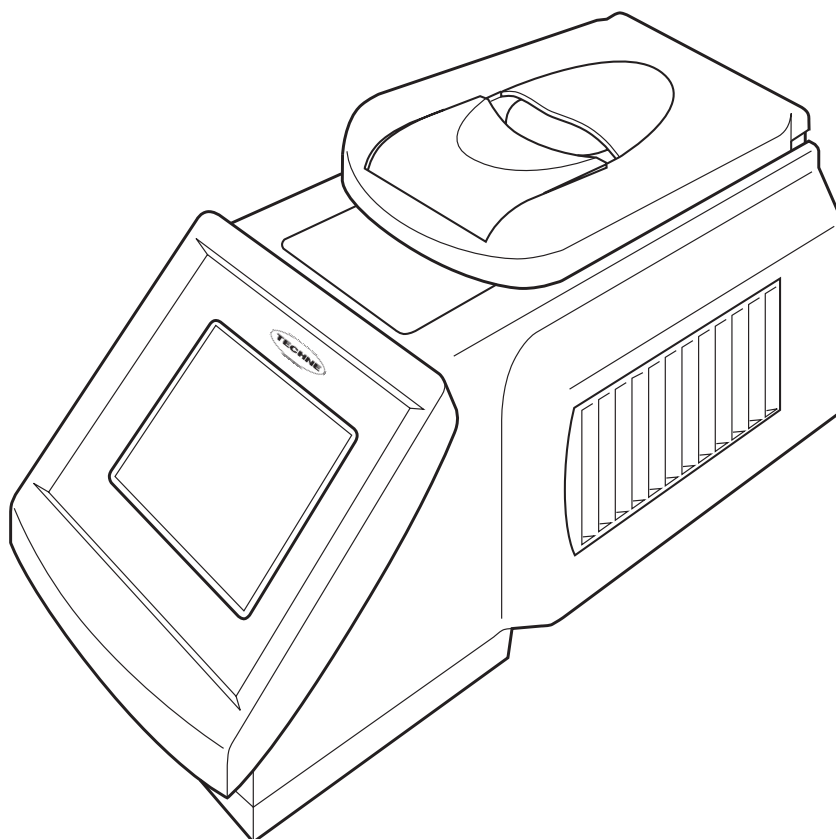¹Aspects of the PCR process are claimed in USA Patent Nos. 4,683,195; 4,683,202, and 4,965,188. Use of the TC-512 in such processes does not convey a licence to practice the processes themselves.

SAFETY AND INSTALLATION

OPERATOR SAFETY

Please read this manual carefully before using the thermal cycler. If the equipment is not used in the manner described in this manual the protection provided by the equipment might be impaired.

All Bibby Scientific Ltd instruments are designed to conform to international safety requirements and are fitted with an over-temperature cut-out. If a safety problem should be encountered then switch off the unit at the mains socket and remove the plug from the electricity supply.

WARNING

HIGH TEMPERATURES ARE DANGEROUS as they can cause serious burns to operators and ignite combustible material. Users should be aware of the following potential hazards:

- USE CARE AND WEAR PROTECTIVE GLOVES TO PROTECT HANDS
- DO NOT use combustible substances near hot objects
- DO NOT operate the instrument in the vicinity of inflammable liquids or gases
- DO NOT place any liquid directly into the instrument, always use the appropriate plastic consumables

After heating samples, remember that parts of the instrument, namely the lid and block and any associated accessories may be very hot.

UNPACKING

When unpacking please ensure that the following have been removed from the packaging:

- Thermal cycler
- Mains cable
- Guarantee card
- Decontamination certificate

The user is advised to keep the original packaging in case the instrument ever needs to be returned for service or repair. Bibby Scientific Ltd accepts no responsibility for damage incurred unless the unit is correctly packed and transported in its original packaging.

Please note that the screen of the thermal cycler is easily damaged by sharp objects such as pens, pencils and fingernails. This type of damage will be considered as misuse and invalidate the guarantee for this component.

INSTALLATION

The instrument should be carried using both hands. Never move or carry the instrument when in use or connected to the mains electricity supply.

1. All Bibby Scientific Ltd instruments are supplied with a power cable; this may be integral or plug-in.
2. Before connecting the instrument to the mains electricity supply, check the voltage against the

rating plate (located on the back of the unit). Ensure that the voltage selector switch (located above the On/Off switch) is set to the appropriate voltage for the local supply. Please note that the unit must be earthed to ensure proper electrical safety. Connect the mains cable to a suitable plug according to the table below.

Connections	220V-240V, 50/60Hz Supply	110V-120V, 50/60Hz Supply
Live	brown	black
Neutral	blue	white
Earth	green/yellow	green

3. UK ONLY: The fused plug supplied with the mains cable is fitted with a 10 Amp fuse to protect the instrument and the user.
4. Units showing 230V, 50/60Hz on the rating plate also operate between 210 and 260V; units with 120V, 50/60Hz also operate between 100 and 130V and units marked with 100V will operate between 90 and 110V. In all cases the heating rate will degrade by approximately 8% at the extremes of the voltage range.
5. Plug the mains cable into the socket on the back of the instrument.
6. Place the unit on a suitable flat bench or in a fume cupboard if required, ensuring that the air inlet vents on the underside are free from obstruction. Position the instrument with a minimum distance all round of 200mm from walls or other items and between each unit if multiple units are being used to permit efficient air flow for each instrument.
7. Switch on the instrument:
 - I mains switch On
 - O mains switch Off

REPLACEMENT CABLE

Should the mains lead need replacement, a cable of 1mm² of harmonized code H05VV-F connected to an IEC320 plug should be used. **IF IN DOUBT CONSULT A QUALIFIED ELECTRICIAN.**

WORKING CONDITIONS

The thermal cycler is designed to operate under the following conditions:

- Indoor use
- Ambient temperature range +5°C to +40°C
- Altitude to 2000m
- Mains supply fluctuations not exceeding 10%
- Over voltage category II IEC 60364-4-443
- Pollution degree 2

Note: The control specifications are quoted at an ambient temperature of 20°C. The specification may deteriorate outside an ambient temperature of 10°C to 30°C.

The instrument has been tested for radio frequency interference and is certified under EN61326.

GUARANTEE

The instrument is guaranteed against any defect in material or workmanship for the period specified on the enclosed guarantee card. This period is effective from the date of purchase; within this period all defective parts will be replaced free of charge provided that the defect is not the result of an accident, misuse or negligence. The screen of the thermal cyclers can be damaged by sharp objects such as pens, pencils and fingernails. Damage of this sort will be considered misuse and will invalidate the guarantee for this component.

Servicing under this guarantee should be obtained from the supplier of the instrument.

This manual has been prepared for the convenience of Bibby Scientific Ltd's customers and nothing in this manual shall be taken as a warranty, condition or representation concerning the description, merchantability, fitness for purpose or otherwise of the unit or components.

Notwithstanding the description and specification(s) of the instruments contained in the operator's manual, Bibby Scientific Ltd reserves the right to make such changes as it sees fit to the instruments or to any of the components.

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
INSTRUMENT CONNECTIONS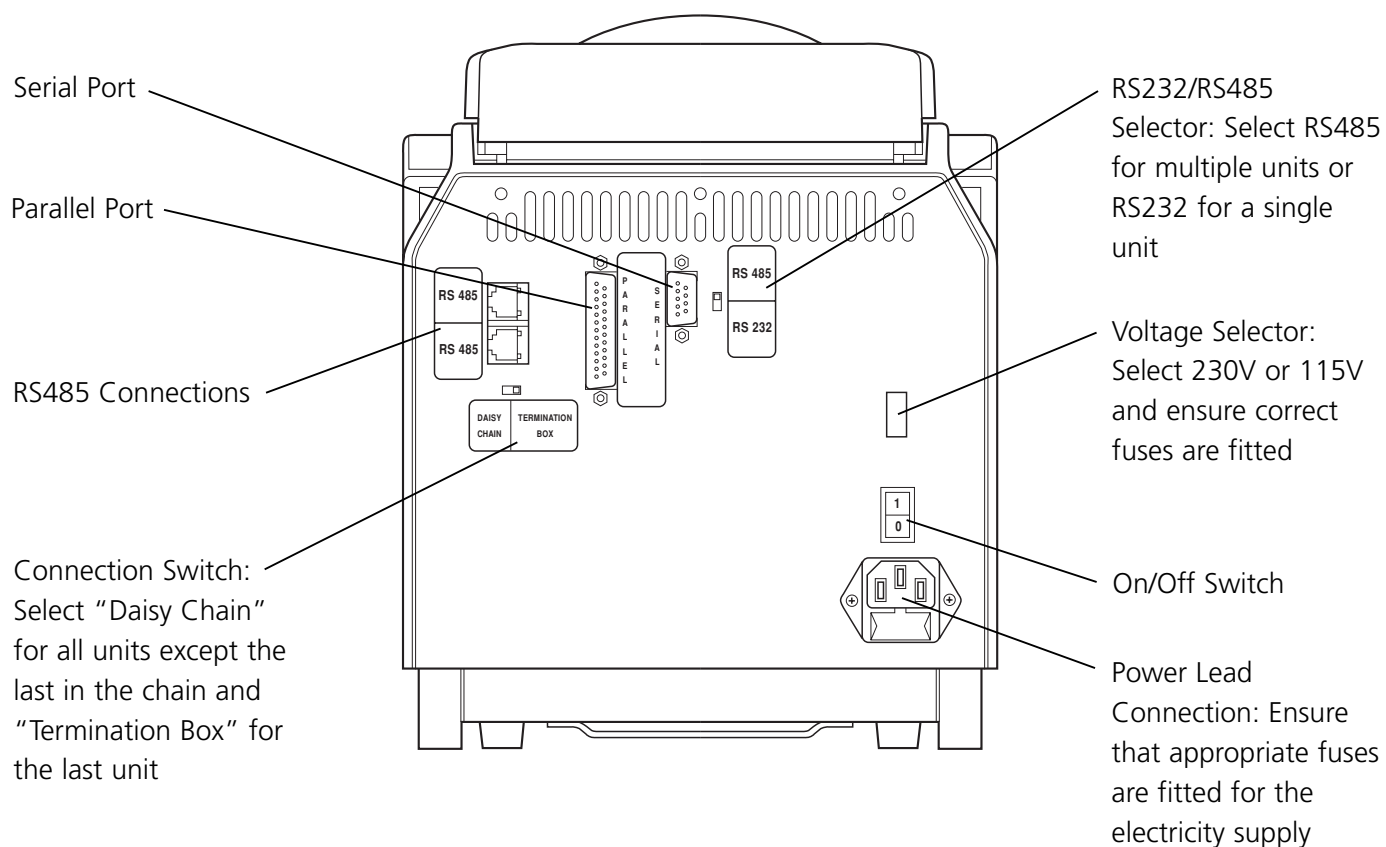

CONTACT INFORMATION

For technical, sales or servicing information, contact your local Bibby Scientific Ltd dealer or:

UK and Rest of World

Bibby Scientific Ltd
Beacon Road, Stone,
Staffordshire, ST15 0SA, UK
Tel: +44 (0)1785 812121
Fax: +44 (0)1785 813748
e-mail: techne@bibby-scientific.com
www.bibby-scientific.com

France

Bibby Scientific France SAS
ZI du Rocher Vert - BP 79,
77793 Nemours Cedex,
France
Tel: +33 1 64 45 13 13
Fax: +33 1 64 45 13 00
e-mail: bsf@bibby-scientific.fr
www.bibby-scientific.com

Middle East

Techne Middle East Ltd.
PO Box 27887, Engomi 2433,
Nicosia,
Cyprus
Tel: +357 22 660427
Fax: +357 22 660356
e-mail: sales@techneme.com

North and South America

Techne Inc.
3 Terri Lane, Suite 10,
Burlington, NJ, 08016, USA
Toll free: 800-225-9243
Tel: +1 609-589-3560
Fax: +1 609-589-2571
e-mail: labproducts@techneusa.com
www.techneusa.com

Italy

Bibby Scientific Italia Srl
Via Alcide de Gasperi 56,
20077 Riozzo di Cerro al Lambro,
Milano, Italia
Tel: +39 (0)2 98230679
Fax: +39 (0)2 98230211
e-mail: marketing@bibby-scientific.it
www.bibby-scientific.it

Far East

Bibby Scientific (Asia) Ltd
Room 607, Yen Sheng Centre,
64 Hoi Yuen Road, Kwun Tong,
Kowloon, Hong Kong
Tel: +852 3583 1581
Fax: +852 3583 1580
e-mail: bibby@bibby-scientificasia.com
www.bibby-scientificasia.com

TC-512 SPECIFICATION

Technical Specification at an ambient temperature of 20°C

Temperature

Gradient Range	20°C to 70°C
Maximum Gradient	30°C
Minimum Gradient	1°C
Number of temperature probes	4
Block temperature range	4°C to 99°C
Typical heating rate	3.0°C/sec
Typical cooling rate	1.3°C/sec
Block uniformity (without gradient)	< $\pm 0.3^{\circ}\text{C}$
Column uniformity with a 10°C gradient	< $\pm 0.3^{\circ}\text{C}$
Control accuracy	< $\pm 0.3^{\circ}\text{C}$
Gradient accuracy	< $\pm 0.3^{\circ}\text{C}$
Temperature set point adjustment	$\pm 0.1^{\circ}\text{C}$

Heated Lid

The heated lid is only operational if the block temperature is set above 35°C	
Selectable heated lid temperature	100°C to 115°C
Heated lid enable/disable	Yes
Warm up time from ambient	4 minutes
Over-temperature cut-out	Yes
Regulated heated lid pressure	Yes

Programs

Password protected	Yes
Incremental/decremental hold time	Yes
Incremental/decremental temperature	Yes
Variable programmable ramp rate steps	$\pm 0.1^{\circ}\text{C}$
Programmable ramp rate	0.1°C
Gradient calculator	Yes
Maximum programs stored internally	84
Maximum number of segments per loop	>60
Maximum number of loops per program	30
Maximum programs stored externally	500
Program naming capacity	Alpha numeric
Program interface	Touch screen
Pause facility	Yes
End of program alarm	Yes (can be disabled)
Auto restart on power failure	Yes
Run completion time	Yes

Miscellaneous

Height x Width x Length	260mm x 220mm x 420mm
Weight	12.5kg
Power consumption	620W (including the heated lid)
Parallel port	Yes
Serial port RS232/RS485 adaptable	Yes
Connection to PC	Yes

TC-512 OPERATION AND MAINTENANCE


The TC-512 has many features which increase the flexibility of the instrument and improve the ease of use. The following features are described in more detail in this chapter:

- Gradient
- Adjustable heated lid
- Memory card
- Adjustable screen
- PC connections

This chapter also includes useful information relating to the type of consumables which can be used in the thermal cyclers.

GRADIENT

A gradient of temperatures across the block can be useful in optimising the annealing conditions for primers; a gradient can be set around a temperature in any step of a program. The set temperature is the temperature required in the middle columns of the block and the gradient is the variation at the two extremes of the block; the left hand columns being the coolest and the right hand columns the hottest.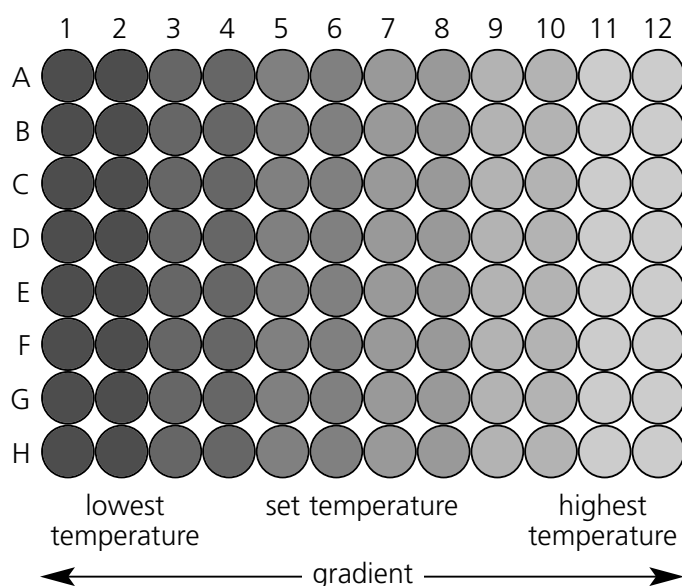

*Schematic diagram
of the 96 well block*

The maximum temperature gradient which can be set is 30°C; the lowest temperature is 20°C and the hottest temperature is 70°C. Examples are given in the table below:

Lowest temperature (°C)	Set temperature (°C)	Highest temperature (°C)	Gradient (°C)
20	35	50	30
40	55	70	30
45	55	65	20
50	55	60	10


It is important to remember that the greater the gradient, the poorer the uniformity across the block, therefore fine-tuning may be required to establish the optimum temperature. While the gradient is mainly linear there are small variations at the extremes of the gradient due to a loss of heat from the hottest part of the block and a gain of heat from the adjacent column on the coldest. This is the reason why the first column does not reach the lowest ideal temperature of the gradient and the last column cannot reach the highest temperature.

A gradient calculator is included in the software which can be used to display more accurately the temperatures of the block for any given gradient. The gradient calculator is located in the System User main menu.

ADJUSTABLE HEATED LID


The heated lid of the TC-512 can be switched on or off depending on the user's requirements. Details for editing the default lid parameters and those of an individual program can be found on pages 41 and 34 respectively.

The lid is unlocked by lifting the back of the handle as shown in the diagram below, this handle can then be used to lift the lid. To close, simply lower the lid onto the block containing the samples and press the handle shut.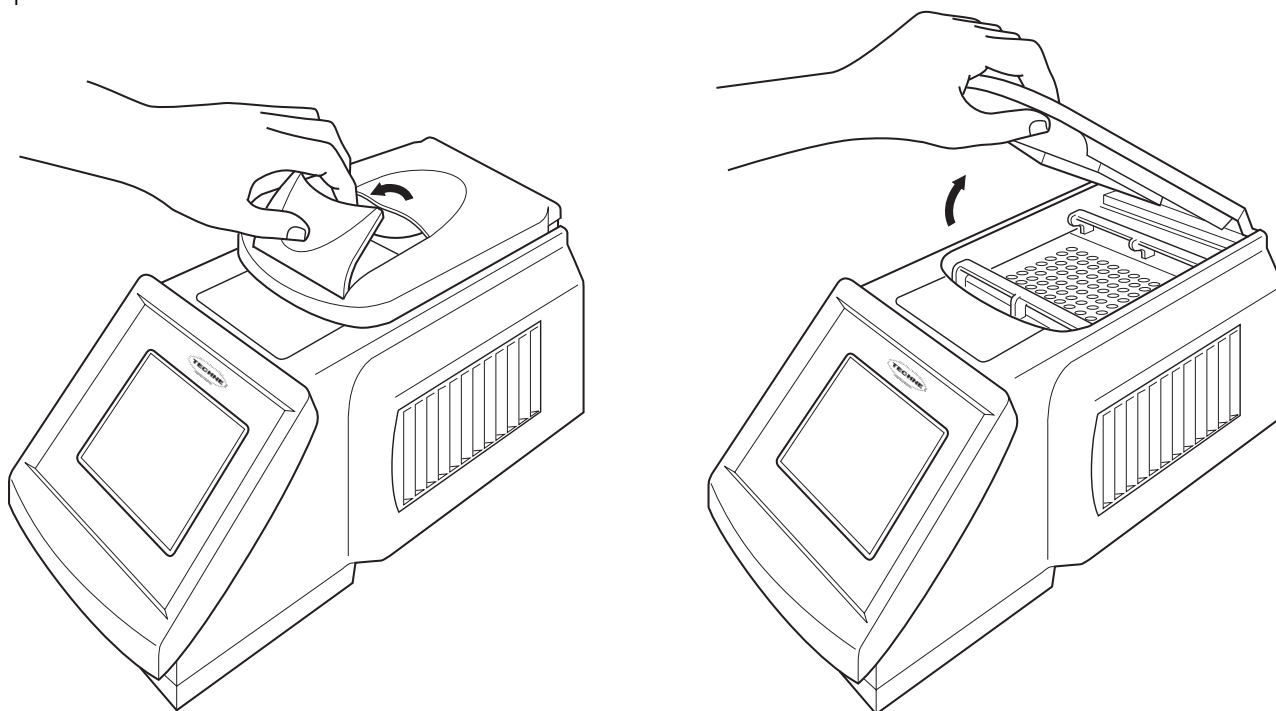

The pressure of the heated lid must be adjusted for each consumable type used in the thermal cycler, to ensure that sufficient pressure is present to prevent any evaporation of samples.

Prior to setting the lid pressure, obtain the lowest pressure by rotating the orange disc anticlockwise. Then shut the lid and latch it. To obtain the correct pressure adjust with the lid in the down position by gently rotating the orange disc clockwise until it is possible to just feel the pressure being applied. Finally, rotate the orange disc a further quarter of a turn; the lid is now at the correct pressure for use with the consumable. The pressure does not need to be adjusted again, unless a different thermal block or consumable is used.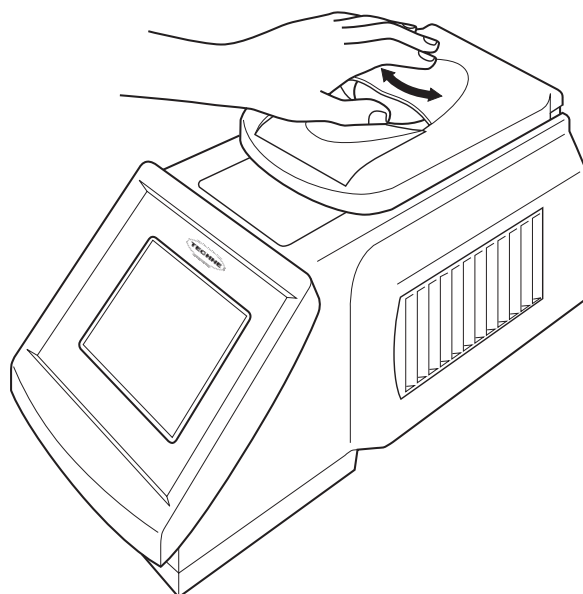


A heated lid mat (Part No. 6104978) is supplied with the TC-512; this is for use with 96-well PCR plates sealed with an adhesive seal. The rubber mat should be placed in top of the sealing film to ensure good thermal contact with the heated lid. Do not over-tighten the lid as this will cause the silicone rubber membrane to deform and will reduce the effectiveness of the heated lid.

MEMORY CARD (not included, must be ordered separately)

The memory card has several functions such as storing programs and temperature logs of completed programs which have been run. It also enables upgrades of the control software to be carried out by either copying software from a new unit to an older one or by obtaining a copy of the software from Bibby Scientific Ltd. The memory card can store up to 500 three segment programs; the number of programs is reduced as the number of segments in each program is increased, to a maximum of 1500 segments in a single program.

The memory card locates in the slot under the silver lid on the front of the instrument. To insert the memory card first lift the heated lid and then the cover and push the memory card into the slot until the button on the left rises. The main label on the memory card should be facing towards the front when the card is positioned correctly.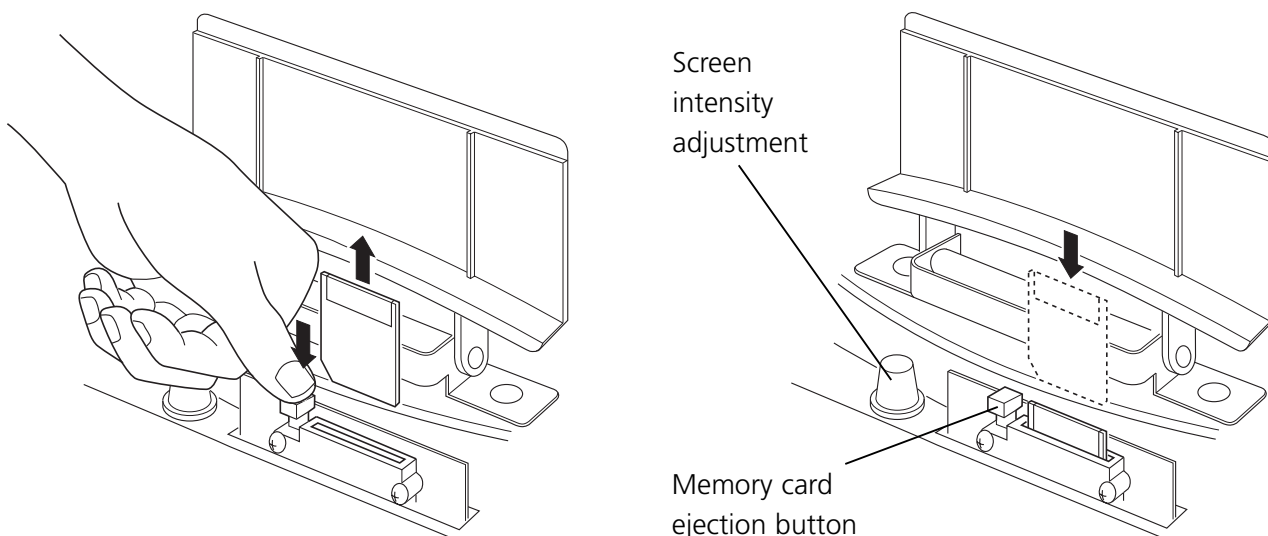

ADJUSTABLE SCREEN

The brightness of the screen can be adjusted according to the light conditions in the laboratory.

The brightness control is located under the front cover of the instrument. After lifting the front cover the small black knob on the left hand side is rotated clockwise to make the screen brighter and anticlockwise to make it darker.

CONSUMABLES


Bibby Scientific Ltd does not recommend any specific tube or reaction vessel other than those described in this manual. The tubes must be able to withstand a pressure of 1 atmosphere at 100°C in addition to the temperatures being used in the thermal cycler block. The tube must not deform to the point of fracture.

We recommend using reaction volumes between 10µl and 50µl. To test the tube, put 25µl of water in each of 5 tubes and subject them to a typical program. At the end of the program measure the volume using a pipette. A loss of more than 1-2µl indicates sample evaporation and a vapour leak. The volume lost and the change in reagent concentrations determines the minimum volume that can be used. Typical volume losses of 1µl in 30 cycles allow the use of samples of 10µl or less.


PC CONNECTIONS

It is possible to connect either a single thermal cycler or a series of instruments to a computer. The options are:

1. To connect a single thermal cycler to a computer part number FGEN232 is required which contains a connecting cable, software disk and instructions.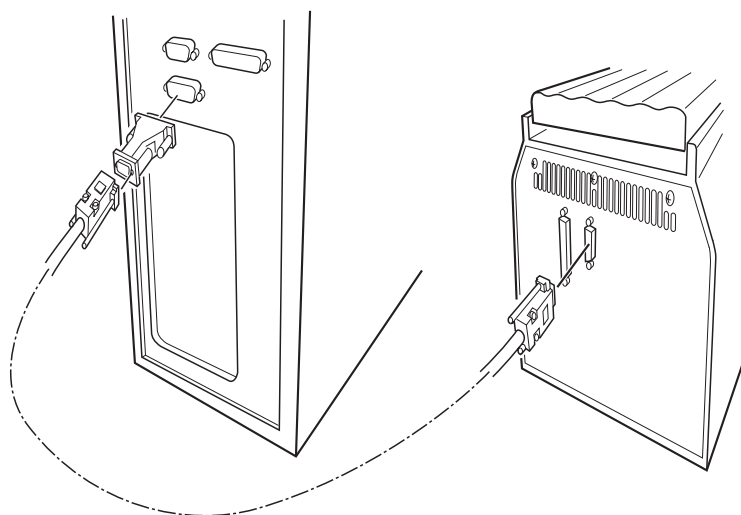

2. To connect several units via a daisy chain to a single computer part number FTGEN485 is required which contains an RS485 set. Included are the connector for the computer, 1m cable for connecting the first instrument to the computer, software disk and instructions. For each unit in the chain a connecting cable (Part number 6103557) is required.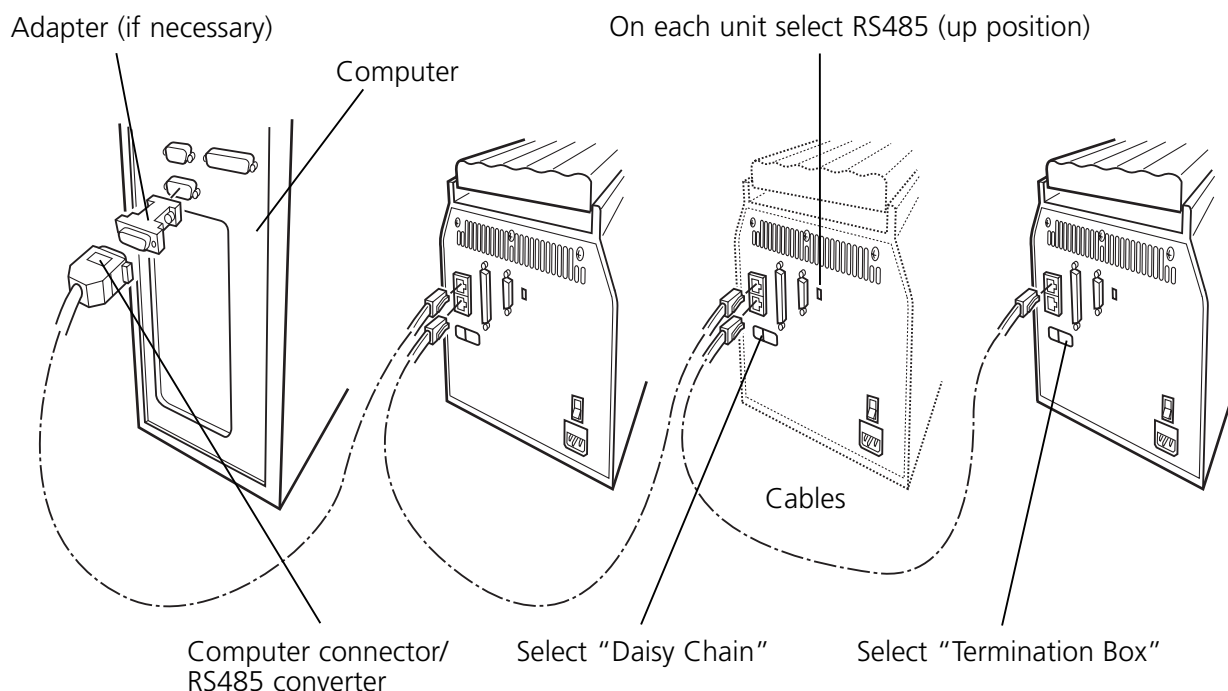

CLEANING THE TC-512

Prior to cleaning the thermal cycler it is important to disconnect the unit from the power supply and if it has been in use allow both the block and heated lid to cool to below 50°C prior to attempting to clean it.

The peltier block, in particular the sample wells should be cleaned regularly to ensure that the maximum heat transfer from the block to the sample occurs. The block should be cleaned if a spillage occurs before the instrument is used again. Use a cloth or cotton bud dipped in 50% propan-2-ol solution (diluted with water) to clean the block ensuring that no deposits are left in the wells of the block.

The case and screen of the thermal cycler can be cleaned using water, ethanol, hexane or 50% methanol (diluted with water). Neither the case nor screen should be immersed in the liquid. Solvents such as acetone should not be used and neither should abrasive cleaners.

If the thermal cycler is being used with any radioactive substances, the peltier block which is made of aluminum, should be cleaned with an agent which is suitable for nonferrous metals, such as Neutracon (from Decon Labs Ltd). However, remember that other parts of the unit are made of ferrous materials and may be damaged by spillage onto them. Neutracon should be used according to the manufacturer's instructions.

REMOVING THE BLOCK


The thermal cycling block can be removed for detailed cleaning or replacement.

First ensure that the unit is disconnected from the power supply and has been allowed to cool if the unit was previously in use. The two quick release handles located at the front and back of the block are used to unlock the block. The padlock symbols indicate the direction in which to slide the handles:

open to unlock and closed to lock. When the block is unlocked it will simply lift out of the unit. Never lift or carry the block by one quick release handle, always use both handles or support the block from underneath.

When refitting the block, slide the quick release handles in the opposite direction to lock the block into the thermal cycler. The software will register and calibrate the unit for the new block.

It may be necessary to edit the heated lid parameters for a new application; for example if the in-situ PCR block is fitted for use with hybridisation slides then it is necessary to switch the heated lid OFF as excess heat can cause evaporation of samples.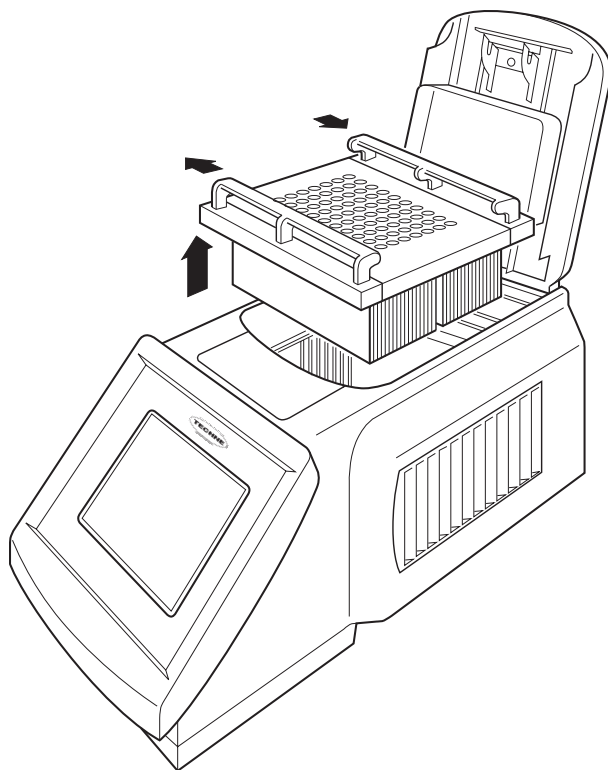

Removal of the TC-512 thermal block

PROGRAMMING THE TC-512

START UP MENU

When the TC-512 is first switched on the screen shown below is displayed and will remain until the unit has finished performing a series of self tests. This screen displays information such as the serial number of the instrument, the type of block being used and the version of software running the thermal cycler.

Bibby Scientific TC-512 TEST UNIT	
Serial number	xxxxxx/xx
Software version	v.vv
Hardware version	h.hh
Block ID	FSG - 60x0.5ml
Last system test	dd/mm/yyyy
Comms address	01
Copyright protected 2003	
Performing self tests	
0 1 2 3	

Bibby Scientific TC-512 TEST UNIT	
Serial number	xxxxxx/xx
Software version	v.vv
Hardware version	h.hh
Block ID	FSG - 60x0.5ml
Last system test	dd/mm/yyyy
Comms address	01
Copyright protected 2003	
System User	System Admin

After the self tests have been completed the screen opposite is shown unless the unit is powered up by auto-restart. If there is a program running and a power failure occurs, the auto-restart function, if enabled, will immediately restart the program once power is restored. The instrument does not perform the self tests at this point. Once the program is complete and has been stopped then the self tests will be initiated.

The System User section of the software allows the user to program the thermal cycler for his/her experiments and view or print the data logs. The System Admin contains the settings for the instrument itself.

THE SYSTEM USER MENU

In this section of the software the user is able to write new programs and edit existing ones. Additionally the program defaults for parameters such as the heated lid and final extension steps can be altered. The gradient calculator and data log files can also be accessed through their sub-menus.

The top screenshot shows the 'System User' menu. It has a header 'Bibby Scientific TC-512 TEST UNIT'. Below it is a table of system information:

Serial number	xxxxxx/xx
Software version	v.vv
Hardware version	h.hh
Block ID	FSG - 60x0.5ml
Last system test	dd/mm/yyyy
Comms address	01
Copyright protected 2003	

At the bottom are two buttons: 'System User' (highlighted) and 'System Admin'.

The bottom screenshot shows the main menu. It has the same header 'Bibby Scientific TC-512 TEST UNIT'. Below it are five buttons: 'Programs', 'Program Defaults', 'Data Logs', 'Gradient Calculator', and 'Exit'.

To access the programming software press the "System User" button on the start-up screen. The System User menu will then appear.

The options within the System User menu include:

Programs: The main functions are reading, writing, editing and running of new or previously stored programs.

Program Defaults: Here the user is able to edit the default settings for the heated lid, initial denaturation, final extension, final hold and the hot start functions, which will then automatically be inserted into any future programs.

Data Logs: Allows the user access to the temperature log files from the previously run programs; the logs can be printed if required. This function is only accessible if a SmartMedia memory card is inserted into the appropriate slot (see page 24).

Gradient Calculator: Calculates the temperature for each column of the block for a user-specified temperature and gradient.

Exit: Returns to the start-up menu.

WRITING AND EDITING PROGRAMS

There are three options available for creating programs on the TC-512:

1. Copy and edit one of the pre-written templates.
2. Edit a pre-written template and save it using a different name.
3. Create a new program with user-specific parameters.

There are 5 pre-written templates. These templates cannot be altered; however, they can be copied, given a new name and then edited to save time when writing a new program.

The Ice Bucket and Ligation 15 Degrees programs are simple temperature holds, with the former having an infinite hold at 4°C and the latter at 15°C. In both of these programs the heated lid is switched off and the initial denaturation, hot start and final extension functions disabled.

The 2 Step and 3 Step PCR templates along with the RT PCR template have the following default parameters:

Parameter	2 Step PCR	3 Step PCR	RT PCR
Heated Lid	105°C	105°C	105°C
Heated lid before program	On	On	On
Pause before program	Off	Off	Off
Initial denaturation	05m00s 94.0°C	05m00s 94.0°C	05m00s 94.0°C
Hot start	Disabled	Disabled	Disabled
Final extension	05m00s 72°C	05m00s 72°C	07m00s 72°C
Final hold	10°C	10°C	10°C

The thermal cycling conditions are different for each of these three templates:

	2 Step PCR	3 Step PCR	RT PCR
Cycles			1
Segment			40m00s 45.0°C
Segment			05m00s 94.0°C
Cycles	30	30	40
Segment	00m30s 92.0°C	00m30s 92.0°C	01m00s 94.0°C
Segment	01m00s 68.0°C	00m30s 55.0°C	00m50s 55.0°C
Segment		00m30s 72.0°C	01m00s 72.0°C

The RT PCR template is for those experiments where the starting material is RNA rather than DNA. RNA requires reverse transcription by a reverse transcriptase enzyme into complementary DNA (cDNA) prior to the amplification by the DNA polymerase in PCR. The extra 45°C step is for the reverse transcription and the 94°C step is to denature the reverse transcriptase before the PCR.

Bibby Scientific
TC-512 TEST UNIT

Programs

Program Defaults

Data Logs

Gradient Calculator

Exit

OPTION 1: COPYING AN EXISTING TEMPLATE

To access the program templates select "Programs" in the System Users menu.

Select Program
14526 bytes of memory free

b 2 STEP PCR TEMPLATE
b 3 STEP PCR TEMPLATE
b ICE BUCKET
b LITIGATION 15 DEGREES
b RT PCR TEMPLATE

▼ ▲

New

Exit

The cursor arrows can be used to scroll up and down the stored programs. An individual program is opened by pressing on the individual program box (shown opposite).

The programs are pre-fixed by b, i or r. The b corresponds to the pre-programmed templates, i for those stored on the internal memory and r for remote and those stored on the SmartMedia memory card. Selecting "New" will open an empty program. See page 34 for more details.

Once the template is open select "Copy". It is also possible to run or print the program by pressing the appropriate button.

Program 3 STEP PCR TEMPLATE

Last Modified	01/11/1998	00:00:00
Heated lid	105.0°C	
Heated lid before program	On	
Pause before program	Off	
Initial denaturation	05m00s	94.0°C
Hot Start	Disabled	
Number cycles	30	

▼ ▲

Copy

Delete

Edit

Run

Print

Exit

Enter the new program name

Q	W	E	R	T	Y	U	I	O	P
A	S	D	F	G	H	J	K	L	
Num	Z	X	C	V	B	N	M	Enter	
Cap	←	Space				→	BS	Cancel	

It is then necessary to enter a new program name. There are several text features within the text screen. Pressing the following buttons allows access to various options:

Num: Numerical keys

Cap: Capital and lowercase letters

BS: Backspace, removes the previous character

Arrow cursors: Move the cursor backwards and forwards

Space: Introduces a space.

When finished press "Enter" to return to the Select Program menu.

Select Program
14526 bytes of memory free

b ICE BUCKET
b LITIGATION 15 DEGREES
b RT PCR TEMPLATE
i B ACTIN
I GAPDH

▼▲

New Exit

Select the copied program.

Program B ACTIN

Last Modified	15/05/2006	11:29:17
Heated lid	105.0°C	
Heated lid before program	On	
Pause before program	Off	
Initial denaturation	05m00s	94.0°C
Hot Start	Disabled	
Number cycles	30	

▼▲

Copy Delete Edit

Run Print Exit

The program cannot be changed at this point so press "Edit" to make any changes required.

The program can be copied, deleted, printed or run by selecting the appropriate option.

Editing B ACTIN

Initial denaturation	05m00s	94.0°C
Hot Start	Disabled	
Number cycles	30	
Seg	MAX°C/s	00m30s 92.0°C
Seg	MAX°C/s	00m30s 55.0°C
Seg	MAX°C/s	00m30s 72.0°C
Next step		

▼▲

Ins Del

PWord Save As Save Exit

All of the parameters can be edited simply by selecting each line in turn. The cursors can be used to navigate through the program as all the parameters cannot be viewed at the same time. To save any changes select "Save" and a prompt will ask:

Save

Yes No

and exit back to the program. For details of all the features and parameters see pages 34.

Select Program
14526 bytes of memory free

b 2 STEP PCR TEMPLATE
b 3 STEP PCR TEMPLATE
b ICE BUCKET
b LITIGATION 15 DEGREES
b RT PCR TEMPLATE

▼ ▲

New Exit

OPTION 2: EDITING AND SAVING AN EXISTING TEMPLATE

Select the template from the program menu.

Program 3 STEP PCR TEMPLATE

Last Modified	01/11/1998	00:00:00
Heated lid	105.0°C	
Heated lid before program	On	
Pause before program	Off	
Initial denaturation	05m00s	94.0°C
Hot Start	Disabled	
Number cycles	30	

▼ ▲

Copy Delete Edit

Run Print Exit

Select "Edit"

Editing 3 STEP PCR TEMPLATE

Initial denaturation	05m00s	94.0°C
Hot Start	Disabled	
Number cycles	30	
Seg	MAX°C/s	00m30s 92.0°C
Seg	MAX°C/s	00m30s 55.0°C
Seg	MAX°C/s	00m30s 72.0°C
Next step		

▼ ▲

Ins Del

PWord Save As Exit

All of the parameters can be edited simply by selecting each line in turn. The cursors can be used to navigate through the program as all the parameters cannot be viewed at the same time. Once all the changes have been made, select "Save As"; there is no option to save. For details of all the features and parameters see page 34.

Enter the new program name

Q	W	E	R	T	Y	U	I	O	P
A	S	D	F	G	H	J	K	L	
Num	Z	X	C	V	B	N	M	Enter	
Cap	←	Space				→	BS	Cancel	

Enter a new program name. There are several text features within the text screen. Pressing the following buttons allows access to various options:

Num: Numerical keys

Cap: Capital and lowercase letters

BS: Backspace, removes the previous character

Arrow cursors: Move the cursor backwards and forwards

Space: Introduces a space.

When finished press "Enter" to return to the program. Select "Run" to start the program.

Select Program
14526 bytes of memory free

b 2 STEP PCR TEMPLATE
b 3 STEP PCR TEMPLATE
b ICE BUCKET
b LITIGATION 15 DEGREES
b RT PCR TEMPLATE

▼ ▲

New Exit

Editing B ACTIN

Initial denaturation	Disabled
Hot Start	Disabled
Next Step	
Final extension	Disabled
Final hold	Disabled

▼ ▲

Ins Del

PWord Save As Save Exit

OPTION 3: CREATING A NEW PROGRAM

The final option is to create a completely new program by selecting "New" in the Select Program menu. This section also contains details of the individual parameters and functions.

To create a new program, select "New" in the Select Program menu.

Type in a program name; for details see the previous page.

The cursor icons can be used to scroll up and down to access more parameters. Scrolling up the menu displays the default parameters:

- Heated lid
- Pause before the program
- Initial denaturation
- Hot start
- Final extension
- Final hold

The default parameters are defined in the System user menu and information about the programming of these parameters can be found on page 41. All of the default parameters can be set uniquely within an individual program.

The keys in the editing menu are as follows:

- "Del" Delete will remove only the highlighted step. The default parameters cannot be deleted.
- "Ins" Insert will add a next step if no "Next step" is present.
- "Save As" Allows the program to be saved with an alternative name.
- "Save" Saves the program.
- "PWord" Adds a password which protects the program so other users cannot change it.
- "Exit" This returns the screen to the Select Program menu.

Next Step

The Next step line is inserted automatically between the Hot Start and Final extension. Select "Next Step" to open the Edit program function menu. If Next Step is not displayed then selecting "Ins" will add a Next Step below the highlighted line.

Edit program function						
Next step						
Segment	7	8	9	←	→	
Number Cycles	4	5	6	Continue		
Inc Dec	1	2	3	Cancel		
Pause	0	.	BS	OK		

The next step enables one of the following to be added to the program:

Segment: Enables the temperature, time and heating rate to be inserted into the segment. Additionally, a gradient across the block can be added.

Number Cycles: Up to 99 cycles can be added

Inc Dec: Allows increasing and decreasing temperatures and times to be set. See page 36 for further details.

Pause: Adds a pause as a step. User intervention will then be required before the program can move on to the next step.

INSERTING A SEGMENT

Selecting "Segment" inserts a temperature step of 94°C for 30s at the maximum ramp rate.

Edit program function						
Segment 94°C ▽ 0°C 00m30s MAX 0°C/s						
Segment	7	8	9	←	→	
Number Cycles	4	5	6	Continue		
Inc Dec	1	2	3	Cancel		
Pause	0	.	BS	OK		

Segment 94°C ▽ 0°C 00m30s MAX 0°C/s			
-------------------------------------	--	--	--

If required change the temperature using the numerical keypad. Then use the → to move across the screen to the gradient option ▽ 0°C.

The gradient on the TC-512 can be set from 1 to 30°C across the temperature range 20 to 70°C. Outside this range the accuracy of the different columns cannot be guaranteed. If a gradient greater than 30°C or outside the temperature range is set then the software will automatically set it to the maximum, 30°C or within the permitted range. If temperatures are set outside 20 to 70°C range then the following error message is displayed:

WARNING! Gradient has been adjusted to be within cycler limits Yes
--

Further information relating to the gradient function can be found on page 22. Use the → to move across the screen to set the time.

Edit program function									
Segment		94.0°C		▢ 0°C		00m30s		MAX 0°C/s	
Segment		h	7	8	9	←	→		
Number Cycles		m	4	5	6	Continue			
Inc Dec		s	1	2	3	Cancel			
Pause			0	.	BS	OK			

The time can be set in hours (h), minutes (m) or seconds (s). Type in the time required followed by the unit. Use the → to move across the screen to the ramp rate.

Edit program function									
Segment		95.0°C		0°C		00m15s		MAX 0°C/s	
Segment		Max	7	8	9	←	→		
Number Cycles			4	5	6	Continue			
Inc Dec			1	2	3	Cancel			
Pause			0	.	BS	OK			

The ramp rate is automatically set to the maximum 3°C per second. A number less than 2°C/s can be inserted. Selecting "Max" will reset the ramp to the maximum.

Selecting "OK" leaves the Edit program function menu and returns to the program.

Selecting "Continue" adds another new step to the program.

If a segment or the Inc/Dec function is inserted then the software will automatically assign 1 cycle to this stage or loop, as it is referred to in the software.


Edit program function									
Number cycles		01							
Segment			7	8	9	←	→		
Number Cycles			4	5	6	Continue			
Inc Dec			1	2	3	Cancel			
Pause			0	.	BS	OK			

INSERTING CYCLE NUMBERS

Selecting "Number Cycles" allows the number of cycles to be increased up to 99. This function allows the user to easily program the 2 or 3 step section of the PCR which usually involves between 20 and 45 repeats of 2 or 3 temperature steps. The program will automatically repeat in the programmed sequence each of the temperature steps for the required number of cycles.

INSERTING INCREMENTAL OR DECREMENTAL TIME OR TEMPERATURE


Under normal circumstances, the Hold Temperature of all steps is constant. However, it is possible to automatically increase or decrease the temperature of a specified step of a programme. Touchdown amplification of DNA requires a higher temperature during the initial cycles to ensure that only the specific piece of DNA is amplified and not non-specific products. A decreasing, increasingly less stringent annealing temperature is then used for the later cycles of the PCR to increase the yield of the product.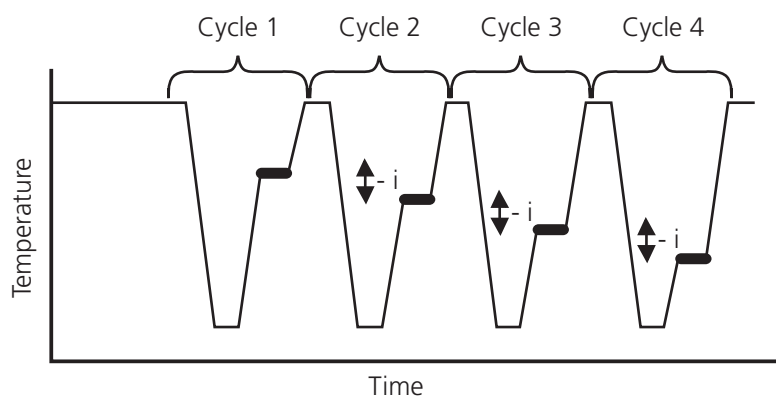

Decremental temperature for Touchdown DNA amplification

The Hold Temperature of the incremented/decremented step is the defined Hold Temperature plus the summation of the increments/decrements. The first cycle is never altered, only subsequent cycles. If decrement is selected, the Hold Temperature is prevented from falling below 4°C. If increment is selected the Hold Temperature is prevented from rising above 99°C.

Similarly, under normal circumstances, the Hold Time of all steps is constant. However, it is possible to automatically increment or decrement the duration of a specified step of a programme. When amplifying large fragments of DNA it is necessary to ensure that the extension is fully complete after each cycle, so the extension time is increased with each cycle to compensate for the increased product quantity and decreasing concentrations of reagents.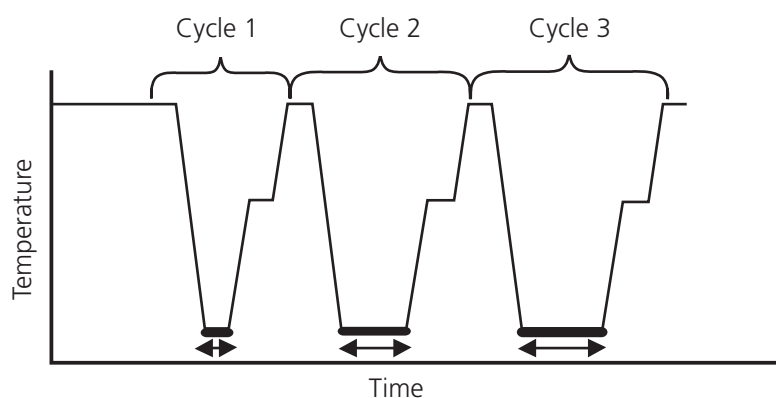

Incremental time for amplification of long DNA products

The Hold Time of the incremented/ decremented step is the defined Hold Time plus the summation of the increments/decrements. The first cycle is never altered, only subsequent cycles. If you select a decrement, the Hold Time is prevented from falling below one second.

Edit program function					
PAUSE					
Segment	7	8	9	←	→
Number Cycles	4	5	6	Continue	
Inc Dec	1	2	3	Cancel	
Pause	0	.	BS	OK	

INSERTING A PAUSE

Selecting "Pause" adds a stop to the program, which when running requires user intervention before the program can continue.

Editing		B ACTIN	
Heated lid		105.0°C	
Heated lid before program		On	
Pause before program		Off	
Initial denaturation	10m00s	75.0°C	
Hot Start		95.0°C	
Next step			
Final extention	07m00s	72.0°C	
▼▲	Ins	Del	
PWord	Save As	Save	Exit

SAVING CHANGES TO A PROGRAM

To change a saved program, select the program and when opened select "Edit" to change the parameters.

Select "Save" to save the changes under the existing program name or "Save As" to save the changes as a new program.

RUNNING A PROGRAM

To run any previously saved program simply follow these steps:

Bibby Scientific TC-512 TEST UNIT			
Programs		Program Defaults	
Data Logs	Gradient Calculator	Exit	

Select "Programs" from the main System User menu.

Select Program
14526 bytes of memory free

b 2 STEP PCR TEMPLATE
b 3 STEP PCR TEMPLATE
b ICE BUCKET
b LITIGATION 15 DEGREES
b RT PCR TEMPLATE

▼ ▲

New Exit

Select the program required, using the cursor keys to view all the available programs.

Program 3 STEP PCR TEMPLATE

Last Modified	01/11/1998	00:00:00
Heated lid	105.0°C	
Heated lid	On	Off
Pause be	Disabled	
Initial de		
Hot Star		
Number		

Do you wish to log this Program?

Yes No

▼ ▲

Copy Delete Edit

Run Print Exit

Press "Run".

"If a SmartMedia card is inserted in the cycler then the software will ask if the Data Log needs to be saved.

If "Yes" is selected then the data log file will require a file name to be inserted.

Program B ACTIN

Data Log None

Waiting Heated Lid 03m40s


Stop Pause

The run screen is then displayed; if the heated lid is on then the screen will show the message "Waiting Heated Lid" and the time until the lid will be ready. It is possible to stop or pause the program.


If "Stop" is pressed then it is necessary to confirm that that the program is to be stopped by pressing "Stop" again.

If pressing "Stop" was an error then pressing "Continue" will enable the program to continue from where it was stopped.


If "Pause" was selected then press "Unpause" to restart the program. The pause function might be required for a manual hot start.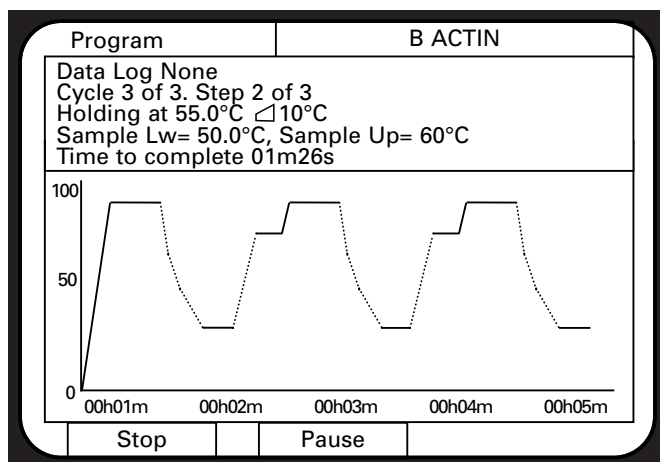


As the program continues the temperatures are displayed on the screen, as shown opposite.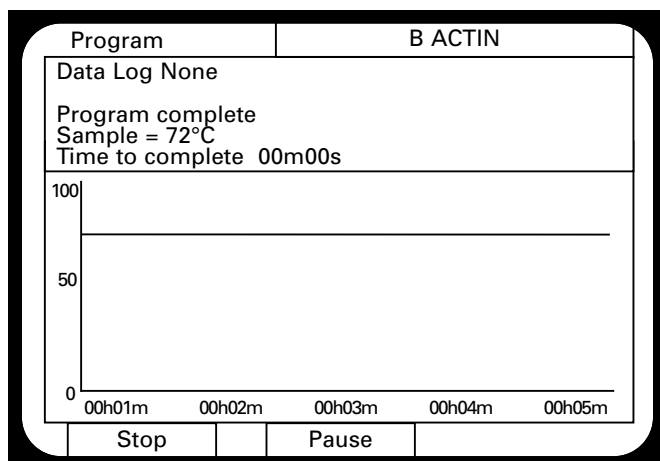

This screen is displayed when the run is complete. If there is a final hold in the program press "Stop" to end the program.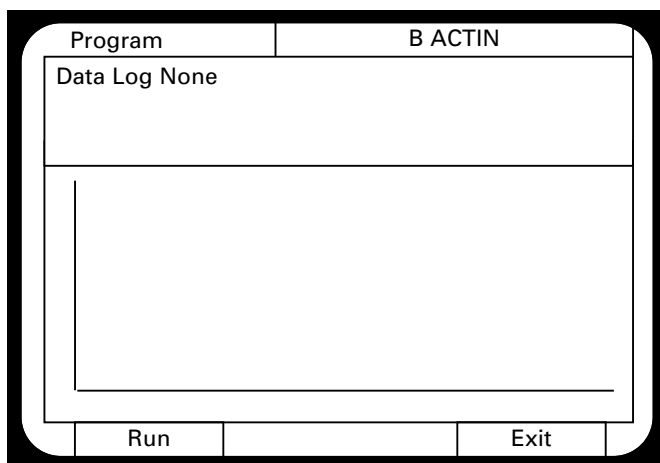

Selecting "Run" at this point will start the program from the beginning again and "Exit" closes the run screen and displays the selected program again.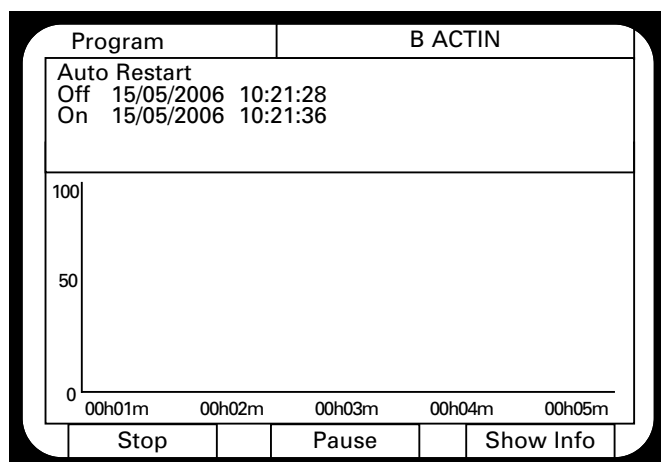



This screen is displayed if there is a power loss during the running of the program and the auto-restart function is switched on. The display shows the times at which the power failed and resumed.

The "Show Info" button takes the screen back to the program information and display.

When the program has finished or been stopped then the instrument will perform its normal start-up tests.

PROGRAM DEFAULTS

It is possible to change the pre-defined parameters which are used for every program by editing the Program Defaults. This information can be found in the System User menu.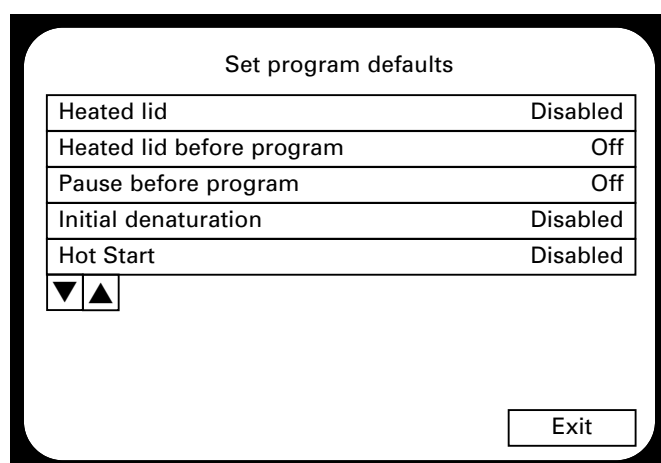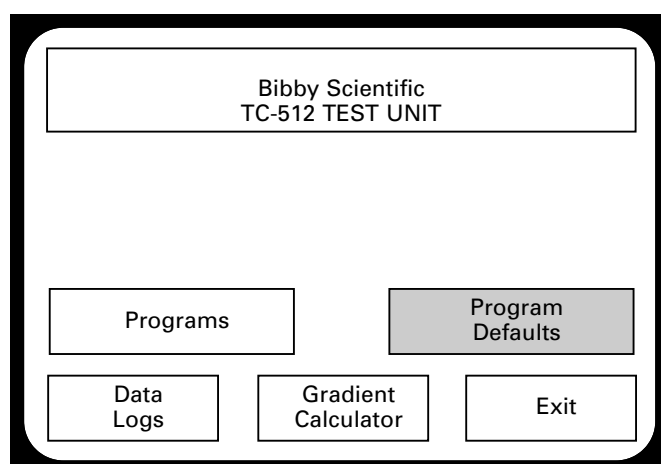

To access all the parameters it is necessary to move up and down the menu using the cursor arrows. The parameters which can be edited are:

Heated lid: Select the "Heated Lid" and then in the temperature editing window it is possible to enable/disable the heated lid. If "Enabled" is pressed the text changes from xxxxx to allow a temperature to be typed in. The lid can be programmed between 100 and 115°C.

Heated lid before program: Selecting this option simply moves between On and Off. This option ensures that the heated lid is at the correct temperature before starting the actual program.

Pause before program: This allows the insertion of a pause in the experiment before the experiment starts. After the pause there must be user intervention before the program can move on to the next step of the program.

Edit initial denaturation setting

Initial denaturation		05m00s			94°C	
h	7	8	9	←	→	
m	4	5	6	Enable		
s	1	2	3	Cancel		
	0	.	BS	OK		

Initial denaturation: Many experimental protocols require a high temperature incubation to remove any secondary structure of the nucleic acid template.

Selecting "Enable" allows a temperature and time to be added. The cursors are used to move between the time and the temperature.

Edit initial denaturation setting

Initial denaturation		xxxxxx			xxxxx°C	
h	7	8	9	←	→	
m	4	5	6	Enable		
s	1	2	3	Cancel		
	0	.	BS	OK		

On the left of the numbers there are 3 buttons for setting the time: h for hours, m for minutes and s for seconds. The correct button should be pressed after the temperature has been typed.

Hot start: Many standard enzymes used in molecular biology techniques can hybridise to and extend non-specific products, resulting in multiple bands on an agarose gel, a greatly reduced yield or no amplification of the specific product. The 'Hot Start' programming step is used to pause the instrument at a specific temperature, typically greater than 70°C, after the initial template denaturation. The reason is to allow the manual addition of unmodified *Taq* DNA polymerase which may lose activity if added during the initial 5 minute denaturation. Heat-activated *Taq* or Hot Start enzymes do not require this step.

Final Extension: Many PCR experiments require a final step at the end of the cycling phase to allow the polymerase to complete the synthesis of the DNA strands. By selecting "Enable" it is possible to put a time and temperature for the step. A typical example is 72°C for 7 minutes.

Final Hold: Frequently the samples are required for further analysis such as agarose gel electrophoresis and the samples are better stored at a temperature lower than ambient. Selecting "Enable" allows the samples to be held indefinitely between 4 and 99°C. This step requires the user to stop the program manually on the instrument. Samples are normally held between 4 and 10°C.

DATA LOGS

The data log contains information relating to the performance of the thermal cycler, which may need to be recorded as part of the experimental procedure, for example, Good Laboratory Practice. The log file contains the temperatures and times of the steps throughout the run, in addition to information relating to the instrument such as serial number and last test date.

Bibby Scientific
TC-512 TEST UNIT

Programs Program Defaults

Data Logs Gradient Calculator Exit

To access the log files select "Data Logs" in the main System User menu.

Set program defaults

B ACTIN

▼ ▲

Print data log Delete data log Exit

All of the programs for which a data log has been generated are displayed. The cursors can be used to access further files. To select an individual log simply press it to select. There is not an option to view the log file on the actual thermal cycler as there is too much information to display. It is possible to either print it or delete it.

PRINTING A DATA LOG

To print the data log it is necessary to have a printer connected directly to the thermal cycler. The printer must be switched on and set on-line before the "Print data log" button is pressed. The printed profile is 80 columns wide and uses only standard characters in the ASCII range, 1 to 127, so all common dot matrix, pin-fed, continuous stationery printers can be used. For example one printer which could be used is the Epson dot matrix fan-form LX300+.

Example of the Data Log

TC-512 program				30/11/98		14:27:25	
				System Configuration			
				System Version:		0.49 1.00	
				Unit Serial No:		88192/01	
				Last System Test:		19/11/1988 09:34:24	
Program: DEPT PROGRAM							
Date last edited: 12/10/1988 17:46:11							
Data log file: B ACTIN							
Event	Date & Time	Func	Cycle	Step	Temp	Set point	Ramp Rate
Program	30/11/98 14:27:26	01	01				
Ramping	30/11/98 14:27:26	01	01	01	4.0°C	4.0°C	- MAX °C/s
Holding	30/11/98 14:27:51	02	01	01	6.3°C	4°C	
Ramping	30/11/98 14:37:52	02	01	01	4.0°C	95.0°C	+ MAX °C/s
Holding	30/11/98 14:38:29	02	01	01	88.1°C	95.0°C	
Ramping	30/11/98 14:38:59	02	01	02	95.0°C	55.0°C	- MAX °C/s
Holding	30/11/98 14:39:26	02	01	02	59.2°C	55.0°C	
Ramping	30/11/98 14:39:56	02	01	03	55.0°C	72.0°C	+ MAX °C/s
Holding	30/11/98 14:40:05	02	01	03	65.5°C	72.0°C	
Ramping	30/11/98 14:41:05	02	02	01	72.0°C	95.0°C	+ MAX °C/s
Holding	30/11/98 14:41:16	02	02	01	88.5°C	95.0°C	
	Signed:						

The data log contains information relating to the events which occurred during the program. The recorded events and descriptions are shown below:

Start: Start of program event is generated when the run button is first pressed.

Stop: Created by pressing the "Stop" button to stop a run prematurely.

Start of program: Created whenever a new program starts.

Start of cycle: Created when the current program starts a new cycle.

Start of segment: Created when a new segment (step) starts.


Start of hold time: Created when the segment reaches the hold window.

End of program: Created when the program finishes.

Pause on: Created when the "Pause" button is first pressed to pause a program.


Pause off: Created when the "Unpause" button is subsequently pressed to end a pause.

Typical graph
from a data log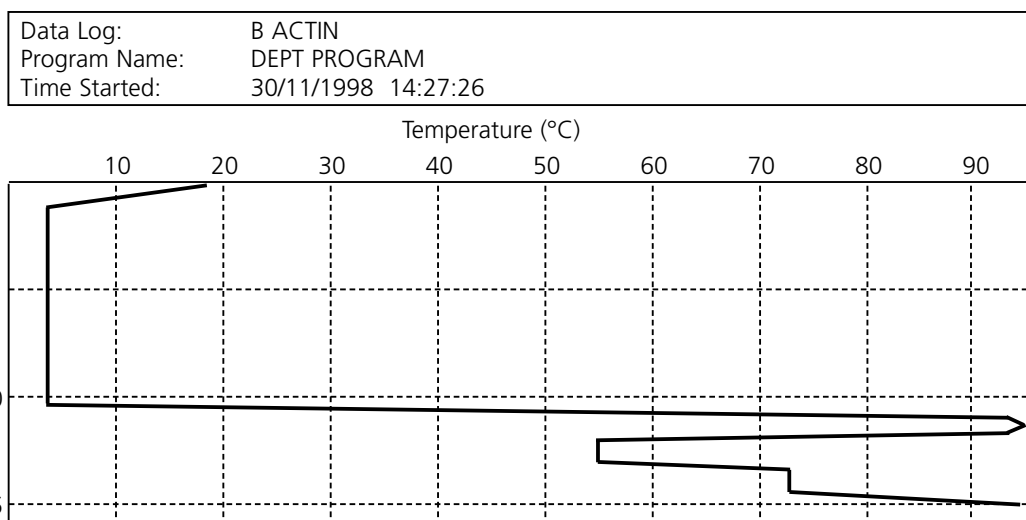


GRADIENT CALCULATOR

The gradient is set within the program (see page 35 for further details), however, it is not possible to view the actual temperatures of each column in the program. The gradient calculator calculates the temperatures for each of the 12 columns (24 for the 384 well block) and allows the user to determine the optimal amplification conditions for the assay.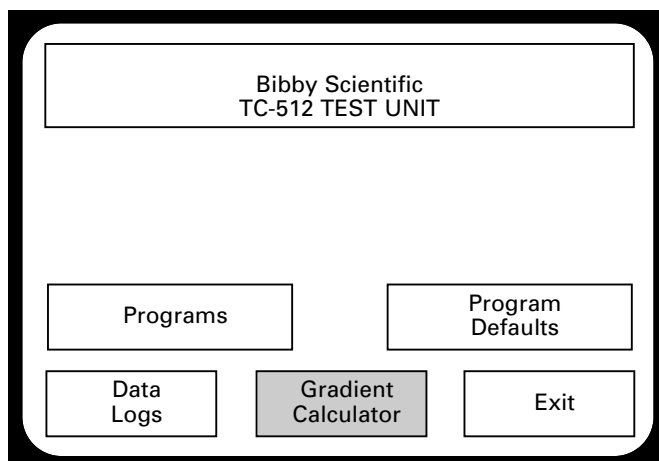

The Gradient Calculator is found in the System User menu.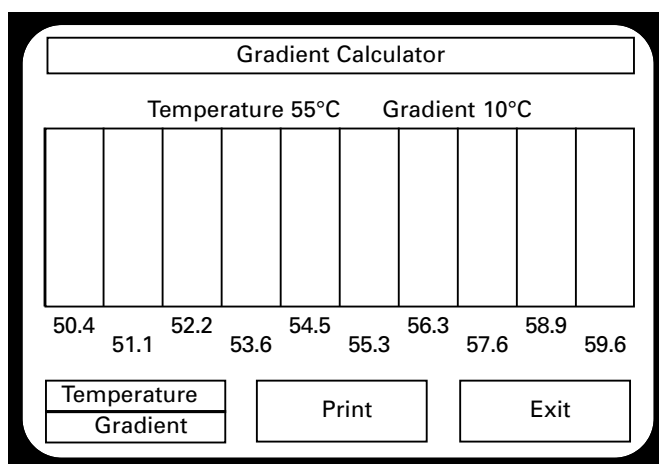

Selecting "Temperature" enables the temperature of the middle of the gradient to be set. The gradient can be set between 20 and 70°C.

Selecting "Gradient" allows the range of the gradient to be set. The minimum is 1°C and the maximum is 30°C.

An example for the 60 x 0.5ml block is shown. The gradient is 10°C with the middle of the block set to 55°C, so the block will be approximately 50°C on the left and 60°C on the right and 55°C in the middle of the block.

It is possible to print the column and gradient information by selecting "Print". Selecting "Exit" returns the screen to the System User menu.

SYSTEM ADMINISTRATION

This section of the software contains settings for the instrument, rather than the actual programs. From various sub-menus it is possible to change passwords for programs and the instrument, test the instrument, delete current programs stored on the memory card and completely reset the instrument.

Bibby Scientific TC-512 TEST UNIT	
Serial number	xxxxxx/xx
Software version	v.vv
Hardware version	h.hh
Block ID	FSG - 60x0.5ml
Last system test	dd/mm/yyyy
Comms address	01
Copyright protected 2003	
System User	System Admin

To access the System Admin simply press the "System Admin" button on the start-up screen.

Enter password																

Q	W	E	R	T	Y	U	I	O	P							
		A	S	D	F	G	H	J	K	L						
Num	Z	X	C	V	B	N	M	Enter								
Cap	←	Space						→	BS	Cancel						

The software will then ask for a password. Type in the default password TECHNE using the touch screen as shown opposite.

To exit the password screen press "Cancel".

Pressing "Enter" takes the user through to the system administration menu.

Bibby Scientific TC-512 TEST UNIT		
		Program Passwords
Sys Admin Password	System Test	Controller Reset
Card Utils	System Config.	Exit

The System Administration Menu.

The following functions are located within the System Administration menu:

Program Passwords: Enables an administrator to change the password for any program.

Sys Admin Password: Allows the administrator to change the administrator password.

System Test: Initiates a simple diagnostic test.

Controller Reset: Allows the user to restore the factory settings.

Card Utilities: Using this function SmartMedia cards can be formatted or all files can be deleted.

System Configuration: This sub-menu allows the user to change the format of dates, times and temperature units. Language, auto restart settings and the comms address can also be altered. Additionally information relating to the heated lid and cycle count of the block is stored in this section of the software.

Select Program

b 2 STEP PCR TEMPLATE
b 3 STEP PCR TEMPLATE
b ICE BUCKET
b LITIGATION 15 DEGREES
b RT PCR TEMPLATE

▼ ▲

Change PWord Exit

PROGRAM PASSWORDS

Selecting the "Program Passwords" function allows the user to change the passwords for any password protected program. First touch the screen to select the program for which the password needs to be changed and then select "Change PWord". To change the password simply type the new password by touching the letters/numbers and then confirm by pressing "Enter". It is then necessary to confirm the new password by re-typing it.

New system administrator password

Q	W	E	R	T	Y	U	I	O	P
A	S	D	F	G	H	J	K	L	
Num	Z	X	C	V	B	N	M	Enter	
Cap	←	Space	→	BS	Cancel				

SYSTEM ADMINISTRATION PASSWORD

Selecting "Sys Admin Password" enables the user to change the initial administrator password.

The screen shown opposite is then displayed. To change the password simply type the new password by touching the letters/numbers and then confirm by pressing "Enter". It is then necessary to confirm the new password by re-typing it.

Bibby Scientific
TC-512 TEST UNIT

RTC OK

Exit

SYSTEM TEST

This is a diagnostic test to ensure that the Real Time Clock settings are correct.

Bibby Scientific
TC-512 TEST UNIT

*** Controller Reset ***
WARNING!!
This will reset the unit to
system defaults. All stored
files will be lost.

Calibration data is retained

Confirm
Reset

Exit

CONTROLLER RESET

If "Confirm Reset" is pressed the instrument default settings will be restored.

WARNING! All files and programs which have been stored on the internal memory will be lost.

Bibby Scientific
TC-512 TEST UNIT

SYSTEM CARD GENERATOR
CARD ERASE

Insert a SmartMedia memory card and press
required function
ALL DATA WILL BE ERASED

System Card

Erase
Card


Exit

CARD UTILITIES

This function has 2 options;

1. To erase any files or programs from an existing SmartMedia memory card.
2. To format a new card for use on the TC-512. In either case all existing data will be deleted from the card.

Selecting the "Erase Card" option will erase all the files and programs from the inserted SmartMedia memory card, leaving an empty card ready for storing new programs.
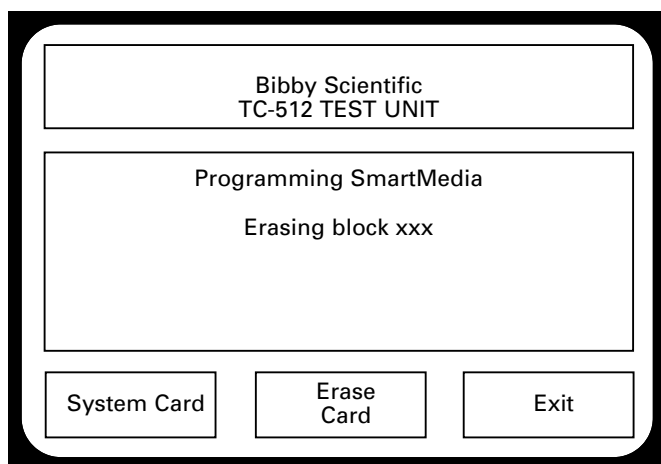


To format a new SmartMedia card;

1. Open and lift the heated lid of the thermal cyclor and then lift the silver lid on the front of the unit.
2. Remove any memory card already present in the slot by pressing the button on the left-hand side of the slot.
3. Then insert the new memory card as described on page 24.
4. Once the card is inserted correctly then press "System Card" on the Card Utilities screen.

The screen will then display an "Erasing block xxxx" message.


The erasing block message will then change to: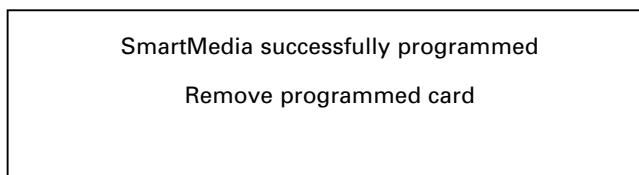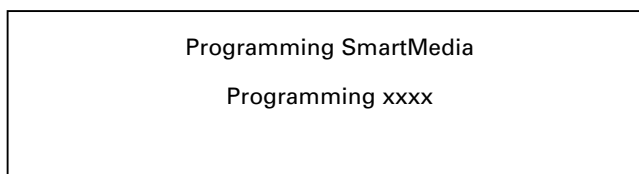

Remove the formatted card and press "Exit".

SYSTEM CONFIGURATION

This contains all the settings for programming the instrument such as units of temperature, date format, language, and date and time settings. Other important features are enabling the auto restart in case of a power failure and setting the comms address of the instrument if multiple units are linked to a computer. This part of the software also contains information relating to the cycle count for the block and the heated lid.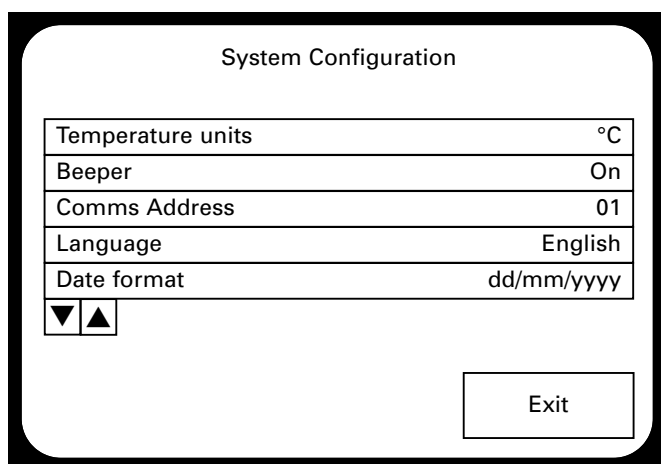

The cursor buttons can be used to scroll up and down the menu to access additional parameters.

A parameter is selected by touching the box on the screen. Then, depending on the parameter the screen will either toggle between the options or display the numerical interface over the page.

Today's Date

Enter date 05/01/2006

7	8	9
4	5	6
1	2	3
0	.	BS

←→

Cancel

OK

To enter the comms address, correct date or time type in the correct numbers and press "OK". "BS" (which is short for backspace) and the cursors can be used to delete and move across the characters.

The parameters which are editable in the systems configuration menu are:

Temperature units: This allows the user to change the units used in the thermal cycling programming of the TC-512. There are 2 options: degrees Centigrade (°C) or degrees Fahrenheit (°F).

Beeper: If required the instrument can be programmed to beep at the end of the program and also when buttons are pressed on the touch screen. This can be enabled or disabled.

Comms Address: This enables the user to specify a unique identification number for each instrument if multiple instruments are connected together and controlled via a computer. It is possible to select a communications address between 01 and 64.

Language: There are 5 languages for the user to select from: English, Français, Español, Deutsch and Italiano.

Date Format: This allows the user to change the date format from dd/mm/yyyy to mm/dd/yyyy.

Date: The date can be corrected by the user by using the numerical interface shown above.

Time: The time can be corrected by the user by using the numerical interface shown in above.

Auto restart: This allows a program to be re-started in the event of a power failure when the instrument is in the middle of running a program. The options are Enabled or Disabled.

Baud Rate: This is the speed of communication between the thermal cycler and a computer. The rate can be selected from the range of 300 to 38400 depending on individual requirements.

FREQUENTLY ASKED QUESTIONS

Q1 What is "Hot start"?

- A1** The Hot Start step is used to pause the machine at a specific temperature, typically around 70°C, after the initial template denaturation. The reason is to allow the manual addition of unmodified *Taq* DNA polymerase which may lose activity if added during the initial 5min denaturation. Heat-activated *Taq* or Hot Start enzymes do not require this step.

Q2 What is the incremental function? (See page 36).

- A2** Incremental timing and temperature are used to increase or decrease either the time or temperature incrementally over the number of cycles in a stage.

Incrementation of extension time is used with 'Long PCR' which is when large template fragments are to be amplified (e.g. 27kb lambda DNA, 40kb genomic DNA).

Decremental temperature is used for protocols such as 'Touchdown PCR' where the first cycle starts with a high annealing temperature and over the number of cycles there is a gradual decrease in the temperature. This ensures that only the specific product is amplified.

Q3 What is a gradient thermal cycler?

- A3** Gradient blocks enable a particular temperature step in a protocol to be altered so that the temperature varies across the block. By specifying a temperature and the gradient to be applied, each column will achieve a different temperature, with the set temperature in the middle block, the lowest on the left and the highest on the right. The gradient is the total difference across the block, for example if the set temperature was 60°C and the gradient 10°C then the temperatures would range from 55°C to 65°C from left to right.

Q4 Why is a temperature gradient required?

- A4** The annealing temperature of the specific primers used in DNA amplification often requires optimisation. Instead of running multiple experiments with different annealing temperatures the gradient thermal cycler allows the testing of 12 different annealing temperatures simultaneously.

If you require further technical or applicational assistance please contact Bibby Scientific Ltd at: technehelp@bibby-scientific.com.

For servicing information please contact: service@bibby-scientific.com.

We are continually striving to improve our thermal cyclers and software. If you have any comments and suggestions on how we can do things better please send them to us at: techne@bibby-scientific.com

ADDITIONAL INFORMATION

Brief fault finding notes and a list of replacement parts are given in this section.

Note: that this equipment should only be dismantled by properly trained personnel. Removing the outer cover exposes potentially lethal mains voltages. There are no user serviceable parts within this equipment.

FAULT FINDING

Should you have any problems with your TC-512 which cannot be easily remedied, you should contact your supplier and return the unit if necessary. Please include details of the fault observed and remember to return the unit in its original packing. Bibby Scientific Ltd accept no responsibility for damage to units which are not properly packed for shipping: if in doubt, contact your supplier, giving the full serial number of the unit and software version number (shown when the unit is first switched on).

FUSES

If neither the power light nor display on the front panel is lit, one of the two fuses may have blown. Check that there is no external cause, such as a faulty plug or lead. Check both fuses and replace the faulty fuse with a new one of the correct value (fuse values are given on the label next to the power inlet). Note that fuses should only be replaced by a qualified electrician.

The holder for the two fuses is built into the mains input socket. First remove the power cable and then gently prise the fuse drawer open with a flat-bladed screwdriver or similar tool.

Each fuse can be removed by using the screwdriver as a lever.

Exchange the faulty fuse in the fuse holder for a working fuse of the correct value. Finally, replace the fuse drawer in the fuse compartment and push the drawer shut.

Fuses which blow repeatedly indicate a serious fault and you should return the unit to your supplier for repair.

ACCESSORIES

Product no	Description	Quantity
FTUB05TW	0.5ml thin wall microtubes	Pack of 1000 tubes with caps
FTUB02TW	0.2ml thin wall microtubes	Pack of 1000 tubes with caps
FMITUBC	0.2ml thin wall microtubes	250 strips of 8 tubes
FMICAPC	0.2ml microtube caps	250 strips of 8 caps
F96T02	96-well flat top, non-skirted plates	Pack of 25 plates
F96PLHS	96-well half skirted plates	Pack of 25 plates
F96SEAL	96-well sealing mat (silicon)	Pack of 50 mats
FMIP384	384-well skirted plates	Pack of 50 plates
FHSEAL	Heat sealing film	Pack of 100

Product Number	Description	Quantity
FTGSM1	Smart Media card, 5V	1
FTGSM2	Smart Media card, 3V	1
FHSEALSD	Heat sealer	1
FHSPA96	Heat sealer plate adapter for 96-well plates	1
FHSPA384	Heat sealer plate adapter for 384-well plates	1
FTC51BHD	TC-512 block for 96 x 0.2ml microtubes or 96-well plate	1
FTC51BFD	TC-512 block for 96 x 0.2ml microtubes or 96-well fully skirted plate	1
FTC51B5D	TC-512 block 60 x 0.5ml microtubes	1
FTC51BID	Flat plate block for in-situ (non gradient)	1
FTCFB384	TC-512 block for 384-well fully skirted plate	1

PC ACCESSORIES

Product Number	Description	Quantity
FGEN232	Cable and software to connect 1 cyclor to PC (RS232 only)	1
FGEN485D	Power pack, 230V, 485 converter, cable and software to connect 2+ cyclers to PC	1
FGENONE	One cyclor cable extension set	1
FGENTWO	Two cyclor cable set	1
FGENFOUR	Four cyclor cable set	1
FGENTEN	Ten cyclor cable set	1

REPLACEMENT PARTS

Product Number	Description	Quantity Required
FCABRTUK	Mains cable and plug UK 230V	1
FCABRTEU	Mains cable and Schuko plug 230V	1
FCABRTUS	Mains cable and US 3-pin plug 120V	1
FCABRTUS	Mains cable and US 3-pin plug 100V	1
6500342	Fuse T6.3A, 230V units	2
6008445	Fuse T8A, 120V units	2
6500324	Fuse T10A, 100V units	2
6104172	Non-volatile Memory Module	1

**Bibby Scientific Ltd**

Beacon Road, Stone,
Staffordshire, ST15 0SA
United Kingdom
Tel: +44 (0)1785 812121
Fax: +44 (0)1785 813748
email: techne@bibby-scientific.com
www.bibby-scientific.com

Bibby Scientific France SAS

ZI du Rocher Vert - BP 79,
77793 Nemours Cedex,
France
Tel: +33 1 64 45 13 13
Fax: +33 1 64 45 13 00
e-mail: bsf@bibby-scientific.fr
www.bibby-scientific.com

Techne Inc.

3 Terri Lane, Suite 10,
Burlington, NJ, 08016, USA
Toll free: 800-225-9243
Tel: +1 609-589-3560
Fax: +1 609-589-2571
email: labproducts@techneusa.com
www.techneusa.com

Bibby Scientific Italia Srl

Via Alcide de Gasperi 56,
20077 Riozzo di Cerro al Lambro,
Milano, Italia
Tel: +39 (0)2 98230679
Fax: +39 (0)2 98230211
e-mail: marketing@bibby-scientific.it
www.bibby-scientific.it

Bibby Scientific (Asia) Ltd

Room 607, Yen Sheng Centre,
64 Hoi Yuen Road, Kwun Tong,
Kowloon, Hong Kong
Tel: +852 3583 1581
Fax: +852 3583 1580
e-mail bibby@bibby-scientificasia.com
www.bibby-scientificasia.com