



Synthesis of a primer test sequence using StratoSpheres™ DNA

Andrew Coffey
Varian, Inc.

Introduction

The synthesis of primer sequences for PCR is perhaps the most commonly used application for oligonucleotide synthesis. 5' ATA CCG ATT AAG CGA AGT TT 3' is a commonly used 20-mer test sequence for evaluating synthesis supports or automated instrumentation. The StratoSpheres DNA dT support matches the performance of the leading commercial macroporous PS/DVB.

Oligonucleotide Synthesis

The oligonucleotides were prepared on 0.2 μ M scale using an Applied Biosystems 392 DNA/RNA Synthesizer and standard chemistry. Following synthesis, the oligonucleotides were

cleaved using ammonium hydroxide. Deprotection of the side chain protecting groups (A, C and G) was accomplished by heating to 55 °C overnight as required. The "DMT on" oligonucleotides were diluted with water in order to give an on-scale response during HPLC analysis.

HPLC Analysis

Column: Pellicular SAX, 4 x 250 mm


Eluent: A = 25 mM Tris-HCl, 0.5% ACN, pH 8.0; B = 25 mM Tris-HCl, 0.8 M Ammonium Chloride, 0.5% ACN, pH 8.0

Gradient: 0 - 100% B in 26 min

Flow Rate: 1.5 mL/min

Temp: 60 °C

Detector: UV 260 nm

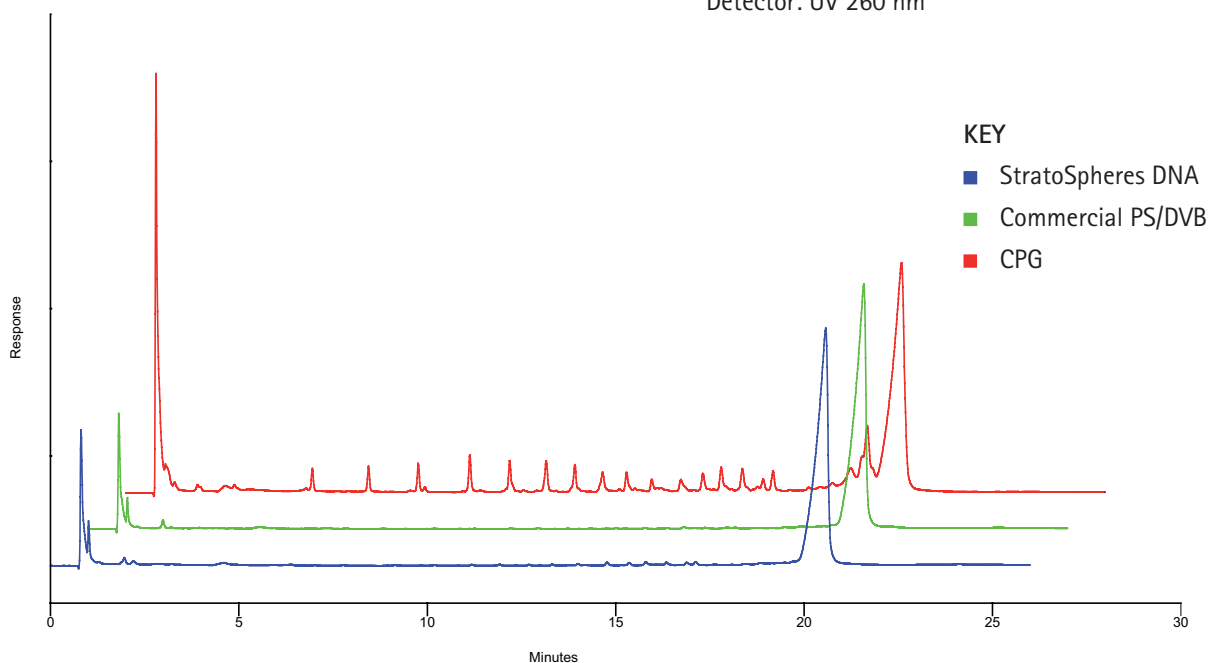

J. Shanagar (2005) Purification of a synthetic oligonucleotide by anion exchange chromatography: method optimisation and scale-up. *J. Biochem. Biophys. Methods*, 64, 216-225.

These data represent typical results.

For further information, contact your local Varian Sales Office.