



# Agilent 1260 Infinity Standard Degasser

User Manual



**Agilent Technologies**

# Notices

© Agilent Technologies, Inc. 2006-2008, 2010

No part of this manual may be reproduced in any form or by any means (including electronic storage and retrieval or translation into a foreign language) without prior agreement and written consent from Agilent Technologies, Inc. as governed by United States and international copyright laws.

## Manual Part Number

G1322-90012

## Edition

06/10

Printed in Germany

Agilent Technologies  
Hewlett-Packard-Strasse 8  
76337 Waldbronn

**This product may be used as a component of an in vitro diagnostic system if the system is registered with the appropriate authorities and complies with the relevant regulations. Otherwise, it is intended only for general laboratory use.**

## Warranty

**The material contained in this document is provided “as is,” and is subject to being changed, without notice, in future editions. Further, to the maximum extent permitted by applicable law, Agilent disclaims all warranties, either express or implied, with regard to this manual and any information contained herein, including but not limited to the implied warranties of merchantability and fitness for a particular purpose. Agilent shall not be liable for errors or for incidental or consequential damages in connection with the furnishing, use, or performance of this document or of any information contained herein. Should Agilent and the user have a separate written agreement with warranty terms covering the material in this document that conflict with these terms, the warranty terms in the separate agreement shall control.**

## Technology Licenses

The hardware and/or software described in this document are furnished under a license and may be used or copied only in accordance with the terms of such license.

## Restricted Rights Legend

If software is for use in the performance of a U.S. Government prime contract or subcontract, Software is delivered and licensed as “Commercial computer software” as defined in DFAR 252.227-7014 (June 1995), or as a “commercial item” as defined in FAR 2.101(a) or as “Restricted computer software” as defined in FAR 52.227-19 (June 1987) or any equivalent agency regulation or contract clause. Use, duplication or disclosure of Software is subject to Agilent Technologies’ standard commercial license terms, and non-DOD Departments and Agencies of the U.S. Government will

receive no greater than Restricted Rights as defined in FAR 52.227-19(c)(1-2) (June 1987). U.S. Government users will receive no greater than Limited Rights as defined in FAR 52.227-14 (June 1987) or DFAR 252.227-7015 (b)(2) (November 1995), as applicable in any technical data.

## Safety Notices

### CAUTION

A **CAUTION** notice denotes a hazard. It calls attention to an operating procedure, practice, or the like that, if not correctly performed or adhered to, could result in damage to the product or loss of important data. Do not proceed beyond a **CAUTION** notice until the indicated conditions are fully understood and met.

### WARNING

A **WARNING** notice denotes a hazard. It calls attention to an operating procedure, practice, or the like that, if not correctly performed or adhered to, could result in personal injury or death. Do not proceed beyond a **WARNING** notice until the indicated conditions are fully understood and met.

# Contents

<b>1</b>	<b>Introduction</b>	<b>5</b>
	Introduction to the Vacuum Degasser	6
	How the Electronics Work	7
<b>2</b>	<b>Site Requirements and Specifications</b>	<b>11</b>
	Site Requirements	12
	Physical Specifications	15
	Performance Specifications	16
<b>3</b>	<b>Installing the Standard Degasser</b>	<b>17</b>
	Unpacking the Standard Degasser	18
	Optimizing the Stack Configuration	20
	Installing the Standard Degasser	23
	Flow Connections to the Standard Degasser	26
	Operational Hints for the Standard Degasser	30
	Transporting the Standard Degasser	33
<b>4</b>	<b>Using the Vacuum Degasser</b>	<b>35</b>
	When to Use a Vacuum Degasser?	36
	Solvent Information	37
	Prevent Blocking of Solvent Filters	38
	Operation Modes of the Vacuum Degasser	40
<b>5</b>	<b>Troubleshooting and Diagnostics</b>	<b>43</b>
	Overview of the Degasser's Indicators	44
	Status Indicators	45
<b>6</b>	<b>Maintenance</b>	<b>47</b>
	Warnings and Cautions	48
	Introduction to Maintenance	50

<b>7</b>	<b>Parts and Materials for Maintenance</b>	<b>59</b>
	Cover Parts	60
	Power and Status Light Pipes	61
	Accessory Kit Contents	62
<b>8</b>	<b>Identifying Cables</b>	<b>63</b>
	Cable Overview	64
<b>9</b>	<b>Appendix</b>	<b>71</b>
	General Safety Information	72
	The Waste Electrical and Electronic Equipment (WEEE) Directive (2002/96/EC)	75
	Radio Interference	76
	Sound Emission	77
	Agilent Technologies on Internet	78



# 1

## Introduction

Introduction to the Vacuum Degasser 6


How the Electronics Work 7



## Introduction to the Vacuum Degasser

The Agilent 1200 Series vacuum degasser, model G1322A, comprises a 4-channel vacuum container, including 4 tubular plastic membranes, and a vacuum pump. When the vacuum degasser is switched on, the control circuit turns on the vacuum pump which generates a partial vacuum in the vacuum container. The pressure is measured by a pressure sensor. The vacuum degasser maintains the partial vacuum by turning on and off the vacuum pump depending on the signal from the pressure sensor.


The LC pump draws the solvents from their bottles through the special tubular plastic membranes of the vacuum container. As the solvents pass through the vacuum tubes any dissolved gas in the solvents permeates through the membranes into the vacuum container. The solvents will be almost completely degassed when leaving the outlets of the vacuum degasser.

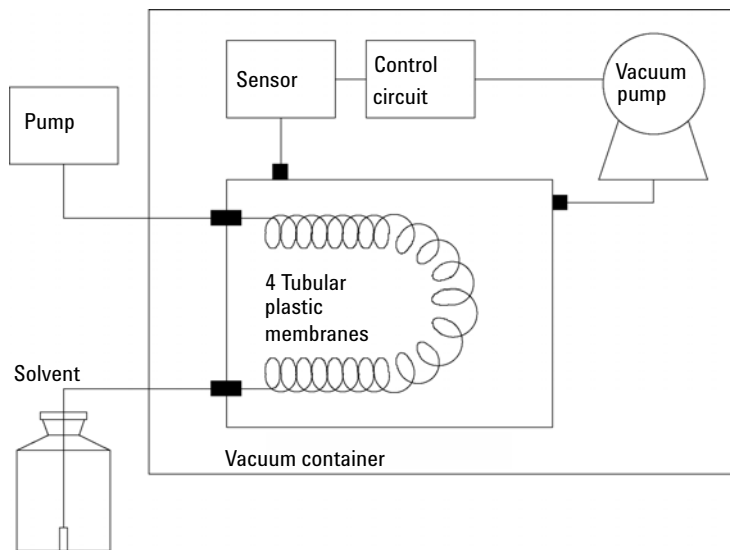

**Figure 1** Overview (only one of the four solvent channels is shown)

## How the Electronics Work

The vacuum degasser has two different normal operation modes and a continuous mode. In operation mode 1 the vacuum degasser works around a defined set point (115 Torr). Due to environmental conditions it is possible that the vacuum degasser cannot reach the pre-defined set point. Under this condition the operation mode 2 becomes active and the vacuum pump is activated in defined time intervals (vacuum level 115 to 190 Torr). In case of a malfunction of the vacuum degasser (vacuum level above 190 Torr) the instrument is turned into the error mode.

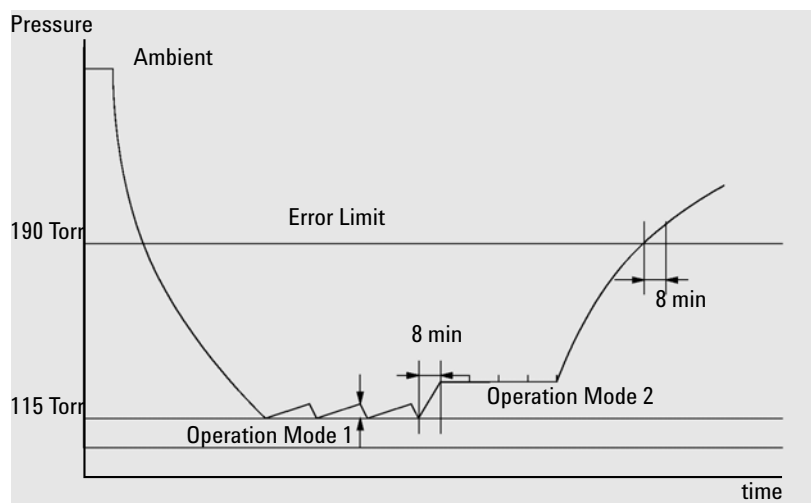

**Figure 2** Operation Modes of the Vacuum Degasser

The main function of the vacuum degasser control assembly is to control the vacuum pump and to check the vacuum in the vacuum container.

The power section of the control assembly comprises a switching power supply that generates + 24 V from line voltage. The + 24 V is used to drive the vacuum pump and the solenoid valve. The electronic control circuit uses + 12 V which is generated from the + 24 V.

The pressure sensor is connected to the vacuum chamber and checks for the correct vacuum in the system.

The amplifier and comparator circuit determines the working range of the vacuum that has to be built up. When the vacuum degasser is turned on and the vacuum in the chamber is not within working range (above error limit of 190 Torr), the amplifier and comparator circuit sends a signal to the vacuum pump driver and the timers of the vacuum pump (timer 1) and the solenoid valve (timer 2).

The vacuum pump is turned on immediately while the solenoid valve closing is delayed by 15 seconds. This time delay allows the vacuum pump to start without load before it is connected to the vacuum chamber. The status indicator turns yellow when the pump is activated. The status lamp is turned off as soon as the vacuum is below the error limit.

When the vacuum in the chamber reaches its operation mode 1 (approximately 115 Torr) the amplifier and comparator circuit turns off the solenoid valve and the vacuum pump is turned off by a timer with a delay of 15 seconds.

As soon as the pressure sensor detects that the limit of the operation mode 1 has been exceeded (e.g. when dissolved gas from the solvent diffused into the vacuum chamber) the vacuum pump is started again as described before.

The pressure signal is available at the auxiliary output. It allows to monitor the vacuum system. The upper limit of operation mode 1 is 600 mV. Values below 600 mV on the pressure output indicate sufficient vacuum in the chamber. If the 600 mV are exceeded the vacuum pump will be started to keep the vacuum within its working limit.

The amplifier and comparator circuit also activates the timer 3 when the vacuum in the vacuum chambers is below operation mode 1. The timer 3 is reset when operation mode 1 is reached within a maximum of 8 minutes. If operation mode 1 is not reached and the time (8 minutes) of timer 3 elapsed, the timer mode (operation mode 2) is activated. In this mode the vacuum pump is automatically turned on every 2 minutes for a time frame of 30 seconds. Timers 1 and 2 are activated as described earlier.

The error monitor continuously checks the error limit of the degasser (190 Torr). If the error limit is exceeded (for example, leak in chamber), the error timer is activated and the yellow status indicator lamp is turned on. The vacuum pump is turned on continuously. If the vacuum pump cannot reach either of its two operation modes within 8 minutes (error timer limit) the error driver is activated. The error driver will deactivate the vacuum pump and solenoid valve. The status lamp turns red and the error output on the remote connector will be activated.




The error output at the remote connector provides a *contact closure* (potential free open collector maximum load 35 V DC/50 mA) as long as the error condition is active. The error condition is set (closed) when the status light shows the error condition (red).

The continuous mode overwrites all other operation modes of the degasser. When activated (switch SW1 on the electronic board or via the auxiliary cable) the vacuum pump is forced into continuous mode and is turned on as long as the degasser is switched ON.

# 1 Introduction

## How the Electronics Work

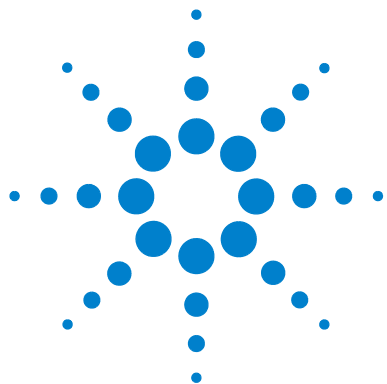

## 2

# Site Requirements and Specifications

Site Requirements 12

Physical Specifications 15

Performance Specifications 16

This chapter provides information on environmental requirements, physical and performance specifications.



## Site Requirements

A suitable environment is important to ensure optimal performance of the instrument.

### Power Considerations

The module power supply has wide ranging capability. It accepts any line voltage in the range described in [Table 1](#) on page 15. Consequently there is no voltage selector in the rear of the module. There are also no externally accessible fuses, because automatic electronic fuses are implemented in the power supply.

#### WARNING

**Hazard of electrical shock or damage of your instrumentation can result, if the devices are connected to a line voltage higher than specified.**

→ Connect your instrument to the specified line voltage only.

---

#### WARNING

**Module is partially energized when switched off, as long as the power cord is plugged in.**

**Repair work at the module can lead to personal injuries, e.g. electrical shock, when the cover is opened and the module is connected to power.**

→ Always unplug the power cable before opening the cover.

→ Do not connect the power cable to the instrument while the covers are removed.

---

#### CAUTION

Unaccessible power plug.

In case of emergency it must be possible to disconnect the instrument from the power line at any time.

→ Make sure the power connector of the instrument can be easily reached and unplugged.

→ Provide sufficient space behind the power socket of the instrument to unplug the cable.

---

## Power Cords

Different power cords are offered as options with the module. The female end of all power cords is identical. It plugs into the power-input socket at the rear. The male end of each power cord is different and designed to match the wall socket of a particular country or region.

### **WARNING**

#### **Absence of ground connection or use of unspecified power cord**

**The absence of ground connection or the use of unspecified power cord can lead to electric shock or short circuit.**

- Never operate your instrumentation from a power outlet that has no ground connection.
  - Never use a power cord other than the Agilent Technologies power cord designed for your region.
- 

### **WARNING**

#### **Use of unsupplied cables**

**Using cables not supplied by Agilent Technologies can lead to damage of the electronic components or personal injury.**

- Never use cables other than the ones supplied by Agilent Technologies to ensure proper functionality and compliance with safety or EMC regulations.
- 

### **WARNING**

#### **Unintended use of supplied power cords**

**Using power cords for unintended purposes can lead to personal injury or damage of electronic equipment.**

- Never use the power cords that Agilent Technologies supplies with this instrument for any other equipment.
-

## Bench Space

The module dimensions and weight (see [Table 1](#) on page 15) allow you to place the module on almost any desk or laboratory bench. It needs an additional 2.5 cm (1.0 inches) of space on either side and approximately 8 cm (3.1 inches) in the rear for air circulation and electric connections.

If the bench should carry an Agilent system, make sure that the bench is designed to bear the weight of all modules.

The module should be operated in a horizontal position.

## Condensation

### CAUTION

Condensation within the module

Condensation will damage the system electronics.

- Do not store, ship or use your module under conditions where temperature fluctuations could cause condensation within the module.
  - If your module was shipped in cold weather, leave it in its box and allow it to warm slowly to room temperature to avoid condensation.
-

# Physical Specifications

**Table 1** Physical Specifications

Type	Specification	Comments
Weight	7 kg (16 lbs)	
Dimensions (height × width × depth)	80 × 345 × 435 mm (3.1 × 13.5 × 17 inches)	
Line voltage	100 – 120 / 220 – 240 VAC	Wide-ranging capability
Line frequency	50 or 60 Hz, ± 5 %	
Power consumption	250 VA / 210 W / 717 BTU	Maximum
Ambient operating temperature	0–55 °C (32–131 °F) 1	
Ambient non-operating temperature	-40 – 70 °C (-40 – 158 °F)	
Humidity	< 95 %, at 25 – 40 °C (77 – 104 °F)	Non-condensing
Operating Altitude	Up to 2000 m (6562 ft)	
Non-operating altitude	Up to 4600 m (15091 ft)	For storing the module
Safety standards: IEC, CSA, UL	Installation Category II, Pollution Degree 2	For indoor use only. Research Use Only. Not for use in Diagnostic Procedures.

<sup>1</sup> This temperature range represents the technical specifications for this instrument. The mentioned temperatures may not be suitable for all applications and all types of solvents.

## NOTE


The Agilent 1260 Infinity Standard Degasser has been tested for evaporation of solvents into the atmosphere by an independent institute with approved methods. The tests were performed with methanol (BIA Nr. 7810) and acetonitrile (NIOSH, Nr. 1606). Evaporation of these solvents into the atmosphere when operating the degasser was below the limits of detection.

## Performance Specifications

**Table 2**    Performance Specifications Agilent 1260 Infinity Standard Degasser

Type	Specification
Maximum flow rate	10 mL/min per channel
Number of channels	4
Internal volume per channel	Typically 12 mL per channel
Materials in contact with solvent	PTFE, PEEK
pH range	1 – 14
Analog output (AUX)	For pressure monitoring, range 0 – 3 V



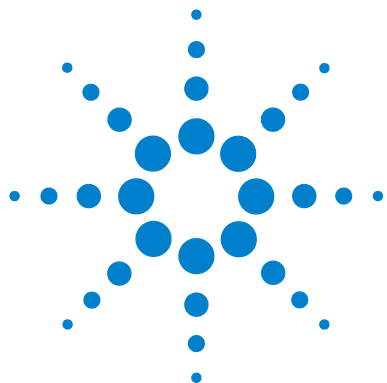

## 3 Installing the Standard Degasser

Unpacking the Standard Degasser	18
Damaged Packaging	18
Delivery Checklist	19
Accessory Kit Contents	19
Optimizing the Stack Configuration	20
Installing the Standard Degasser	23
Flow Connections to the Standard Degasser	26
Operational Hints for the Standard Degasser	30
Priming the Degasser	30
Priming with a Syringe	30
Priming with the Pump	32
Transporting the Standard Degasser	33

This chapter provides information on unpacking, checking on completeness, stack considerations and installation of the module.



## Unpacking the Standard Degasser

### Damaged Packaging

If the delivery packaging shows signs of external damage, please call your Agilent Technologies sales and service office immediately. Inform your service representative that the instrument may have been damaged during shipment.

#### CAUTION

"Defective on arrival" problems

If there are signs of damage, please do not attempt to install the module. Inspection by Agilent is required to evaluate if the instrument is in good condition or damaged.

- Notify your Agilent sales and service office about the damage.
  - An Agilent service representative will inspect the instrument at your site and initiate appropriate actions.
-

## Delivery Checklist

Ensure all parts and materials have been delivered with the degasser. The delivery checklist is shown below. To aid in parts identification, please see [“Parts and Materials for Maintenance”](#) on page 59. Please report missing or damaged parts to your local Agilent Technologies Sales and Service Office.

**Table 3** Standard Degasser Delivery Checklist

Description	Quantity
Standard degasser	1
Power cable	1
Remote cable	1
Auxiliary cable	As ordered
User manual	1
Accessory kit ( <a href="#">“Accessory Kit Contents”</a> on page 19)	1


## Accessory Kit Contents

Accessory Kit (p/n G1322-68705)

p/n	Description
5062-8534	Syringe Reorder number (pack of 10)
9301-1337	Syringe adapter
0100-1710	Mounting Tool for Tubing Connections
5062-2463	Corrugated waste tube (reorder pack), 5 m
G1322-67300	Kit of 4 solvent tubes for connection degasser to MCGV (Quaternary Pump) including labels

## Optimizing the Stack Configuration


If your standard degasser is part of a system, you can ensure optimum performance by limiting the configuration of the system stack to the following configuration. This configuration optimizes the system flow path, ensuring minimum delay volume (from point of solvent mixing to head of column) and dead volume (from point of injection to detector outlet).

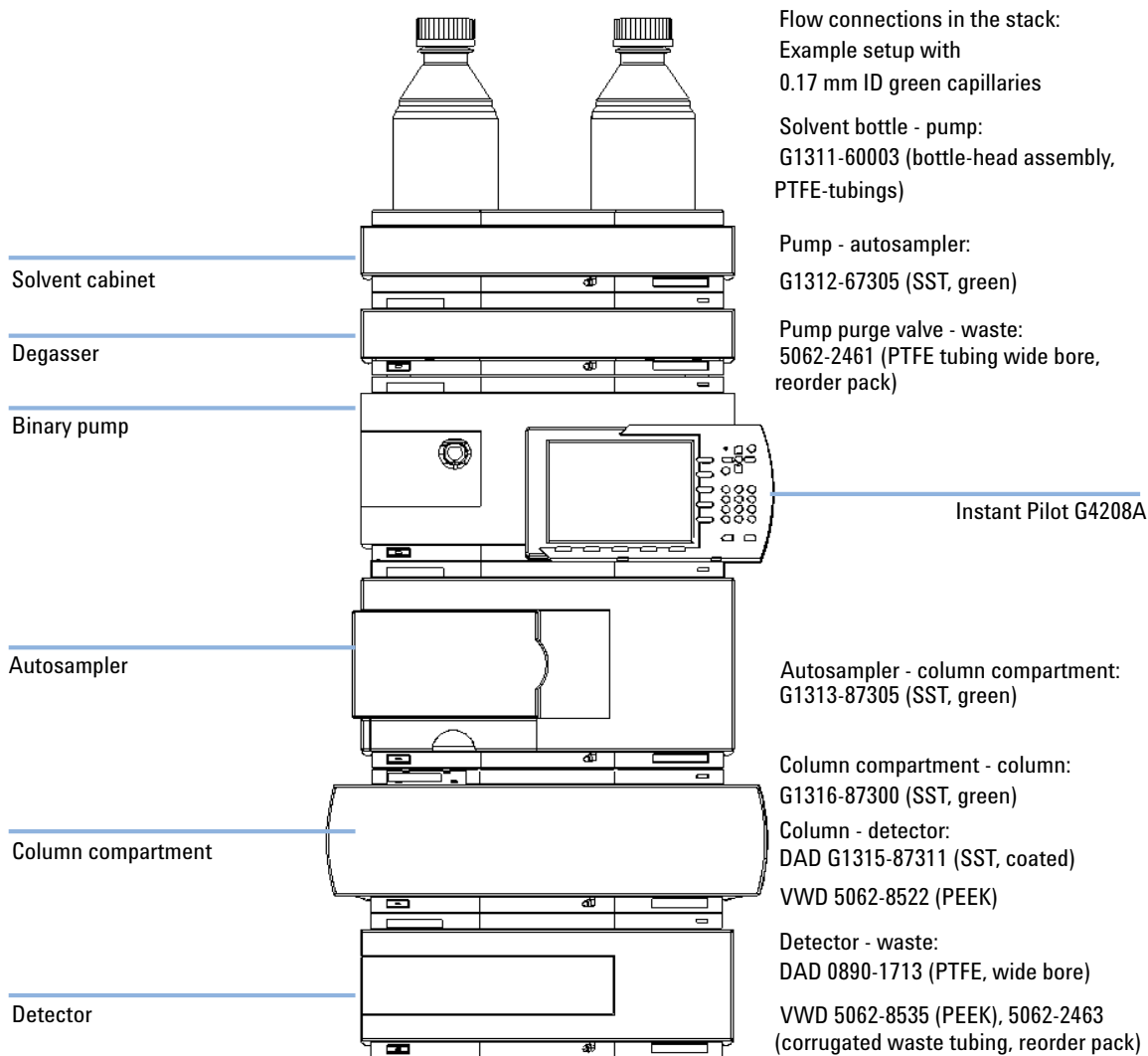

**Figure 3** Recommended Stack Configuration (Front View)

### 3 Installing the Standard Degasser

#### Optimizing the Stack Configuration

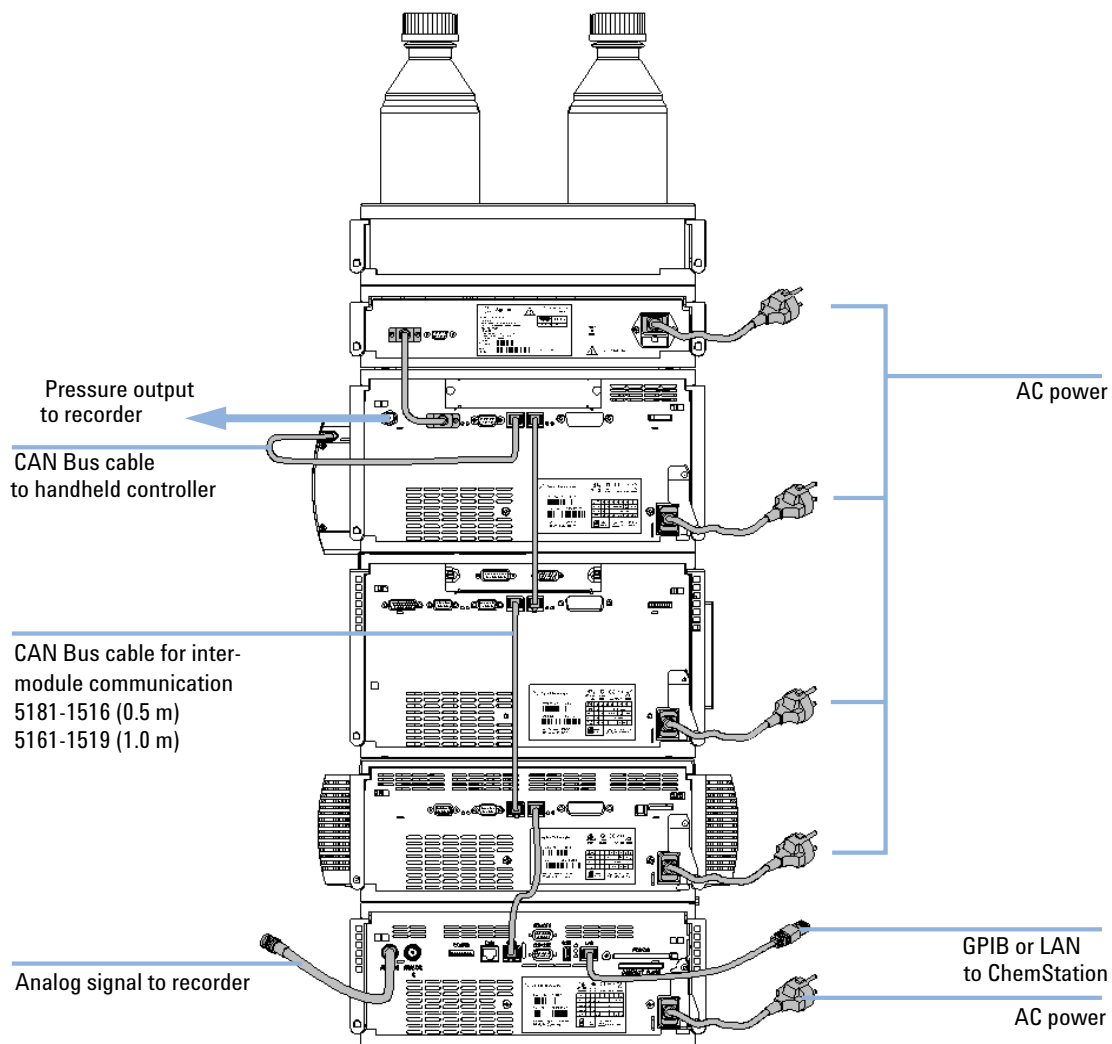

**Figure 4** Recommended Stack Configuration (Rear View)

# Installing the Standard Degasser

Parts required	#	Description
	1	Standard degasser
	1	Power cord
	1	Interface cable as ordered

See [“Cable Overview”](#) on page 64.

Preparations	
	<ul style="list-style-type: none"> <li>• Locate bench space</li> <li>• Provide power connections</li> <li>• Unpack the module</li> </ul>

## CAUTION

"Defective on arrival" problems

If there are signs of damage, please do not attempt to install the module. Inspection by Agilent is required to evaluate if the instrument is in good condition or damaged.

- Notify your Agilent sales and service office about the damage.
- An Agilent service representative will inspect the instrument at your site and initiate appropriate actions.

- 1 Place the module on the bench in a horizontal position.
- 2 Ensure the power switch on the front of the module is OFF (switch stands out).
- 3 Connect the power cable to the power connector at the rear of the module.

### 3 Installing the Standard Degasser


#### Installing the Standard Degasser

- 4 Connect the interface cable to the degasser. The interface cable (remote cable) is a one way connection to send a not-ready signal from the degasser to the other modules to shut down the whole system after an error condition of the degasser.

#### NOTE

In an Agilent 1260 Infinity system, the individual modules are connected through a CAN cable. The Agilent 1260 Infinity Standard Degasser is an exception. The standard degasser can be connected via the APG remote connector to the other modules of the stack. The AUX output allows to monitor the vacuum pressure in the degasser chamber. A G4208A Instant Pilot can be connected to the CAN bus at any of the modules in the system except for the degasser. The chromatography data system can be connected to the system through a LAN cable at any of the modules (except for the degasser), preferably at the detector. For more information about connecting the Instant Pilot or a data system refer to the respective user manual.


Power socket

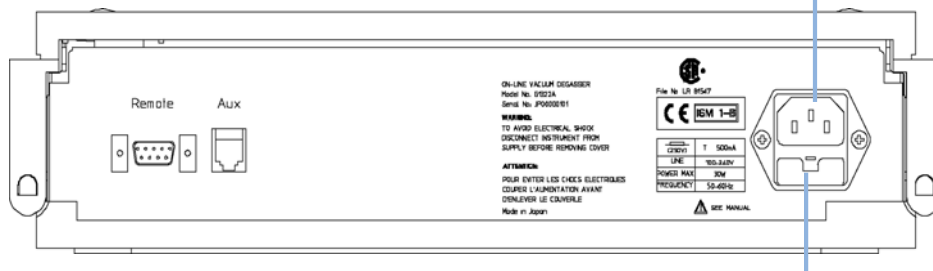

Fuse holder

**Figure 5** Rear of the Standard Degasser



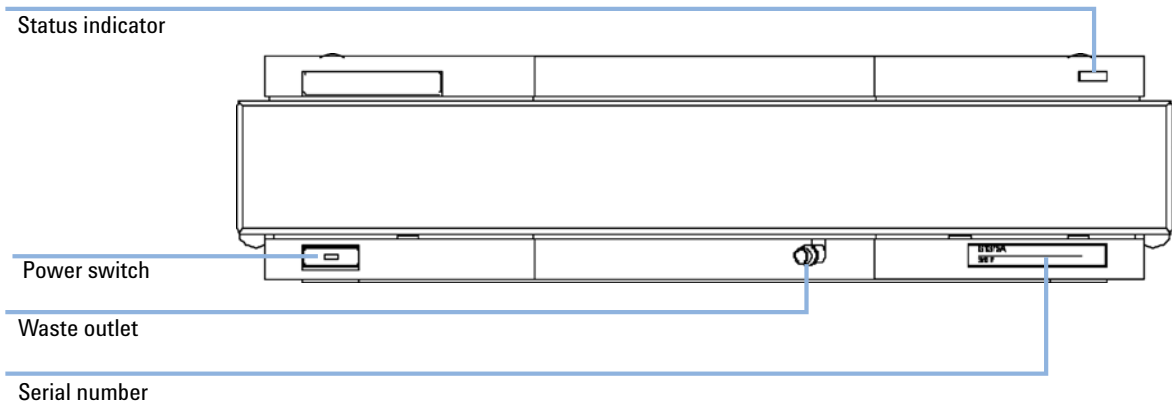

**Figure 6** Front of the Standard Degasser

- 5 Press in the power switch to turn on the vacuum degasser.

**NOTE**

The power switch stays pressed in and a green indicator lamp in the power switch is ON when the vacuum degasser is turned ON. When the line power switch stands out and the green light is OFF, the vacuum degasser is turned OFF.

## Flow Connections to the Standard Degasser

Parts required	#	Description
	1	Solvent cabinet including solvent bottles (filled with solvent) and bottle head assemblies
	1	Standard degasser
	1	Solvent outlet tubes
	1	Syringe with adapter

**Preparations**

- Install the degasser


### WARNING

**Toxic, flammable and hazardous solvents, samples and reagents**

**The handling of solvents, samples and reagents can hold health and safety risks.**


- When working with these substances observe appropriate safety procedures (for example by wearing goggles, safety gloves and protective clothing) as described in the material handling and safety data sheet supplied by the vendor and follow good laboratory practice.
  - The amount of substances should be reduced to the minimal volume required for the analysis.
  - Do not operate the instrument in an explosive atmosphere.
-

- 1 Place solvent cabinet with the bottle(s) on top of the degasser.
- 2 Remove the front cover by pressing the snap fasteners on both sides.

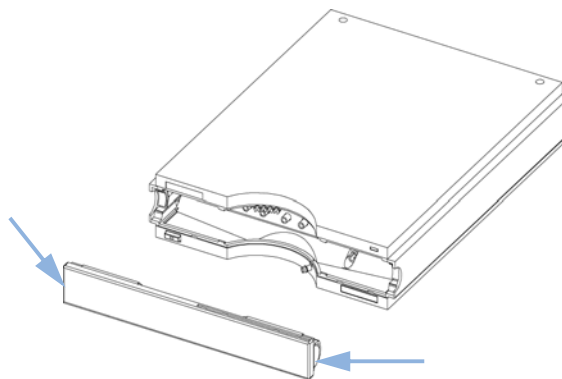

**Figure 7** Removing the Front Cover

- 3 If the degasser is not used with an Agilent 1260 Infinity pump, connect the waste tube from the accessory kit to the waste outlet and place into your waste system.
- 4 Put the bottle head assemblies into solvent bottles containing your mobile phase.
- 5 Connect the solvent tubes from the bottle head assemblies to the inlet connectors A to D (typically the left connection of the channel) of the degasser. Use the mounting tool shown in [Figure 8](#) on page 27 to fix the tube screw. Fix the tubes in the clips of the degasser.

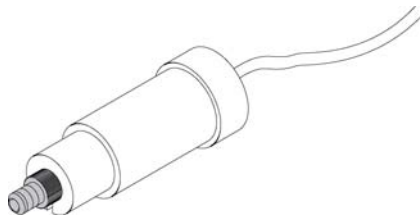

**Figure 8** Mounting Tool

### 3 Installing the Standard Degasser


#### Flow Connections to the Standard Degasser

- 6 Connect the outlet tubes to the output ports (typically right connection of the channel) of the degasser.
- 7 Prime the degasser before first use (see [“Priming the Degasser”](#) on page 30).

#### NOTE

Atmospheric gases can diffuse through the tubing and dissolve in the mobile phase solvents. For best chromatographic results, keep the length of tubing between the degasser and the pump as short as possible.

---

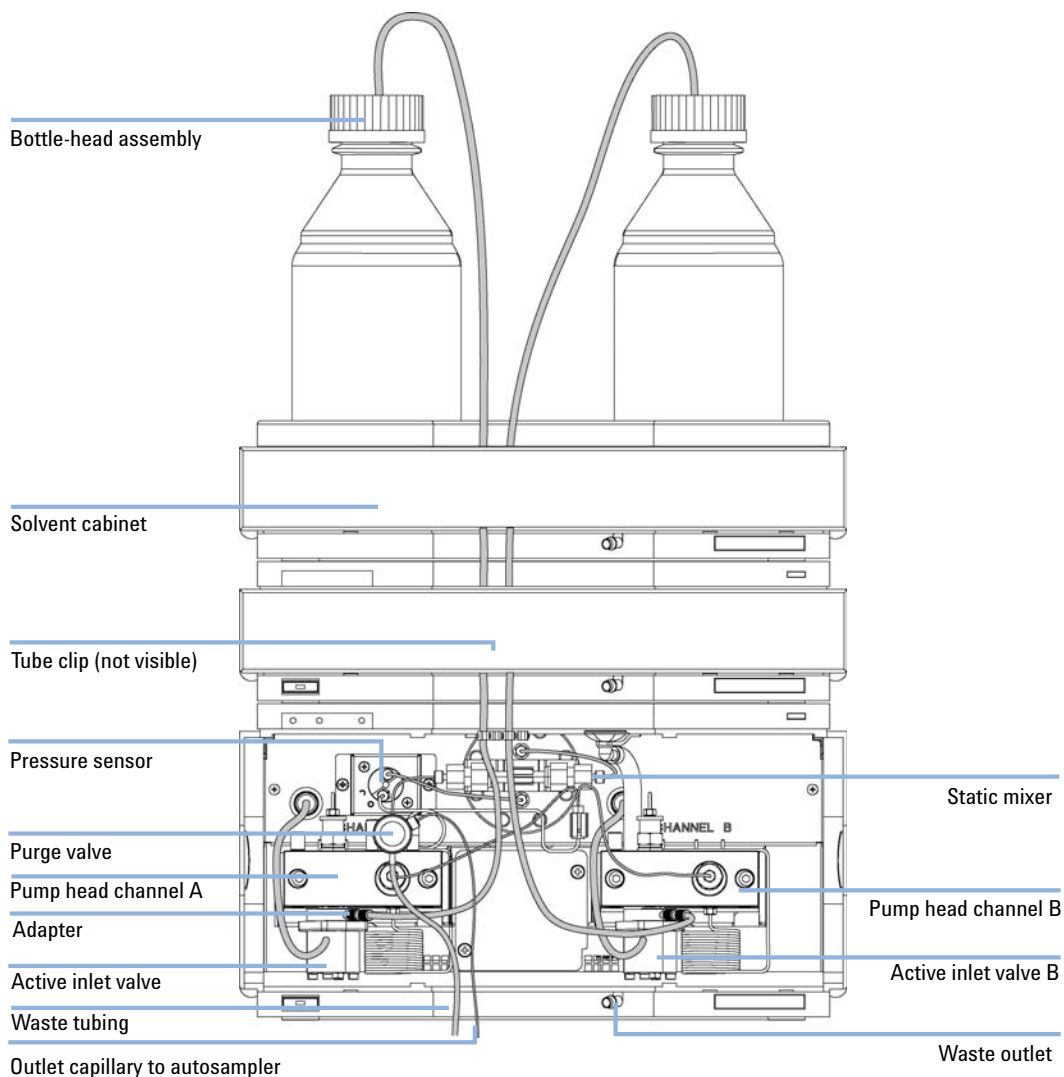

**Figure 9** Flow Connections to the Standard Vacuum Degasser

## Operational Hints for the Standard Degasser

### Priming the Degasser

The standard degasser can be primed either by drawing solvent through the degasser with a syringe or by pumping with the connected pump.

Priming the standard degasser with a syringe is recommended, when:

- standard degasser is used for the first time, or vacuum tubes are empty, or
- changing to solvents that are immiscible with the solvent currently in the vacuum tubes.

Priming the standard degasser by using the pump at high flow rate (3 – 5 mL/min) is recommended, when:

- pumping system was turned off for a length of time (for example, overnight) and if volatile solvent mixtures are used, or
- if solvents have been changed.

### Priming with a Syringe

Before using a new degasser or new tubings for the first time:

#### **WARNING**

**Toxic, flammable and hazardous solvents, samples and reagents**

**The handling of solvents, samples and reagents can hold health and safety risks.**

- When working with these substances observe appropriate safety procedures (for example by wearing goggles, safety gloves and protective clothing) as described in the material handling and safety data sheet supplied by the vendor and follow good laboratory practice.
  - The amount of substances should be reduced to the minimal volume required for the analysis.
  - Do not operate the instrument in an explosive atmosphere.
-

- 1** Prime all tubings with at least 30 ml of iso-propanol no matter whether the channels will be used with organic mobile phase or with water.  
If you are changing to a solvent that is immiscible with the solvent currently in the tubing continue as follows:
- 2** Replace the current solvent with iso-propanol, if current solvent is organic or with water, if current solvent is an inorganic buffer or contains salt.
- 3** Disconnect solvent outlet tube of the channel that is supposed to be primed from your pump.
- 4** Connect syringe adapter to solvent outlet tube.
- 5** Push syringe adapter onto syringe.
- 6** Pull syringe plunger to draw at least 30 ml of solvent through degasser and tubing.
- 7** Replace the priming solvent with the new solvent of your choice.
- 8** Pull syringe plunger to draw at least 30 ml of solvent through degasser and tubing.
- 9** Disconnect syringe adapter from solvent tube.
- 10** Connect solvent tube to your pump.
- 11** Repeat step 3 on page 31 to step 10 on page 31 for the other solvent channels.

**NOTE**

When priming the standard degasser with a syringe the solvent is drawn through the degasser tubes very quickly. The solvent at the degasser outlet will therefore not be fully degassed. Pump for approximately 10 min with your selected flow rate before starting any application. This will allow the standard degasser to properly degas the solvent in the degasser tubes.

**NOTE**

The pump should never be used for priming empty tubings (never let the pump run dry). Use the syringe to draw enough solvent for completely filling the tubings to the pump inlet before continuing to prime with the pump.

## **Priming with the Pump**

When the pumping system has been turned off for a certain time (for example, overnight) oxygen will rediffuse into the solvent channels between the standard degasser and the pump. Solvents containing volatile ingredients will slightly lose these, if left in the degasser without flow for a prolonged period of time. Therefore priming of the standard degasser and the pumping system is required before starting an application.

- 1**   Open the purge valve of your pump and set flow rate to 5 ml/min.
- 2**   Flush the standard degasser and all tubes with at least 30 mL of solvent.
- 3**   Set flow to required value of your application and close the purge valve.
- 4**   Pump for approximately 10 minutes before starting your application.
- 5**   Repeat step 1 on page 32 through step 4 on page 32 with other channels, where needed.



## Transporting the Standard Degasser

### WARNING

#### Solvents leaking out

**Solvents remaining in the solvent channels may leak out during transport. This can possibly cause personal damage.**


→ Drain any remaining solvents from the degassing channels before transporting the degasser.

---

- 1 Disconnect the solvent tubes from solvent inlets from front panel.
- 2 Disconnect one of the solvent tubes from your pump.
- 3 Connect syringe adapter to solvent tube of this solvent channel.
- 4 Push syringe adapter onto syringe.
- 5 Pull syringe plunger to draw solvent out of degasser and tubing. Continue to draw solvent into syringe until the solvent channel is completely empty.
- 6 Repeat step 2 on page 33 through step 5 on page 33 for the remaining solvent channels.

### **3 Installing the Standard Degasser**

#### **Transporting the Standard Degasser**

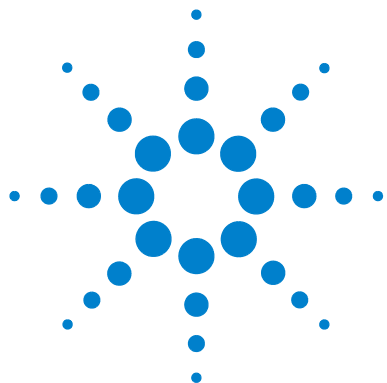

## 4 Using the Vacuum Degasser

When to Use a Vacuum Degasser?	36
Solvent Information	37
Prevent Blocking of Solvent Filters	38
Operation Modes of the Vacuum Degasser	40
Normal Operation Mode 1	41
Normal Operation Mode 2 (Timing Mode)	41
Continuous Mode	42
Error Mode	42



## When to Use a Vacuum Degasser?

### WARNING

#### *Unspecified Conditions*

**Operating the instrumentation under conditions other than their intended use might result in a potential safety hazard or might damage the instrumentation.**

→ Never operate your instrumentation under conditions other than specified by the vendor.

---

Pumps that mix the solvent on the low pressure side need degassing and must be equipped with a degasser or alternative degassing systems (for example, helium). Isocratic pumps and high-pressure mixing pumps do not always require degassing. However for the following conditions the standard degasser is also recommended for an isocratic or a high pressure mixing pump:

- if your detector is used with maximum sensitivity in the low UV wavelength range,
- if your application requires optimum injection precision, or
- if your application requires highest retention time reproducibility (mandatory at flow rates below 0.5 mL/min),
- if your sample or detection is sensitive to dissolved oxygen in the mobile phase (degradation).


Generally a degasser should be used when negative effects due to dissolved gas in the mobile phase exceed the limits that are acceptable for the user. Negative effects that can be caused by dissolved gas are:

- Unstable flow due to unstable pumping conditions. This may result in a high ripple (unstable pressure at constant flow and with constant mobile phase composition) or high standard deviations of peak retention times and peak areas especially at low flow rates.
- Baseline noise on detectors that are sensitive to changes in the refractive index (e.g. RI detector or UV detector in the low UV range, both at maximum sensitivity).
- Sample degradation.
- Fluorescence quenching due to dissolved oxygen.
- Baseline drift in electrochemical detectors due to dissolved oxygen especially in reduction mode.

## Solvent Information

Always filter solvents through 0.4 µm filters, small particles can permanently block the capillaries. Avoid the use of the following steel-corrosive solvents:

- Solutions of alkali halides and their respective acids (for example, lithium iodide, potassium chloride, and so on).
- High concentrations of inorganic acids like sulfuric acid, especially at higher temperatures (replace, if your chromatography method allows, by phosphoric acid or phosphate buffer which are less corrosive against stainless steel).
- Halogenated solvents or mixtures which form radicals and/or acids, for example:

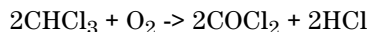

This reaction, in which stainless steel probably acts as a catalyst, occurs quickly with dried chloroform if the drying process removes the stabilizing alcohol.

- Chromatographic grade ethers, which can contain peroxides (for example, THF, dioxane, di-isopropylether). Such ethers should be filtered through dry aluminium oxide which adsorbs the peroxides.
- Mixtures of carbon tetrachloride with 2-propanol or THF dissolve stainless steel.

## Prevent Blocking of Solvent Filters

Contaminated solvents or algae growth in the solvent bottle will reduce the lifetime of the solvent filter and will influence the performance of the pump. This is especially true for aqueous solvents or phosphate buffers (pH 4 – 7). The following suggestions will prolong lifetime of the solvent filter and will maintain the performance of the pump.

- Use sterile, if possible amber solvent bottles to slow down algae growth.
- Filter solvents through filters or membranes that remove algae.
- Exchange solvents every two days or refilter.
- If the application permits add 0.0001 – 0.001 M sodium azide to the solvent.
- Place a layer of argon on top of your solvent.
- Avoid exposure of the solvent bottles to direct sunlight.

### Prevent Blocking of Solvent Filters

The solvent filters are on the low-pressure side of the pumping system. A blocked filter therefore does not affect the pressure readings of the pump. The pressure readings cannot be used to identify blocked filters. If the solvent cabinet is placed on top of the standard degasser the filter condition can be checked in the following way:

Remove the tubing at the inlet port of the standard degasser. If the filter is in good condition the solvent will freely drip out of the solvent tube (due to hydrostatic pressure). If the solvent filter is partly blocked no solvent or only very little solvent will drip out of the solvent tube.

### **Cleaning the Solvent Filters**

- Remove the blocked solvent filter from the bottle-head assembly and place it in a beaker with concentrated nitric acid (35 %) for one hour.
- Thoroughly flush the filter with bidistilled water (remove all nitric acid).
- Replace the filter.

#### **NOTE**

Never use the system without solvent filter installed.

---

# Operation Modes of the Vacuum Degasser

The vacuum degasser allows various operation modes. Operation mode 1 and 2 are the normal operation modes of the degasser:

- In operation mode 1 the vacuum degasser works at 115 Torr.
- In operation mode 2 the vacuum degasser works in the range between 115 to 190 Torr.
- The continuous mode can be selected for highest degassing efficiency of the degasser. In this mode the vacuum level is below 115 Torr.
- The error mode is activated in case the degasser cannot achieve a vacuum level of 190 Torr.

(1 Torr =  $1.33 \times 10^{-3}$  bar)

**Table 4**    Operation Modes of the Vacuum Degasser

	Operation Mode 1	Operation Mode 2	Continuous Mode	Error Mode
Pressure Range [Torr]	115	115 -190	Below 115	Above 190
DC Voltage Readings [mV] (rough values)	Lower Limit: 590 - 600 Upper Limit: 600 - 610	600 - 800	Below 600	Above 800
Operation mode	Pump switched on and off at upper and lower limit during operation	Pump switched on every 2 minutes for 30 seconds	Pump permanently running	8 Minutes of continuous pumping before shutdown
Status Indicator	OFF - below 800 mV YELLOW - above 800 mV	OFF - below 800 mV YELLOW - above 800 mV	OFF - below 800 mV YELLOW - above 800 mV	YELLOW
Failure Actions	Switch to Operation Mode 2	Switch to Error Mode	Switch to Error Mode	Shut down module, switch status indicator to RED

**NOTE**

The voltage readings given for the various modes are approximate values. All values are set in the factory according to the vacuum level. Depending on the vacuum sensor batch variations this will result in different voltage readings. Regardless of the readings of the various stages no adjustment should be performed on the degasser electronics.



**NOTE**

See service manual (troubleshooting and diagnostics) for setting up test meter connections to the degasser for reading the pressure sensor output.

## Normal Operation Mode 1

When the degasser is turned ON, the vacuum pump runs and is connected to ambient through the solenoid valve. The solenoid valve activates about 15 seconds after turning ON the degasser (you can hear it click). The vacuum pump then begins to pull a vacuum on the vacuum chamber. The voltage measurements begin to decrease rapidly. The yellow status LED turns off when the vacuum level reaches 190 Torr (DC voltage readings around 800 mV). The normal operation mode vacuum level (115 Torr) is achieved when the DC voltage measures approximately 590 to 600 mV. After achieving the normal operation mode vacuum level, the solenoid valve turns off. The vacuum pump continues to run for a few seconds, then it turns OFF. When the DC voltage measurement rises back to approximately 600 to 610 mV, the turn on process begins again. If the vacuum level of normal operation mode one cannot be achieved within 8 minutes the instrument turns into normal operation mode 2.

## Normal Operation Mode 2 (Timing Mode)

Under certain operational conditions (large amount of dissolved gas in mobile phase, high flow rates) the 115 Torr trigger level for operation mode 1 cannot be reached. The standard degasser automatically turns into operation mode 2. Normal operation mode 2 is a fixed timing mode. Every 2 min the degasser is turned ON for 30 s. This assures a pressure level in the range of 115 – 190 Torr. The pressure signal measured with the auxiliary cable is in the range of 600 – 800 mV. If the vacuum level of normal operation mode two cannot be achieved within 8 min the instrument turns into error mode.

## Continuous Mode

The continuous mode is activated either by switching SW1 on the main board to 1 (removal of cover is required, see “[Removing and Refitting the Top Cover](#)” on page 52, for identifying SW1, see *Service Manual, Repairs, Overview of Internal Parts*), or by connecting pin 1 (white cable) and pin 3 (green cable) of the auxiliary cable to each other. When turning ON the vacuum degasser the vacuum pump will run continuously. This will establish a vacuum level that is below the trigger level (600 mV / 115 Torr) of the normal operation mode 1. If the vacuum level of normal operation mode 2 cannot be achieved within 8 minutes the instrument turns into error mode.

### NOTE


When set to continuous mode the life time of the vacuum pump will be significantly reduced.

### When to use Continuous Mode

In continuous mode the vacuum pump runs continuously thus generating the highest degassing efficiency of the degasser. This mode is only recommended for very sensitive applications (e.g. RI detection).

## Error Mode

The error level for the standard degasser is 190 Torr (approximately 800 mV). This level cannot be achieved when there is a failure in the degasser (for example, leaks, etc.). When the error level is exceeded the yellow status indicator lamp is turned on and the vacuum pump runs continuously. If the degasser cannot reach one of the normal operation modes within 8 min the status indicator turns red and the vacuum pump is turned OFF.

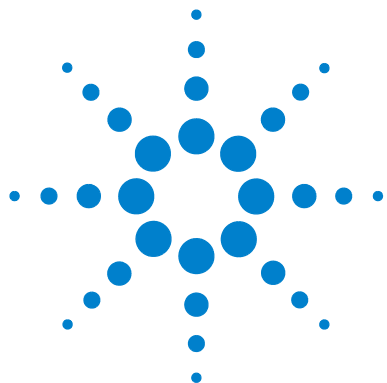

## 5 Troubleshooting and Diagnostics

Overview of the Degasser's Indicators 44

Status Indicators 45

Power Supply Lamp 45

Instrument Status Indicator 46



## Overview of the Degasser's Indicators

### Status Indicators


The standard degasser is provided with two status indicators which indicate the operational state (ready, busy, and error states) of the standard degasser. The status indicators provide a quick visual check of the operation of the standard degasser (see [“Status Indicators”](#) on page 45).

### Hardware Symptoms

A red status lamp at the vacuum degasser indicates a problem with the vacuum system or with the electronic control. The vacuum degasser generates an error output on the remote lines.

## Status Indicators

Two status lamps are located on the front of the standard degasser. The left lamp indicates the power supply status, the right lamp indicates the standard degasser status.

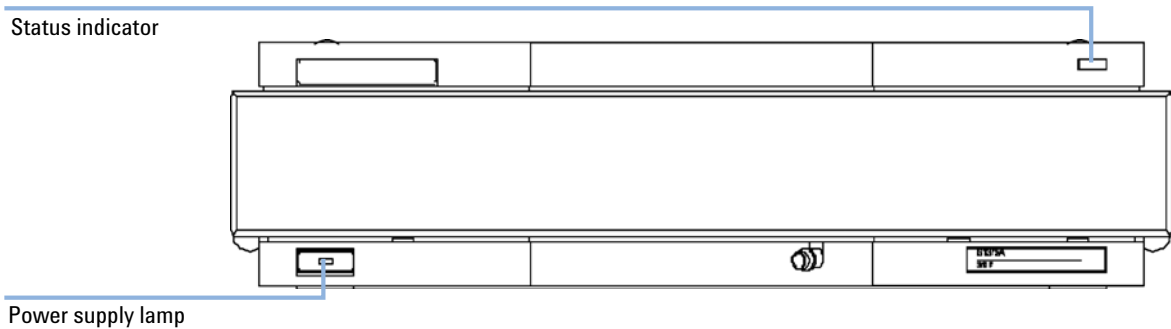

**Figure 10** Location of Status Lamps

## Power Supply Lamp

The power supply lamp is integrated into the main power switch. When the lamp is illuminated (green), the power is ON.

## Instrument Status Indicator

The instrument status indicator indicates one of three possible instrument conditions.

- When the lamp is OFF, the standard degasser is in *ready* condition (only if the power supply lamp is ON, otherwise, the instrument is switched OFF, or there is a defect in the power supply). A ready condition exists when there is sufficient vacuum in the standard degasser.
- A *busy* condition is indicated ,when the lamp is yellow. A busy condition exists when the vacuum pump of the standard degasser is working to generate or maintain vacuum in the chambers. This is the case when the standard degasser is turned ON for the first time or the pressure rises above its limit during normal operation.
- An *error* condition is indicated, when the lamp is red. An error condition exists when the standard degasser detects an internal defect that does not allow the vacuum to be built up in a certain time frame (approximately 8 minutes).

### CAUTION

#### Internal Leak or Electronic Failure

If the error LED is on, there is either an internal leak in the vacuum system or an electronic failure. In case of an internal leak it is possible that solvent may enter the vacuum chamber and solvent may leak into the waste drain.

- To prevent any damage of the standard degasser, switch off the degasser and remove the solvent bottles from the solvent cabinet to stop any gravity-caused flow of solvent into the vacuum chamber.



## 6 Maintenance

Warnings and Cautions	48
Introduction to Maintenance	50
Simple Repairs - Maintenance	50
Using the ESD Strap	50
Cleaning the Instrument	51
Removing and Refitting the Top Cover	52
Assembling the Main Cover	55
Exchanging the Fuses of the Power Inline Filter	56
Exchanging the Status Light Pipe	58

This chapter describes the maintenance of the module.



## Warnings and Cautions

### WARNING

#### Toxic, flammable and hazardous solvents, samples and reagents

**The handling of solvents, samples and reagents can hold health and safety risks.**

- When working with these substances observe appropriate safety procedures (for example by wearing goggles, safety gloves and protective clothing) as described in the material handling and safety data sheet supplied by the vendor and follow good laboratory practice.
  - The amount of substances should be reduced to the minimal volume required for the analysis.
  - Do not operate the instrument in an explosive atmosphere.
- 

### WARNING

#### Electrical shock

**Repair work at the module can lead to personal injuries, e.g. shock hazard, when the cover is opened.**

- Do not remove the metal top cover of the module. No serviceable parts inside.
  - Only certified persons are authorized to carry out repairs inside the module.
- 

### WARNING

#### Personal injury or damage to the product

**Agilent is not responsible for any damages caused, in whole or in part, by improper use of the products, unauthorized alterations, adjustments or modifications to the products, failure to comply with procedures in Agilent product user guides, or use of the products in violation of applicable laws, rules or regulations.**

- Use your Agilent products only in the manner described in the Agilent product user guides.
-



**CAUTION**

Safety standards for external equipment

- If you connect external equipment to the instrument, make sure that you only use accessory units tested and approved according to the safety standards appropriate for the type of external equipment.
-

## Introduction to Maintenance


### Simple Repairs - Maintenance

The standard degasser is designed for easy repair. The most frequent repairs such as exchanging power fuses and status light pipes can be performed by the user, but require opening the main cover of the standard degasser. These repairs are described in this chapter.

### Using the ESD Strap

Electronic boards are sensitive to electronic discharge (ESD). In order to prevent damage, always use an ESD strap supplied in the standard accessory kit when handling electronic boards and components.

- 1 Unwrap the first two folds of the band and wrap the exposed adhesive side firmly around your wrist.
- 2 Unroll the rest of the band and peel the liner from the copper foil at the opposite end.
- 3 Attach the copper foil to a convenient and exposed electrical ground.

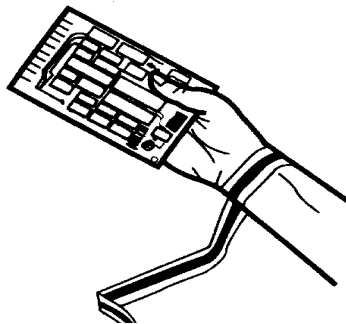

**Figure 11**    Using the ESD Strap

## Cleaning the Instrument

### **WARNING**

**Liquid dripping into the electronic compartment of your module.**

**Liquid in the module electronics can cause shock hazard and damage the module.**

- Do not use an excessively damp cloth during cleaning.
  - Drain all solvent lines before opening any fittings.
- 

The vacuum degasser case should be kept clean. Cleaning should be done with a soft cloth slightly dampened with water or a solution of water and a mild detergent. Do not use an excessively damp cloth that liquid can drip into the vacuum degasser.

## Removing and Refitting the Top Cover

**Tools required**            Screwdriver Pozidriv #1

- Preparations**
- Switch OFF the degasser at the main power switch
  - Disconnect the power cable and remote cable
  - Disconnect all solvent tubes from the ports of the degasser
  - Remove solvent cabinet from the degasser
  - Remove degasser from the stack.


### **WARNING**

**Toxic, flammable and hazardous solvents, samples and reagents**


**The handling of solvents, samples and reagents can hold health and safety risks.**

- When working with these substances observe appropriate safety procedures (for example by wearing goggles, safety gloves and protective clothing) as described in the material handling and safety data sheet supplied by the vendor and follow good laboratory practice.
- The amount of substances should be reduced to the minimal volume required for the analysis.
- Do not operate the instrument in an explosive atmosphere.


**1** Remove the front panel.

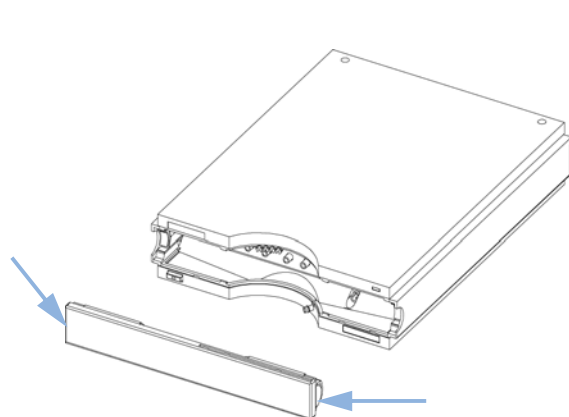


**2** Unclip the clips on the top cover .

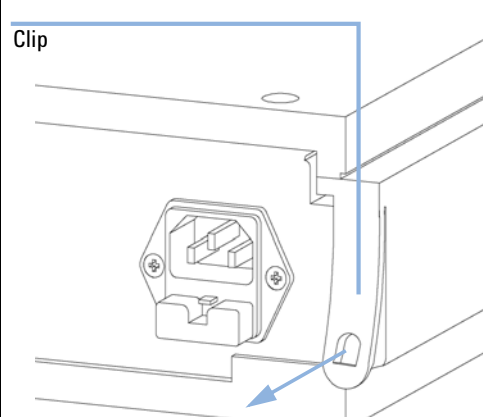


**3** Lift the cover away.

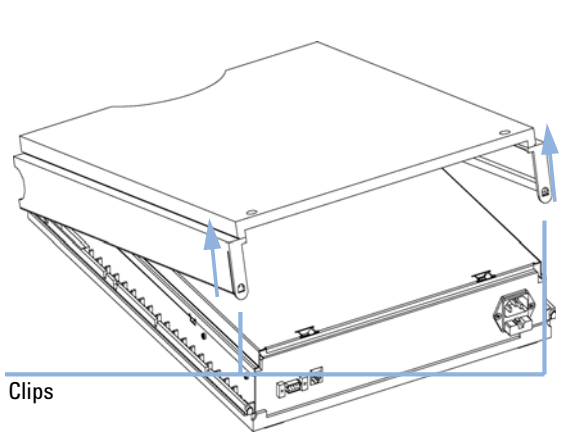


**4** Unscrew the screws on the top plate and remove the plate by lifting its back and then sliding the plate to the front.

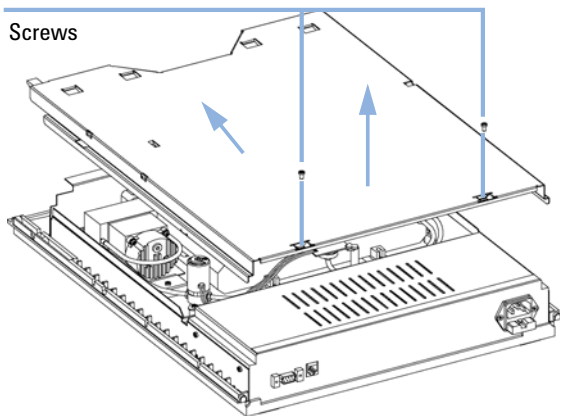

**5** Place the metal cover onto the housing. Lower the front of the plate first, then slide panel to the back. Assure the four metal tabs of the panel slide underneath the Z-plane. Fix the two holding screws.

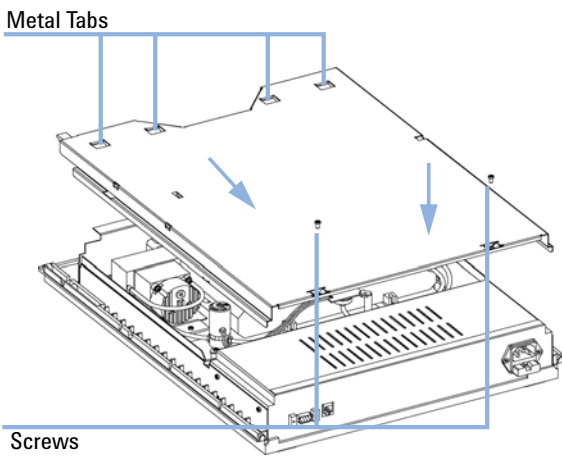


**6** Replace the top cover. Ensure the clips are seated correctly.

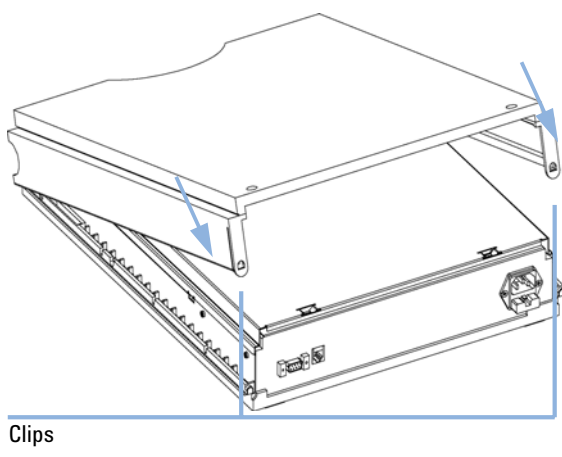

## 6 Maintenance

### Introduction to Maintenance

**7** Reinstall the front cover.

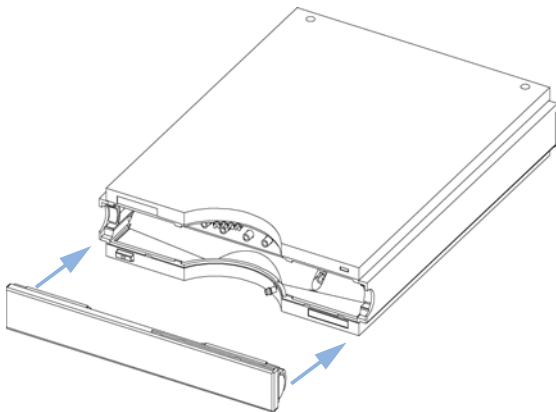

**8** Reinstall the degasser in your system stack and connect the cables and capillaries and turn on the degasser.

## Assembling the Main Cover

When	• If cover is broken.		
Parts required	#	p/n	Description
	1	5065-9989	Cover kit (includes base, top, left and right)

### CAUTION



Wrong assembly

In case you insert the left or right side in the opposite position, you may not be able to remove the side from the top part.

→ Take care not to mix up left and right side.

### NOTE

The cover kit contains all parts, but it is not assembled.
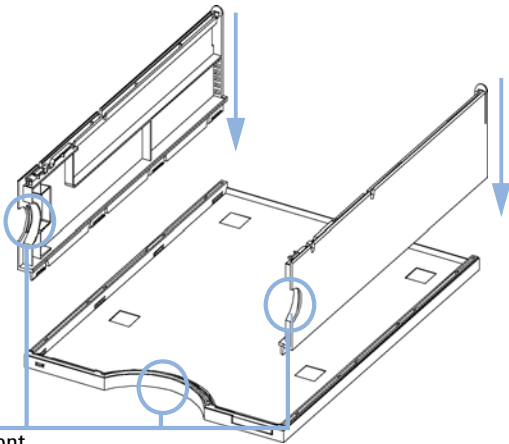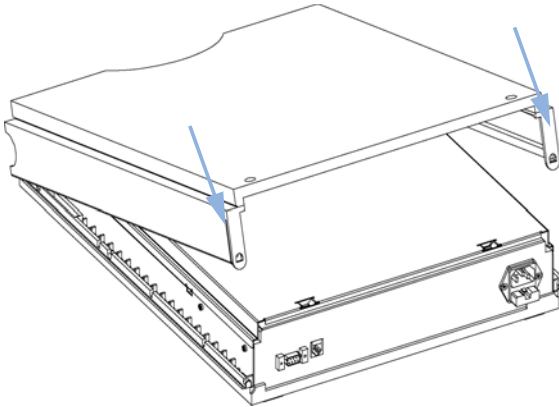
<p><b>1</b> Place the top part on the bench and insert the left and right side into the top part.</p>  <p>Front</p>	<p><b>2</b> Replace the cover.</p> 
<p><b>Next Steps:</b></p> <p><b>3</b> Replace the degasser in the stack and reconnect the cables and capillaries.</p> <p><b>4</b> Turn ON the degasser.</p>	

# Exchanging the Fuses of the Power Inline Filter

When	• If the standard degasser appears dead		
Tools required	• Test meter (if available)		
Parts required	#	p/n	Description
	1	2110-0458	Fuse: 250V, T 500 mA , compatible to all supported line voltages

**WARNING**

*Use of unsupplied cables*

**The use of cables which haven't been supplied by Agilent Technologies can lead to damage of the electronic components or personal injury.**

→ ?Never use cables other than the ones supplied by Agilent Technologies to ensure proper functionality and compliance with safety or EMC regulations.


**NOTE**

The Agilent 1260 Infinity Standard Degasser power supply has wide-ranging capability (see [Table 1](#) on page 15). It accepts any line voltage in the range mentioned below. Consequently there is no voltage selector in the rear of the standard degasser. There are two externally accessible fuses, that protect the power supply. These fuses are identical for all accepted line voltages.

- 1 Switch OFF the power switch at the front of the instrument.
- 2 Remove the power cable from the power connector at the rear of the instrument.



- 3 Press down the clip of the fuse holder and pull out of the power socket.

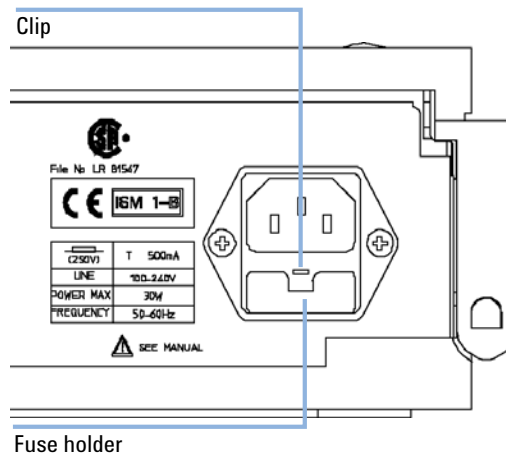



**Figure 12** Fuse Holder and Clip

- 4 Remove the fuses from the fuse holders.
- 5 Ensure the fuse wires inside the fuses are not broken. If a test meter is available, check the resistance of each fuse. A good fuse shows a low resistance (approximately 0 Ohm).
- 6 If a fuse is defective (wire broken or high resistance), insert a new fuse.
- 7 Reinsert the fuse holders and the power cable.
- 8 Switch ON the power switch.

# Exchanging the Status Light Pipe

<b>When</b>	• If part is broken		
<b>Tools required</b>	Screwdriver Pozidriv #1		
<b>Parts required</b>	<b>#</b>	<b>p/n</b>	<b>Description</b>
	1	5041-8384	Status light pipe
<b>Preparations</b>	• Remove the front cover and top cover, see <a href="#">“Removing and Refitting the Top Cover”</a> on page 52.		
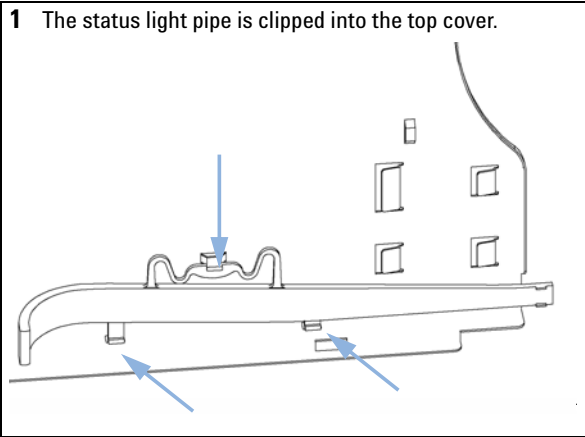
<p><b>1</b> The status light pipe is clipped into the top cover.</p> 	<p><b>2</b> Replace the top cover, see <a href="#">“Removing and Refitting the Top Cover”</a> on page 52.</p>
---------------------------------------------------------------------------------------------------------------------------------------------------------	---------------------------------------------------------------------------------------------------------------

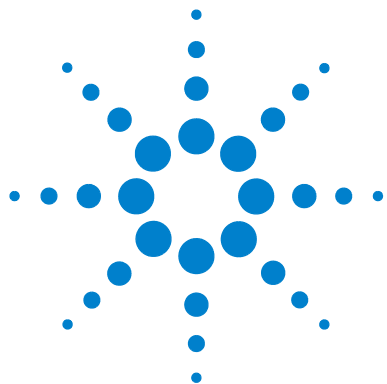

## 7

# Parts and Materials for Maintenance

Cover Parts 60


Power and Status Light Pipes 61

Accessory Kit Contents 62



## Cover Parts


Item	p/n	Description
1	5065-9989	Cover kit (includes base, top, left and right)
2	5065-9990	Front cover
3	5042-8901	Name plate
4	5041-8387	Tube clip

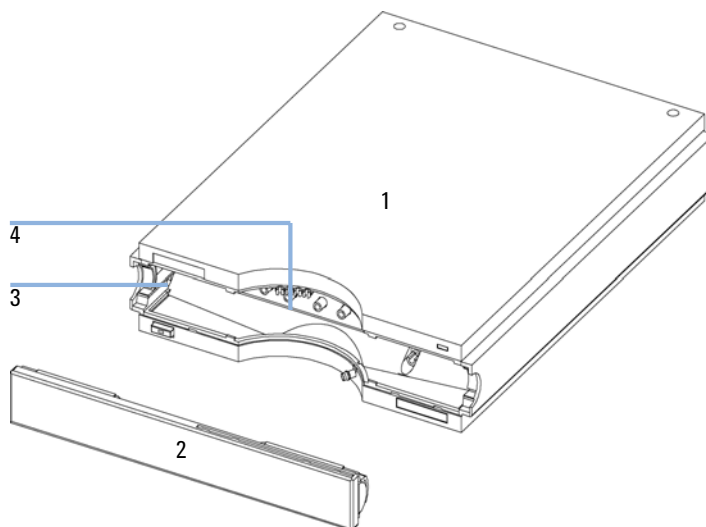

**Figure 13**   Cover Parts

## Power and Status Light Pipes

Item	p/n	Description
1	5041-8383	Power switch coupler
2	5041-8382	Power switch light pipe
3	5041-8381	Power switch button
4	5041-8384	Status light pipe

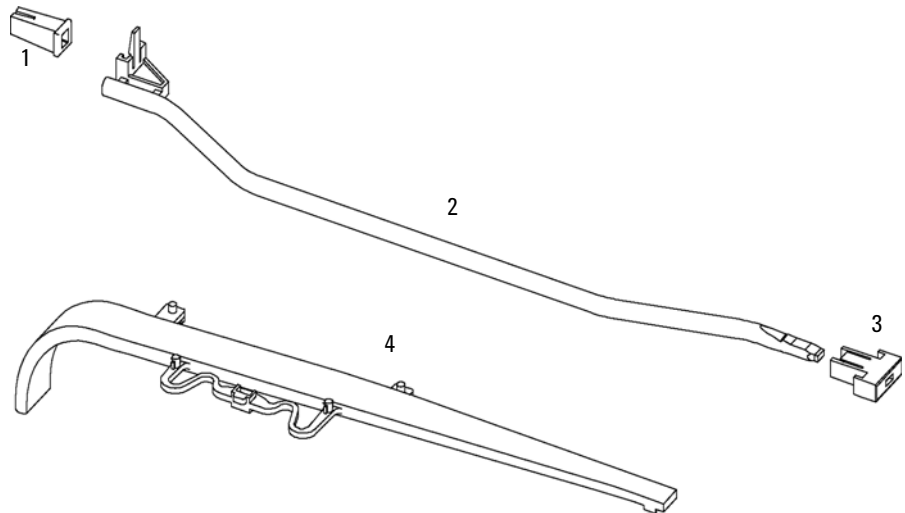

**Figure 14** Power and Status Light Pipes

## Accessory Kit Contents

Accessory Kit (p/n G1322-68705)

<b>p/n</b>	<b>Description</b>
5062-8534	Syringe Reorder number (pack of 10)
9301-1337	Syringe adapter
0100-1710	Mounting Tool for Tubing Connections
5062-2463	Corrugated waste tube (reorder pack), 5 m
G1322-67300	Kit of 4 solvent tubes for connection degasser to MCGV (Quaternary Pump) including labels



## 8 Identifying Cables

Cable Overview	64
Remote Cables	66
Auxiliary Cable	68



## Cable Overview

**NOTE**

Never use cables other than the ones supplied by Agilent Technologies to ensure proper functionality and compliance with safety or EMC regulations.

### Analog cables

p/n	Description
35900-60750	Agilent module to 3394/6 integrators
35900-60750	Agilent 35900A A/D converter
01046-60105	Analog cable (BNC to general purpose, spade lugs)

### Remote cables

p/n	Description
03394-60600	Agilent module to 3396A Series I integrators  3396 Series II / 3395A integrator, see details in section <a href="#">“Remote Cables”</a> on page 66
03396-61010	Agilent module to 3396 Series III / 3395B integrators
5061-3378	Agilent module to Agilent 35900 A/D converters (or HP 1050/1046A/1049A)
01046-60201	Agilent module to general purpose

### BCD cables

p/n	Description
03396-60560	Agilent module to 3396 integrators
G1351-81600	Agilent module to general purpose



### CAN cables

p/n	Description
5181-1516	CAN cable, Agilent module to module, 0.5 m
5181-1519	CAN cable, Agilent module to module, 1 m

### LAN cables

p/n	Description
5023-0203	Cross-over network cable, shielded, 3 m (for point to point connection)
5023-0202	Twisted pair network cable, shielded, 7 m (for point to point connection)


### External Contact Cable

p/n	Description
G1103-61611	External contact cable - Agilent module interface board to general purposes

### RS-232 cables


p/n	Description
G1530-60600	RS-232 cable, 2 m
RS232-61600	RS-232 cable, 2.5 m Instrument to PC, 9-to-9 pin (female). This cable has special pin-out, and is not compatible with connecting printers and plotters. It's also called "Null Modem Cable" with full handshaking where the wiring is made between pins 1-1, 2-3, 3-2, 4-6, 5-5, 6-4, 7-8, 8-7, 9-9.
5181-1561	RS-232 cable, 8 m

# Remote Cables

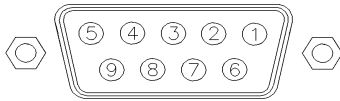

One end of these cables provides a Agilent Technologies APG (Analytical Products Group) remote connector to be connected to Agilent modules. The other end depends on the instrument to be connected to.


## Agilent Module to 3396A Integrators

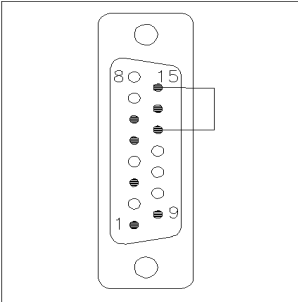p/n 03394-60600	Pin 3394	Pin Agilent module	Signal Name	Active (TTL)
	9	1 - White	Digital ground	
	NC	2 - Brown	Prepare run	Low
	3	3 - Gray	Start	Low
	NC	4 - Blue	Shut down	Low
	NC	5 - Pink	Not connected	
	NC	6 - Yellow	Power on	High
	5,14	7 - Red	Ready	High
	1	8 - Green	Stop	Low
	NC	9 - Black	Start request	Low
	13, 15		Not connected	


## Agilent Module to 3396 Series II / 3395A Integrators

Use the cable Agilent module to 3396A Series I integrators (p/n 03394-60600) and cut pin #5 on the integrator side. Otherwise the integrator prints START; not ready.


## Agilent Module to 3396 Series III / 3395B Integrators

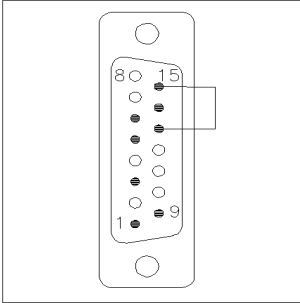p/n 03396-61010	Pin 33XX	Pin Agilent module	Signal Name	Active (TTL)
	9	1 - White	Digital ground	
	NC	2 - Brown	Prepare run	Low
	3	3 - Gray	Start	Low
	NC	4 - Blue	Shut down	Low
	NC	5 - Pink	Not connected	
	NC	6 - Yellow	Power on	High
	14	7 - Red	Ready	High
	4	8 - Green	Stop	Low
	NC	9 - Black	Start request	Low
	13, 15		Not connected	

## Agilent Module to Agilent 35900 A/D Converters

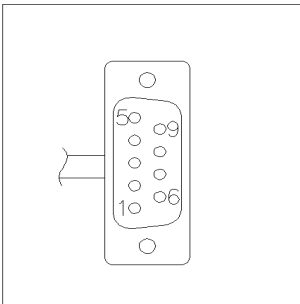p/n 5061-3378	Pin 35900 A/D	Pin Agilent module	Signal Name	Active (TTL)
	1 - White	1 - White	Digital ground	
	2 - Brown	2 - Brown	Prepare run	Low
	3 - Gray	3 - Gray	Start	Low
	4 - Blue	4 - Blue	Shut down	Low
	5 - Pink	5 - Pink	Not connected	
	6 - Yellow	6 - Yellow	Power on	High
	7 - Red	7 - Red	Ready	High
	8 - Green	8 - Green	Stop	Low
	9 - Black	9 - Black	Start request	Low

Agilent Module to General Purpose

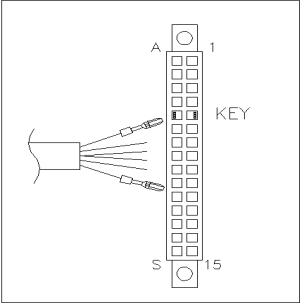
p/n 01046-60201	Pin Universal	Pin Agilent module	Signal Name	Active (TTL)
		1 - White	Digital ground	
		2 - Brown	Prepare run	Low
		3 - Gray	Start	Low
		4 - Blue	Shut down	Low
		5 - Pink	Not connected	
		6 - Yellow	Power on	High
		7 - Red	Ready	High
		8 - Green	Stop	Low
		9 - Black	Start request	Low

Auxiliary Cable



One end of this cable provides a modular plug to be connected to the Agilent vacuum degasser. The other end is for general purpose.

Agilent Vacuum Degasser to general purposes

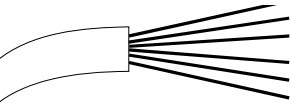p/n G1322-81600	Color	Pin Agilent 1100	Signal Name
	White	1	Ground
	Brown	2	Pressure signal
	Green	3	
	Yellow	4	
	Grey	5	DC + 5 V IN
	Pink	6	Vent

## **8 Identifying Cables**

### **Cable Overview**



## 9 Appendix

General Safety Information [72](#)

The Waste Electrical and Electronic Equipment (WEEE) Directive  
(2002/96/EC) [75](#)

Radio Interference [76](#)

Sound Emission [77](#)

Agilent Technologies on Internet [78](#)



## General Safety Information

The following general safety precautions must be observed during all phases of operation, service, and repair of this instrument. Failure to comply with these precautions or with specific warnings elsewhere in this manual violates safety standards of design, manufacture, and intended use of the instrument. Agilent Technologies assumes no liability for the customer's failure to comply with these requirements.

### **WARNING**

**Ensure the proper usage of the equipment.**

**The protection provided by the equipment may be impaired.**

→ The operator of this instrument is advised to use the equipment in a manner as specified in this manual.

---

## Safety Standards

This is a Safety Class I instrument (provided with terminal for protective earthing) and has been manufactured and tested according to international safety standards.

## Operation

Before applying power, comply with the installation section. Additionally the following must be observed.

Do not remove instrument covers when operating. Before the instrument is switched on, all protective earth terminals, extension cords, auto-transformers, and devices connected to it must be connected to a protective earth via a ground socket. Any interruption of the protective earth grounding will cause a potential shock hazard that could result in serious personal injury. Whenever it is likely that the protection has been impaired, the instrument must be made inoperative and be secured against any intended operation.



Make sure that only fuses with the required rated current and of the specified type (normal blow, time delay, and so on) are used for replacement. The use of repaired fuses and the short-circuiting of fuse holders must be avoided.

Some adjustments described in the manual, are made with power supplied to the instrument, and protective covers removed. Energy available at many points may, if contacted, result in personal injury.

Any adjustment, maintenance, and repair of the opened instrument under voltage should be avoided whenever possible. When inevitable, this has to be carried out by a skilled person who is aware of the hazard involved. Do not attempt internal service or adjustment unless another person, capable of rendering first aid and resuscitation, is present. Do not replace components with power cable connected.

Do not operate the instrument in the presence of flammable gases or fumes. Operation of any electrical instrument in such an environment constitutes a definite safety hazard.






Do not install substitute parts or make any unauthorized modification to the instrument.

Capacitors inside the instrument may still be charged, even though the instrument has been disconnected from its source of supply. Dangerous voltages, capable of causing serious personal injury, are present in this instrument. Use extreme caution when handling, testing and adjusting.

When working with solvents please observe appropriate safety procedures (e.g. goggles, safety gloves and protective clothing) as described in the material handling and safety data sheet by the solvent vendor, especially when toxic or hazardous solvents are used.

# Safety Symbols

**Table 5** Safety Symbols

Symbol	Description
	The apparatus is marked with this symbol when the user should refer to the instruction manual in order to protect risk of harm to the operator and to protect the apparatus against damage.
	Indicates dangerous voltages.
	Indicates a protected ground terminal.
	Indicates eye damage may result from directly viewing the light produced by the deuterium lamp used in this product.
	The apparatus is marked with this symbol when hot surfaces are available and the user should not touch it when heated up.

## WARNING

### A WARNING

alerts you to situations that could cause physical injury or death.

- Do not proceed beyond a warning until you have fully understood and met the indicated conditions.

## CAUTION

### A CAUTION

alerts you to situations that could cause loss of data, or damage of equipment.

- Do not proceed beyond a caution until you have fully understood and met the indicated conditions.

## The Waste Electrical and Electronic Equipment (WEEE) Directive (2002/96/EC)

### Abstract

The Waste Electrical and Electronic Equipment (WEEE) Directive (2002/96/EC), adopted by EU Commission on 13 February 2003, is introducing producer responsibility on all electric and electronic appliances starting with 13 August 2005.

#### NOTE

This product complies with the WEEE Directive (2002/96/EC) marking requirements. The affixed label indicates that you must not discard this electrical/electronic product in domestic household waste.

Product Category:

With reference to the equipment types in the WEEE Directive Annex I, this product is classed as a Monitoring and Control Instrumentation product.

---



#### NOTE

Do not dispose off in domestic household waste

To return unwanted products, contact your local Agilent office, or see [www.agilent.com](http://www.agilent.com) for more information.

---

## Radio Interference

Cables supplied by Agilent Technologies are screened to provide optimized protection against radio interference. All cables are in compliance with safety or EMC regulations.

### **Test and Measurement**

If test and measurement equipment is operated with unscreened cables, or used for measurements on open set-ups, the user has to assure that under operating conditions the radio interference limits are still met within the premises.

## Sound Emission

### Manufacturer's Declaration

This statement is provided to comply with the requirements of the German Sound Emission Directive of 18 January 1991.

This product has a sound pressure emission (at the operator position) < 70 dB.

- Sound Pressure  $L_p < 70$  dB (A)
- At Operator Position
- Normal Operation
- According to ISO 7779:1988/EN 27779/1991 (Type Test)

## **Agilent Technologies on Internet**

For the latest information on products and services visit our worldwide web site on the Internet at:

<http://www.agilent.com>

Select Products/Chemical Analysis

It will provide also the latest firmware of the modules for download.

# Index

## A

Agilent  
     on internet 78  
 algae growth 38  
 ambient operating temperature 15  
 ambient non-operating temperature 15  
 analog output 16, 16  
 APG remote connector 24  
 AUX output 24  
 auxiliary  
     cable 68

## B

bench space 14  
 busy condition 46

## C

cable  
     auxiliary 19, 68  
     CAN 24  
     power 19, 52  
     remote 19, 66  
 cables  
     analog 64  
     BCD 64  
     CAN 65  
     external contact 65  
     LAN 65  
     overview 64  
     remote 64  
     RS-232 65  
 CAN cable 24  
 changing solvents 30  
 cleaning the instrument 51

condensation 14  
 control circuit 6  
 cover 60

## D

dead volume 20  
 defect on arrival 18  
 delay volume 20  
 delivery checklist 19  
 dimensions 15

## E

electronic waste 75  
 electronics  
     how does it work 7  
 error condition 46  
 exchanging  
     status light pipe 58

## F

flow connections 26, 26  
 flow rate 16  
 frequency range 15  
 front of the vacuum degasser 25  
 fuse 56

## H

highest injection precision 36  
 highest retention time reproducibility 36  
 humidity 15

## I

information

    on sound emission 77  
 installation  
     bench space 14  
 internal leak 46  
 internet 78  
 introduction to the degasser 6

## L

line frequency 15  
 line voltage 15

## M

main cover, assembling 55  
 materials in contact with solvent 16  
 maximum flow rate 16  
 maximum sensitivity 36  
 mounting tool 27, 27

## N

non-operating altitude 15  
 non-operating temperature 15  
 number of channels 16

## O

operating altitude 15  
 operating temperature 15  
 operational hints 30

## P

packaging  
     damaged 18  
 parts  
     cover 60

## Index

- damaged 19
- light pipes 61
- missing 19
- performance specifications 16
- pH range 16
- physical specifications 15
- power light pipe 61
- power supply lamp 45
- power cable 52
- power consideration 12
- power consumption 15
- power cords 13
- power supply 56
- power switch 23
- pressure sensor readings 24
- pressure sensor 6
- priming
  - with a pump 30, 32
  - with a syringe 30
- principle of operation 6

## R

- radio interference 76
- ready condition 46
- rear of the vacuum degasser 24
- remote
  - cable 66

## S

- safety class I 72
- safety
  - general information 72
  - standards 15
  - symbols 74
- screwdriver pozidriv #1 52, 58
- site requirements
  - power cords 13
- snap fastener 27

- solvent cabinet 52
- solvent filters
  - cleaning 39
  - prevent blocking 38
- solvent information 40
- sound emission 77
- specification
  - performance specifications 16
  - physical 15
- stack configuration 20
- status light pipe 58, 61
- status indicator 45
- syringe adapter 31, 33

## T

- tool 27
  - mounting tool 27
- tools
  - screwdriver pozidriv #1 52, 58
- top cover, removing and refitting 52
- transportation 33

## V

- vacuum pump 6
- voltage range 15
- voltage selector 56
- volume per channel 16

## W

- waste
  - electrical and electronic equipment 75
- WEEE directive 75
- weight 15
- when to use a degasser 42





## **In This Book**

This manual contains user information about the Agilent 1260 Infinity Standard Degasser. The manual describes the following:

- introduction,
- site requirements and specifications,
- installing the degasser,
- using the degasser,
- troubleshooting and diagnostics,
- maintenance,
- parts and materials for maintenance,
- identifying cables,
- appendix.

© Agilent Technologies 2006-2008, 2010

Printed in Germany  
06/10



G1322-90012



**Agilent Technologies**