

Agilent 4200 FlexScan FTIR

Operation Manual



Agilent Technologies

Notices

© Agilent Technologies, Inc. 2010, 2011

No part of this manual may be reproduced in any form or by any means (including electronic storage and retrieval or translation into a foreign language) without prior agreement and written consent from Agilent Technologies, Inc. as governed by United States and international copyright laws.

Manual Part Number

0024-401

Edition

Second edition, April 2011 Printed in USA Agilent Technologies, Inc.

Errata Statement

NOTICE: This document contains references to A2 Technologies. Please note that A2 Technologies is now part of Agilent Technologies. For more information, go to www.agilent.com/chem

Warranty

The material contained in this document is provided "as is," and is subject to being changed, without notice, in future editions. Further, to the maximum extent permitted by applicable law, Agilent disclaims all warranties, either express or implied, with regard to this manual and any information contained herein, including but not limited to the implied warranties of merchantability and fitness for a particular purpose. Agilent shall not be liable for errors or for incidental or consequential damages in connection with the furnishing, use, or performance of this document or of any information contained herein. Should Agilent and the user have a separate written agreement with warranty terms covering the material in this document that conflict with these terms, the warranty terms in the separate agreement shall control.

Technology Licenses

The hardware and/or software described in this document are furnished under a license and may be used or copied only in accordance with the terms of such license.

Restricted Rights Legend

If software is for use in the performance of a U.S. Government prime contract or subcontract, Software is delivered and licensed as "Commercial computer software" as defined in DFAR 252.227-7014 (June 1995), or as a "commercial item" as defined in FAR 2.101(a) or as "Restricted computer software" as defined in FAR 52.227-19 (June 1987) or any equivalent agency regulation or contract clause. Use, duplication or disclosure of Software is subject to Agilent Technologies' standard commercial license terms, and non-DOD Departments and Agencies of the U.S. Government will receive no greater than Restricted Rights as defined in FAR 52.227-19(c)(1-2) (June 1987). U.S. Government users will receive no greater than Limited Rights as defined in FAR 52.227-14 (June 1987) or DFAR 252.227-7015 (b)(2) (November 1995), as applicable in any technical data.

Safety Notices

CAUTION

A **CAUTION** notice denotes a hazard. It calls attention to an operating procedure, practice, or the like that, if not correctly performed or adhered to, could result in damage to the product or loss of important data. Do not proceed beyond a **CAUTION** notice until the indicated conditions are fully understood and met.

WARNING

A WARNING notice denotes a hazard. It calls attention to an operating procedure, practice, or the like that, if not correctly performed or adhered to, could result in personal injury or death. Do not proceed beyond a WARNING notice until the indicated conditions are fully understood and met.

Contents

Contents

1.	Introduction	5
2.	Getting Started	9
	Unpacking	9
	Pa	cking list 10
	Component	ts 11
	Installing th	ne battery12
	Power swit	ch LED status 13
	Starting the	e software 14
	Ch	arging the handheld PC 15
	Log	gging on to MicroLab Mobile for the first time 15
	Performanc	e verification 16
	Dia	agnostic values 17
	Per	rformance validation 19
3.	Analyzing Sam	ples 23
	External ref	flectance, grazing angle and diffuse reflectance 23
	Re	flectance standards and gain settings 24
	Ste	eps 25
	Co	llecting a background spectrum 25
	Co	llecting a sample spectrum 26
	Internal ref	lectance (ATR) 27
	Ste	eps 28
	Cle	eaning the sample probe 28
	Co	llecting a background spectrum 29

Contents

Collecting a sample spectrum30Using the solid sample ATR press31

- 4. Maintenance 35
- 5. Spare Parts 37
- 6. Specifications 39

Technical 39

Environmental conditions 40

Handheld PC data station 40



The Agilent 4200 FlexScan FTIR instrument is a portable, handheld, battery-operated analyzer designed to measure a variety of solid and liquid samples in locations difficult or impossible to analyze in a laboratory. It can be used for quantitative or qualitative analysis of materials. The system is small, lightweight and packaged in a weather-resistant enclosure suitable for outdoor use. It comes standard with internal battery power for operation of up to 4 hours.

The 4200 FlexScan FTIR can be safely used in operating temperatures ranging from 0 to 50 degrees Celsius (32 to 122 degrees Fahrenheit) and in humidity environments as high as 95% for extended periods of time.





Fire Hazard

4200 FlexScan FTIR systems are NOT intrinsically safe. Use the system only in atmospheres that have been tested for flammable materials. If the equipment is used in a manner not specified in this manual, the protection provided by the equipment may be impaired.

CAUTION

Although the 4200 FlexScan FTIR system is designed to withstand significant shock and vibration found in a field environment, use caution when handling and avoid dropping the system when it is out of the shipping case. Dropping the system may result in equipment failure.

The 4200 FlexScan FTIR uses technology known as Fourier transform infrared (FTIR) spectroscopy, which is the current state-of-the-art technique for molecular compound identification and quantification. FTIR uses an infrared (IR) light source to pass through the sample and onto a detector, which precisely measures the amount of light absorbed by the sample. This absorbance creates a unique spectral fingerprint that is used to identify the molecular structure of the sample and determine the exact quantity of a particular compound in a mixture. In addition to the capability to measure samples onsite that in the past were typically measured in the lab, this system is a non-destructive tool (NDT) that allows the analysis to be performed without damaging the sample.

The heart of the Agilent FTIR spectrometer system is a patented Michelson interferometer design mounted on shock-dampening mechanisms to protect the components from shock and vibration in the field. This proprietary design is the key to successfully making FTIR technology compact, lightweight, rugged and field-portable.

The 4200 FlexScan FTIR system can be configured for a number of different sampling configurations to accommodate the analysis of a range of liquids, solids, powders, pastes and gels. Information on the optimal use of the different sampling configurations of the 4200 FlexScan system is listed in the following categories:

- Agilent 4200 FlexScan FTIR External Reflectance System: For the chemical analysis of solid or hard surface samples, especially thickness measurement of thin coatings and films on reflective surfaces. Typical applications include thickness measurement, analysis of anodization of metals, and measurement of surface cleanliness. The External Reflectance System uses a 45 degree specular reflectance probe or a grazing angle specular reflectance probe. The 45 degree probe is optimal for analysis of thin films greater than 1 micron in thickness. Specular reflectance is an ideal method for the analysis of surface contamination.
- Agilent 4200 FlexScan FTIR Grazing Angle System: The grazing angle sample interface works in a similar manner to the external reflectance sample interface; however, the angle of incidence is 82 degrees. The increased angle of incidence increases. The grazing angle probe is optimal for analysis of ultra thin films less than 1 micron in thickness. The grazing angle probe should be used when very low level contamination needs to be analyzed (nanometer thickness levels).
- Agilent 4200 FlexScan FTIR Internal Reflectance System: For the chemical analysis of liquids, powders, pastes and gels. The Internal Reflectance System uses a diamond optical sampling probe that incorporates the attenuated total reflectance (ATR) technique. This method is optimal for the identification of a variety of samples and has the advantage that no sample preparation is required to obtain the proper IR absorbance levels. The diamond ATR probe for the 4200 FlexScan FTIR is a single reflection ATR. The single reflection ATR sampling technique is ideal for identification of thick or highly-absorbing samples where small IR pathlengths are required (typically 2 micron depth of penetration).

The 4200 FlexScan FTIR is equipped with a handheld PC (Socket PDA) as standard and can be interfaced to the optional laptop computer for development of new methods and for detailed viewing and manipulating the spectra. The 4200 FlexScan software user interface is intuitive, simple to use and requires no specialized technical training. The spectral display is represented with easy to read status bars and report formats, which increase ease of use on the smaller handheld PC display. With the touch of a button, the system provides valuable information about the identity and amount of chemical substances present in a material.



2. Getting Started

Unpacking 9 Components 11 Installing the battery 12 Power switch LED status 13 Starting the software 14 Performance verification 16

Unpacking

To unpack your spectrometer system:

1 After receiving the Agilent 4200 FlexScan FTIR delivery, do not immediately open the shipping container. Instead, place the shipping case in a room-temperature environment and allow several hours for the contents of the container to reach the room ambient temperature. This is to avoid unnecessary condensation on the components prior to the initial setup and installation process.

Your 4200 FlexScan FTIR has been packed to ensure that it arrives at your facility ready to go to work. Custom foam inserts protect the instrument from damage during travel. Whenever your instrument is shipped via common carrier, it needs to be repacked in the shipping case. Using this case, your 4200 FlexScan can be removed and re-packed in the case in a matter of seconds, while maintaining instrument security.

Getting Started

NOTE

2 The shipping enclosure should contain the standard items listed below as well as any optional items ordered. Inspect the enclosure closely to make sure all items have been removed from the shipment packaging. Also ensure that all the items in the packing list below have arrived undamaged and are in good working condition. Contact Agilent immediately if any items are found to be missing or damaged in shipment.

Keep all original packing material for storing, shipping and transporting the system in the future.

3 Remove the 4200 FlexScan FTIR system from its shipping case and place it on its side on a flat, stable surface. The instrument must be kept away from hot surfaces and any sources of electromagnetic interference.

Packing list

Plastic shipping case

- One Agilent 4200 FlexScan FTIR spectrometer system
- One Agilent 4200 FlexScan FTIR mirrored reference cap
- One Agilent 4200 FlexScan FTIR diffuse reference cap
- One Agilent 4200 FlexScan FTIR polystyrene frequency check standard
- Two Agilent 4200 FlexScan FTIR Li-ion batteries
- One Agilent 4200 FlexScan FTIR Operation Manual
- One Agilent MicroLab Software Operation Manual
- One Socket handheld PC system (PDA)
- One Agilent 4200 FlexScan FTIR wrist strap
- One shipping container

Cardboard box

- One handheld PC USB cable with power connector
- One battery charger
- One power supply for battery charger
- One AC line cord
- One installation disk (needed as backup copy only)
- One Agilent 4200 FlexScan FTIR USB cable

Components

The 4200 FlexScan FTIR is designed to measure samples in a handheld configuration. Sample measurement can be started using the handle-mounted trigger or the handheld PC touch screen.



Figure 1. Agilent 4200 FlexScan FTIR

Agilent 4200 FlexScan FTIR Operation Manual

Installing the battery

The 4200 FlexScan FTIR comes with two lithium-ion batteries, which are fully charged. Each battery gives approximately 4 hours of run time. A battery charger is also included; refer to the battery charger manual for information on its operation. There is an LCD display on the front of the battery indicating the level of charge. Five (5) bars indicate a full charge.

The 4200 FlexScan FTIR has a door on the front of the handle. The door on the front gives access to the system battery. The door must remain closed and locked while the system is in use. The 4200 FlexScan should arrive with one battery in the shipping case.

To install the battery:

- 1 Release the battery door tab on the top of the electrical housing (see Figure 1).
- 2 Slide the battery into the compartment with the tab of the battery facing downward and towards the front side of the 4200 FlexScan FTIR. The V-grove on the top of the battery interfaces with the tab on the top of the battery compartment.
- **3** Slide the battery towards the back of the compartment until you feel the battery engage.
- **4** Continue to slide another eighth of an inch until completely engaged. If the battery is not fully engaged, the system will not operate properly.
- **5** Close the battery compartment door.

Depending on temperature and age of the battery, the system should generally operate for a period of four hours on battery power.

To remove the battery from the compartment, open the door and pull on the black plastic tab to slide the battery out of the 4200 FlexScan FTIR housing.



Electrical Shock Hazard Avoid shock hazard. Never allow liquid to run into the 4200 FlexScan FTIR battery compartment.

Power switch LED status

The power button contains a two-colored LED. The color displayed on the LED can indicate the state of the system. The table below lists the colors of the LED and the status of the instrument.

Table 1. Power switch LED status

Status	LED color	Action (% duty cycle)
System OFF	No LED illumination	N/A
System OFF	Red	100%
System startup	Red / green	Green 0.5 s on / red 0.5 s on
System ON	Green	100%
Low battery	Green	50% (0.5 s on / 0.5 s off)
Critically low battery	Red	50% (0.5 s on / 0.5 s off)
Dead battery	No LED illumination	N/A
Firmware update	Red / green	Green blinks rapidly twice then red blinks rapidly twice

CAUTION

The power button is a momentary switch to avoid accidental power off. In order to turn the system on or off, hold the button down for two (2) seconds.

NOTE

The 4200 FlexScan FTIR contains a lithium-ion rechargeable battery pack. Agilent has fully charged the 4200 FlexScan system prior to shipment. The 4200 FlexScan battery life under continuous use is approximately 4 hours.

Starting the software

The 4200 FlexScan FTIR systems are supplied with the Agilent MicroLab Mobile software, which has already been installed on your handheld PC.

To start the software once the 4200 FlexScan FTIR has been powered up:

1 Turn on the handheld PC data station by depressing the power button on the top right corner of the data station (see Figure 2).



Figure 2. Data station showing power button location

2 The system is provided with a shortcut to the MicroLab Mobile software in the Start menu. Upon initial power-up of the handheld PC, look in the computer display window and tap Start. Then tap the MicroLab Mobile icon. This will launch the MicroLab Mobile software application.

NOTE

If the you see this message upon loading your software:

Instrument failed to initialize. Instrument returned error code: -2. The application will exit; then power-cycle the instrument and restart the application.

the spectrometer is not connected to the computer. When using the handheld PC with MicroLab Mobile, verify that the instrument displays a green light on the power button and verify that the 'Bluetooth' communications of the handheld PC are turned 'On'. The instrument must be connected via Bluetooth for the MicroLab Mobile software to launch.

Charging the handheld PC

To connect the handheld PC to an AC power source for charging the battery:

- **1** Connect the USB cable to the handheld PC.
- **2** Connect the computer AC power cable to the 'Y' connector of the USB cable.
- 3 Connect the handheld PC power cable to the AC power source.

Alternatively, if purchased, place the PC in the optional handheld charging cradle. For more information on the operation of the handheld PC, see the supplied Socket User Manual.

Logging on to MicroLab Mobile for the first time

As mentioned above, the 4200 FlexScan FTIR systems are supplied with the Agilent MicroLab Mobile software already installed on your system. (A software disk for the Agilent MicroLab PC software is also included. This allows use of the 4200 FlexScan with a desktop or notebook computer.)

In addition, the handheld PC is provided with a shortcut to the MicroLab Mobile software already installed in the Start menu.

To start the software:

1 Tap **Start** then tap the **MicroLab Mobile** icon.

Upon initial launch of the system, the MicroLab Mobile software automatically loads and a MicroLab application splash screen appears briefly showing software version, copyright and trademark information.

The software will launch and display a logon screen.

NOTE If the 4200 FlexScan FTIR instrument has not been turned on first, you will receive an 'Instrument Failed to Initialize' warning message on the handheld PC. If this happens, tap OK, power-cycle the 4100 ExoScan instrument, and restart the MicroLab Mobile application.

2 For detailed instructions on initial software logon and user management, refer to the MicroLab Software Operation Manual.

For additional instructions on adding a user to the system and other initial software setup procedures, refer to the MicroLab Software Operation Manual.

CAUTION

If a 2.4 GHz cordless phone or a cellular telephone is used near the 4200 FlexScan FTIR, you may experience problems with the wireless communication and spectroscopic performance. Try turning off the phone or moving further away from the instrument and attempt the application again.

Performance verification

The 4200 FlexScan FTIR spectrometer systems have been thoroughly tested at the factory, so no alignment steps are required. However, it is recommended to run the performance test initially to ensure the instrument is running properly. To run and interpret the performance test, refer to the MicroLab Software Operation Manual. The performance test measures the energy level of the instrument (based on the interferogram voltage or height). A successful test will be indicated by a green circle at the top of the software screen. You are now ready to analyze a sample. A yellow or red circle indicates the instrument is operating outside of the factory-defined parameters. These parameters are listed below as well as on the Advanced Features software page.

NOTE

If the circle at the top of the software screen is yellow or red, contact Agilent technical support for assistance.

Like any measurement device, it is important to verify that the 4200 FlexScan FTIR is performing properly before using the system to make critical measurements. The 4200 FlexScan provides both diagnostic values and performance validation tests to demonstrate the system's performance. The diagnostic values provide an easy to understand, quick assessment of the instrument's function. If the instrument is not functioning properly, one of the diagnostic values will be out of specification. Typically, the instrument will not allow data collection if the diagnostic values are out of specification. Performance validation tests are longer, more involved tests that measure how well the instrument is working. The tests listed under 'Performance validation' check the instrument's sensitivity (performance), stability and frequency precision (laser calibration).

Each industry has different requirements for instrument verification. In general, the diagnostic values should be checked on a daily or weekly basis. The instrument will not collect data if the diagnostic values are grossly out of specification, but it is good practice to verify that the instrument is working properly. Performance validation should be run on a quarterly or bi-annual basis. Highly-regulated industries may require the performance validation to be run monthly, depending on the use of the instrument. The performance validation verifies the key aspects to the instrument's ability to measure good data. If the performance validation tests are within specification, the instrument should provide data that works well with methods developed for the 4200 FlexScan. As with any instrument, however, results can always be verified by running a known sample with the sample specific method.

Diagnostic values

The following values can also be viewed on the Diagnostics page in the MicroLab Mobile or MicroLab PC software. These values can be used to determine if the instrument is working properly. They can be used on a daily basis to determine that the instrument is operating as intended. In each case, the Optimal value indicates that the system is running at its intended performance level. The Marginal value indicates that the instrument is still functioning, but at a lower performance level. The Critical value indicates that the system is not working correctly. Contact Agilent technical support for assistance if there is a problem with your instrument.

Getting Started

NOTE

When using the external reflectance or grazing angle sample interface, the appropriate reflectance cap must be in place to make these measurements. The diffuse gold reference cap should be used with the 45° reflectance probe when measuring carbon black or other low reflectance samples. The gold mirror reference cap should be used with the 45° reflectance probe when measuring thin films or highly reflecting samples. The gold mirror reference cap should always be used with the grazing angle sample probe.

Table 2. Diagnostic values

-

Value	Optimal (green)	Marginal (yellow)	Critical (red)	Comment
Energy	29,000–21,000	>29,000 or	> 31,000 or	Indicates the overall
(centerburst)		<21,000	< 18,000	alignment and proper gain
				adjustment for the system.
Battery life	AC or	30–5 min	<5 min	
	>30 min			
Source	1.9 A	>2.2 or	>2.5 or	Indicates a problem with
		<1.6	<1.0	the source control voltage
				or that the source has
				burnt out. Both voltage and
				amperage are displayed;
				however, amperage
				provides a sufficient
				diagnostic.
Laser signal	12,000–4,000	>12,000 or	>15,000 or	Can verify gross alignment
		<4,000	< 3,000	errors even if the
				reflectance cap is not in
D	05 44	05		place.
Detector	35–44	<35	<30	Indicates a problem with
temperature		or >44	or >48	the cooling circuitry or an
				ambient temperature over
	10 75	. 75	. 00	the specified range.
UPU 	10—/5	>/5	>UU<	indicates the ambient
temperature				temperature over the
				specified range.

Performance validation

The MicroLab PC and MicroLab Mobile software have three tests appropriate for performance validation of the 4200 FlexScan FTIR. All of these tests can be accessed from the Advanced Features, System Check page of the MicroLab PC and MicroLab Mobile software. The 4200 FlexScan system should be warmed up for at least 30 minutes before conducting any of these tests. These tests can be run on a quarterly, bi-annual or annual basis to determine that the instrument is performing within its specification.

Performance (signal-to-noise) test

This test measures the signal-to-noise level at two regions of the IR spectrum: 2500 cm⁻¹ and 1000 cm⁻¹. The signal-to-noise is defined as the reciprocal of the root mean squared (RMS) noise in the defined region for a blank sample measured with a blank background. Both the background and sample are measured at 4 cm⁻¹ resolution with a one minute collection time. This test takes 2 minutes per test. The user can specify a number of tests to be conducted. For 45 degree external reflectance and grazing angle reflectance measurements, the appropriate reflective reference cap should be in place for the entire test. No reference cap is needed when using the ATR sample interface. At least 10 tests should be collected in order to obtain an accurate picture of the performance. The table below gives the minimum performance for each sample interface.

Sample interface	Specification @ 2500 cm ⁻¹	Specification @ 1000 cm ⁻¹
ATR	2500	6000
External reflectance	1500	2800
Grazing angle	1500	2800

 Table 3.
 Minimum performance for the Signal-to-Noise test

Stability test

This test measures the short-term stability at two regions of the spectrum: 3000 cm⁻¹ and 1000 cm⁻¹. Stability is a measurement of the baseline differences observed over the selected time period. The test measures a background at the beginning, then one sample every minute for the duration of the test as specified in the 'Number of minutes' field in the software. For 45 degree external reflectance and grazing angle reflectance measurements, the appropriate reflectance reference cap should be in place for the entire test. No reference cap is needed when using the ATR sample interface. The test results are expressed as %Transmittance (difference versus 100%) of maximum deviation during the Stability test. The table below gives the minimum performance for this test.

Table 4. Minimum performance for the Stability test

Time	Specification @ 3000 cm ⁻¹	Specification @ 1000 cm ⁻¹
30 min	2%	1%

Laser Frequency Calibration test

The Laser Frequency Calibration test measures the frequency (Xaxis) accuracy. The test is conducted by measuring a spectrum of a polystyrene film. Absorbance frequencies from this spectrum are compared to frequencies set by NIST SRM 1921 polystyrene film. The software enables administrative-level users to use the test results to set the laser calibration. Laser calibrations should only be set after consultation with an Agilent technical support engineer. For this test, a background is measured first. For 45 degree external reflectance and grazing angle reflectance sample interfaces, the reflective reference cap is required for background measurements. After the background collection, a spectrum of polystyrene film is measured as instructed by the software. When using the 45 degree external reflectance and grazing angle reflectance interfaces, the polystyrene sample should be placed between the reference cap and the sample interface. For the ATR sample interface, the reference cap is not required for a background, but the polystyrene test sample must be pressed firmly against the diamond sample interface during the sample collection. The specification is listed below for each sample interface. The difference between any of the listed frequencies and the standard value should be no greater than the specification.

Sample interface	Specification
ATR	<2 cm ⁻¹
External reflectance	<1 cm ⁻¹
Grazing angle	<1 cm ⁻¹

 Table 5.
 Minimum performance for the Laser Frequency Calibration test

Getting Started

This page is intentionally left blank.



3. Analyzing Samples

External reflectance, grazing angle and diffuse 23 Internal reflectance (ATR) 27

The Agilent 4200 FlexScan FTIR instrument is provided with a sample probe interface as specified on your order. This section covers instructions for how to analyze a sample with all sampling probes offered on the 4200 FlexScan.

External reflectance, grazing angle and diffuse reflectance

The 4200 FlexScan FTIR instrument can be provided with three different reflectance interfaces: external (or specular) reflectance, grazing angle reflectance or diffuse reflectance. Each of these sample interfaces relies on light reflecting from the sample surface. When coming in contact with a sample, the infrared (IR) light passes through the top surface of the sample before being reflected back out of the sample and into the detector of the 4200 FlexScan system. This small and consistent pathlength provides the advantage of no sample preparation required to get good measurement results on a variety of samples. The key to obtaining good results with external reflectance is making sure the tip of the probe makes stable contact with the sample surface and does not move during the measurement process. The front of the sample interface defines the IR light focus or each of the reflectance sample interfaces.

The 45 degree probe is optimal for analysis of thin films greater than 1 micron in thickness. The grazing angle probe is optimal for analysis of ultra thin films less than 1 micron in thickness. Diffuse reflectance is specifically designed for measurement of rough surfaces that scatter rather than reflect light. Specular reflectance is an ideal method for the analysis of surface contamination. The grazing angle probe should be used when very low level contamination needs to be analyzed (nanometer thickness levels).

Each reflectance sample interface requires the use of a reference sample for taking the background measurement prior to analyzing the sample. The reference sample is incorporated into the 4200 FlexScan FTIR by means of a metal disk mounted into a cap that snaps onto the tip of the reflectance probe. The 4200 FlexScan system offers two reference caps designed for this purpose – a diffuse gold reference cap with a rough surface finish and a stainless steel mirror reference cap with a very smooth surface finish.

The diffuse gold reference cap should be used with the 45 degree reflectance probe when measuring low-reflectance samples. The mirror reference cap should be used with the 45 degree reflectance probe when measuring thin films or highly-reflecting samples. The mirror reference cap should always be used with the grazing angle sample probe. The diffuse cap should always be used with the diffuse reflectance sample interface.

CAUTION

Ensure that the sample makes solid contact with the entire surface area of the external reflectance probe tip and is not held at an angle against the sample, and that no movement occurs between the sample and the probe during the measurement.

Reflectance standards and gain settings

The 4200 FlexScan FTIR is designed to measure a variety of sample types. In order to measure the best quality spectra, it is important to use the correct reflectance standard and to appropriately set the detector gain. Detector gain is an electronic setting that maximizes the detector response from the available light. The detector gain must be set such that both the reference material and the sample are within the linear range of the detector.

The response of the detector is reported on the MicroLab PC and MicroLab Mobile software on the Diagnostics page, available from the Advanced Features page; it is labeled Energy. The energy value has a range from 1000 to 30,000.

First, select the reference standard that is most similar to the sample; the diffuse standard should be used for samples that scatter light, while the mirror standard should be used for samples that reflect light. With the appropriate reference in place, observe the energy value. The value should be close to 26,000 with the reference standard in place. If the energy value is below 23,000 or above 29,000 counts, click the Gain Adjust button.

The Gain Adjustment screen shows the Energy value at the top and has up/down arrows to adjust the gain. With the reference in place, adjust the gain to bring the energy to 26,000 counts. Once the appropriate gain is set, it can either be saved as the System Default or Method gain. The System Default gain is used for all methods, while the Method gain is used only for the currently selected method.

Once the gain has been set, it is useful to check the energy from a sample. The sample must have energy greater than 1000. Also, the sample should have energy less than that of the reference.

Steps

The steps involved in performing an external reflectance analysis are:

- **1** Collect a background spectrum using the reflectance cap.
- 2 Collect a sample spectrum.

Collecting a background spectrum

To ensure accuracy of the measurement, it is recommended that the system is configured in the software to collect a background spectrum before every sample is analyzed. This provides a baseline profile of the system conditions using the gold reference cap. By collecting a background automatically before each sample measurement, negative effects of changes in the environment can be avoided.

To collect a background spectrum:

- 1 From the software, choose a method and click **Start** to begin the measurement.
- 2 The software will collect a background automatically if the method selected was set up as recommended in the paragraph above. The software may prompt you to install the reference cap. Install the appropriate gold reflectance reference cap tightly against the tip of the sample probe. The reference cap has a black nylon strap with Velcro that can be used for convenience and security. To install, slide the nylon strap into the loop located on the front of the 4200 FlexScan FTIR near the sample interface. Overlap the strap around the loop and secure using the Velcro at the end of the strap.
- **3** For detailed software instructions for collecting a background, refer to the MicroLab Software Operation Manual.

Collecting a sample spectrum

To collect a sample spectrum:

- 1 Once the background is finished collecting, the software will instruct you to put the probe in contact with the sample.
- **NOTE** It is recommended to use the provided strap to secure the 4200 FlexScan FTIR to the user and avoid damage due to dropping the system. To install the strap, hook both ends (2 spring clips) of strap to the rod on the lower back panel of 4200 FlexScan (just above the handle). Place the strap around the shoulder or neck.
 - 2 The software will now display a graph showing the intensity (absorbance) of the sample measurement. This screen is used to ensure that the sample has been correctly placed at the sampling interface. For 4200 FlexScan FTIR applications using reflectance, the value should be between 0.01 and 2 AU. Use the real-time data viewer in the software to ensure good contact between the sample and the probe, and then choose to collect the data using the 4200 FlexScan FTIR software.

- **3** Click the **Next** button on the software screen to proceed with the analysis. Clicking the Home button will stop the measurement and return you to the Home screen.
- **4** For detailed software instructions for measuring a sample, refer to the MicroLab Software Operation Manual.
- **5** After completing the sample measurement, immediately clean the sample from the accessory using the instructions provided above.
- **6** For detailed software instructions on reviewing results and handling sample data, refer to the MicroLab Software Operation Manual.

Internal reflectance (ATR)

The 4200 FlexScan FTIR instrument can be provided with a dedicated sampling probe interface called the attenuated total reflectance (ATR) probe. The ATR takes advantage of the physical properties of light when encountering two materials with differences in index of refraction. When coming in contact with a sample having a different index of refraction, the infrared (IR) light creates an evanescent wave that provides a very small and specific depth of penetration into the sample before reflecting back into the detector of the 4200 FlexScan FTIR system. This small and consistent pathlength provides the advantage of no sample preparation required to get good measurement results on a variety of samples. The key to obtaining good results with an ATR is making good contact between the sample and ATR crystal of the sampling device. The ATR technique can be used for analysis of liquids, pastes, powders and even some solid samples.

The Agilent ATR uses a type IIa synthetic diamond crystal as the interface between the sample and the infrared energy. The diamond provides the advantage of extreme hardness as well as chemical resistivity. Diamond can accept samples with a pH range from 1 to 14, which means hard or abrasive samples and even strong acids can be safely analyzed. The Agilent 4200 FlexScan FTIR ATR sampling probe is a single reflection system. The diamond is the world's most durable substance and the ZnSe monolithic substrate is used as a focusing lens to provide the maximum IR signal.

The single reflection ATR is most suitable for higher absorbing samples such as rubber, polymers, paints and fibers. Powder and solid samples are best measured on the single reflection ATR as well due to the optional sample press device that can apply high pressure to powder and solid samples to ensure good contact with the diamond sampling surface. The single reflection ATR is also a good choice when the amount of available sample is limited. The single reflection diamond is 2 mm in diameter with a 200 μ m active area and provides an approximate 2 micron depth of penetration for IR energy at 1700 cm⁻¹. The single reflection ATR element protrudes slightly above the probe's metal mounting plate.

CAUTION

Although diamond is a very hard material, the ATR window is relatively thin and can crack under extreme pressure conditions. Make sure the sample makes contact with the entire surface area of the diamond and not just a point within the diamond. Avoid the use of the sample press on samples that may be sharp or pointed.

Steps

The steps involved in performing an ATR analysis are:

- 1 Clean the ATR sample mounting window.
- 2 Collect a background spectrum.
- **3** Collect a sample spectrum.

Cleaning the sample probe

Before using the 4200 FlexScan FTIR ATR system on a new sample, it is important as a first step to ensure the sampling surface is clean. Kimwipes or a suitable substitute (such as cotton swabs) should be used to clean the sensor and sampling area. Clean the sample interface with a suitable solvent such as acetone, methanol, ethanol or isopropyl alcohol. Visually inspect the probe tip after cleaning to make sure there is no film, haze or other residue present. CAUTION

Do not break the spectrometer seal and attempt to clean interior surfaces. Breaking the seal will void the warranty.

Collecting a background spectrum

To ensure accuracy of the measurement, it is recommended that the system is configured in the software to collect a background spectrum before every sample is analyzed. This provides a baseline profile of the system conditions with no sample loaded on the instrument. By collecting a background automatically before each sample measurement, negative effects of changes in the environment can be avoided.

To collect a background spectrum:

1 From the software, choose a method and click **Start** to begin the measurement.

The software will collect a background automatically if the method selected was set up as recommended in the paragraph above.

The diamond should not be in contact with the sample or any substance other than the air in the atmosphere during the background measurement.

- **2** For detailed software instructions for collecting a background, refer to the MicroLab Software Operation Manual.
- **CAUTION** To ensure an accurate background spectrum is collected, perform a visual inspection of the diamond ATR sample mounting surface, looking for any haze or film present on the diamond from measuring the previous sample. If a film is observed, repeat the cleaning procedure above until the diamond surface is clear of any residue.

Collecting a sample spectrum

To collect a sample spectrum using the ATR probe:

1 Place a small amount of material to be measured against the diamond ATR crystal. The crystal is the clear, circular shaped material held in place by the surrounding metal disk at the end of the probe.

For liquid samples, simply immerse the tip of the sample interface in the sample. Small samples can be measured by placing a drop of sample on the center of the diamond interface.

Solid samples require establishing good contact between the diamond sampling interface and the sample. Press the diamond firmly into the sample. Tilt the instrument to ensure that the diamond surface is flat to the sample.

In the case of a volatile sample, larger amounts of sample can be applied without concern of leakage or damage to the instrument. However, using the smallest amount of sample possible will ease in the cleaning process.

In all cases, please be sure that the sample covers the entire surface area of the diamond crystal.

If the sample is a powder or solid, the optional sample press may be needed to ensure good contact with the sample.

- 2 Use the real-time data viewer in the software to ensure good contact between the sample and the diamond, and then choose to collect the data using the 4200 FlexScan FTIR software.
- **3** Click the **NEXT** button on the software screen to proceed with the analysis.
- **4** For detailed software instructions for measuring a sample, refer to the MicroLab Software Operation Manual.
- 5 After completing the sample measurement, immediately clean the sample from the accessory using the instructions provided above. It is important to make sure the sample mounting window is free of any residue from the previous sample.

- **6** For detailed software instructions on reviewing results and handling sample data, refer to the MicroLab Software Operation Manual.
- CAUTION The sample mounting window and accessory window are made from type IIa synthetic diamond, which is extremely chemically-resistant. However, it is still possible the diamond or metal mounting plate can be damaged by extreme samples. Only measure samples with a pH between 1 and 14. Do not leave extremely acidic samples on the metal mounting plate for an extended time.

CAUTION

Although diamond is a very hard material, the ATR window is relatively thin and can crack under extreme pressure conditions. Ensure that the sample makes contact with the entire surface area of the diamond and not just a point within the diamond. Avoid using the sample press on samples that may be sharp or pointed.

Using the solid sample ATR press

The 4200 FlexScan FTIR ATR system is offered with an optional sample press attachment that is removable when not required. The sample press enables improved measurement of solid or powder samples. The press ensures that the sample is in good contact with the diamond ATR sample interface, which ensures the best possible results for these sample types. The press is designed to deliver 15 pounds of force to the sample or over 2500 pounds per square inch of pressure. It is not for use with liquids, gels or pastes, as the press may come directly in contact with the IR energy and alter the results. These types of samples make good contact without the use of a press.

Installing

For initial installation of the ATR sample press:

- 1 Clean any remaining sample residue off the diamond sample interface and the stainless steel sampling tube.
- **2** Hold the 4200 FlexScan FTIR with the diamond sample interface pointing up.
- **3** Open the press arm.
- **4** Slide the ring of the press onto the stainless steel sampling tube.
- 5 Rotate the press in the clockwise direction while pressing down until the ring locks in place. The press arm will be located at 12:00, 4:00 or 8:00. There should be an audible click when the ring locks into place.
- **CAUTION** Installation of the solid sample press with the press arm in the downward position can damage the diamond sample interface. Always install the press in the open position.

Removing

To remove the ATR sample press:

- 1 Hold the 4200 FlexScan FTIR with the diamond sample interface pointing up.
- **2** Open the sample press arm so that the press tip is not in contact with the diamond sample surface.
- **3** Twist the press ring in a counterclockwise direction approximately one-quarter of a full rotation (¹/₄ turn).
- **4** Slide the press off the stainless steel sampling tube.

Measuring a sample

To collect a sample spectrum using the ATR press:

- 1 From the software, choose a method and click **Start** to begin the measurement.
- **2** Once the background is finished collecting, the software will instruct you to put the diamond in contact with the sample.

Place a small amount of material to be measured against the diamond ATR crystal. The crystal is the clear, circular shaped material held in place by the surrounding metal disk at the end of the probe.

For liquid samples, simply immerse the tip of the sample interface in the sample. Small samples can be measured by placing a drop of sample on the center of the diamond interface.

Solid samples require establishing good contact between the diamond sampling interface and the sample. Firmly press the diamond into the sample. Tilt the instrument to ensure that the diamond surface is flat to the sample.

In the case of a volatile sample, larger amounts of sample can be applied without concern of leakage or damage to the instrument. However, using the smallest amount of sample possible will ease in the cleaning process.

In all cases, ensure that the sample covers the entire surface area of the diamond crystal.

- **3** Use the real-time data viewer in the software to ensure good contact between the sample and the diamond, and then choose to collect the data using the 4200 FlexScan FTIR software.
- 4 Click the **NEXT** button on the software screen to proceed with the analysis.
- **5** For detailed software instructions for measuring a sample, refer to the MicroLab Software Operation Manual.
- 6 After completing the sample measurement, immediately clean the sample from the accessory using the instructions provided above. It is important to make sure the sample mounting window is free of any residue from the previous sample.
- 7 For detailed software instructions on reviewing results and handling sample data, refer to the MicroLab Software Operation Manual.

Analyzing Samples

This page is intentionally left blank.



When the outside of the Agilent 4200 FlexScan FTIR system needs cleaning, disconnect the USB cable and ensure that the battery compartment is closed. Use a damp, soft cloth and a mild detergent to wipe down the outside of the unit. Do not completely immerse the 4200 FlexScan in water. Although the main optical housing is sealed, water can penetrate the electrical compartment.

WARNING

Hazard



User servicing of parts and components beyond the procedures outlined in this manual may pose a safety risk to you, damage your instrument, and invalidate the instrument warranty. Perform only those procedures described in this manual. For all other technical problems, contact your authorized Agilent representative.

WARNING Hazard



The 4200 FlexScan FTIR was designed to operate in remote locations. However, analyzing certain materials and some maintenance procedures may be hazardous if the following guidelines are not practiced. Maintenance

This page is intentionally left blank.



Part number	Description
0024 - 001	Agilent 4200 FlexScan FTIR system with external reflectance
0024 - 002	Agilent 4200 FlexScan FTIR system with diffuse reflectance
0024 - 003	Agilent 4200 FlexScan FTIR system with grazing angle
0024 - 004	Agilent 4200 FlexScan FTIR system with diamond ATR
0023 - 300 -	Diffuse gold reflectance reference cap for diffuse reflectance
0023-301	Polystyrene frequency reference cap for diffuse reflectance
0023 - 304	Diffuse gold reflectance reference cap for external reflectance
0023 - 305	Polystyrene frequency reference cap for external reflectance
0023 - 307	Mirrored stainless steel reflectance reference cap for external reflectance
0023 - 520	ATR sample press
800 - 2002	Handheld PC (PDA)
800 - 2003	Handheld PC docking station (includes USB and power cables)
800 - 2004	Replacement handheld PC battery
800 - 2005	Replacement lithium-ion battery
430 - 0018	USB cable

For sampling accessories, upgrades, service contracts, repair and refurbishment services, contact Agilent.

Spare Parts

This page is intentionally left blank.



Technical

- **Interferometer geometry:** High-throughput Michelson interferometer with fixed and moving flat mirrors
- Standard beamsplitter: Zinc selenide
- Maximum spectral resolution: 4 cm⁻¹
- Laser: Low-powered solid state
- Source: Wire-wound element
- Spectral range: 4000 to 650 cm⁻¹
- **Sample interface:** External reflectance (45°), grazing angle (82°), diffuse reflectance (normal incidence), or single reflection diamond internal reflectance.
- Detector: 1.3 mm diameter, thermoelectrically-cooled dTGS
- Optical compartment: 5.5 x 4.75 x 3.9 in (14 x 12.1 x 9.9 cm);
 3.2 lb (1.45 kg)
- Electrical compartment: 7.5 x 4 x 2.5 in (19 x 10.2 x 6.4 cm); 2.5 lb (1.16 kg)
- Cord: 6.5 ft (1.98 m) extended, 20 in (7.8 cm) coiled;
 2.1 lb (0.95 kg)
- **Battery:** Li-ion, 10.8 V; 4400 mAh

Specifications

Environmental conditions

- **Temperature:** 0 to 50 °C; 32 to 122 °F (during use)
- **Temperature:** Minus 30 to 60 °C; -22 to 140 °F (storage)
- **Humidity:** Up to 90%, non-condensing

Handheld PC data station

- **Processor:** 520 MHz Intel[®] PXA270
- Memory: 64 MB low power RAM
- Hard drive: 128 MB NAND flash