

Agilent 6100 Series Quadrupole LC/MS Systems

Quick Start Guide

Where to find more information	1
Configuration Diagrams	4
ChemStation Views	6
Basic Operation	20
Operating Tips	27
Troubleshooting Tips	29

This manual applies to ChemStation Software (version B.04.02 SP1 and later until superseded) and Agilent 6100 Series LC/MS instruments.

Where to find more information

Online Help Use online Help for in-depth information beyond what is shown in this *Quick Start Guide*. Display online Help in any of the following ways:

- Click the **Help** button in the toolbar.
- Select **Help Topics** from the **Help** menu.
- Click the **Help** button on most dialog boxes to show task-specific help.



Agilent Technologies

User Guides Besides this Quick Start Guide, the following guides are shipped with your system and are also available in PDF format on the installation disk.

- *Agilent 6100 Series Quadrupole LC/MS Systems Familiarization Guide*
- *Agilent 6100 Series Quadrupole LC/MS Systems Concepts Guide*
- *Agilent 6100 Series Quadrupole LC/MS Systems Installation Guide*
- *Agilent 6100 Series Quadrupole LC/MS Systems Maintenance Guide*

What's New in B.04.02 SP1?

- Support for new G6120B, G6130B, and G6150B Single Quadrupole LC/MS systems.
- Support for Agilent Jet Stream Technology on the 6150.
- Support for Microsoft Windows XP Professional 32-Bit with Service Pack 3.
- Support for Microsoft Windows Vista Business 32-Bit with Service Pack 2.
- Easy Sequence, a new interface to plan, generate and queue sequences.

What's New in B.04.02?

New ChemStation Software Features


- Easy Sequence: A new easy-to-use user interface for sequence planning, generation and queueing.
- Full support for all modules in the Agilent 1200 LC series, including the Agilent 1200 Rapid Resolution SL LC System and the new Agilent 1290 Infinity LC System, which consists of the following hardware modules.
 - Agilent G4220A 1290 Binary Pump
 - Agilent G4226A 1290 High Performance Autosampler
 - Agilent G1316C 1290 Thermostatted Column Compartment
 - Agilent G4212A 1290 DAD
 - Agilent G4227A Infinity FlexCube
- A new user interface for LC instrument control in the Method and Run Control view with:

**For Agilent 1100
/1200 LC
Systems users**


- Flexible sizing of the Instrument Control panel to make best use of the available desktop space.
- Show/hide functionality to reduce the display to the most essential information.
- Graphical representation of multiple devices of the same kind (for example, two pumps).
- Graphical representation of all valves.
- Direct access to method parameters and all important functions for each module.
- Status display for the individual modules and resulting overall instrument status.
- Tooltips display most important information on specific functions.
- Tabular display of the sample information for the current run or sequence line.
- Direct access to Data Analysis parameters and settings for OpenLAB ECM (with G2189BA ChemStation OpenLAB option only).

Configuration Diagrams

1100 or 1200 Series LC

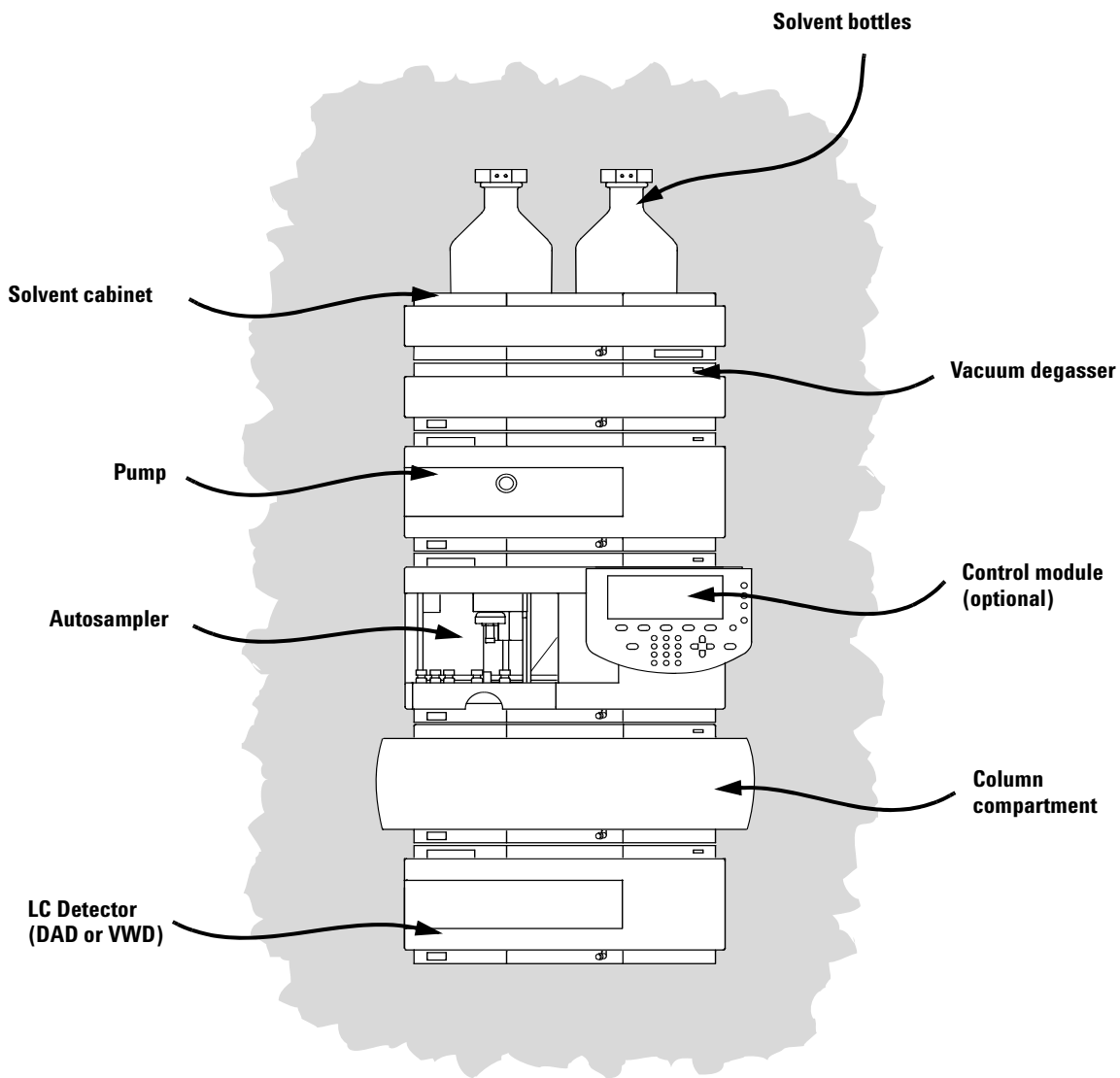


LC Cabling

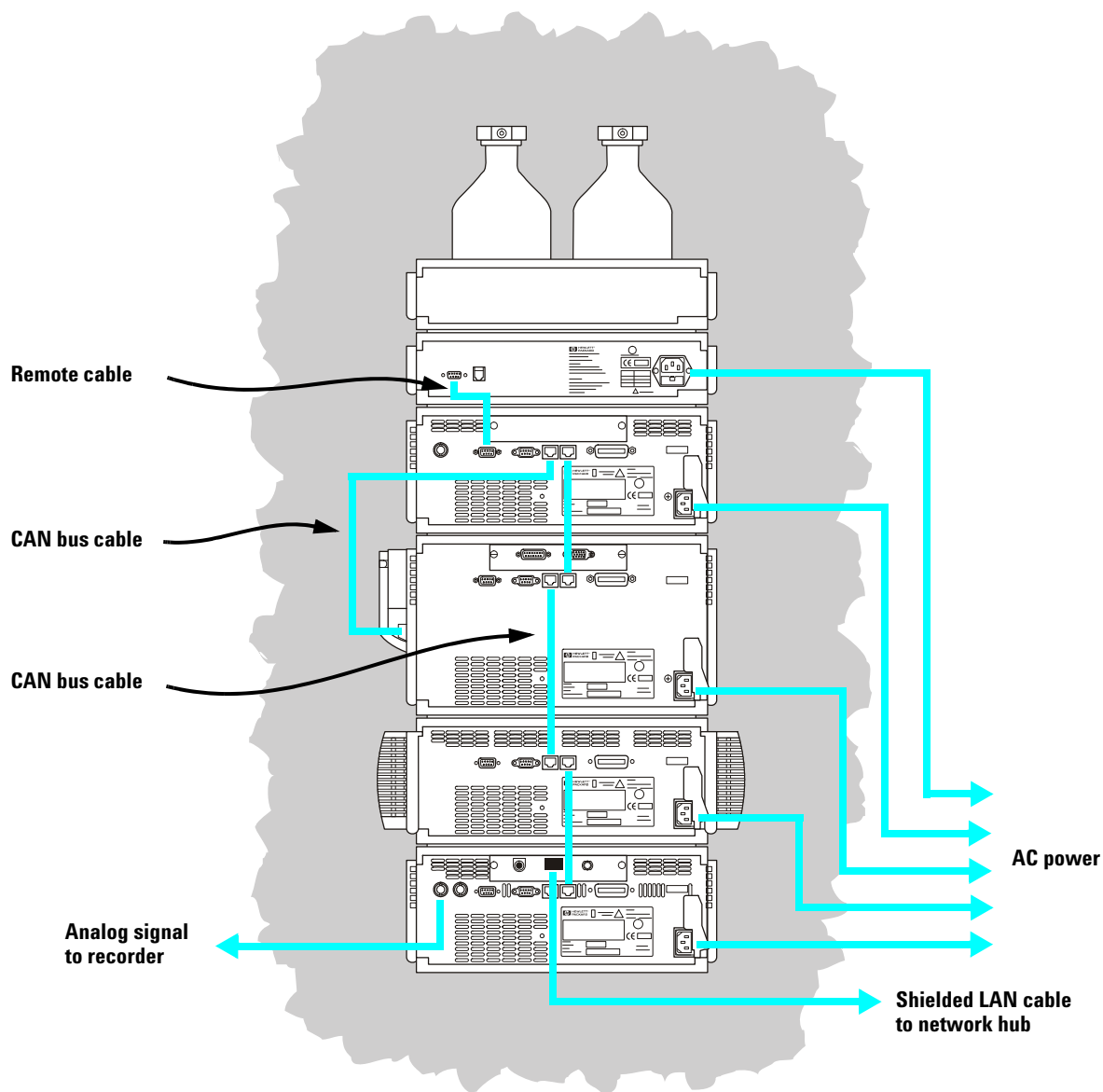

ChemStation Views

The LC/MS Quad ChemStation software consists of several views, which are described below. Switch to a different view by clicking one of the view buttons in the lower left corner of the screen (shown below) or from the **View** menu.

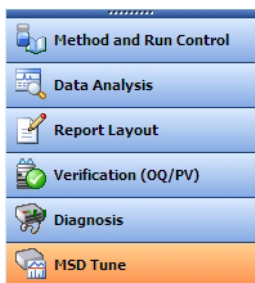

Method and Run Control View Allows you to set up methods and adjust instrument parameters to inject samples and acquire data one sample at a time or in automated sequences. See the next page for more information.

Data Analysis View Allows you to perform various data evaluation tasks on chromatograms and spectra, such as integration, quantitation, checking peak purity, deconvolution, and report generation. See [page 11](#) for more information.


Report Layout View Allows you to design custom report templates to use with the data generated by your ChemStation. See [page 14](#) for more information.

Verification (OQ/PV) View For use by Agilent service engineers to determine if your system is operating in a predictable manner. This is useful to show Good Laboratory Practice (GLP) compliance and may be required by some government agencies. See [page 16](#) for more information.

Diagnosis View Allows you to run tests to diagnose instrument problems and access information on how to resolve these problems. The early maintenance feedback (EMF) feature can notify you when it is time to perform system maintenance before a problem occurs. See [page 18](#) for more information.

MSD Tune View In this view, you can calibrate your LC/MS quad automatically. You can also set MS parameters manually for specific types of molecules. See [page 19](#) for more information.

Method and Run Control View

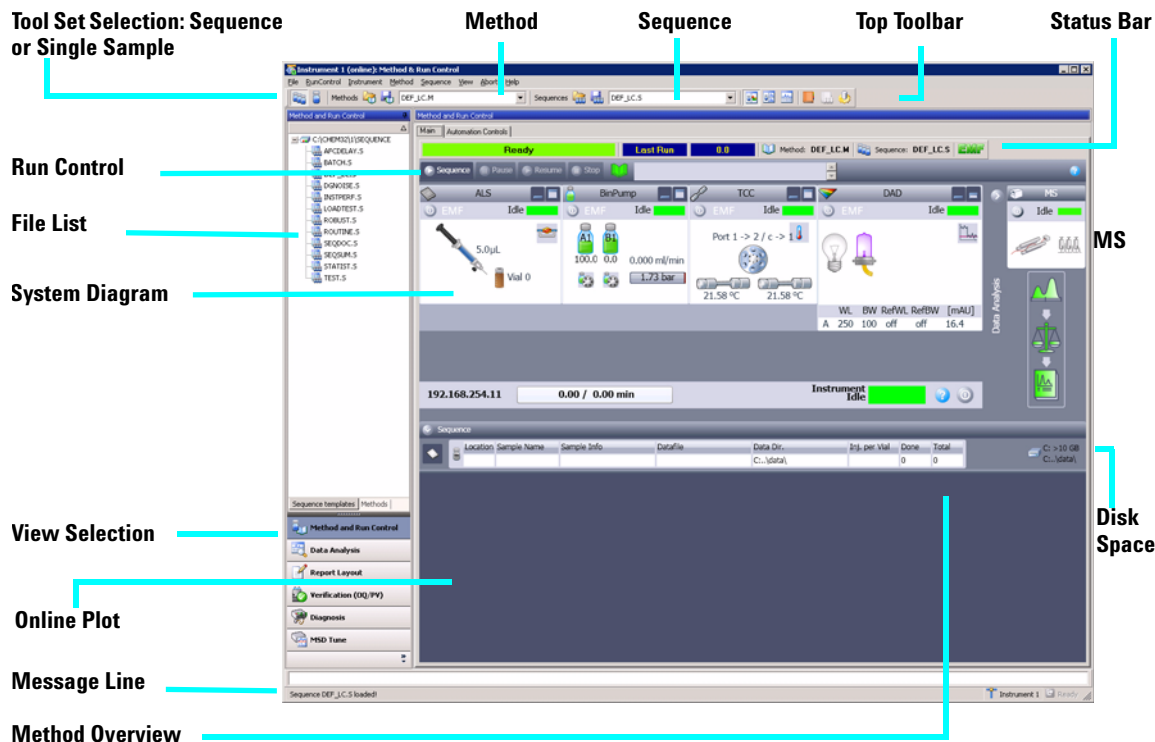

NOTE


The user interface for Method and Run Control can be configured in two ways during installation. See *Agilent 6100 Series Single Quad LC/MS System Installation Guide* for information on installation and configuration.

If your instrument is configured with the Classic user interface, then see ChemStation online Help and the following manuals on the ChemStation disk for information and instructions using the Classic interface:

- *Agilent 6100 Series Single Quad LC/MS Systems Quick Start Guide*
- *Agilent 6100 Series Quadrupole LC/MS Systems Familiarization Guide*
- *Agilent 6100 Series Quadrupole LC/MS Systems Concepts Guide*

Status and Run Bars: Main Tab

Right-click the Method icon to see the following commands: Run Time Checklist, Method Information, Edit Entire Method, Method Audit Trail, Print Method, and Help.

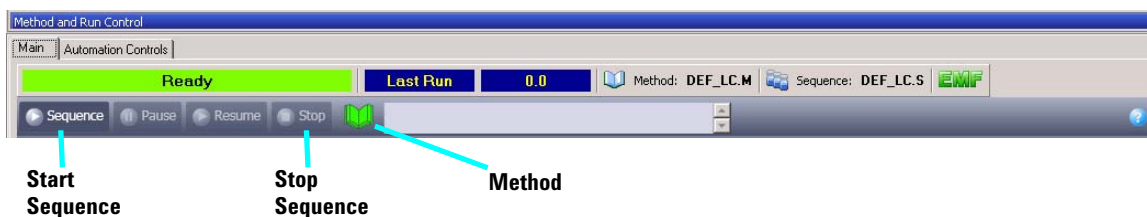

System Diagram Each icon of the system diagram represents one component or module of your system. Click on an icon if you want to edit the method parameters or go to the online Help for that particular component.

Injector**Pump & Solvents****Column****Detectors**

**Integration
Calibration
Reporting**

Single Sample Tool Set

This toolset allows you to work on methods and run an analysis for a single sample. Move the cursor over a button to view a description of it.



Click the Single Sample icon to display this toolset.



Sequence Tool Set

This toolset allows you to work on sequences and run automated analyses of multiple samples. Move the cursor over a button to view a description of it.



Click the Sequence icon to display this toolset.




Flow Injection Analysis (FIA)

This toolset allows you to inject multiple samples directly into the detector, bypassing the chromatographic column. The results are sent to a single data file. FIA can be used for method development or for applications that do not require chromatography.



To enable FIA from the system diagram:

- Click the FIA icon.
- Select Edit FIA Series.
- Mark the Enable FIA Series in Method option and edit the FIA Table.

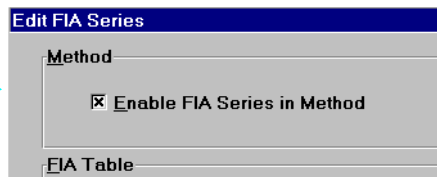

Data Analysis View

Navigation Tools **Cursor Tools**

File List

Tool Set Selection

Signal View Selection

Display Graphics Tools

DAD Signal

LC/MS quad Signal

Message Line

The various tool sets available in Data Analysis View are shown on the following pages. You can view a description of a button by moving the cursor over it.

Common Tool Set This toolset is always present in Data Analysis view.




Graphics Tool Set Allows you to manipulate the graphic display.



Integration Tool Set Allows you to perform integration and reporting tasks on a chromatogram.



Calibration Tool Set Allows you to perform calibration tasks for quantitation.

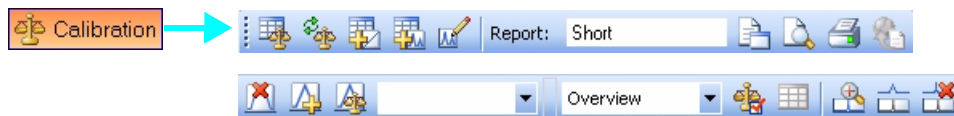


Signal Tool Set Allows you to graphically work with the UV or MS signal.




Purify Tool Set Allows you to graphically work with your purification data.




Spectral Tool Set Allows you to perform spectral evaluation tasks.

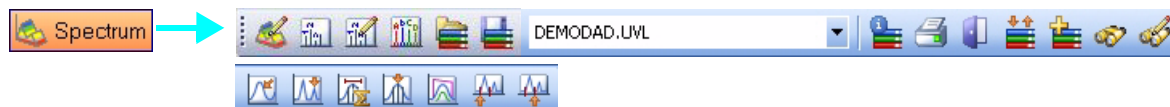

Deconvolution Tool Set Allows you to transform mass spectra from multiply-charged ions into a calculated molecular weight.

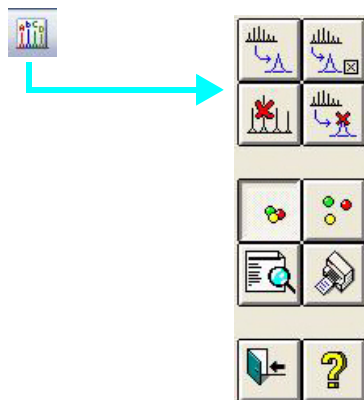

Report Layout View

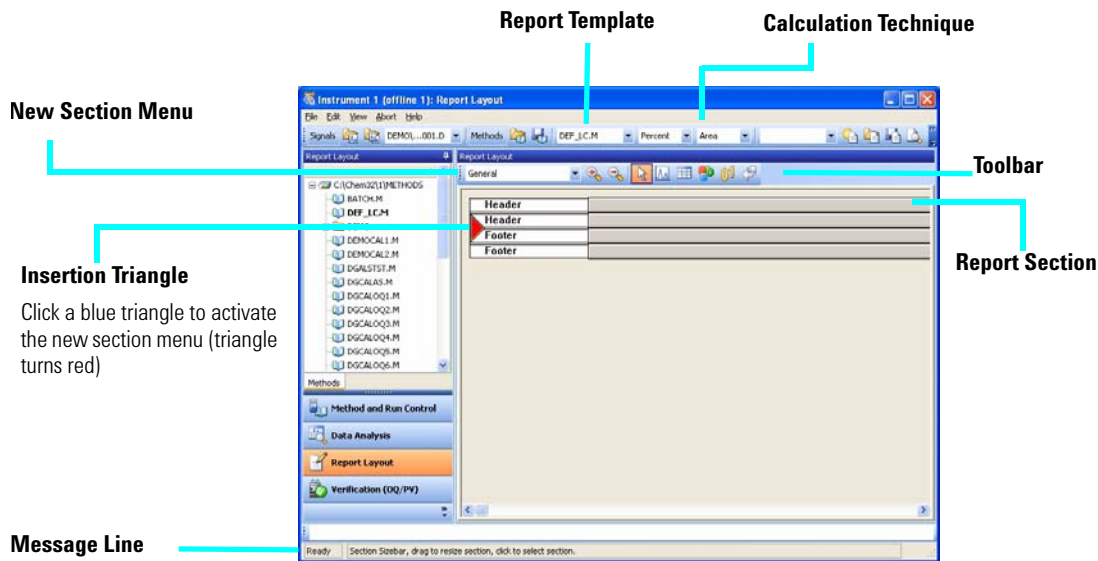

Report Layout

Tool Set This toolset is present in the Report Layout view. Move the cursor over a button to view a description of it.




Tips

- Edit or create a report layout as described in the online Help.
- To test your report layout, select a calculation technique from the drop-down list box to define how the results should be calculated.
- Load a data file. The results are loaded into the report template using the selected calculation.
- Select **File / Add to Report Styles** to add your completed report template to the list of available report styles. Now you can use your customized report within a method.


Sample Report Layouts (in \CHEM32\REPSTYLE)

Template	Description
SIMPLE.FRP	Includes a chromatogram and calibration curves per calibrated compounds.
NESTED.FRP	Shows how to nest subsections.
SPECTRA.FRP	Shows how to include spectra.
PURITY.FRP	Shows how to include peak purity data.
ESTD.FRP	Uses the same elements as the standard ESTD reports so you can customize your external standard reports.
AREAPCT.FRP	Uses the same elements as the standard area percent report so you can customize your area percent reports.
LIBRARY.FRP	A customized report that uses library search results. You must have a method loaded that specifies the spectra library.

Print Preview button

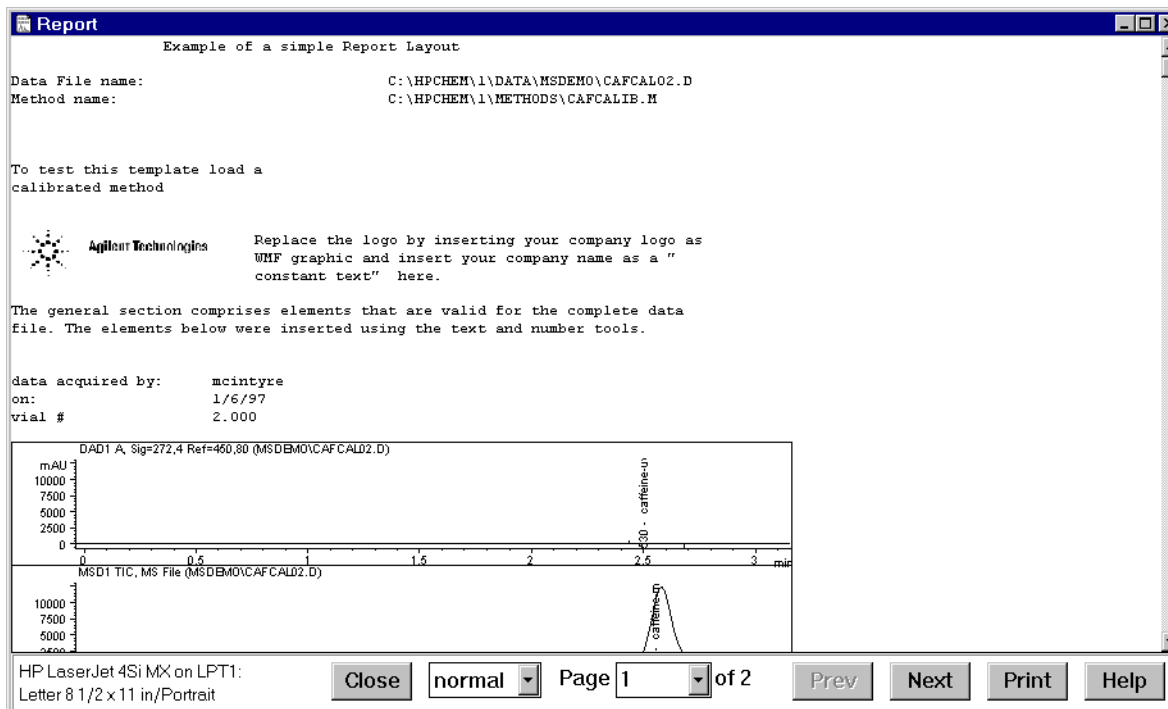

Verification (OQ/PV) View

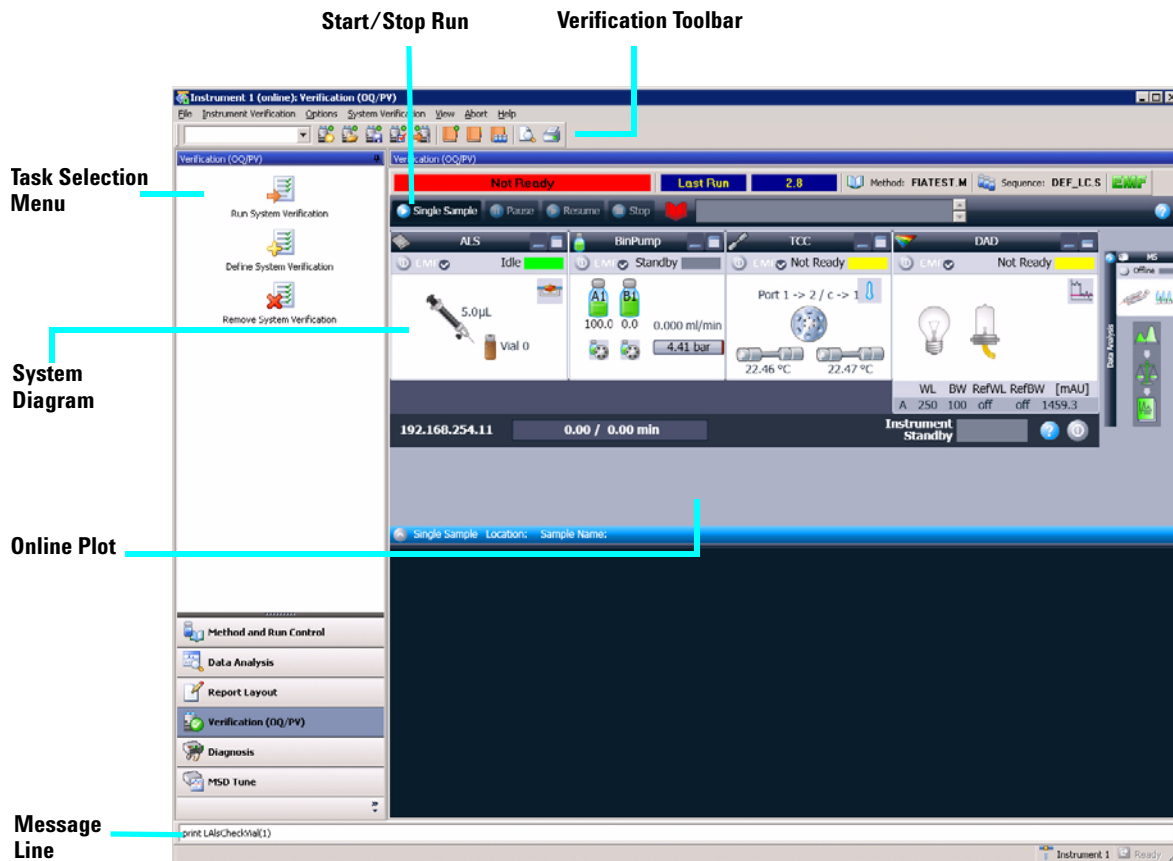

The Verification (OQ/PV) view lets Agilent service engineers test whether your analytical instruments and the ChemStation software are operating correctly according to predefined performance criteria.

Operation Qualification (OQ) Operation Qualification is the documented verification that the equipment-related system or subsystem performs as intended throughout representative or anticipated operating ranges.

Performance Verification (PV) Performance Verification is the documented verification that the process and or the total process-related system performs as intended throughout all anticipated operating ranges.

System Diagram Each icon of the system diagram represents one component or module of your system.

Verification Toolbar The Verification toolbar is displayed when you select **Show Top Toolbar** from the **View** menu. Move the cursor over a button to view a description of it.




Available Tests The following verification tests are supplied with your ChemStation software for use by an Agilent service engineer. Refer to the online Help for more information on these tests.

- VWD Wavelength Accuracy
- Intensity
- Holmium
- Temperature Accuracy
- Noise, Flow, Temperature
- DAD Wavelength Accuracy
- Injector Precision*
- Detector Linearity/Carry-over*
- Injector Linearity*
- Gradient Composition


* These tests can be used to verify LC/MS quad performance.

Diagnosis View

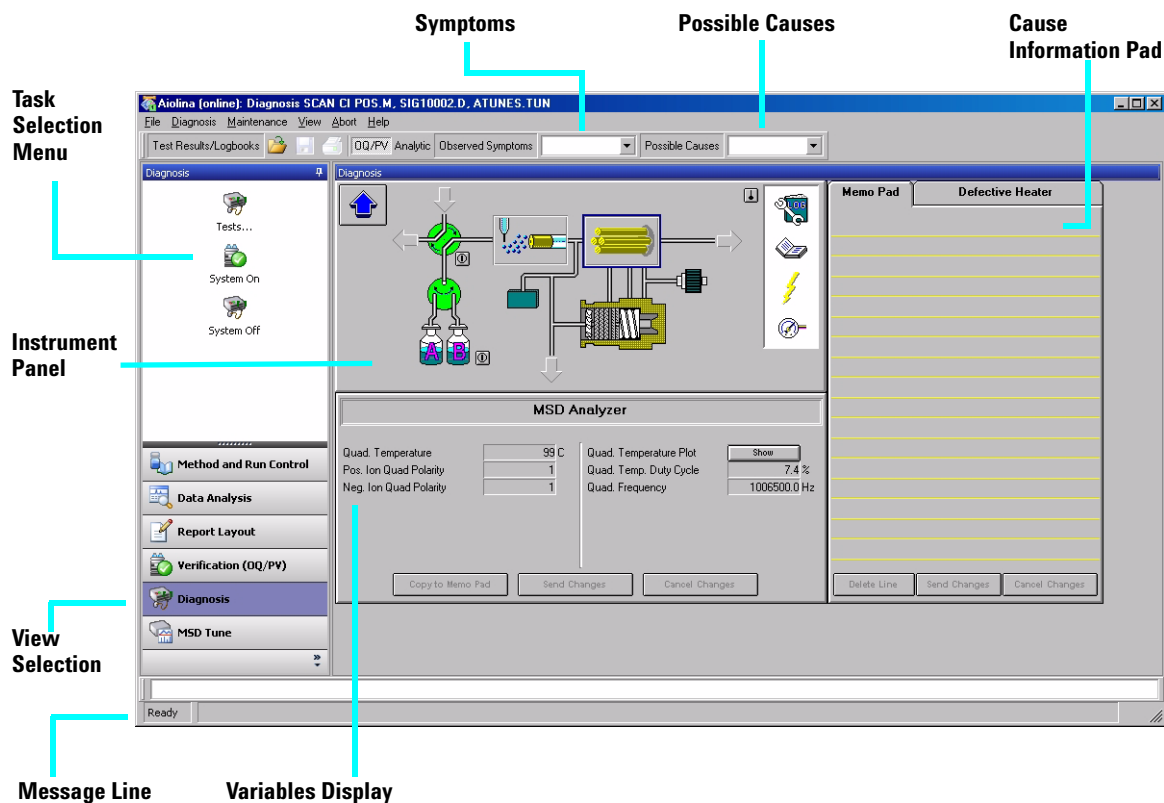


This view allows you to carry out diagnostic and maintenance activities on your system.

Diagnosis Toolbar The Diagnosis toolbar is displayed when you select **Show Top Toolbar** from the **View** menu. Move the cursor over button to view a description of it.

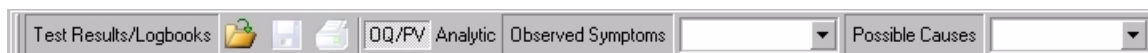

MSD Tune View

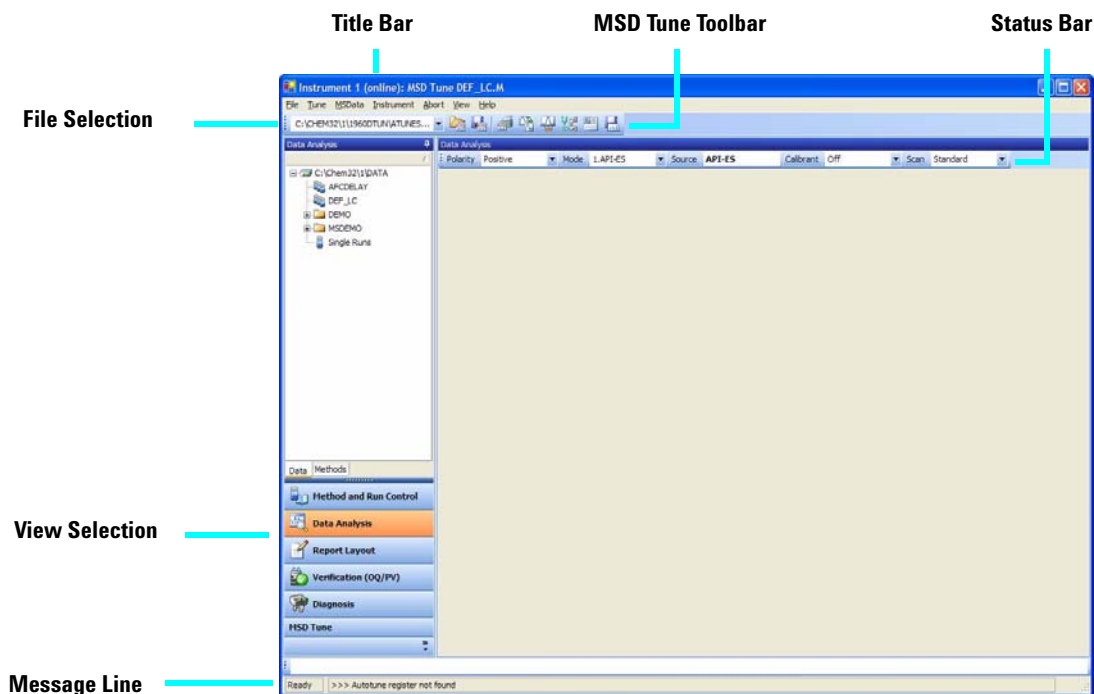

MSD Tune Toolset

This toolset is present in MSD Tune view.



Load an LC/MS quad tune file.



Calibrate the mass axis.



Save the current tune file.



Change spray chamber parameters.



Generate a profile and scan report.



Edit LC/MS quad parameters for manual tuning.





Autotune the LC/MS quad and print a tune report.



Acquire multiple scans and send results to data file.

Basic Operation

Start up and Shut Down the System

Task	Instruction	Comment
To start up the system	<ol style="list-style-type: none"> 1 From the Method and Run Control view, click the  button on the system status bar. When you move mouse over the button, it turns green if system is currently idle. 2 Set up the LC conditions (pump, column heater, and detector). 3 Set up the LC/MS quad spray chamber conditions as described in the MSD Spray Chamber dialog box topic in the online Help. <p>Allow 15 minutes for the system to warm up.</p>	<p>This procedure assumes the system is under vacuum and the LC and LC/MS quad are properly connected. It also assumes the liquid flow path is properly set up.</p>
To put the system in Standby Mode	<ol style="list-style-type: none"> 1 Flush the system with pure mobile phase to ensure the flow path is clear of buffers. This avoids plugging the nebulizer. A plugged nebulizer can cause a high back pressure, which can damage the LC flow cell. It is good practice to flush the flow path (including the column) for 5-10 minutes with a mobile phase without buffers such as 50:50 water/acetonitrile. 2 Click the  button on the system status bar. All modules are set to standby mode (indicated by the gray color in the system diagram). 	<p>Put the system in standby mode overnight or whenever you won't be analyzing samples for an extended time. The standby state for the LC/MS quad leaves the nebulizer and drying gas on at low flow.</p> <p>Standby for the LC/MS quad is:</p> <ul style="list-style-type: none"> • 20 psi for the nebulizer, • 3 L/min for drying gas, • 300°C for drying gas temperature, • 325°C for the APCI vaporizer, if present, and • the MS stream selection valve is set LC to Waste.

Tune the LC/MS quad

When the LC/MS quad is used as a detector for the LC, a mass spectrum is associated with each data point in the LC chromatogram. To obtain high quality, accurate mass spectra, the LC/MS quad must be optimized to:

- Maximize sensitivity
- Maintain acceptable resolution
- Ensure accurate mass assignment

Tuning is the process of adjusting LC/MS quad parameters to achieve these goals. After the LC/MS quad parameters have been optimized, they must be saved in a tune file (.tun). This tune file is then specified in the method that is used to acquire data for your samples.

Frequent tuning, automated or manual, is not required. Once tuned, the LC/MS quad is very stable. Tuning should generally be needed no more often than monthly, weekly at most.

Wait ~12 hours after pumpdown before tuning or operating your LC/MS quad. The analyzer takes about 12 hours to reach thermal equilibrium. Tune files that are created or data that is acquired before the LC/MS quad is at thermal equilibrium may have incorrect mass assignments and other inaccuracies.

Task	Instructions	Comments
To use Autotune	<div><div>1</div><div>From the MSD Tune view, select Autotune from the Tune menu or click the Autotune toolbar button.</div></div> <div><div>2</div><div>Review the tune report that is printed automatically when tuning is completed.</div></div>	Use autotune for automated adjustment of the LC/MS quad performance.

Task	Instructions	Comments
To use Check Tune	<ol style="list-style-type: none"> 1 Select Check Tune from the Tune menu. 2 Review the Check Tune report. Adjustments are suggested for values that are outside of acceptable ranges. 	Use Check Tune to quickly determine whether the LC/MS quad is correctly tuned. It performs a single profile scan of the tune masses and compares the peak widths and mass axes with target values.
To use Manual Tune	<ol style="list-style-type: none"> 1 Select Manual Tune from the Tune menu. 2 Set the desired mass resolution (adjusting width gain and width offset). 3 Calibrate the mass axis (adjusting mass gain and mass offset). 4 Optimize the ion transmission through the source ion optics (fragmentor, skimmer(s), lens 1, lens 2, octopole peak and octopole knee). 5 Adjust the signal strength (setting iris and adjusting the multiplier gain). <p>Note that fragmentor and gain are method parameters. The fragmentor affects ion transmission and fragmentation. For more information, see the online Help.</p>	Use Manual Tune when you want to: <ul style="list-style-type: none"> • Achieve maximum sensitivity by sacrificing some resolution • Tune specifically for the very low end (<150 amu) of the mass range • Tune with a compound other than the standard calibrants

Acquire LC/MS Data

Modes of Acquisition

There are three modes of acquiring data:






- Running a method for a single sample
- Running a sequence for multiple samples
- Running an FIA series

Note the following about acquiring data:

- All three acquisition modes require an appropriate method.
- Samples may be injected either manually or with an ALS.
- A run must always be started from the software.
- An FIA method cannot be used in a sequence but multiple FIA methods can be run sequentially.

To edit a method and start a run




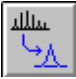

Once you know the acquisition mode you want to use, you need to set up an appropriate method. Methods are set up in the Method and Run Control view.

Instructions	Comments
1 Activate the single sample toolset by clicking the  button on the toolbar.	
2 Load a method by selecting Method / Load Method or by clicking the  button on the toolbar and selecting a method from the list.	
3 Click the  button to begin editing the method. You can also begin by selecting Edit Entire Method from the Method menu or by right-clicking the  icon on the Run Control Bar.	
4 Select the method sections you want to edit (select all sections to become familiar with the method parameters that are available) and click OK.	A series of dialog boxes is displayed, which allows you to set up your method and instrument parameters.
5 Add any method comments you want to appear on your reports.	Click the Help button on any of the method dialog boxes for descriptive information on the items available in each.
6 Set up the instrument parameters, such as pump, injector, DAD or VWD, column thermostat, MS quad signals, and MS quad spray chamber.	
7 Set up the Data Analysis parameters, such as signal details, integration events, report parameters, instrument curves, calibration curves, calibration table, and ion parameters.	
8 Complete the Run Time Checklist.	
9 Save the method using a different name. Select Method / Save Method As or click the  button on the toolbar.	Once you are familiar with the options that are available, you can use the system diagram menus for quick access to particular method parameters, rather than using the edit entire method process.
10 When you are ready to begin a run, click the Start button.	

Use Deconvolution (optional - requires additional license)

Deconvolution is a process that transforms mass spectra from multiply-charged ions into a calculated molecular weight.

To deconvolute a mass spectrum

Instructions	Comments
1 In the Data Analysis view, load an electrospray MS data file. The TIC is displayed.	
2 Generate the MS spectrum that you want to deconvolute (select the Spectral task tool set and use the spectrum selection tools to pick the spectrum of interest).	Note that these operations are performed on data acquired in full scan mode.
3 Click the Enter Deconvolution Tool button to set up the deconvolution display area.	
4 When the spectrum is displayed, click the Label Ions button to locate masses to be used in deconvolution.	
5 Examine the ions that were found. Very noisy data or data with unresolved regions often need special settings to find ions.	You can use the Ion Label Options button to optimize the set of ions found.
6 Click the Edit Deconvolution Parameters button to change any of the parameters.	
7 Click the Run Deconvolution button to begin the deconvolution process.	
8 Select specific components from the Component list to look at individual components or groups of components.	Optionally, click the Preview Deconvolution Report button or click the Print Report button.
9 Optionally, select components from the Component List box and then click the Delete Components button.	
10 Repeat the process.	This action removes the peaks for the selected components from the original spectrum. You can then continue with step 4 above, using this new spectrum.



Deconvolution Report

The Deconvolution report contains summary information about each component selected, along with detailed information on how each peak in a component contributes to the component's molecular weight. In the first part of the report, the components are ranked by percent relative abundance, which is useful for estimating the percent of impurities.

Note that the actual molecular weight may differ from the computed molecular weight even if the data fit a Gaussian curve perfectly, due to other errors such as errors in mass axis assignment or unresolved chemical impurities.

The Deconvolution software is optional and is part of the G2720BA Bioanalysis Software package.

Operating Tips

- Back up your data and methods **regularly** to avoid loss of data if the files are accidentally overwritten, deleted, or if a hardware problem develops with your disk drive.
- Put the system in standby mode overnight or whenever you won't be analyzing samples for an extended time.
- Verify that the tune file you are using is appropriate for your samples.
- Save Tune reports in an MS Logbook for future reference.
- Regular system maintenance can reduce problems. Keep a maintenance record.
- Use the Maintenance Logbook and EMF features in Diagnosis view to help you keep track of when maintenance is needed and to keep an online maintenance record.
- Flush the sample path and clean the spray chamber, capillary tip, and spray shield daily or at the end of each shift. Check the foreline pump fluid level every week.
- The spray chamber vent hose must be connected to a lab vent that is used **only** for the source (completely separate from the vent hose for the foreline pump). Otherwise, waste products can migrate into the spray chamber vent producing chemical noise.
- Samples need to be filtered. They should be salt and detergent-free if no chromatography is used.
- If a UV detector is available, use it in series with the LC/MS quad. Try to minimize chromatographic peak broadening by using low dispersion tubing.
- To avoid chromatographic band broadening, be sure that all tubing connections are free of dead volume. Use zero dead-volume (ZDV) fittings when possible.
- Use finger-tight fittings on the LC inlet union. Swaged fittings can compress the frit.
- If you are using APCI, the optimal flow rate is 1.0 ml/min. The range is 0.5 to 1.5 ml/min.

- Use the following table as a guide to using SIM, condensed scan, and full scan acquisition modes.

Task	Mode
Acquire electrospray data for samples containing large, multiply-charged analytes.	Full Scan
Analyze a mixture with unknown components (small molecules).	Scan
Analyze a mixture with known components in unknown amounts (quantitate).	Scan or SIM
Identify the presence of a few known compounds at low levels within a mixture.	SIM
Use fast chromatography on SL MS quads without missing peaks.	Fast Scan
Use fast chromatography on 6150B AJS-HT MS quads without missing peaks.	Ultra Fast Scan

Troubleshooting Tips

No peaks

- ✓ Verify that there is spray from the nebulizer.
- ✓ Verify that the capillary voltage is set correctly.
- ✓ Verify that the MS system is tuned correctly.
- ✓ Verify that MS vacuum pressures are within normal ranges.
- ✓ Check the drying gas flow and temperature.
- ✓ Verify that the fragmentor is set correctly.
- ✓ Check if any error message indicates a problem with electronics.
- ✓ Verify that sufficient sample is present in vial.

Poor mass accuracy

- ✓ Recalibrate the mass axis.
- ✓ Verify that the ions used for tuning span the mass range of the sample ions and that they show strong, stable signals.

Low signal

- ✓ Check the solution chemistry. Verify that the solvent you're using is appropriate for your sample. Mixed samples can exhibit signal suppression of one or more components.
- ✓ Verify that the sample is fresh and has been stored correctly.
- ✓ Verify that the MS system is tuned correctly.
- ✓ Check the nebulizer condition.
- ✓ Clean the capillary entrance.
- ✓ Check the capillary for damage and contamination.
- ✓ Verify that the spray chamber parameters are set correctly, such as drying gas heater and nebulizer pressure.

Unstable signal

- ✓ Verify that the drying gas flow and temperature are correct for the solvent flow you are using.
- ✓ Verify that the solvent is thoroughly degassed. Do **not** use ultrasonic degassing with protein samples.
- ✓ Verify that the LC backpressure is steady; this indicates a steady solvent flow.

High spectral noise

- ✓ Use appropriate mass filter values.
- ✓ Check the spray shape. Nebulizer may be damaged or incorrectly set.
- ✓ Verify that drying gas flow and temperature are correct for the solvent flow you are using.
- ✓ Verify that the solvent is thoroughly degassed. Do **not** use ultrasonic degassing with protein samples.
- ✓ Verify that the LC backpressure is steady; this indicates a steady solvent flow.
- ✓ If you are using water as part of the mobile phase, verify that it is de-ionized ($>18\text{M}\Omega$).

Droplets, not spray, exiting the nebulizer

- ✓ Verify that the nebulizing gas pressure is set high enough for the LC flow being used.
- ✓ Check the position of the needle in the nebulizer.
- ✓ Stop the solvent flow and remove the nebulizer assembly. Use a magnifying glass to examine the end of the nebulizer for damage.

No flow

- ✓ Verify that the LC is on and there is sufficient solvent in the correct bottle.
- ✓ Check for LC error messages.
- ✓ Check for blockages. Repair or replace any blocked components.
- ✓ Check for leaks.
- ✓ Verify that the MS stream selector valve is set to **LC to MSD**.

Undesired fragmentation

- ✓ Fragmentor is set too high.
- ✓ Ionization is causing fragmentation (APCI vs. Electrospray).
- ✓ APCI temperature is too high.

Safety Warnings

Intended Purpose

Use of the LC/MS quad in a manner not intended by the manufacturer is prohibited.

Safety Class

The LC/MS quad is a Safety Class 1 instrument and has been designed and tested in accordance with IEC Publication 1010-1, Safety Requirements for Electrical Equipment for Measurement, Control, and Laboratory Use.

Stack Configuration

WARNING

Do not stack more than three LC modules of any kind on top of the LC/MS instrument. Stacking more than three LC modules is potentially unstable and dangerous.

Placing the LC modules on top of the LC/MS quad is also inconvenient. It puts the solvent bottles and some LC controls out of easy reach and requires the LC stack to be disassembled and removed for LC/MS quad maintenance.

WARNING

Use only Nitrogen as the drying and nebulizer gas. Air, oxygen, or other gases, when combined with solvents and high voltages in the spray chamber, can result in an explosion.

WARNING

Do not connect the LC/MS quad to a power source that is not equipped with a protective earth contact. You can create a shock hazard for the operator and can damage the instrument.

WARNING

Do not interrupt the protective conductor inside or outside the LC/MS quad or disconnect the protective earth terminal. You can create a shock hazard for the operator and can damage the instrument.

WARNING

Do not adjust, maintain, or repair the opened instrument while it is connected to a power source. These operations should be performed only by trained persons who are aware of the hazards involved.

WARNING

Before you perform maintenance on the electronics or turbo controllers, turn off the main circuit breaker and disconnect the power cord. The power switch on the front of the LC/MS quad does not completely cut power to the instrument.

www.agilent.com

In this book

This book contains:

- [Where to find more information](#)
- [What's New in B.04.02?](#)
- [Configuration Diagrams](#)
- [ChemStation Views](#)
- [Basic Operation](#)
- [Operating Tips](#)
- [Troubleshooting Tips](#)
- [Safety Warnings](#)

© Agilent Technologies, Inc. 2010

Printed in USA
Fourth Edition, April 2010



G1960-90070



Agilent Technologies