



Analysis of Carbohydrates in Lemonade using HPLC

Angelika
Gratzfeld-Huesgen


Food

Abstract

The following carbohydrates have been analyzed: glucose, galactose, raffinose, fructose, mannitol, sorbitol, lactose, maltose, cellobiose, and sucrose. Food carbohydrates are characterized by a wide range of chemical reactivity and molecular size. Because carbohydrates do not possess chromophores or fluorophores, they cannot be detected with UV-visible or fluorescence techniques. Nowadays, however, refractive index detection can be used to detect concentrations in the low parts per million (ppm) range and above, whereas electrochemical detection is used in the analysis of sugars in the low parts per billion (ppb) range.

Sample preparation

Degassed drinks can be injected directly after filtration. More complex samples require more extensive treatment, such as fat extraction and deproteinization. Sample cleanup to remove less polar impurities can be done through solid-phase extraction on C18 columns.

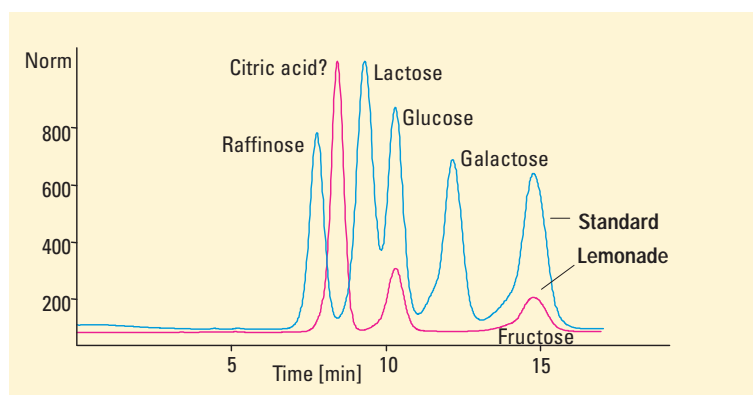

Conditions

Column

300 ~ 7.8 mm Bio-Rad HPXP, 9 μ m

Mobile phase water

Column compartment 80 °C

Flow rate 0.7 ml/min

Detector refractive index

Sample preparation

Samples were directly injected.

Figure 1
Analysis of carbohydrates in lemonade



Agilent Technologies
Innovating the HP Way

Chromatographic conditions

The HPLC method presented here was used to analyze mono-, di-, and trisaccharides as well as sugar alcohols.

HPLC method performance

Limit of detection

<80 ng with S/N = 2

Repeatability of

RT over 10 runs <0.05 %

areas over 10 runs 2 %

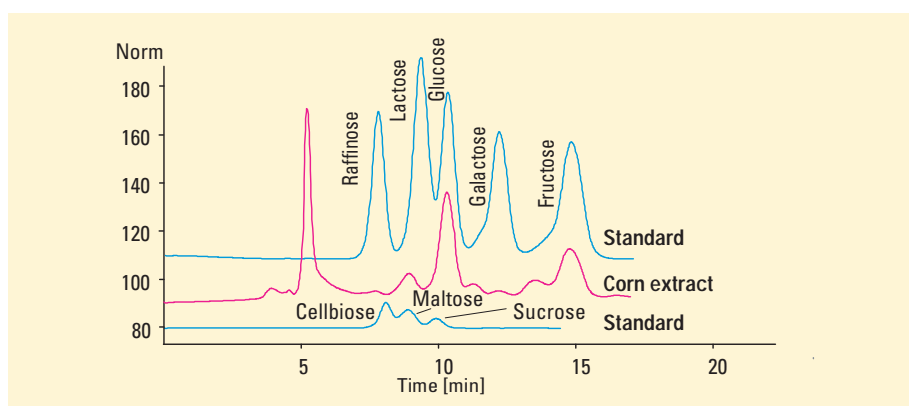


Figure 2
Analysis of carbohydrates in corn extract


References

1. *Official Methods of Analysis, Food Compositions; Additives, Natural Contaminants*, 15th ed; AOAC: Arlington, VA, 1990, Vol. 2; AOAC Official Method 980.13: Fructose, glucose, lactose, maltose, sucrose in milk chocolate; AOAC Official Method 982.14: Glucose, fructose, sucrose, and maltose in presweetened cereals; AOAC Official Method 977.20: Separation of sugars in honey; AOAC Official Method 979.23: Saccharides (major) in corn syrup; AOAC Official Method 983.22: Saccharides (minor) in corn syrup; AOAC Official Method 984.14: Sugars in licorice extracts.

Equipment

Agilent 1100 Series

- degasser
 - isocratic pump
 - autosampler
 - thermostatted column compartment
 - refractive index detector
- Agilent ChemStation + software

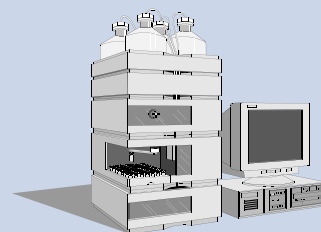

Angelika Gratzfeld-Huesgen is application chemist at Agilent Technologies, Waldbronn, Germany.

For more information on our products and services, visit our worldwide website at <http://www.agilent.com/chem>

© Copyright 1997 Agilent Technologies
Released 09/97
Publication Number 5966-0637E



Agilent Technologies
Innovating the HP Way