


Anorectics (Fen-Phen) Analysis on ZORBAX Bonus-RP

Application
Pharmaceutical
Robert Ricker

The popular but controversial dietary aide, Fen-Phen, is a mixture of two basic drugs. These compounds elute from ZORBAX Bonus-RP with reduced peak tailing compared to that on an alkyl-C8 column.

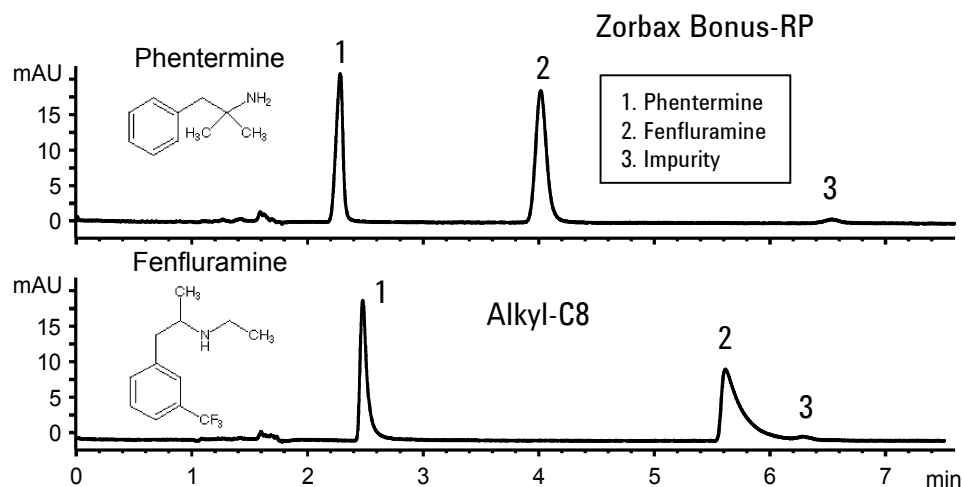

Conditions: HPLC System: Agilent 1100 with quaternary pump
Column: ZORBAX Bonus-RP, 4.6 x 150 mm, Agilent P/N: 883668-901
Mobile Phase: [MeOH: ACN (50:50)] : K₂HPO₄, 25 mM, pH 7.2, (55: 45)
Detector: UV 254 nm; Flow: 1 mL/ min.; Temperature: ambient

Highlights

- ZORBAX Bonus-RP provides unique selectivity and peak shape that significantly improves resolution between peak 2 and an impurity.
- ZORBAX Bonus-RP is a unique polar-linked bonded phase that provides:
 - Good peak shape (improved quantitation and resolution)
 - Alternate selectivity (improved resolution)
 - Long life over a wide pH range (reproducible separations)



Agilent Technologies

*Robert Ricker is an application chemist
based at Agilent Technologies, Wilmington,
Delaware.*

For more information on our products and
services, visit our website at:
www.agilent.com/chem

Copyright© 2002 Agilent Technologies, Inc.
All Rights Reserved. Reproduction,
adaptation or translation without prior
written permission is prohibited, except as
allowed under the copyright laws.

Agilent shall not be liable for errors
contained herein or for incidental or
consequential damages in connection with
the furnishing, performance, or use of this
material.

Information, descriptions, and specifications
in this publication are subject to change
without notice.

Printed in the USA
April 25, 2002
5988-6415EN



Agilent Technologies