

User's guide

Cleaning sequence and reprocessing list audit trails

1) Note that this tool will only work with Galaxie 1.9 SP2b or newer version

2) Copy the files on a USB drive, and make sure the following files are in the same directory:

- CleanSequenceAuditTrail.exe
- ForEachParser.exe
- CleanSequenceAuditTrail.ini
- CleanReprocessingAuditTrail.ini
- Clean.Bat
- orlando_filapi.dll (If you use a newer version than 1.9SP2b, don't add this .dll)

3) In both CleanSequenceAuditTrail.ini and CleanReprocessingAuditTrail.ini, modify the Galaxie Data folder path if needed. The tool will clean all the sequences and reprocessing lists found under the folder indicated (also in children folders).

For example: "C:\Galaxie\Data\Group1\".

4) Launch the Clean.bat. The tool works without any intervention. The windows close automatically when the tool has finished.

5) Note: the tool creates automatically a backup file in the folder indicated in the *.ini file:

e.g.: C:\Galaxie\Data\Group1\example.SEQU.Backup

NOTICE: This document contains references to Varian. Please note that Varian, Inc. is now part of Agilent Technologies. For more information, go to www.agilent.com/chem.



Agilent Technologies