


Agilent CombiZorb Scavengers . . .

For removal of unwanted components from combinatorial chemistry reaction mixtures

Application Note

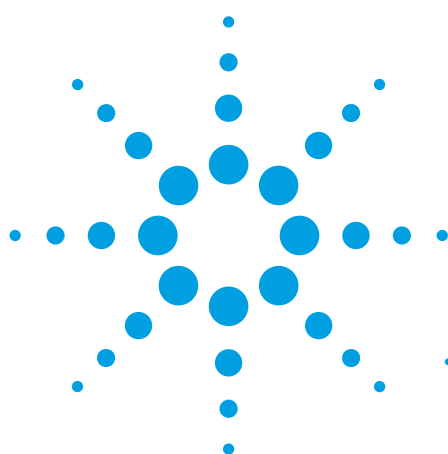

Agilent CombiZorb Scavengers are based on either ultrapure (99.995%) porous silica microspheres or macroporous polystyrene divinylbenzene (MP PS/DVB). CombiZorb Scavengers provide more capacity per unit volume compared to gel-type PS scavengers.

CombiZorb S Scavengers

- **No swelling** - usable in different formats like cartridges, columns, or 96-well plates
- Based on high purity (99.995%) porous silica microspheres with low metal content
 - No interfering reactions or catalysis with metals
 - No contamination from the adsorbent
- **Broad solvent compatibility**
 - Use with polar and non-polar solvents
 - No need to switch solvents between reaction and scavenging steps
- No static charge
- Easy to handle when packing columns or transferring the adsorbent

CombiZorb MP Scavengers

- **Significantly less swelling** (+50%) than standard PS/DVB resins (+500%)
 - Higher, more consistent capacity/volume ratio than resins that swell
 - No need to pre-swell
- **Broad solvent compatibility**
 - Use with polar or non-polar solvents
 - No need to switch solvents between reaction and scavenging steps
- High purity

Agilent also makes several other products for the combinatorial chemistry environment.

- ZORBAX Rapid Resolution columns (15, 30, 50 mm) for fast screening
- CombiHT columns for fast preparative LC in the 15-50 mg range
- The Agilent 1100 LC and MSD instruments for liquid chromatography and mass spectral analysis

Ordering Information

CombiZorb Scavenger	Part No.	Size
S-Triamine	550145-138	10g
	550245-238	100g
S-Monoamine	550145-118	10g
	550245-218	100g
S-Sulfonic Acid	550145-124	10g
	550245-224	100g
S-Tertiary Amine	550145-148	10g
	550245-248	100g
MP-Aldehyde	510100-150	10g
	510200-250	100g
MP-Isocyanate	510100-160	10g
	510200-260	100g
MP-Piperidinomethyl	510100-170	10g
	510200-270	100g
MP-Sulfonyl Chloride	510100-114	10g
	510200-214	100g
MP-Sulfonyl Hydrazide	510100-134	10g
	510200-234	100g
Variety Pack*	550550-901	6-5g bottles

S = silica (based on high purity, 99.995% pure porous silica microspheres)
MP = macroporous polystyrene/divinylbenzene (PS/DVB)

* Variety pack contains a 5 g bottle of each of the six scavengers:
MP-Aldehyde, MP-Isocyanate, S-Triamine, S-Monoamine,
S-Sulfonic Acid and S-Tertiary Amine

Request a quote for special orders and larger quantities of CombiZorb Scavengers.



Agilent Technologies

Innovating the HP Way

CombiZorb Product Information



CombiZorb Scavengers	Structure	Characteristics	Loading Capacity	Applications
CombiZorb S Scavengers				
S-Triamine		Spherical, ultra-pure silica; Density: 1.3 g/ml; Particle Size: 20-80 µm; no swelling	1.3-1.6 meq/g	Scavenging acids, acid chlorides, acid anhydrides, chloroformates, isocyanates, aldehydes, and many other electrophiles
S-Monoamine		Spherical, ultra-pure silica; Density: 1.3 g/ml; Particle Size: 20-80 µm; no swelling	0.6-0.9 mmol/g	Scavenging acids, acid chlorides, acid anhydrides, chloroformates, isocyanates, aldehydes, and many other electrophiles
S-Sulfonic Acid		Spherical, ultra-pure silica; Density: 1.3 g/ml; Particle Size: 20-80 µm; no swelling	0.4-0.6 mmol/g	Scavenging amines and other basic compounds
S-Tertiary Amine		Spherical, ultra-pure silica; Density: 1.3 g/ml; Particle Size: 20-80 µm; no swelling	0.6-0.9 mmol/g	Scavenging acids generated from condensation reactions and may catalyze acylation reactions
CombiZorb MP* Scavengers				
MP-Aldehyde		Spherical, macroporous polystyrene/DVB particle; Density: 0.6 g/ml; Particle Size: 30-100 µm; swelling: 20-30% in THF or CH ₂ Cl ₂	1.2-1.8 mmol/g	Scavenging hydrazines, hydroxylamines, primary amines, organometallic reagents (organolithium and Grignard reagents), and other nucleophiles
MP-Isocyanate		Spherical, macroporous polystyrene/DVB particle; Density: 0.6 g/ml; Particle Size: 30-100 µm; swelling: 20-30% in THF or CH ₂ Cl ₂	0.9-1.3 mmol/g	Scavenging primary and secondary amines, anilines, hydrazines, and other nucleophiles
MP-Piperidinomethyl		Spherical, macroporous polystyrene/DVB particle; Density: 0.6 g/ml; Particle Size: 30-100 µm; swelling: 30-50% in THF or CH ₂ Cl ₂	1.7-2.5 mmol/g	Use as catalyst for acylation reactions in place of tertiary amine and for scavenging acids generated from condensation reactions. ¹
MP-Sulfonyl Chloride		Spherical, macroporous polystyrene/DVB particle; Density: 0.6 g/ml; Particle Size: 30-100 µm; swelling: 30-50% in THF or CH ₂ Cl ₂	1.6-2.0 mmol/g	Scavenging alcohols, phenols, and amines, more widely used as solid phase reagents in "catch and release" applications. ^{2,3}
MP-Sulfonyl Hydrazide		Spherical, macroporous polystyrene/DVB particle; Density: 0.6 g/ml; Particle Size: 30-100 µm; swelling: 30-50% in THF or CH ₂ Cl ₂	1.5-2.5 mmol/g	Scavenging aldehydes and ketones and used as solid phase reagents in some "catch and release" applications. ^{4,5}

*MP = macroporous polystyrene divinylbenzene

References

1. R.J. Booth and J.C. Hodges, *J. Am. Chem. Soc.* 1997, 119, 4882
2. T. Takahashi, S. Ebata, and T. Doi, *Tetrahedron Lett.* 1998, 39, 1369
3. E.W. Baxter, J.K. Rueter, S. Nortey, and A.B. Reitz, *Tetrahedron Lett.* 1998, 39, 979
4. Y. Hu, S. Baudart, and J. Porce, Jr., *J. Org. Chem.* 1999, 64, 1049
5. O. Galliglu and A. Auar, *Eur. Polym. J.* 1989, 25, 313

For the latest information on the complete line of Agilent Technologies columns and supplies for analytical instruments, see our online catalog at www.agilent.com, or contact your local Agilent sales office. For all other areas contact Agilent or your local authorized distributor.

Information, descriptions and specifications in this publication are subject to change without notice.

