

## Effect of Bonded-Phase on Resolution and Retention of Polypeptides

Application **Biochemical** 

Reversed-phase HPLC continues to be the analysis technique of choice for the analysis of proteins and peptides. The mobile phase for these separations generally contains aqueous buffers with ACN and TFA. Selectivity changes between samples are often accomplished with the use of different bonded phases.



## Highlights

- · Differences in bonded-phase selectivity can significantly improve resolution in reversed-phase analysis of peptides and proteins. See change in separation between peaks 2 and 3, peaks 5 and 6, and peaks 8 and 9.
- · Change of temperature, in addition to change of bonded phase, can be used to obtain desired separations.
- · Sterically-protected bonded phases of ZORBAX SB packings are well suited to chromatography at very low pH.

## **Conditions:**

ZORBAX 300 SB-CN, SB-C3, SB-C8, SB-C18 (4.6 x 150 mm, Agilent P/N: 883995-905, 883995-909, 883995-906, 883995-902) Mobile Phase: 15-53% in 20 min. A: 5:95 ACN:Water with 0.10 % TFA (v/v%) B: 95:5 ACN:Water with 0.085% TFA (v/v%) Injection 10µL, 2-6µg protein (in 6M Guanidine HCl pH 7.0), 1.0mL/min., 35° C, Detect. UV (215 nm)



Robert Ricker is an application chemist based at Agilent Technologies, Wilmington, Delaware.

For more information on our products and services, visit our website at: www.agilent.com/chem

Copyright<sup>©</sup> 2002 Agilent Technologies, Inc. All Rights Reserved. Reproduction, adaptation or translation without prior written permission is prohibited, except as allowed under the copyright laws.

Agilent shall not be liable for errors contained herein or for incidental or consequential damages in connection with the furnishing, performance, or use of this material.

Information, descriptions, and specifications in this publication are subject to change without notice.

Printed in the USA April 25, 2002 5988-6310EN



**Agilent Technologies**