

Epoxy Resin Analysis on Agilent PLgel MIXED-E with Gel Permeation Chromatography

Application Note

Materials Testing and Research, Polymers

Authors

Greg Saunders and Ben MacCreath
Agilent Technologies, Inc.
Essex Road
Church Stretton
S46 6AX
UK

Introduction

Epoxy resins are thermosetting copolymers formed by mixing epoxide with polyamine. They are extremely durable after curing, and so are widely used as general-purpose adhesives, coatings and composites.

Analysis of epoxy resins by gel permeation chromatography is straightforward with Agilent PLgel 3 μ m MIXED-E columns, which are ideal for low molecular weight samples that contain oligomeric fractions, as well as polymers, up to 30,000 MW.



Agilent Technologies

Analysis of polysiloxane

In this analysis of epoxy resin, the ultra high efficiency achieved with Agilent PLgel 3 μ m particles gives baseline resolution of the sample components with the lowest MW component eluting within 8 minutes.

Conditions

Columns	Agilent PLgel 3 μ m MIXED-E, 300 \times 7.5 mm (p/n PL1110-6300)
Eluent	THF
Flow rate	1.0 mL/min
Detector	RI
System	Agilent PL-GPC 50

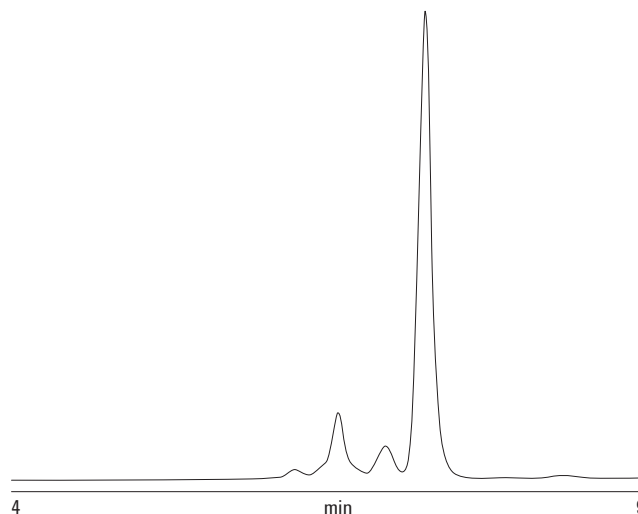


Figure 1 Fast elution of low molecular components of an epoxy resin on an Agilent PLgel 3 μ m MIXED-E column.

For More Information

For more information on our products and services, visit our Web site at www.agilent.com/chem.

www.agilent.com/chem

Agilent shall not be liable for errors contained herein or for incidental or consequential damages in connection with the furnishing, performance, or use of this material.

Information, descriptions, and specifications in this publication are subject to change without notice.

© Agilent Technologies, Inc., 2011

Printed in the USA

April 18, 2011

5990-7980EN



Agilent Technologies