

This chromatogram shows the complete separation of ß-amyloid peptides (ßAP) using Agilent ZORBAX 300 SB-C8 column. The inset box shows the high recovery following the HPLC separation, as determined by direct protein assay of collected peaks. The separation of these peptides was not possible at column temperatures below 80°C.

## Highlights

- Sterically-protected bonded phases (StableBond) permit use of low pH mobile phases, even at high temperature.
- Use of elevated column temperature can dramatically improve separation and recoveries of difficult samples.
- Note the greater retention of ßAP(1-42) than ßAP(1-43).



2.5 μg each peptide in 20 μL 6 M urea / 5% HOAc Flow: 1 mL/min.



Robert Ricker is an application chemist based at Agilent Technologies, Wilmington, Delaware.

For more information on our products and services, visit our website at: www.agilent.com/chem

Copyright<sup>©</sup> 2002 Agilent Technologies, Inc. All Rights Reserved. Reproduction, adaptation or translation without prior written permission is prohibited, except as allowed under the copyright laws.

Agilent shall not be liable for errors contained herein or for incidental or consequential damages in connection with the furnishing, performance, or use of this material.

Information, descriptions, and specifications in this publication are subject to change without notice.

Printed in the USA April 25, 2002 5988-6326EN



**Agilent Technologies**