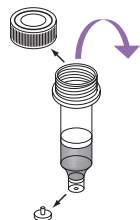
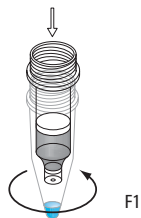


1 Dilute and filter sample

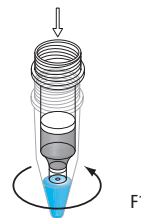
- Dilute 8-10 μ L human serum/plasma sample to 200 μ L with Buffer A.
- **NOTE:** Consult cartridge certificate for true capacity.
- Filter through 0.22 μ m spin filter.

2 Remove buffer

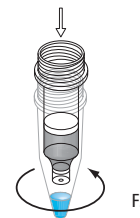
- Remove cartridge cap and plug and remove buffer from top of resin bed with transfer pipette.
- **Never let frits or resin bed run dry.**

3 Apply sample and incubate

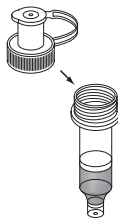
- Add 200 μ L diluted serum/plasma sample. Cap cartridge loosely or leave open.
- Place in 1.5 mL collection tube labeled "Flow-through fraction 1" (F1).
- Centrifuge 1 minute at $100 \times g$.
- Remove cartridge from centrifuge and let sit for 5 minutes at room temperature.

4 Wash and collect flow-through F1

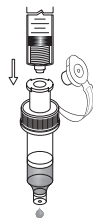
- Add 400 μ L Buffer A.
- Centrifuge 2.5 min at $100 \times g$.
- Collect in F1 tube.

5 Wash and collect flow-through F2

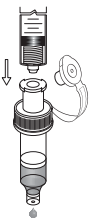
- Place spin cartridge in new collection tube labeled "Flow-through fraction 2" (F2).
- Add 400 μ L Buffer A. Centrifuge 2.5 min at $100 \times g$.
- Collect in F2 tube.

6 Prepare for elution

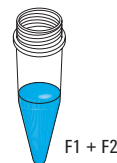
- Remove spin cartridge from F2 tube and attach Luer-Lock adapter tightly to top of cartridge.

7 Elute bound fraction

- Fill 5 mL Luer-Lock plastic syringe with 2.5 mL Buffer B and attach to Luer-Lock adapter.
- Slowly push Buffer B through cartridge to elute bound proteins into new collection tube.
- Save eluate with targeted high-abundance proteins for analysis or discard.

8 Re-equilibrate

- Fill new 5 mL plastic syringe with 5 mL Buffer A and attach to Luer-Lock adapter.
- Slowly push Buffer A through cartridge to re-equilibrate the cartridge for the next sample or store wetted with Buffer A (at 4 °C).
- Re-cap both ends for storage.

9 Analyze F1 + F2

- Fractions F1 and F2 can be analyzed individually or combined.
- Concentrate and analyze these fractions containing low-abundance proteins.

**Agilent Technologies**

For more detailed instructions or information on accessories refer to the Agilent Human 14 Multiple Affinity Removal Spin Cartridge Instruction Guide

Visit www.agilent.com/chem/bioreagents for more products that will help make your proteomics research more efficient:

- Multiple Affinity Removal LC columns for automated serum/plasma processing
- Tryptic digestion kit for digesting gel-isolated proteins
- C18 Cleanup Pipette Tips and Spin Tubes for purifying proteomics samples
- Spin filters and concentrators for cleaning, concentrating, and buffer-exchanging samples
- MALDI-MS matrices and peptide/protein standards for your proteomics studies
- mRP-C18 column for extremely high recovery protein fractionation or desalting



Multiple Affinity Removal LC column



Protein In-Gel Tryptic Digestion Kit



Cleanup C18 Pipette Tips



mRP Column

For More Information

For more information on our products and services, visit our Web site at www.agilent.com/chem/bioreagents.

Or call your local authorized Agilent distributor.

Agilent shall not be liable for errors contained herein or for incidental or consequential damages in connection with the furnishing, performance, or use of this material.

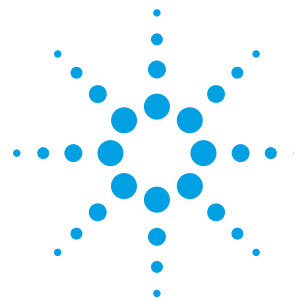
Information, descriptions, and specifications in this publication are subject to change without notice.

© Agilent Technologies, Inc. 2007

Printed in the USA
May 16, 2007
5969-1598



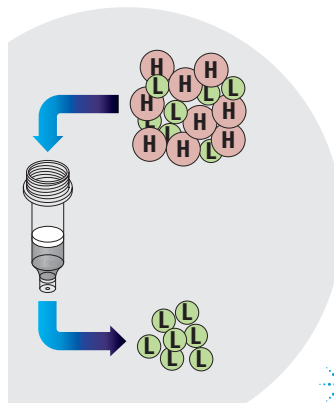
Agilent Technologies



Agilent Human 14 Multiple Affinity Removal Spin Cartridge

Part Number 5188-6560

Quick Reference Guide



Agilent Technologies