

Metanephrines in Urine

Application Note

Bioanalysis

Introduction

An HPLC separation of four metanephrines in a urine sample was accomplished with an Agilent Polaris C18-A column.



Agilent Technologies

Conditions

Column	Agilent Polaris C18-A, 4.6 × 150 mm, 5 µm (p/n A2000150X046)
Sample	Standard and patient urine sample
Sample concentration	Urine standard - normetanephrine 3543 nmol/L, metanephrine 1496 nmol/L, 3-methoxytyramine 1638 nmol/L, internal standard (4-O-methyldopamine)
Mobile phase	0.1 M KH ₂ PO ₄ , pH 6.0, 20 mg/L 1-octanesulfonic acid, sodium salt, 1 g/L sodium EDTA, 6% acetonitrile
Flow rate	1.0 mL/min
Injection volume	5 µL
Temperature	30 °C
Detector	ECD, Antec Decade, +0.63 V Glassy Carbon – Ag/AgCl

For More Information

These data represent typical results. For more information on our products and services, visit our Web site at www.agilent.com/chem.

Peak identification (for both figures)

1. Normetanephrine
2. Metanephrine
3. 3-methoxytyramine (3-MT)
4. 4-O-methyldopamine (IS)

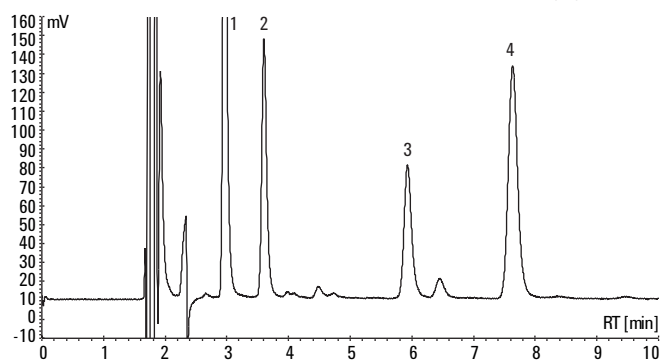


Figure 1. Metanephrines in urine with an Agilent Polaris C18-A column and electrochemical detection on a standard urine sample.

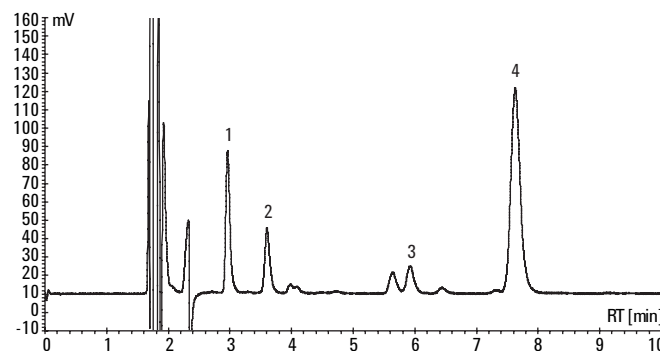


Figure 2. Metanephrines in urine with an Agilent Polaris C18-A column and electrochemical detection on a patient urine sample.

www.agilent.com/chem

Agilent shall not be liable for errors contained herein or for incidental or consequential damages in connection with the furnishing, performance, or use of this material.

Information, descriptions, and specifications in this publication are subject to change without notice. This publication was originally published in 2009.

© Agilent Technologies, Inc., 2011
Published in USA, July 12, 2011
SI-02160



Agilent Technologies