



Monitoring the Purification of Synthetic Amyloid Peptide Ab (1-42) by High-Temperature Reversed-Phase HPLC

Application


Biochemical

Robert Ricker

This example illustrates the use of elevated column temperature to examine the purity of isolated Ab (1-42) peptide. The high column temperature, 80°C, is required to obtain good recovery (>75%) and peak shape for this highly hydrophobic and aggregation-prone peptide. Chromatography at lower temperature results in increasing loss of the peptide (5% recovery at 25°C), and marked band-tailing. This example compares the purity of the crude synthetic peptide in Panel A, to the purified peptide in Panel B. The analysis takes advantage of isocratic separation (no re-equilibration time) using highly efficient 3.5 µm particle diameter, 300Å pore size, Agilent ZORBAX SB-C18 column packing material. The high stability of the StableBond packing permits operation at elevated temperature and low pH, without damage to the column.

Highlights

- *High temperature separation enhances recovery and improves peak shape for this badly behaved peptide.*
- *Highly stable, sterically-protected silane chemistry permits the use of extreme separation conditions, without the consequence of rapid column break down.*
- *Small particle-diameter packing material results in improved resolution.*

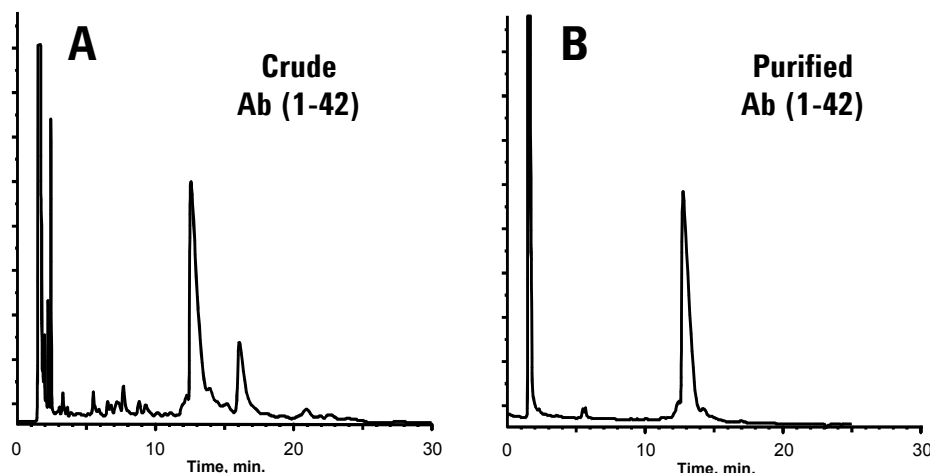

Conditions:
LC: HP1090
Columns: ZORBAX 300SB-C18, 3.5µm, 4.6 x 150mm, Agilent P/N: 863973-902
Mobile Phase: 29% ACN / 0.1% TFA in Water; 80°C
Sample: 20µL, Ab(1-42); Flow: 1.0 mL / min.; UV 210 nm



Agilent Technologies

*Robert Ricker is an application chemist
based at Agilent Technologies, Wilmington,
Delaware.*

For more information on our products and
services, visit our website at:
www.agilent.com/chem

Copyright© 2002 Agilent Technologies, Inc.
All Rights Reserved. Reproduction,
adaptation or translation without prior
written permission is prohibited, except as
allowed under the copyright laws.

Agilent shall not be liable for errors
contained herein or for incidental or
consequential damages in connection with
the furnishing, performance, or use of this
material.

Information, descriptions, and specifications
in this publication are subject to change
without notice.

Printed in the USA
April 25, 2002
5988-6325EN



Agilent Technologies