




# Separation of Retinal (Vitamin A) Isomers: Comparison of Mobile-Phase Composition

Application  
Food Analysis  
Robert Ricker

HPLC separation of vitamin A isomers (retinal isomers) is important in the analysis of various food stuffs, metabolism of visual pigments, liver function, and the corpus luteum. HPLC has traditionally been carried out using 1,4-dioxane. Efforts here were to find a more friendly solvent and a column that provides good chromatographic performance.

## Highlights

- Various solvent combinations can be used in conjunction with ZORBAX Sil to provide adequate resolution of Vitamin A isomers.
- ZORBAX Sil silica provides good peak shape for these retinoids.

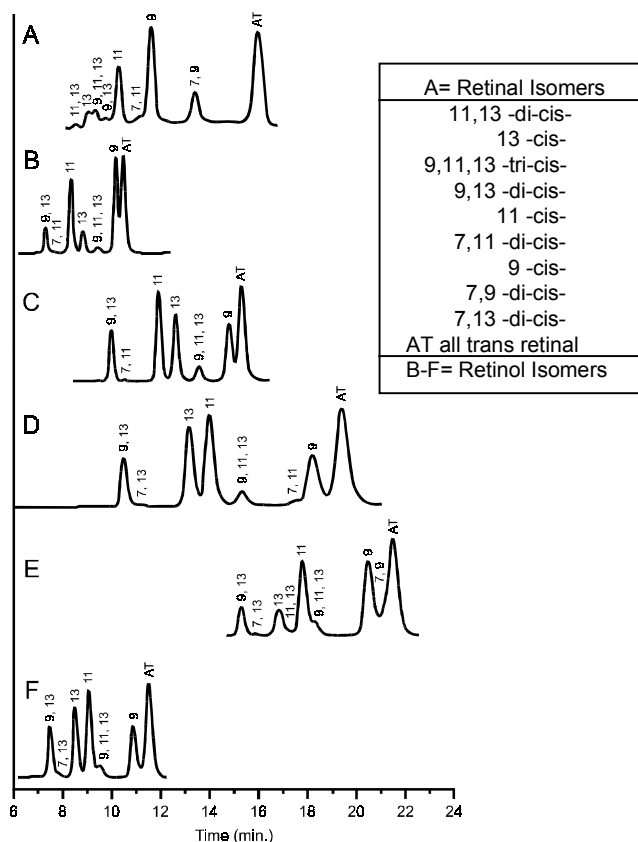

Courtesy of Dr. G. Nöll Physiologisches Inst. -- Justus Liebig Uni. Giessen

Conditions:  
ZORBAX Sil (4.6 x 250 mm) (Agilent P/N: 880952-701)  
Injection volume 20µL, 25°C



Agilent Technologies

Chromatogram	Mobile-Phase	Ratio (V/V)	Flow Rate mL/min	Pressure (bar)	$\lambda$ nm	All Trans at min
A. retinal	n-hexane/t-BME	97:3	2	58	371	15.94
B. retinol	n-hexane/1,4-dioxane	93:7	2	54	320	10.07
C. retinol	n-hexane/1,4-dioxane	94:6	2	67	325	14.90
D. retinol	n-hexane/BME	93:7	4	118	325	19.29
E. retinol	n-heptane/BME	94:6	3	140	325	21.50
F. retinol	n-heptane/BME	93:7	4	187	325	11.25

*Robert Ricker is an application chemist based at Agilent Technologies, Wilmington, Delaware.*

For more information on our products and services, visit our website at:  
[www.agilent.com/chem](http://www.agilent.com/chem)

Copyright© 2002 Agilent Technologies, Inc.  
All Rights Reserved. Reproduction, adaptation or translation without prior written permission is prohibited, except as allowed under the copyright laws.

Agilent shall not be liable for errors contained herein or for incidental or consequential damages in connection with the furnishing, performance, or use of this material.

Information, descriptions, and specifications in this publication are subject to change without notice.

Printed in the USA  
April 25, 2002  
5988-6361EN



**Agilent Technologies**