

## SPE Method for Anabolic Steroids

### Application Brief

**Condition:** Add 5 mL methanol to the cartridge. Apply vacuum and discard the eluant. Repeat with 5 mL water. Do not allow the sorbent to go dry at any point during this step.

**Load:** Attach reservoir to the cartridge using the coupling adaptor. Add 10 mL urine or 4 mL serum to the cartridge. Apply vacuum so that sample takes 30–60 seconds to pass through the cartridge. Serum may be diluted 1/1 with water to aid in handling.

**Rinse:** Add 3 mL 50% methanol to the cartridge. Apply vacuum and discard eluant. Leave vacuum on for 30 seconds after all of the 50% methanol has passed through the cartridge. Centrifuge the cartridge at 1000–1500 rpm for 5 minutes.<sup>1</sup>

**Elution:** Place a collection tube beneath the cartridge. Add 3 mL methanol to the cartridge. Apply vacuum and collect the eluant. Concentrate to dryness using gentle heat (<45°C) and under a stream of nitrogen (transfer to a small vial may be necessary). Dissolve the residue in 100 µL methanol. Inject 1–2 µL into GC.

<sup>1</sup> Centrifuging removes additional water which aids in sample concentration.

### Equipment

- AccuBOND<sup>®</sup> ODS (C18) 6 mL/500 mg cartridge (P/N 188-1356)
- coupling fitting (P/N 5185-5794)
- 25 mL sample reservoir (P/N 700-4007)
- vacuum manifold (P/N 5185-5754, 10-port)  
(P/N 5185-5765, 20-port)

### Reagents

- water (HPLC grade)
- methanol (HPLC grade)



Compound	k	$\bar{x}$	Std. Dev.
Dehydroisoandrosterone (prasterone)	9.55	76	7
5 $\alpha$ -Androstan-17 $\beta$ -ol-3-one (stanolone)	9.85	22	6
19-Nortestosterone (nandrolone)	9.95	70	5
Mesterolone	10.24	90	8
Testosterone	10.37	70	4
1-Dehydrotestosterone (boldenone)	10.51	98	6
17 $\alpha$ -Methyltestosterone	10.55	76	5
1-Dehydro-17 $\alpha$ -methyltestosterone (methandrostenolone)	10.77	76	6
Norethandrolone	10.97	77	5
1-Dehydrotestosterone acetate	11.30	81	4
19-Nortestosterone-17-propionate	11.35	74	9
4-Chlortestosterone-17-acetate (clostebol)	12.47	60	8
Stanozolol	12.76*	46	10
1-Dehydrotestosterone benzoate	15.22	57	17
19-Nortestosterone-17-decanoate	15.48	59	9
1-Dehydrotestosterone undecylenate	16.60	55	12

\*poor peak shape


k = partition ratio (a measure of retention)

$\bar{x}$  = % recovery

## Steroid Extract from Urine

Column: DB-1  
30 m x 0.25 mm I.D.  
P/N: 122-1031  
Film Thickness: 0.1  $\mu$ m  
Carrier: Helium at 40 cm/sec (measured at 55°C)  
Oven: 55°C for 1 min  
55°C – 175°C at 30°C/min  
175°C – 325°C at 10°C/min  
Injector: Splitless; 275°C  
45 sec. purge activation time  
Detector: FID; 300°C


1. Dehydroisoandrosterone (prasterone)  
2. 10-Nortestosterone (nandrolone)  
3. Testosterone  
4. 17 $\alpha$ -Methyltestosterone  
5. 1-Dehydro-17 $\alpha$ -methyltestosterone (methandrostenalone)  
6. Norethandrolone  
7. 1-Dehydrotestosterone acetate  
8. 19-Nortestosterone-17-propionate  
9. 1-Dehydrotestosterone benzoate

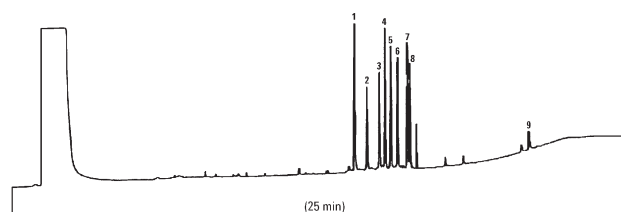

## Steroid Extract from Serum

Column: DB-1  
30 m x 0.25 mm I.D.  
P/N: 122-1031  
Film Thickness: 0.1  $\mu$ m  
Carrier: Helium at 40 cm/sec (measured at 55°C)  
Oven: 55°C for 1 min  
55°C – 175°C at 30°C/min  
175°C – 325°C at 10°C/min  
Injector: Splitless; 275°C  
45 sec. purge activation time  
Detector: FID; 300°C

1. Dehydroisoandrosterone (prasterone)  
2. 10-Nortestosterone (nandrolone)  
3. Testosterone  
4. 17 $\alpha$ -Methyltestosterone  
5. 1-Dehydro-17 $\alpha$ -methyltestosterone (methandrostenolone)  
6. Norethandrolone  
7. 1-Dehydrotestosterone acetate  
8. 19-Nortestosterone-17-propionate

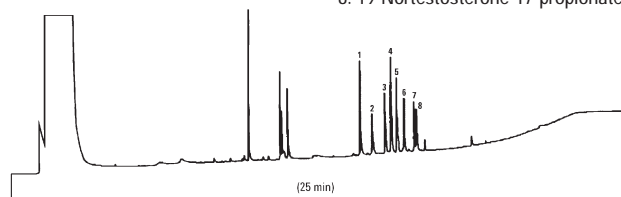

To order or for technical assistance,  
contact your local authorized  
Agilent distributor, or visit our  
website at [www.agilent.com/chem](http://www.agilent.com/chem).

© Agilent Technologies, Inc. 2001

Information, descriptions and specifications in this  
publication are subject to change without notice.

All rights reserved. Reproduction, adaptation,  
or translation without prior written permission  
is prohibited, except as allowed under copyright laws.

Printed in the U.S.A.  
October 25, 2001  
5988-4061EN



Agilent Technologies