

# Studies of Polyvinyl Alcohol under Temperature and Humidity Control

## **Application Note**

Jon Hammerschmidt<sup>1</sup> <sup>1</sup> University of Minnesota

The atomic force microscope (AFM) is capable of imaging insulating materials with atomic resolution, thus making it an excellent tool for studying polymer surfaces. Atomic force microscopy can be used to measure many properties of polymers as well as surface structure, including friction, adhesion, and viscoelasticity. All of these measurements can be done without damaging preparative processes, such as metallic coating or placement in vacuums, which are common with many other analytical techniques. Additionally, AFM measurements can be performed at the molecular level.

Atomic force microscopy was first applied to a polymer surface in 1988 (1); it has subsequently become a standard research tool in the field. This application note will focus on the utility of AFM for investigating various temperature- and humiditydependent properties of polymers.

It is well known that both the environment and the history of a polymer can drastically influence its morphology and its material properties (2). Typically, annealing will drive polymer chains into ordered (crystalline) conformations, depending on interacting chemical functional groups, structure, and stereochemistry of the polymer (2). Not surprisingly, heating a polymer above its glass transition temperature (Tg) gives the polymer chains the mobility to migrate into more energetically favorable orientations. For example, polyvinyl alcohol -(CH2CH0H)n-(PV0H) crystallizes upon heating. For a PVOH sample with only 1-2% water content, annealing at 90°C for up to 10 hours vields approximately 38% crystallinity, whereas heating PVOH to 200 °C for a few minutes produces approximately 60% crystallinity (3).

PVOH is somewhat special in that crystallization can occur in the atactic form. This is attributed to hydrogen atoms and hydroxyl groups (-OH) having roughly similar sizes, which allows the close, regular packing of chains. Strong hydrogen bonding between -OH groups on neighboring chains also tends to hold the molecules in a regular structure despite the lack of stereoregularity (3). For hydrophilic polymers such as PVOH, relative humidity (RH) exerts as strong an influence on the polymer's properties as temperature. When exposed to high humidity, some polymers can have water content of more than 25% (4). Absorption of water causes hydrophilic polymers to swell and elevates the mobility of individual polymer chains. The higher mobility allowed by the solvent greatly affects the polymer's properties. In particular, the glass transition temperature can be lowered, or "plasticized", by more than 60 °C with a solvent (4).

To demonstrate the capabilities of controlled-environment atomic force microscopy, two experiments using a PVOH film were conducted. The film was imaged in situ under precise temperature and humidity control using an Agilent atomic force microscope (AFM) with an environmental chamber and a temperature-controlled sample stage.



#### Temperature

Figure 1 presents lateral force images of a PVOH sample acquired at (a) 22 °C and then at (b) 60 °C. Friction, or lateral force, images were used instead of topography since they often yield better contrast of different morphologies. This is primarily because softer, amorphous regions often produce more dissipative responses than the relatively stiffer, more crystalline regions (5). Humidity was kept constant during imaging, thus resulting in a Tg of approximately 25 °C. Subtle structural changes are especially evident in lateral force images; rounded structures have become longer and thinner, which is characteristic of nanocrystalline regions.

As expected, the polymer started to crystallize slightly when the temperature was elevated. In both images 1(a) and 1(b), the applied force was held constant and the tip-sample adhesive forces were determined to be constant, suggesting that the contact zone did not change appreciably during imaging.

#### Humidity

Figure 2(a) shows topography and friction images of PVOH at 85% RH, where it has approximately 20% water content (4). After exposure to this moist air, the humidity was lowered to 10% RH and the sample was re-imaged. Figure 2(b) reveals large crystalline regions on the surface of the PVOH film at 10% RH. Once again, surface crystallization occurred, but this time crystallization was attributable to extra mobility afforded by the solvent. The swollen film allowed reorganization of polymer chains and then, upon drying, locked a more highly ordered structure in place. The more crystalline film was less dissipative than the amorphous film.

#### Conclusion

The *in situ* control of temperature, humidity, and other environmental parameters is essential when imaging polymers that are sensitive to their surroundings. *In situ* environmental control eliminates many experimental limitations, thus allowing more thorough investigations of polymer morphology and kinetics.



Figure 1. Lateral force images of polyvinyl alcohol acquired at (a) 22 °C and then at (b) 60 °C. Both images are 1.2 μm x 1.2 μm.

Figure 2. Topography (left) and lateral force (right) images of polyvinyl alcohol acquired at (a) 85% RH and then at (b) 10% RH. All images display the same area and are 5 µm x 5 µm.

#### References

- 1. T.R. Albrecht, M.M. Dovek, C.A. Lang, P. Grütter, C.F. Quate, S.N.J. Kuan, C.W. Frank, R.F.W. Pease, J. Appl. Phys., 64, 1178-1184 (1988).
- 2. G. Strobl, The Physics of Polymers: Concepts for Understanding their Structures and Behavior, Springer, Berlin (1996).
- 3. N.G. McCrum, B.E. Read, G. Williams, Anelastic and Dielectric Effects in Polymeric Solids, Wiley, London (1967).
- 4. C.A. Finch, Polyvinyl Alcohol: Properties and Applications, Wiley, London (1973).
- 5. G. Haugstad, W.L. Gladfelter, E.B. Weberg, R.T. Weberg, T.D. Weatherill, Langmuir, 10, 4295-4306 (1994).

### **AFM Instrumentation from Agilent Technologies**

Agilent Technologies offers high-precision, modular AFM solutions for research, industry, and education. Exceptional worldwide support is provided by experienced application scientists and technical service personnel. Agilent's leading-edge R&D laboratories are dedicated to the timely introduction and optimization of innovative and easy-to-use AFM technologies.

#### www.agilent.com/find/afm

#### www.agilent.com

For more information on Agilent Technologies' products, applications or services, please contact your local Agilent office. The complete list is available at:

#### www.agilent.com/find/contactus

Americas	(077) 004 4444
Canada	(8//) 894-4414
Latin America	305 269 7500
United States	(800) 829-4444
Asia Pacific	
Australia	1 800 629 485
China	800 810 0189
Hong Kong	800 938 693
India	1 800 112 929
Japan	81 426 56 7832
Korea	080 769 0800
Malaysia	1 800 888 848
Singapore	1 800 375 8100
Taiwan	0800 047 866
Thailand	1 800 226 008
Europe	
Austria	0820 87 44 11
Belgium	32 (0) 2 404 93 40
Denmark	45 70 13 15 15
Finland	358 (0) 10 855 2100
France	0825 010 700
Germany	01805 24 6333*
	*0.14€/minute
Ireland	1890 924 204
Italy	39 02 92 60 8484
Netherlands	31 (0) 20 547 2111
Spain	34 (91) 631 3300
Sweden	0200-88 22 55
Switzerland (French)	44 (21) 8113811(Opt 2)
Switzerland (German)	0800 80 53 53 (Opt 1)
United Kingdom	44 (0) 7004 666666
Other European Countries:	
www.agilent.com/find/contactus	
Revised: March 23, 2007	

Revised: March 23, 200

Product specifications and descriptions in this document subject to change without notice.

© Agilent Technologies, Inc. 2007 Printed in USA, April 23, 2007 5989-6626EN

