




Surfactants TX-114

Application
Polymers
Robert Ricker

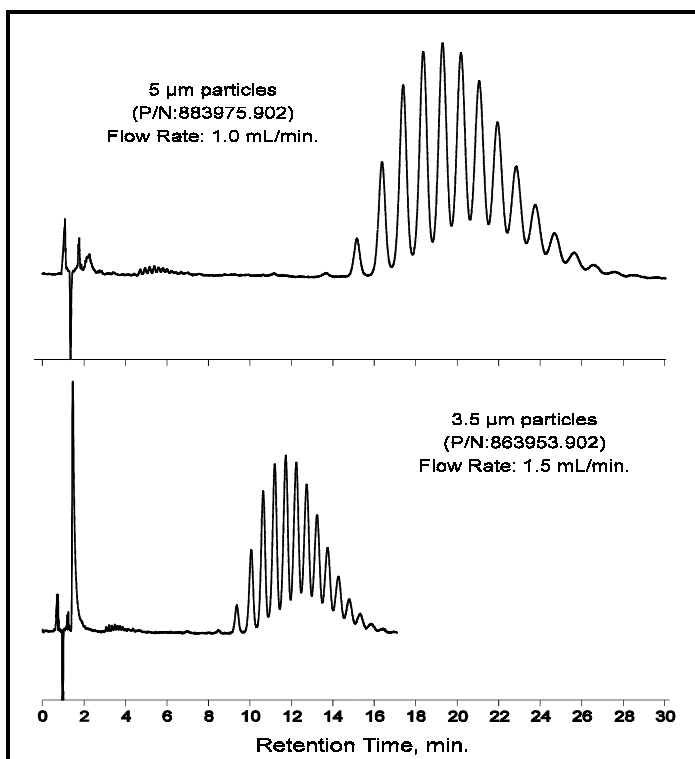

Reproduced from the *Journal of Chromatographic Science*
by permission of Preston Publications, a Division of Preston Industries, Inc.

Conditions:
ZORBAX SB-C18 (4.6 x 150 mm) (Agilent P/N: above)
Mobile Phase: 75% MeOH, 25% H₂O, 50°C, Detec. UV(225 nm)

Highlights

- *StableBond C18 columns are extremely stable at elevated temperatures.*
- *Oligomers of low-molecular-weight polymers can be resolved by reversed-phase techniques.*
- *Smaller packing particles (3.5µm) provide means for faster separations with comparable resolution in half the time.*



Agilent Technologies

*Robert Ricker is an application chemist
based at Agilent Technologies, Wilmington,
Delaware.*

For more information on our products and
services, visit our website at:
www.agilent.com/chem

Copyright© 2002 Agilent Technologies, Inc.
All Rights Reserved. Reproduction,
adaptation or translation without prior
written permission is prohibited, except as
allowed under the copyright laws.

Agilent shall not be liable for errors
contained herein or for incidental or
consequential damages in connection with
the furnishing, performance, or use of this
material.

Information, descriptions, and specifications
in this publication are subject to change
without notice.

Printed in the USA
April 25, 2002
5988-6376EN



Agilent Technologies