NOTICE: Varian, Inc. was acquired by Agilent Technologies in May 2010. This document is provided as a courtesy but is no longer kept current and thus will contain historical references to Varian. For more information, go to **www.agilent.com/chem**.



# Application Note SI-01893

# Synthesis of HIV-1 gp41 using AmphiSpheres™ Resin

Andrew Coffey Varian, Inc.

### Introduction

The 28 residue HIV-1 gp41 ectodomain sequence, H-NEQELLELDKWASLWNWFNITNWLWYIK-NH<sub>2</sub> (de la Torre *et al.*, 2007), is difficult to synthesize. Multiple tryptophan, leucine, asparagine and glutamic acid residues as well as other amino acids requiring side chain protection clearly make this a difficult prospect and would normally require optimization of the synthetic strategy.

## Instrumentation and Conditions

Column: Varian Polaris 5 µm C18-A 150 x 4.6 mm (P/N A2000250X046) Gradient: 10 – 30% MeCN (0.1% TFA), 0 – 30 min Flow rate: 1 mL/min UV detector: 220 nm Injection volume: 20 µL

### Sample Preparation

The peptide was prepared using a peptide synthesizer on a 0.1 mmol scale using 10 fold excess Fmoc-AA-OH (with standard side chain protection as required) and coupling reagent (HBTU) for coupling reactions. Single 9 minute couplings were used throughout. Deprotection reactions (20% piperidine in DMF) were monitored by conductivity meter. Following peptide assembly and cleavage of the final Fmoc group on the synthesizer the peptide was cleaved from the resin by treating with TFA / TIPS / water 95:2.5:2.5 (v/v) for 2-3 h. Cleavage reagents were removed by rotary evaporation followed by trituration of the resultant solid with cold ether. The peptide was then lyophilized from water or water / acetic acid mix.



# **Results and Discussion**

Synthesis of HIV-1 gp41 using AmphiSpheres resin outperformed the synthesis using a commercially available polyether resin (HPLC/MS confirms the major peak as 3653 [M+H]<sup>+</sup>).

### Conclusions

AmphiSpheres 40 RAM successfully synthesized HIV-1 gp41. AmphiSpheres is a significant addition to the StratoSpheres product line from Varian, Inc. AmphiSpheres resins contain hydrophobic polystyrene and hydrophilic polyethyleneglycol components. This imparts a change in the swell characteristics of this amphipathic resin that can help improve synthesis of "difficult" peptide sequences.

### References

de la Torre, B. G. *et al.*, 2007, Int. J. Pept. Prot. Res., **13**, 265-270

These data represent typical results. For further information, contact your local Varian Sales Office.

VARIAN

Varian and the Varian Logo are trademarks of Varian, Inc. <sup>®</sup> Varian, Inc. 2008 Varian, Inc. www.varianinc.com North America: 800.926.3000 – 925.939.2400 Europe: *The Netherlands*: 31.118.67.1000 Asia Pacific: *Australia*: 613.9560.7133 Latin America: *Brazil*: 55.11.3238.0400