

Effect of Mobile-Phase Ionic Strength on Neutral and Strongly Basic Antibodies at pH 7.0

Application


Biochemical

Robert Ricker

Separation of antibodies from antibody conjugants and reaction components is a job well suited to size-exclusion chromatography (SEC). However, mobile-phase ionic strength and pH have a strong influence over proper elution of antibodies and other biomolecules. The chromatograms below show the effect of phosphate buffer concentration on peak shape and retention.

Highlights

- All SEC columns have some electrostatic and hydrophobic characteristics. Therefore, take care to carry out experiments under conditions that promote ideal SEC behavior.
- Strongly basic molecules (right panel) interact with silica-based column packings to a much greater extent than neutral molecules (left panel). Note the complete adsorbance of IgG₃ at low ionic strength; and peak tailing of both antibodies at low and high ionic strength.

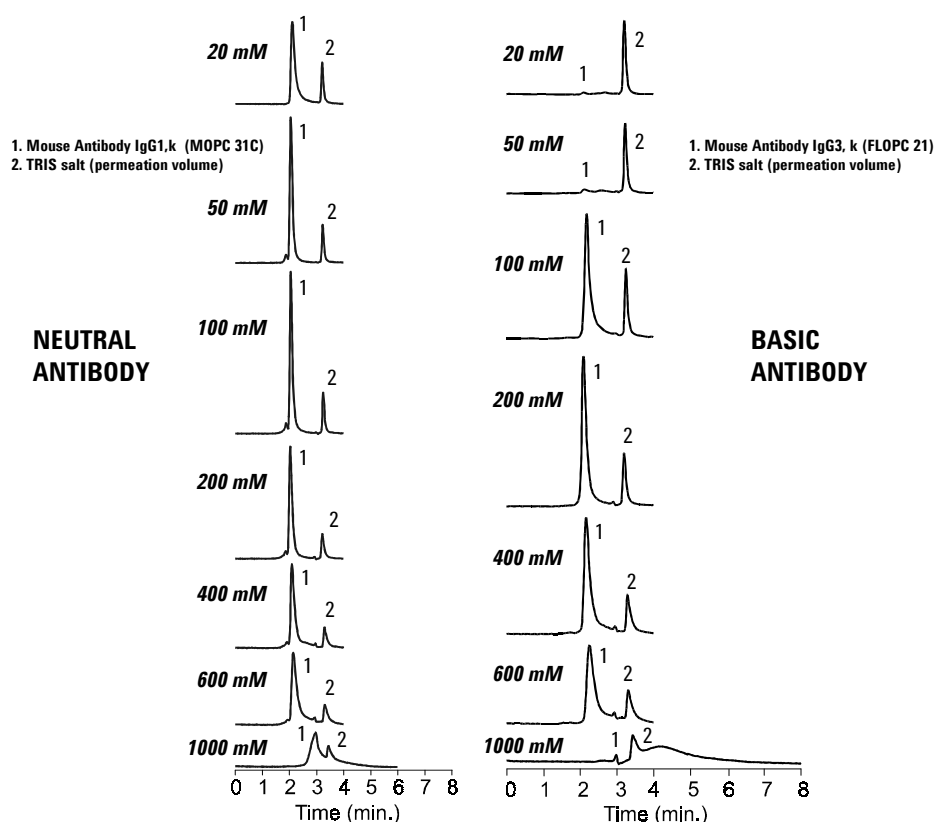

Conditions:
ZORBAX GF-250 (4.6 x 250 mm) (Agilent P/N: 884973-701)
Mobile Phase: Sodium Phosphate, pH 7.0 (concentration as shown)
Injection volume 2µL, 0.96 mL/min, 35°C, Detect. UV (230 nm)



Agilent Technologies

*Robert Ricker is an application chemist
based at Agilent Technologies, Wilmington,
Delaware.*

For more information on our products and
services, visit our website at:
www.agilent.com/chem

Copyright© 2002 Agilent Technologies, Inc.
All Rights Reserved. Reproduction,
adaptation or translation without prior
written permission is prohibited, except as
allowed under the copyright laws.

Agilent shall not be liable for errors
contained herein or for incidental or
consequential damages in connection with
the furnishing, performance, or use of this
material.

Information, descriptions, and specifications
in this publication are subject to change
without notice.

Printed in the USA
April 25, 2002
5988-6315EN



Agilent Technologies