

Using LC/MS/MS 6410 for Analysis of Chloramphenicol, Thiamphenicol, and Florfenicol in Fish Samples

Application Brief

Xiaorong Ran, Tao Bo

Chloramphenicol is a banned compound in the EU. It is a zero tolerance compound, and some methods have already been developed for this analysis. The LC/MS/MS is often used for its greater sensitivity and higher selectivity.

In many cases, not only chloramphenicol but also thiamphenicol and florfenicol are also found. This method uses a simple method for detecting all three compounds within less than 6 minutes. Furthermore, the results gave the good results, showing the performance at the negative mode.

Experimental

LC Conditions

Column	Agilent ZORBAX Eclipse Plus, 2.1 mm × 50 mm, 1.8 μm
Mobile phase	A: Water B: Methanol
Flow rate	0.4 mL/min
Gradient	0–2 min/B 10% to 90%; 2–3 min/B 90%; 3.01/B 10%
Stop time	6 min
Column compartment temperature	45 °C
Injection	5 μL

MSD Condition

Instrument	Agilent 6410A triple quadrupole LC/MS system
Source	ESI –

Highlights

- Using a simple RRLLC method can separate the compounds well within 6 minutes
- Quite high sensitivity in negative mode
- The ISTD method can remove the matrix effect and minimize the sample preparation interference



Agilent Technologies

Sample Preparation

All SPE cartridges are conditioned with 2 mL of water before use.

1. **Honey**, 1 g sample is diluted to 5 mL with water and 25 μ L 10 ppb IS is added. The solution is loaded onto the SPE cartridge and allowed to stand for 5 min. Elution is performed with 10 mL ethyl acetate. The eluate is collected and the solvent is evaporated under a nitrogen stream at 40 °C. The residue is redissolved in 1 mL methanol and put in an ultrasonic bath for 1 min. The solution is filtered, using a syringe filter, before injection. No additional cleanup of the sample solution is performed.
2. **Shrimp**, 1 g of shrimp is defrosted and mixed in a blender. To the 1 g of the mixed shrimp, 3 mL of water and 25 μ L 10 ppb IS are added. The portion is centrifuged for 5 min (8,000 rpm). The supernatant is loaded on the cartridge and allowed to stand for 5 min. Elution is performed with 5 mL ethyl acetate. The eluate is collected and the solvent evaporated under a nitrogen stream at 40 °C. The residue is redissolved in 1 mL methanol and put in an ultrasonic bath for 1 min; the solution is filtered before injection.
3. **Chicken**, 1 g of chicken is defrosted and mixed in a blender. To the 1 g of the mixed chicken, 3 mL of water and 25 μ L 10 ppb IS are added. The portion is centrifuged for 5 min (8,000 rpm). The supernatant is loaded on the cartridge and allowed to stand for 5 min. Elution is performed with 5 mL ethyl acetate. The eluate is collected and the solvent evaporated under a nitrogen stream at 40 °C. The residue is redissolved in 1 mL methanol and put in an ultrasonic bath for 1 min; the solution is filtered before injection.

MRM Setting


Name	Precursor ion	Product ion	Frag (V)	CE (V)	Dwell (ms)
TAP	354.1	185.1*	120	20	60
	354.1	289.9	120	10	60
FF	356	185.1*	120	20	60
	356	335.8	120	5	60
CAP	321.2	152.1*	120	10	60
	321.2	257.1	120	5	60
D5-CAP (ISTD)	326.2	157.2	130	15	60

Results


Name	Linearity (0.5–20 ppb)
TAP	0.994
FF	0.992
CAP	0.994

Sensitivity

1. CAP: 0.5 ppb S/N = 81

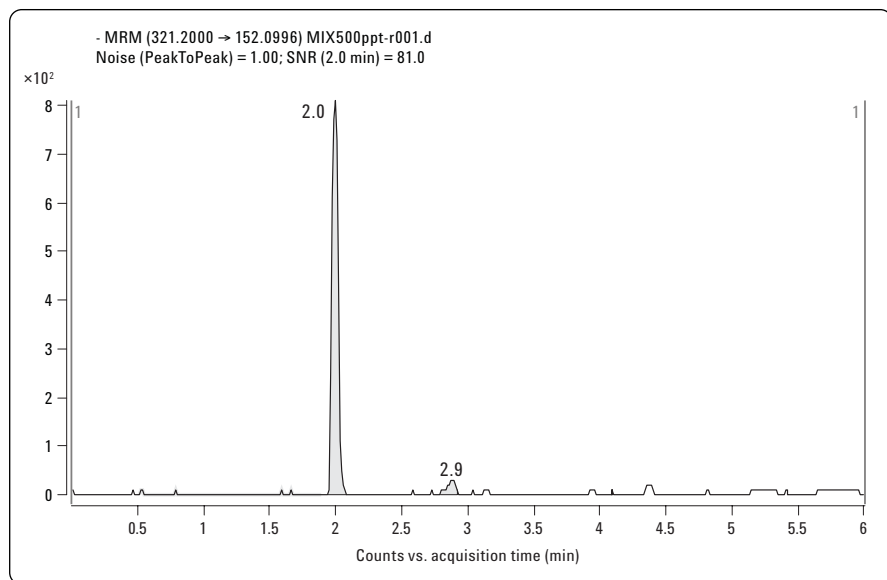

2. FF: 0.5 ppb S/N = 56

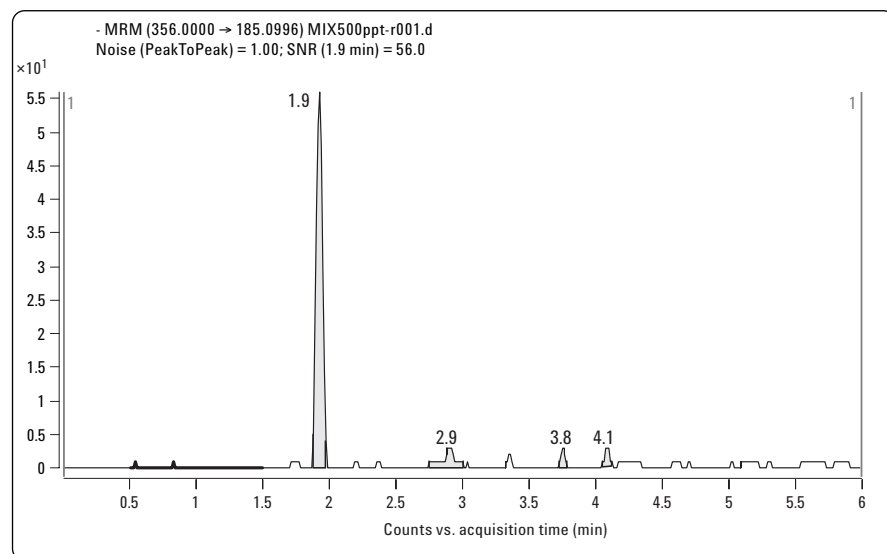


Xiaorong Ran, Tao Bo are application chemist based at Agilent Technologies, Beijing, China

For More Information

For more information on our products and services, visit our Web site at www.agilent.com/chem.

3. TAP 1 ppb, S/N = 12

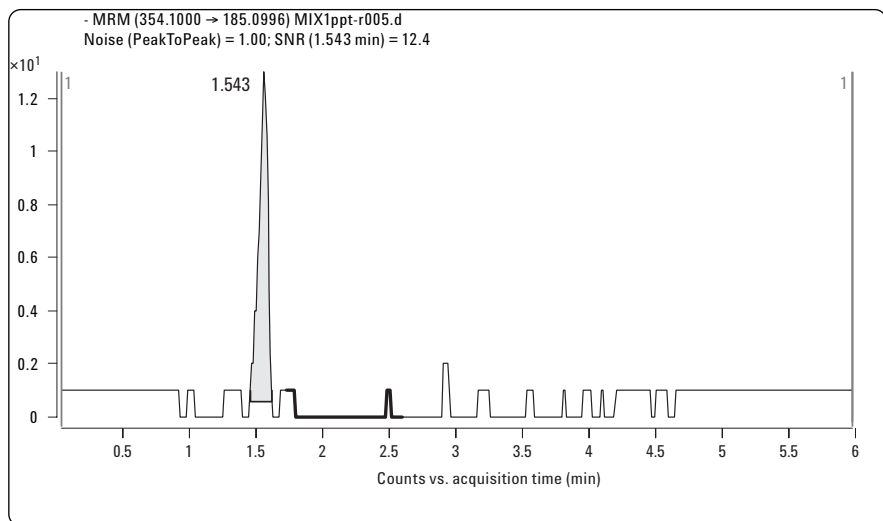

www.agilent.com/chem

Agilent shall not be liable for errors contained herein or for incidental or consequential damages in connection with the furnishing, performance, or use of this material.

Information, descriptions, and specifications in this publication are subject to change without notice.

© Agilent Technologies, Inc., 2009
Printed in the USA
June 15, 2009
5990-3734EN



Agilent Technologies