

Analysis of Fragrance and Allergens using an Agilent J&W FactorFour VF-WAXms GC Column

Application Note

Author

John Oostdijk
Agilent Technologies, Inc.

Introduction

Allergic reactions to perfumes and the individual components responsible for these reactions have been studied since the late 1970s. The fragrance mix used internationally to diagnose fragrance-allergic patients includes eight ingredients: oak moss extract, iso-eugenol, eugenol, cinnamal, hydroxy citronellal, geraniol, cinnamyl alcohol and amyl cinnamal. Analysis of this particular fragrance mix on a VF-WAXms reveals the presence of a number of these common allergens. The MS compatibility of VF-WAXms is clearly evident, especially in the low bleed profile around 250 °C , as shown in Figure 1 on page 2.



Agilent Technologies

Conditions	Peak Identification:	
Technique: GC-MS	1. <i>Linalool</i>	14. <i>Eugenol</i>
Column: VF-WAXms	2. <i>Methyl heptin carbonate</i>	15. <i>Amyl cinnamyl aldehyde</i>
	3. <i>Phenyl acetaldehyde</i>	16. <i>Anisic alcohol</i>
30 m x 0.25 mm x 0.25 µm (part number CP9205)	4. <i>Methyl chavicol</i>	17. <i>Cinnamyl alcohol</i>
Oven: 100 °C to 250 °C with 10 °C / min	5. <i>Methyl octin carbonate</i>	18. <i>Farnesol isomer I + II</i>
Carrier Gas: Helium, 1.0 mL/min	6. <i>Citronellol</i>	19. <i>Farnesol isomer III</i>
Injector: Split 1:30, T = 250 °C	7. <i>Geraniol</i>	20. <i>iso-Eugenol</i>
Detector: GC-MS Iontrap	8. <i>Methyl gamma ionone</i>	21. <i>Hexyl cinnamic aldehyde</i>
Trap: 200 °C	9. <i>Benzyl alcohol</i>	22. <i>Lyral (4,4-isomer)</i>
Manifold: 60 °C	10. <i>Cinnamaldehyde</i>	23. <i>Coumarine</i>
Sample Size: 0.1 µL	11. <i>Hydroxy citronellal</i>	24. <i>Amyl cinnamic alcohol</i>
Sample: Fragrances mixture (500 ppm)	12. <i>Methyl eugenol</i>	25. <i>Benzyl benzoate</i>
	13. <i>Lilial</i>	26. <i>Benzyl salicylate</i>
		27. <i>Benzyl cinnamate</i>

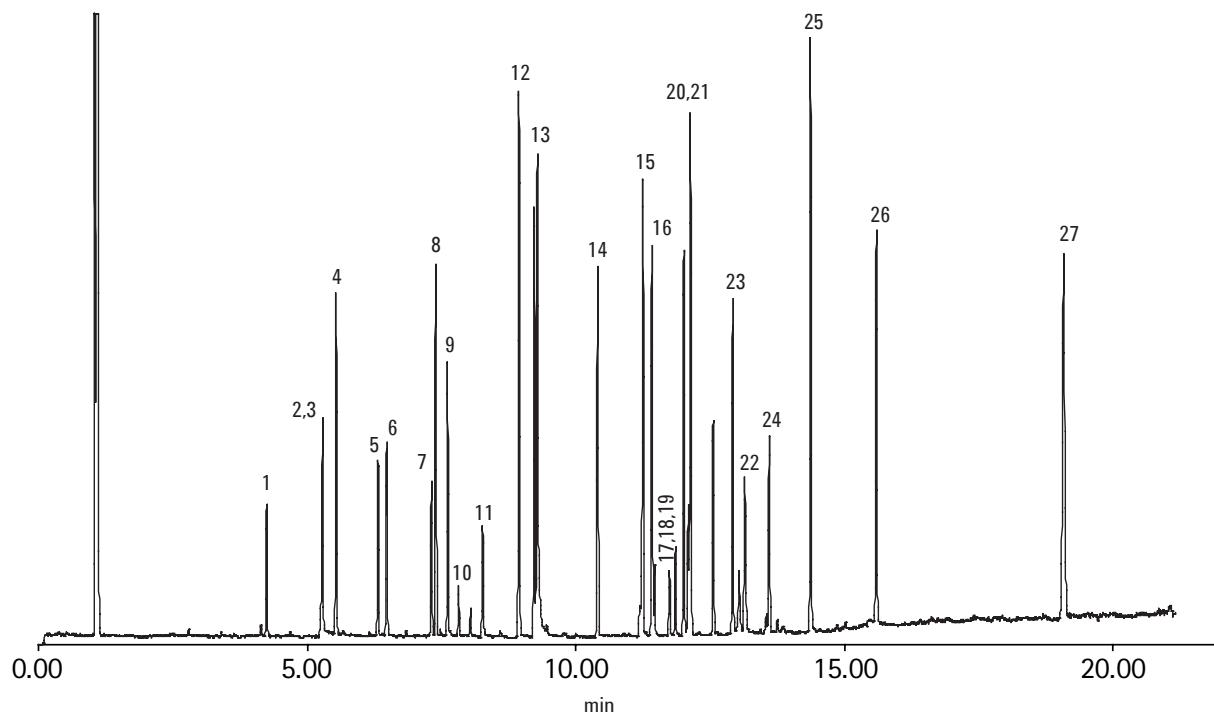


Figure 1. Analysis of fragrance and allergens using VF-WAXms GC columns

www.agilent.com/chem

This information is subject to change without notice.

© Agilent Technologies, Inc. 2010

Published in UK, October 08, 2010

SI-01061



Agilent Technologies