



Analysis of Polyvinyl Pyrrolidone with DMF

Application Note

Materials Testing and Research, Polymers

Authors

Greg Saunders, Ben MacCreath
Agilent Technologies, Inc.

Introduction

Agilent PLgel 10 μ m MIXED-B columns are designed for high MW polymer analysis and demanding eluent conditions such as polyvinyl pyrrolidone in dimethyl formamide.

Polyvinyl pyrrolidone is soluble in both polar organics and water. DMF is used with LiBr to minimize any polyelectrolyte effects. The response in DMF is relatively small using an RI detector. Polyvinyl pyrrolidone can be chromatographed successfully with PLgel MIXED-B columns.



Agilent Technologies

The PLgel MIXED-B spans a wide range of molecular weights, up to 10 million, with a linear calibration curve. It is particularly useful for molecular weight distributions where slightly higher than average MWs are encountered. The 10 μm particle size provides good resolution with relatively low pressures for enhanced lifetimes in demanding conditions.

Conditions

Columns:	2 x PLgel 10 μm MIXED-B, 300 x 7.5 mm (part number PL11110-6100)
Eluent:	DMF + 0.1% LiBr
Flow Rate:	1.0 mL/min
Temperature:	60 °C
Detection:	390-MDS Multi Detector Suite (differential refractive index)

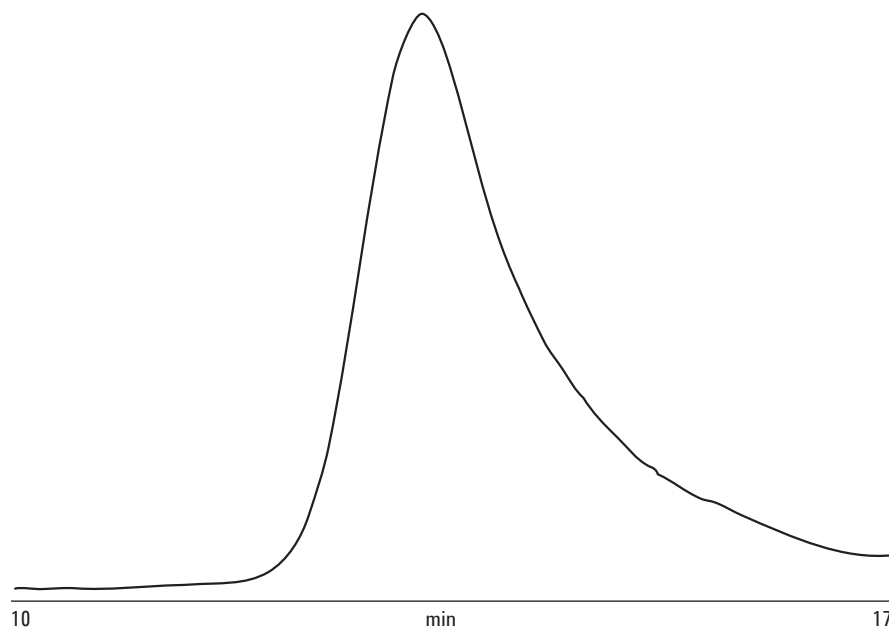


Figure 1. Analysis of polyvinyl pyrrolidone using PLgel 10 μm MIXED-B columns

www.agilent.com/chem

This information is subject to change without notice.

© Agilent Technologies, Inc. 2010

Published in UK, September 29, 2010

SI-1781



Agilent Technologies