

N-terminal Sequencing of 10 pmols of Human Serum Albumin using Routine 3.1 Method on the HP G1005A N-terminal Protein Sequencing System

James Kenny, Hewlett-Packard
California Analytical Division

**HP G1005A
N-terminal Protein
Sequencing System**

AB 96-2

Application Brief

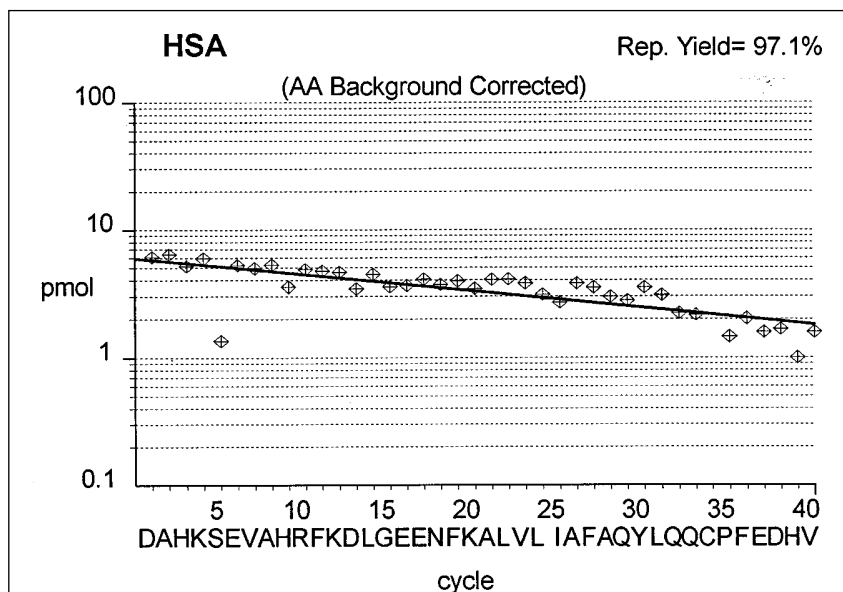


Fig 1. Repetitive yield plot for the sequence analysis of 10 pmols HSA.

Abstract

This Application Brief describes the N-terminal sequence analysis of a 10 pmol of human serum albumin using the HP G1005A N-terminal protein sequence with Routine 3.1 method

Keywords

Sequencing, protein, G1005A

Introduction

The HP G1005A N-terminal Protein Sequencing System performs automated Edman chemistry of peptides and proteins loaded in solution and immobilized on miniature biphasic columns consisting of reverse-phase (RP) sample column mated to a strong anion exchange (SAX) column.^{1,2}

The biphasic support facilitates efficient, simple and rapid sample clean-up and concentration as part of the sample loading procedure. Samples can be directly loaded and retained on the RP column in a variety of buffer matrices (salts, denaturants, detergents), washed free of the contaminants, coupled to the SAX column, and sequenced.

Several performance characteristics of the HP G1005A N-terminal Protein Sequencing System and Routine 3.1 chemistry contribute to the routine sequence analysis of proteins at low picomole level.

Features of the N-terminal Routine 3.1 chemistry are:

- **Extremely efficient Edman-based N-terminal sequencing chemistry.** Typical initial yields for pmol amounts are approximately 50%. Repetitive yields can be greater than 95%, permitting the confident assignment of 40 or more amino acid residues.

- **Reproducible PTH-amino acid retention times and gradient chromatography.** Stable retention times enable chromatogram subtraction or overlay for accurate cycle-to-cycle assignment.
- **Stable and low chemical background.** Confident assignment of cycles yielding less than 1 pmol of PTH-amino acid requires this sufficiently low background.

Methods

The sequence analysis presented here was performed human serum albumin (HSA, 68kDa) protein. A sample aliquot of 10 pmols was diluted in 1 mL of aqueous TFA and directly loaded on a reverse-phase sample column (HP G1073A), then mated to a strong anion exchange (SAX) column (HP G1448A). Prior to sample loading this sequencing column was prepared using Column Prep 3.1/3.5 method. Analysis was carried out using the N-terminal Routine 3.1 sequencer method and PTH_4.M HPLC method on an HP G1005A N-terminal Protein Sequencing System with standard Routine 3.1 reagents and solvents.

Results and Discussion

Shown in Figure 1, the initial yield (cycle 1) of HSA was 6.1 pmol of PTH-Asp, corresponding to an overall repetitive yield of 97.1%.

Representative chromatograms of 10 pmols HSA show the unambiguous assignment of the PTH-amino acid residues and the low chemical background. A PTH-amino acids residues are well-resolved and separated from the Edman chemistry by-products of DPTU (as identified in the figure) and DPU (which elutes after PTH-Leu).

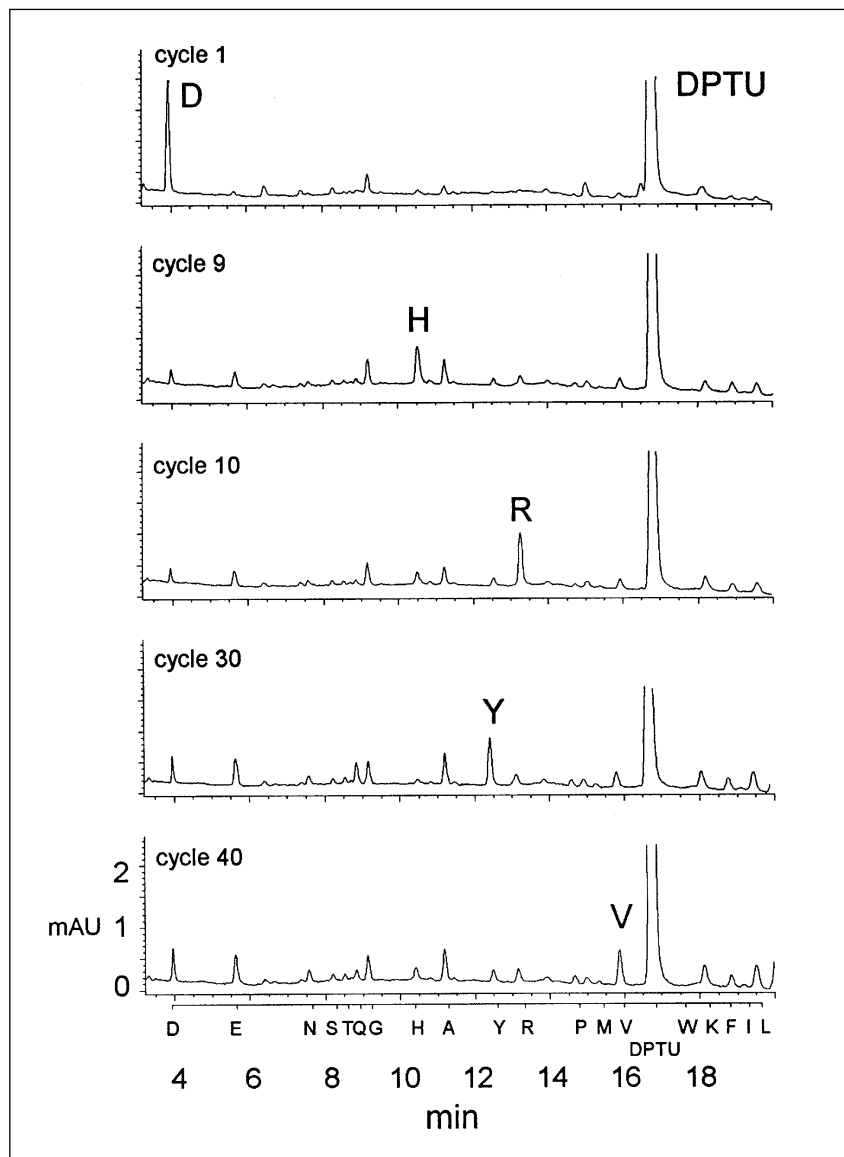


Fig 2. Results of sequence analysis of 10 pmols of HSA using "Routine 3.1" sequencing method.

References

1. Hewlett-Packard Technical Note. "Routine 3.0 Sequencer Methods," TN 95-1.
2. Chad G. Miller, "Adsorptive Biphasic Column technology for Protein Sequence Analysis and Protein Chemical Modification," METHODS: A Companion to Methods in Enzymology, **6**, 315-333 (1994).

Hewlett-Packard shall not be liable for errors contained herein or for incidental or consequential damages in connection with the furnishing, performance or use of this material.

Information, descriptions and specifications in this publication are subject to change without notice.

Copyright © 1996
Hewlett-Packard Company
All rights reserved. Reproduction and adaptation is prohibited.

Printed in the U.S.A. July 1996
(23) 5965-3464E