NOTICE: This document contains references to Varian. Please note that Varian, Inc. is now part of Agilent Technologies. For more information, go to **www.agilent.com/chem.**



QAII Station Operator's Manual

P/N 70-9018 November 2010 Revision K

Limitation of Liability

The information in this document is subject to change without notice. Varian, Inc. makes no warranty of any kind with regard to this material, including, but not limited to, the implied warranties or merchantability and fitness for a particular purpose. Varian, Inc. shall not be liable for errors contained herein or for incidental consequential damages in connection with the furnishing, performance, or use of this material.

All rights are reserved. Reproduction, adaptation, or translation without prior written permission is prohibited, except as allowed under copyright laws.

First Edition (November 2010) **QAII Station Operator's Manual** Part Number 70-9018 Revision K Printed in the United States of America

The following terms are trademarks of Varian, Inc.:

- BenchsaverTM
- BIO-DIS III[®]
- Enhancer Cell[®] Full Flow FiltersTM
- Peak VesselTM
- TruCenterTM
 VanKel[®]

 Practical Solutions[®] QA and QAIITM

VK[®]

Varian, Inc. 13000 Weston Parkway Cary, North Carolina 27513-2250 800.229.1108 919.677.1108 Fax: 919.677.1138 www.varianinc.com



Copyright (c) 1999 - 2010 by Varian, Inc.

Chapter 1	Safety Practices and Hazards 9
	Electrical Hazards 10 Other 10 General 12
	WEEE Directive 13
Chapter 2	Introduction 15
	Base Unit 16 Conventions Used in this Manual 17
Chapter 3	Installation and Setup 19
	Unpacking Your QAII Station 19 QAII Attachments 20 Digital Wobble Gauge 20

Magnetic Tachometer Sensor 21 Magnetic Clip 21 Universal Adapter 22 Electronic Level Sensor 22 Electronic Vibration Sensor 23 Temperature Probe 23 Power Switch Functions 24 Setting up the QAII Station 25

Chapter 4

Operation 27

Placing the Universal Adapter on Apparatus without EaseAlign Centering Rings 28
Operating the QAII Station 29

Main Menu Option 2, Setup 30
Main Menu Option 3, Vibration 30

Running a Test 31
Printing Results 41
Storage Table and Viewing Data 42

Chapter 5

Maintenance and Troubleshooting 45

Periodic Maintenance 45 Calibration 45 Report Center Impact Printer 46 Installing the Cartridge Ribbon 46 Replacing the Paper Roll 47 Toggling Your Printer Online 48 Printer Self Test 49 Printer Configuration 49 Fuse Replacement 51 Troubleshooting 52 QAII Default Values 54

Chapter 6

Service and Warranty 55

Exclusions and Limitations56Obtaining Warranty Service56Warranty Limitations56Exclusive Remedies57

Index 59

Tell Us How We Are Doing 61

This page was intentionally left blank, except for this message.

List of Figures

FIGURE 1. QAI	I Station 16
FIGURE 2. Digi	tal Wobble Gauge 20
FIGURE 3. Mag	gnetic Tachometer Sensor 21
FIGURE 4. Mag	gnetic Clip 21
FIGURE 5. Univ	versal Adapter 22
FIGURE 6. Elec	ctronic Level Sensor 22
FIGURE 7. Elec	ctronic Vibration Sensor 23
FIGURE 8. Tem	perature Probe 23
FIGURE 9. QAI	I Station Rear Panel 25

List of Figures

FIGURE 10.	EaseAlign Mounting Bracket Secured to Dissolution
	Apparatus 28

- FIGURE 11. Placement for RPM Reading 34
- FIGURE 12. Wobble Gauge with Paddle 36
- FIGURE 13. Wobble Gauge with Basket 37
- FIGURE 14. Electronic Level Sensor Placement 38
- FIGURE 15. Vibration Sensor 40

Chapter 1 Safety Practices and Hazards

The QAII Station has been carefully designed so that when used properly you have an accurate, fast, flexible, and safe instrument.

If the equipment is used in a manner not specified by the manufacturer, the protection provided by the equipment may be impaired.

The QAII Station is operated in conjunction with equipment that uses aqueous liquids. Unskilled, improper, or careless use of this instrument can create shock hazards, fire hazards, or other hazards which can cause death, serious injury to personnel, or severe damage to equipment and property.

Information on safety practices is provided with your instrument and operation manuals. Before using your instrument or accessories, you must thoroughly read these safety practices.

Observe all relevant safety practices at all times.

Electrical Hazards

The QAII Station contains electrical circuits, devices, and components operating at dangerous voltages. Contact with these circuits, devices, and components can cause death, serious injury, or painful electric shock.

Panels or covers that are retained by fasteners which require the use of a tool for removal may be opened only by Varian-trained, Varian-qualified, or Varian-authorized service engineers. Consult the manuals or product labels supplied with the QAII Station to determine which parts are operator-accessible.

Application of the wrong supply voltage, connection of the instrument to an incorrectly wired supply outlet, or lack of proper electrical grounding can create a fire hazard or a potentially serious shock hazard and could seriously damage the instrument and any attached ancillary equipment.

Always use a three-wire outlet with ground connection which is adequately rated for the load. The installation must comply with local, state, and federal safety regulations.

Do not connect the instrument to the main power supply until you have made sure that the operating voltage is correctly set for the main power supply in the specific outlet in your laboratory to which the equipment will be connected.

Other

Other specific warnings and cautions appear in the manuals where appropriate and detail the specific hazard, describe how to avoid it, and specify the possible consequences of not heading the warning or caution.

Warning

A 'Warning' message appears in the manual when failure to observe instructions or precautions could result in death or injury. Symbols depicting the nature of the specific hazard are also placed alongside warnings.

Revision K, 11/10 P/N 70-9018	QAII Station	Page 11 Safety Practices and Hazards
F/N /0-5010		

These symbols are also used on warning labels attached to the instrument. When you see one of these symbols, you must refer to the relevant operation or service manual for the correct procedure referred to by that warning label.

The meaning of the symbols that appear alongside warnings in this manual are as follows:



Electrical shock



Caution Refer to accompanying documents

Read all warnings and cautions carefully and observe them at all times.

Caution

A 'Caution' message appears in the manual when failure to observe instructions could result in damage to equipment (Varian supplied and / or other associated equipment).



A 'Note' appears in the manual to give advice or information.



Varian, Inc.

Page 12 Safety Practices and Haz	QAII Station ards Operator's Manual	Revision K, 11/10 P/N 70-9018
Information Symb	ools	
I	Switches main power on	
0	Switches main power off	
\sim	Indicates single-phase alternating current	
CE	Indicates the product complies with the re or more European Union (EU) directives	quirements of one
C Manufactor US	Indicates specific equipment meets conse standards of safety to provide assurance, OSHA, that these products are safe for us for North America	nsus-based required by e in the workplace
	Indicates that this product must not be dis unsorted municipal waste (see "WEEE Dir page 13)	posed of as rective" on

General

CE Compliant Products

The QAII Station has been designed to comply with the requirements of the Electromagnetic Compatibility (EMC) Directive and the Low Voltage Directive (LVD) of the EU.

Varian, Inc. has confirmed that each product complies with the relevant directives by testing a prototype against the prescribed European Norm (EN) standards.

Revision K, 11/10	QAII Station	Page 13
P/N 70-9018	Operator's Manual	Safety Practices and Hazards

Proof that a product complies with the directives is indicated by:

- the CE marking appearing on the rear of the product.
- the documentation package that accompanies the product containing a copy of the declaration of conformity. This declaration is the legal declaration by Varian, Inc. that the product complies with the directives and also shows the EN standards to which the product was tested to demonstrate compliance. The declaration of conformity is signed by the representative of the manufacturing plant.

cTUVus - U.S. and Canadian Product Approvals

The QAII Station has been designed to comply with North American safety requirements.

This product has been tested and certified for the North American market by TUV Rheinland of North America, Inc. The TUVus mark signifies that this product has been tested to U.S. standards and certified for the U.S. market. The cTUV mark signifies that this product has been tested to Canadian standards and certified for the Canadian market. When the two marks are coupled, the cTUVus mark signifies that this product has been tested to standards and certified for both markets.

WEEE Directive

All Varian products that are subject to the WEEE directive shipped after August 13, 2005 are compliant with the WEEE marking requirements. Such products are marked with the "crossed out wheelie bin" WEEE symbol shown on page 12 in accordance with European Standard EN 50419.

This symbol on the product or on its packaging indicates that this product must not be disposed of as unsorted municipal waste. The separate collection and recycling of your waste equipment at the time of disposal will help to conserve natural resources and ensure that it is recycled in a manner that protects human health and the environment.

For more information on collection, reuse, and recycling systems, please contact your local/regional waste administration, your local distributor, or Varian, Inc.

Page 14	QAII Station	Revision K, 11/10
Safety Practices and Hazards	Operator's Manual	P/N 70-9018

This page was intentionally left blank, except for this message.

Chapter 2 Introduction

The QAII Station is designed to help you easily check the critical physical parameters specified in the USP General Chapters <711> Dissolution and <724> Drug Release.

The QAII Station tests and documents the parameters of the following variables:

- shaft wobble (also known as runout)
- instrument level
- vessel temperature
- spindle speed
- vibration

Regulatory authorities emphasize these factors since they can lead to major variances in your dissolution results. Test data from up to 30 dissolution apparatus can be stored in the QAII Station's non-volatile memory. The data can be printed via the Report Center Printer. The QAII Station records the serial numbers of the apparatus in addition to test values.

The built-in printer provides a detailed hard-copy calibration report listing all parameter values for each vessel position. If you use Varian serialized paddles or basket shafts, you can enter the shaft serial numbers for each dissolution apparatus in your lab and have these numbers print on the report as well.

Page 16	QAII Station	Revision K, 11/10
Introduction	Operator's Manual	P/N 70-9018

Base Unit

The base unit contains the keypad, display screen and Report Center Printer. This unit stores information for up to 30 apparatus. The five remote sensors plug into the keyed jacks on the rear panel. The base unit has a rechargeable battery for portable use and its own AC power supply for printing or charging the battery. There is a power switch on the left side panel that controls the battery power to the unit and there is a power switch on the rear panel that controls AC power to the unit. You must operate under AC power in order to print via the Report Center Printer. A certified calibration shaft is included and guaranteed to be within USP specifications.

FIGURE 1. QAII Station





Caution

Panels or covers that are retained by fasteners which require the use of a tool for removal may be opened only by Varian-trained, Varian-qualified, or Varian-authorized service engineers.

QAII Station Operator's Manual Page 17 Introduction

Conventions Used in this Manual

- Items you are asked to press are in bold. For example, "press **TEST** on the keypad" or "Press **ENTER**".
- Key sequences you are asked to press appear like this: **MENU** > 2.



Note Remember to return the warranty card supplied with this manual. Completing and returning the card ensures your right to protection under the terms and conditions of your warranty. It also enables us to better assist you in the event of any problems. Additionally, it guarantees you will be informed of any issues that arise concerning your equipment, such as upgrades, retrofits, or regulatory changes.

Page 18	QAII Station	Revision K, 11/10
Introduction	Operator's Manual	P/N 70-9018

This page was intentionally left blank, except for this message.

Chapter 3 Installation and Setup

Unpacking Your QAII Station

Follow these steps to safely unpack your QAII Station:

- Step 1. Open each carton and check the contents for damage which may have occurred during shipping. Shipping damage rarely occurs, but if it does contact both the carrier who delivered the instruments and the Dissolution Systems Service Department. Though claims for damage should be filed with the carrier, we can help you file a claim.
- Step 2. Carefully remove the QAII Station base unit, digital wobble gauge and magnetic tachometer sensor (attached to an EaseAlign mounting bracket and universal adapter), magnetic clip, electronic level sensor, precision temperature probe, electronic vibration sensor, calibration shaft, and all the cables from the shipping carton.
- Step 3. Remove as much cushioning material and tape as possible.

Page 20	QAII Station	Revision K, 11/10
Installation and Setup	Operator's Manual	P/N 70-9018

Step 4. Place the base unit and accessories on a clear, dry, level section of the bench top close to the dissolution apparatus. The preferred placement of the QAII Station is on the right side of the dissolution apparatus.



QAII Attachments

The QAII comes with attachments necessary to measure and document individual spindle speeds, shaft wobble, shaft perpendicularity, temperature, instrument level, and vibration.

Digital Wobble Gauge

Mounted on a self-aligning bracket (the EaseAlign mounting bracket), the wobble gauge has a built-in display screen and indicates wobble in terms of inches or millimeters. A power switch preserves the life of the lithium battery. The ORIGIN button makes it easy to zero the gauge prior to starting the spindles. A cable transmits the gauge reading to the QAII Station base unit.



FIGURE 2. Digital Wobble Gauge

Varian, Inc.

Revision K, 11/10	QAII Station	Page 21
P/N 70-9018	Operator's Manual	Installation and Setup

Magnetic Tachometer Sensor

The magnetic tachometer sensor is also mounted on the EaseAlign mounting bracket. The sensor detects the rotations per minute of the paddle or basket shaft with the use of a magnetic clip.

FIGURE 3. Magnetic Tachometer Sensor



Magnetic Clip

The magnetic clip is used in conjunction with the magnetic tachometer sensor. Each time the magnet passes the sensor, one rotation is counted.

FIGURE 4. Magnetic Clip



Page 22	QAII Station	Revision K, 11/10
Installation and Setup	Operator's Manual	P/N 70-9018

Universal Adapter

The supplied universal adapter allows the QAII Station to be used in conjunction with any Varian dissolution apparatus, including older models, or with any non-Varian dissolution apparatus. The adapter provides locating pegs to secure the EaseAlign mounting bracket with wobble gauge and tachometer sensor to the dissolution apparatus vessel.

FIGURE 5. Universal Adapter



Electronic Level Sensor

The electronic level sensor allows you to check the level of the vessel plate and drive unit as well as the perpendicularity of each shaft from side-to-side and front-to-back, simultaneously. The sensor has a range of -9.9 to 9.9 degrees.

FIGURE 6. Electronic Level Sensor



P/N 70-9018	QAII Station Operator's Manual	Page 23 Installation and Setup
		-

Electronic Vibration Sensor

The electronic vibration sensor can be placed either on the vessel plate or on the drive unit in the upright position to obtain readings. The sensor can be used to measure vibration on the X, Y, and Z axes individually.

FIGURE 7. Electronic Vibration Sensor



Temperature Probe

The temperature probe allows you to measure the temperature of the medium in the vessel. The temperature probe can also be used to measure the temperature of the water in the water bath or any other aqueous-based solutions.

FIGURE 8. Temperature Probe





Page 24	QAII Station	Revision K, 11/10
Installation and Setup	Operator's Manual	P/N 70-9018

Power Switch Functions

WarningImage: Description of the systemImage: Description of the systemImage

The QAII Station has a main power switch and a battery power / charge switch.

The main power switch is located on the rear panel. See Figure 9, "QAII Station Rear Panel," on page 25. With the switch in the ON position, the QAII Station is on and you can charge the battery and print.

The battery power / charge switch is on the left side panel. When AC power is not present, such as when the power cord is not attached or when the main switch is off, the battery power /charge switch allows battery power for data collection and monitoring. If the AC power is on, the battery power / charge switch is used to charge the battery only.

It is recommended that you fully discharge the battery before the next charge to extend the battery life. When the power is weak, leave the battery switch on overnight without AC power. This drains the remaining power so that it can be recharged. Each charge period takes five to six hours. Note that the internal printer can be operated only with AC power.

Setting up the QAII Station

FIGURE 9. QAII Station Rear Panel





Ensure the AC power setting is at the correct voltage for your power supply. The power setting is indicated on the line fuse holder located next to the on/off switch on the instrument back panel. To change the voltage, see "Fuse Replacement" on page 51.

The electrical connection at the back of the QAII is the primary disconnect for the instrument.

- Step 1. Ensure the power switches on the rear panel and left side panel are in the off position.
- Step 2. Connect the cables between the remote sensors and their corresponding jacks on the QAII Station rear panel. Each cable has a different kind of connector making it impossible to plug it into the wrong jack.
- Step 3. Connect the AC power cord between the receptacle on the rear panel and an appropriate AC power receptacle.



Page 26	QAII Station	Revision K, 11/10
Installation and Setup	Operator's Manual	P/N 70-9018

This page was intentionally left blank, except for this message.

Chapter 4 **Operation**

The QAII Station can be used in conjunction with any Varian dissolution apparatus, including older models, or with any non-Varian dissolution apparatus.

For dissolution apparatus that do not have EaseAlign centering rings, the QAII Station comes with a universal adapter. This adapter mounts to the dissolution apparatus vessel in the same manner as the EaseAlign centering rings. Once this adapter is in place, the procedures for using the QAII Station are the same for all dissolution apparatus. The purpose of the adapter is to provide the locating pegs to secure the EaseAlign mounting bracket to the dissolution apparatus (see "locating pegs" in Figure 10 on page 28).

Page 28	QAII Station	Revision K, 11/10
Operation	Operator's Manual	P/N 70-9018



FIGURE 10. EaseAlign Mounting Bracket Secured to Dissolution Apparatus

Placing the Universal Adapter on Apparatus without EaseAlign Centering Rings

	Note Before beginning operation, the paddles or basket shafts must be set to their proper operating heights.
Step 1.	Raise the dissolution apparatus drive unit.
Step 2.	Position the universal adapter over the vessel and push down gently so the adapter is held in place by the short legs that rest against the inside of the vessel. The bracket should fit without having to be forced and it should sit flat on the dissolution apparatus vessel plate.

Revision K, 11/10	QAII Station	Page 29
P/N 70-9018	Operator's Manual	Operation

Operating the QAII Station

- Step 1. Ensure the power switch is in the OFF position.
- Step 2. Ensure all accessories are connected to the QAII Station rear panel. See "Setting up the QAII Station" on page 25.
- Step 3. Turn on the QAII Station. The monitor screen displays.

RPM 000.00_	TEMP -1.5
WOB .0000	VIB 0.000
LV-X -6.8	LV-Y -6.8
09/30/04	15:35:30

Step 4. Press **MENU** to display the QA Main Menu screen.

QA MAIN MENU	
1 SET CLOCK	2 SETUP
3 VIBRATION	4 TEMP
09/30/04	15:37:37

Following is a description of the QA Main Menu screen options:

Option	Response
1 Set Clock	The QA Main Menu screen displays the date and time of day (in 24-hour format). For example: 14:55 indicates 2:55 pm; 09:05 indicates 9:05 am.
	To change or correct the date or time of day, select option 1, Set Clock. The date displays. Press ENTER to accept the current date or enter the new date and press ENTER . The time displays. Press ENTER to accept the current time of day or enter the correct time in 24-hour format and press ENTER . The QA Main Menu screen displays.
2 Setup	See "Main Menu Option 2, Setup" on page 30.
3 Vibration	See "Main Menu Option 3, Vibration" on page 30.
4 Temp	No action required.

Page 30	QAII Station	Revision K, 11/10
Operation	Operator's Manual	P/N 70-9018

Main Menu Option 2, Setup

Select option **2**, Setup, from the QA Main Menu screen. The System Set Up screen displays.

SYSTEM SET U	JP	
1 POSITION	2 METRIC	
3 SCREEN SAVE	R TIME	
4 EUROPE DATE	STYLE	

Following is a description of the System Set Up screen options:

Option	Response
1 Position	Select option 1 , Position, to select the number of positions being tested on the dissolution apparatus. Select option 1 for six or option 2 for eight. The System Set Up screen displays.
2 Metric	Select option 2 , Metric, to select the vibration unit of measure. Select option 1 for inch or option 2 for metric. The System Set Up screen displays.
3 Screen Saver Time	Select option 3 , Screen Saver Time, to enter the number of minutes (0 - 9) the back lighting for the display screen remains on. Enter the desired duration and press ENTER . The System Set Up screen displays.
4 Europe Date Style	Select option 4 , Europe Date Style, to select the date format as American or European. Select option 1 , America (mm/dd/yy), or option 2 , Europe (dd/mm/yy). The System Set Up screen displays.

Press **MENU** to return to the QA Main Menu screen.

Main Menu Option 3, Vibration

The monitor screen displays the g value only. To display all other values associated with the vibration reading, select option 3, Vibration, from the QA Main Menu screen. Ensure

Revision K, 11/10	QAII Station	Page 31
P/N 70-9018	Operator's Manual	Operation

the vibration sensor is in place prior to selecting this option. See "Measuring and Recording Vibration" on page 40.

Select option **3**, Vibration, from the QA Main Menu screen. The Vibration Analysis screen displays:

VIBR ANALYSIS g 0.000 FRQ 000.00Hz VELOCITY .0000 IN/S DISPLACEMENT 0.000MIL

- "g" represents acceleration expressed in mm/sec² or in/sec².
- "FRQ" represents frequency expressed in Hz.
- "Velocity" is expressed in mm/sec or in/sec.
- "Displacement" is expressed in mm or mils.

Press **MENU** to return to the QA Main Menu screen.

Running a Test

The QAII Station is programmed to portably test up to 30 different dissolution apparatus.

To run a test, complete the procedures through "Measuring and Recording Shaft Verticality" on page 38 for each position, then complete "Measuring and Recording Vessel Plate (Table) Level" on page 39 and "Measuring and Recording Vibration" on page 40.



Page 32	QAII Station	Revision K, 11/10
Operation	Operator's Manual	P/N 70-9018

Entering Storage and Serial Numbers

3 Exit

			Note	
		Press MENU fro temperature, w	om the QA Main Menu so obble, vibration, and lev	creen to view the RPM, el data.
Step 1.	From numb	the monitor scre er displays.	een, press TEST . The mo	ost recently stored storage
Step 2.	Press numb previo	e ENTER to acce per (1 to 30) and pously entered.	ept the storage number of press ENTER . A serial r	or key in a new storage number displays if one was
Step 3.	Press nine (s ENTER to acce digits, and press	pt the serial number or left the serial number or left ENTER . The following s	key in a new number, up to screen displays:
		TEST ON POS	ITION 1	
		1 TEST	2 SKIP NEXT	
		3 EXIT		
		09/30/04	16:11:14	
Step 4.	Selec	t from the follow	ing options on the Test o	on Position 1 screen:
Option		Response		
1 Test		Use this option to run a test on the selected position. See "Entering Shaft Identification Numbers" on page 33.		
2 Skip Nex	t	Select option 2 to skip the position number listed and move to the next position number. The previously stored values for all of the parameters print for any skipped position number.		

Select option 3 to return to the monitor screen.

Revision K, 11/10	QAII Station	Page 33
P/N 70-9018	Operator's Manual	Operation

Entering Shaft Identification Numbers

- Step 1. Select option **1**, Test. A shaft identification number displays if one was previously entered.
- Step 2. Press **ENTER** to accept the current identification number or enter a new number, up to six digits, and press **ENTER**.
- Step 3. Continue to "Measuring and Recording RPM" below.

Measuring and Recording RPM



Immediately after entering the shaft identification number, the parameters being measured display on the screen.



- Step 1. Raise the drive unit to remove the EaseAlign centering rings, if applicable, from the vessels to be tested.
- Step 2. Lower the drive unit until there is a distance of approximately one centimeter from the top of the paddle blade or the bottom rim of the basket to the top of the vessel or adapter.
- Step 3. Place the EaseAlign mounting bracket with the magnetic tachometer sensor and digital wobble gauge on the first vessel position to be tested. The locating holes in the self-aligning mounting bracket slip over the locating pegs (see "locating pegs" in Figure 10 on page 28).

Page 34	QAII Station	Revision K, 11/10
Operation	Operator's Manual	P/N 70-9018

- Step 4. Press down until the mounting bracket is level and firmly seated.
- Step 5. The tip of the wobble gauge pointer should be positioned against the paddle shaft or the bottom rim of the basket.
- Step 6. Place the magnetic clip on the paddle or basket shaft being checked (see Figure 11, "Placement for RPM Reading," below). To avoid damage to the paddle or basket shaft, place the clip at or near the top of the shaft, just below the spindle chuck, then slide it down into its operating position in front of the tachometer sensor.

FIGURE 11. Placement for RPM Reading



- Step 7. Set the desired RPM on the dissolution apparatus.
- Step 8. Turn on the spindles.
- Step 9. On the QAII, press **ENTER** to store the RPM or press **CLEAR** to skip the RPM.



Step 10. Continue to "Measuring and Recording Temperature" on page 35.

Revision K, 11/10	QAII Station	Page 35
P/N 70-9018	Operator's Manual	Operation

Measuring and Recording Temperature

Complete the following steps to record the temperature:

- Step 1. Ensure the temperature probe is in the vessel being measured. Be careful not to entangle the temperature probe with the rotating spindles.
- Step 2. Press **ENTER** to store the temperature or press **CLEAR** to skip the temperature.

 Note
If you press CLEAR, the previously stored value prints for the temperature.

Step 3. Continue to "Measuring and Recording Wobble" below.

Measuring and Recording Wobble

Note
For highest accuracy, runout readings should be recorded with shafts turning at 50 RPM or less.
The shaft or basket wobble value, as measured by the wobble gauge in either millimeters or inches, is the absolute difference between high and low values. After zeroing the gauge, the reading on the display screen may be negative or positive since there is no way for the unit to determine whether you placed the gauge pointer on the shaft high point, low point or somewhere in between. For example, if the reading ranges between - 0.05 mm and + 0.1 mm, the QAII Station displays the runout value as 0.15 mm.

Complete the following steps to record the shaft or basket wobble:

- Step 1. Turn off the spindles.
- Step 2. Press **ON/OFF** on the wobble gauge. The display screen activates.

QAII Station	Revision K, 11/10
Operator's Manual	P/N 70-9018
	QAII Station Operator's Manual

- Step 3. Press **in/mm** on the gauge key to select inches or millimeters as the measurement unit.
- Step 4. Press and hold briefly **ORIGIN** to auto zero the gauge. The screen displays 0.0000 in or 0.00 mm.



Step 5. Ensure the wobble gauge is properly positioned with the tip of the wobble gauge pointer positioned against the paddle shaft or the bottom rim of the basket. See Figure 12, "Wobble Gauge with Paddle," below and Figure 13, "Wobble Gauge with Basket," on page 37.



Note

To avoid damage to the EaseAlign mounting bracket when using Apparatus 2 (paddles), check the clearance between the paddle blade and the bottom of the mounting bracket before rotating the shafts.

FIGURE 12. Wobble Gauge with Paddle



Revision K, 11/10	QAII Station	Page 37
P/N 70-9018	Operator's Manual	Operation



Note

When using Apparatus 1 (baskets), it is necessary to reposition the runout gauge pointer to the bottom rim of the basket before obtaining a reading. Take extra care to avoid damage to the runout gauge pointer. When repositioning the drive unit for proper alignment with the lower rim of the basket, use a thin ruler to hold the pointer safely away from your fingers.

FIGURE 13. Wobble Gauge with Basket



- Step 6. Turn on the spindles.
- Step 7. On the QAII, press **ENTER** to store the wobble value or press **CLEAR** to skip the wobble.



Step 8. Continue to "Measuring and Recording Shaft Verticality" on page 38.

Page 38	QAII Station	Revision K, 11/10
Operation	Operator's Manual	P/N 70-9018

Measuring and Recording Shaft Verticality

This apparatus is used in conjunction with a centering tool to verify proper shaft alignment. Complete the following steps to record the shaft verticality:

- Step 1. Turn off the spindles.
- Step 2. Hold the level sensor flush against the shaft. The perpendicular groove of the sensor should be placed high up on the shaft, approximately 1 inch or 25.4 mm below the bottom of the dissolution apparatus drive unit. This sensor measures in degrees the verticality or perpendicularity of the shaft from side-to-side and front-to-back simultaneously.

FIGURE 14. Electronic Level Sensor Placement



Step 3. On the QAII, press **ENTER** to store the shaft verticality or press **CLEAR** to skip the shaft verticality. Immediately after storing or skipping the shaft verticality, the Test on Position 2 screen displays prompting you to test position 2.



Revision K, 11/10	QAII Station	Page 39
P/N 70-9018	Operator's Manual	Operation

Step 4.	Select option 1 , Test, to repeat the above procedures beginning with step 2 under "Entering Shaft Identification Numbers" on page 33 for each remaining position.
	to the next position number.
	Select option 3 , Exit, to return to the monitor screen.
Step 5.	After testing the last position, continue to "Measuring and Recording

Measuring and Recording Vessel Plate (Table) Level

Vessel Plate (Table) Level" below.

Complete the following steps to record the vessel table level:

- Step 1. Place the level sensor in the upright position on the vessel plate. The sensor measure in degrees the level from side-to-side and front-to-back simultaneously.
- Step 2. Press **ENTER** to store the vessel table level or press **CLEAR** to skip the level.



Step 3. Continue to "Measuring and Recording Vibration" on page 40.

Page 40	QAII Station	Revision K, 11/10
Operation	Operator's Manual	P/N 70-9018

Measuring and Recording Vibration

Complete the following steps to record the vibration:

Step 1. Place the vibration sensor on the vessel plate. The vibration sensor can be placed either on the vessel plate as shown below or on the drive unit to obtain readings.

FIGURE 15. Vibration Sensor



Step 2. On the dissolution apparatus, set the shafts to rotate at the desired speed and activate the spindles.



Step 3. Position the sensing arrow in the direction of the axis to be measured.



Varian, Inc.

Revision K, 11/10 P/N 70-9018	QAII Station Operator's Manual	Page 41 Operation
Step 4.	Press ENTER to store vibration on the X axis or press the X-axis vibration.	CLEAR to skip
Step 5.	Press ENTER to store vibration on the Y axis or press the Y-axis vibration.	CLEAR to skip
Step 6.	Press ENTER to store vibration on the Z axis or press the Z-axis vibration. The monitor screen displays.	CLEAR to skip

Printing Results

To print results, the QAII Station must be plugged into AC power with the main (rear panel) power switch in the ON position.

- Step 1. From the monitor screen, press **PRINT** on the keypad.
- Step 2. Enter the storage number (1 30) of the desired stored data and press **ENTER**. The following screen displays:

TEST REPORT	
1 PRINT	2 SEND DATA
3 COM ID	4 BAUD RATE
09/30/04	08:40:09

Step 3. Select option **1** to print a full report.

The report prints out the following data:

- storage number
- apparatus serial number
- date and time of test
- position (vessel/shaft position)
- shaft ID (serial number)

- actual RPM
- wobble (in or mm)
- temp (vessel temp in °C)
- shaft level (X and Y measured in degrees)

Page 42	QAII Station	Revision K, 11/10
Operation	Operator's Manual	P/N 70-9018

Table level and vibration data display at the end of the report:

- table level (X and Y measured in degrees)
- vibration (g)

- frequency (Hz)
- velocity (mm/sec or in/sec)
- displacement (mm or mil)

Note: Vibration, frequency, velocity, and displacement are expressed for the three dimensions of X, Y, and Z.

Storage Table and Viewing Data

Complete the following steps to manually review the collected data:

- Step 1. From the Test Report screen, press **MENU**. The monitor screen displays.
- Step 2. Press **VIEW DATA**. The Storage View screen displays.

STORAGE VIEW	
1 STORAGE MAP	
2 STORAGE DATA	
03/31/00	08:40:09

- Step 3. Select option 1, Storage Map, to view dissolution apparatus serial numbers stored in the QAII Station memory. To scroll through the list, press ENTER to go to the next entry or CLEAR to return to the previous entry.
- Step 4. Press **MENU** to return to the Storage View screen.
- Step 5. Select option **2**, Storage Data, to review the previous results of an individual apparatus.

Revision K, 11/10	QAII Station	Page 43
P/N 70-9018	Operator's Manual	Operation

- Step 6. Enter the storage number of the apparatus that you would like to review and press **ENTER**. The dissolution apparatus data displays. To scroll through the data, press **ENTER** to go to the next entry or **CLEAR** to return to the previous entry.
- Step 7. Press **MENU** to return to the Storage View screen.
- Step 8. Press **MENU** twice to return to the QA Main Menu screen.

Page 44	QAII Station	Revision K, 11/10
Operation	Operator's Manual	P/N 70-9018

This page was intentionally left blank, except for this message.

Chapter 5 Maintenance and Troubleshooting

Periodic Maintenance



Warning

The QAII contains electrical circuits, devices, and components operating at dangerous voltages. Contact with these circuits, devices, and components can cause death, serious injury, or painful electric shock.

Periodic maintenance needs may vary depending on frequency of instrument usage.

Calibration

We recommend the QAII Station be calibrated at six-month intervals. However, based on your own SOPs, this can be extended for up to a year. Call the Dissolution Systems Service Department for a quote, packing instructions and a Return Authorization Number before sending your unit for re-calibration.

QAII Station Operator's Manual

Report Center Impact Printer

The following is helpful information for using your impact printer.

Installing the Cartridge Ribbon

If the printer is used infrequently, the print impression sometimes becomes weak because the ribbon dries out. If the printed material is difficult to read and you suspect this is the cause of the problem, advance to a new section of the ribbon by pressing the printer toggle switch to the *Paper feed* position. If the printing is still faint, replace the cartridge.

To install the cartridge, complete the following steps:

- Step 1. Toggle the printer off line by pressing the printer toggle switch to the OnLine / Off Line position. When the printer is off line, the Ready LED does not illuminate.
- Step 2. Four small grooves are embossed on the printer cover. Gently push on these grooves to tilt the cover. When the printer cover is tilted up, you can lift it off completely.
- Step 3. Push down on the right side of the ribbon cartridge (marked PUSH) and remove the old cartridge.
- Step 4. Install the new cartridge. If there is already paper in the printer, hold the cartridge between your thumb and index finger, slide it over the paper and into the printer compartment. Ensure the paper is between the ribbon cartridge and the ink ribbon. Ensure the ink cartridge is inserted firmly to prevent weak or irregular printing. The cartridge must be properly seated and aligned for the best printing.
- Step 5. Turn the cartridge knob (marked by an arrow) clockwise to stretch the ribbon taut.
- Step 6. Replace the cover.

Page 4	QAII Station	Revision K, 11/10
Maintenance and Troubleshootin	Operator's Manual	P/N 70-9018

- Step 7. Toggle the printer online by pressing the printer toggle switch to the OnLine / Off Line position. The Ready LED illuminates.
- Step 8. Replace the paper if necessary.

If you get ribbon ink on the printer's plastic cover, remove it immediately. Once dried, it is difficult to remove.

Replacing the Paper Roll

- Step 1. Toggle the printer off line by pressing the printer toggle switch to the OnLine / Off Line position. When the printer is off line, the Ready LED does not illuminate.
- Step 2. Grasp the paper roll cover firmly by the grooves on the side and the front edge. Pull outward to remove the cover.
- Step 3. Press the printer toggle switch to *Paper feed* to advance the paper approximately one inch beyond the paper cutter.
- Step 4. Using scissors, cut the paper feeding to the printer and remove the paper roll.
- Step 5. Pull the remaining paper through the printer mechanism. *Pull the paper from the front (paper cutter side)*. Pulling the paper out of the back of the printer will damage the print mechanism.
- Step 6. Unroll several inches of paper on the new roll.
- Step 7. If it is jagged, cut a straight edge on the paper roll to facilitate the entry of the paper into the printer.
- Step 8. Slide the paper through the slot connecting the paper compartment and the printer compartment. You can slide it in approximately 1/4 inch before it stops.

Page 48	QAII Station	Revision K, 11/10
Maintenance and Troubleshooting	Operator's Manual	P/N 70-9018

Step 9. While holding the paper in place, press the printer toggle switch to the *Paper feed* position and hold until approximately one inch of paper has emerged from the top of the printer.



- Step 10. Release the printer toggle switch.
- Step 11. Turn the paper roll to take up any slack in the paper feeding to the printer.
- Step 12. Place the paper roll into the paper compartment.
- Step 13. Replace the paper roll cover. If the cover is difficult to remove or replace, the left and right edges can be trimmed or shaved with a utility knife allowing the cover to slide easier.
- Step 14. Toggle the printer online by pressing the printer toggle switch to the OnLine / Off Line position. The Ready LED illuminates.

Toggling Your Printer Online

Complete these steps to toggle your printer online:

- Step 1. Toggle the printer online by pressing the printer toggle switch to the OnLine / Off Line position. When the printer is off line, the Ready LED does not illuminate.
- Step 2. Release the switch and it returns to the center position. The Ready LED illuminates and a READY message prints if the PRINT READY command has not been turned off. See "Printer Configuration" on page 49 for

Revision K, 11/10	QAII Station	Page 49
P/N 70-9018	Operator's Manual	Maintenance and Troubleshooting

instructions on turning on and off the PRINT READY command. When you first turn on the instrument, it prints a READY message to assure you that the built-in microprocessor is operating properly.

When you turn off the printer, wait at least three seconds before turning it on again.

Printer Self Test

You can test the print head and ribbon only *after* inserting paper. Do not attempt to print without paper. Follow these steps to perform a printer self test:

- Step 1. Turn off the QAII Station.
- Step 2. Press and hold the printer toggle switch in the *Paper feed* position.
- Step 3. Turn on the QAII Station.
- Step 4. Hold the printer toggle switch until printing begins. The printer prints a list of the current configuration settings and performs a continuous print test.
- Step 5. Press the printer toggle switch to the *OnLine / Off Line* position to stop the printing operation.
- Step 6. The printer is ready to resume normal operation.

Printer Configuration



Step 1. Turn off the QAII Station.

Page 50	QAII Station	Revision K, 11/10
Maintenance and Troubleshooting	Operator's Manual	P/N 70-9018

- Step 2. Press and hold the printer toggle switch in the *OnLine / Off Line* position while turning on the instrument. Hold the printer toggle switch in the *OnLine / Off Line* position for six seconds after the instrument is turned on, then release the switch.
- Step 3. The printer should print: *** SETUP MENU *** and CONFIGURE... [NEXT/OK]. If this message does not print, repeat steps 1 through 3.
- Step 4. The printer toggle switch is used to complete the configuration. Pressing the left side of the printer toggle switch selects NEXT to advance to the next menu item. Pressing the right side of the printer toggle switch selects OK to accept what is stated on this line of the menu item. Each time the switch is pressed, another part of the menu prints. Allow the printer to finish printing before pressing the switch again. See the table of commands below.

(Note		
	The printout is easier to read if the printer cover is removed.		
	*** SETUP MENU***		
	CONFIGURE	[NEXT/OK]	Press NEXT to avoid configuration
	CUSTOM	[NEXT/OK]	Press OK to enter custom mode
	CUSTOM MENU		
	PRINT CUSTOM SETUP	[NEXT/OK]	Press NEXT
	AUTO SEQ = NO	[NEXT/OK]	Press OK
	ZERO = Ø	[NEXT/OK]	Press OK
	POUND SIGN = #	[NEXT/OK]	Press OK
	_(UNDERSCORE)	[NEXT/OK]	Press OK
	ONLINE/OFFLINE = YES	[NEXT/OK]	Press OK
	EXT CH SET = NO	[NEXT/OK]	Press OK
	PRINT READY = YES	[NEXT/OK]	Press NEXT
	PRINT READY = NO	[NEXT/OK]	Press OK
	READY		

Your printer is now configured correctly.

Fuse Replacement

WarningImage: Description of the problem of the

The fuse compartment is located beside the power cord receptacle on the QAII rear panel. See Figure 9, "QAII Station Rear Panel," on page 25.

- Step 1. To check or replace the fuse, remove the power cord from the QAII.
- Step 2. A release tab is located on the right side of the fuse compartment. Insert a slotted screwdriver under the tab. A slight application of pressure forward releases the compartment door.
- Step 3. The fuses are located in the removable holders marked with an arrow on the top of each. The QAII uses one 500 mAmp, metric (5 x 20 mm) standard fuse for each holder.
- Step 4. Replace the fuse in the holder and insert the fuse holder into the fuse compartment with the arrows pointing toward the bottom of the compartment.
- Step 5. The holder is designed for multiple voltages. The voltage displays through a window in the fuse compartment door. To change the voltage, gently pull the wheel to remove it from the holder. Rotate the wheel and snap it back in place displaying the correct voltage.
- Step 6. Push the fuse compartment door closed. It snaps into place.
- Step 7. Replace the power cord.

Troubleshooting

The Dissolution Systems Service Department can assist you if you experience problems or have questions concerning your QAII. Many problems can be traced to simple sources and are easily solved.

Following is a troubleshooting guide which may help you. The Dissolution Systems Service Department can be reached at 800.229.1108 (inside the US) or 919.677.1108 (outside the US). Optionally, you can send a fax to 919.677.1138. You can also e-mail the Dissolution Systems Service Department at dissolution.service@varianinc.com.

Symptom	Probable Cause	Suggested Solution	
The unit does not respond when the main power switch	There is a blown fuse.	Replace the fuse. See "Fuse Replacement" on page 51.	
is turned on.	The unit is not plugged in.	Check the rear panel AC power cord receptacle. Plug in the power cord.	
The unit does not respond when the battery switch is turned on.	The battery is dead.	Recharge the battery. See "Power Switch Functions" on page 24.	
The Report Center Printer does not function.	The printer is disabled.	Ensure the Report Center Printer is online and the Ready LED is illuminated. See "Toggling Your Printer Online" on page 48.	
	THE QAII is not plugged in or the main power switch is not turned on.	Plug the instrument into a wall outlet and ensure the main (rear panel) power switch is in the ON position.	
The Report Center Printer is active but nothing prints.	The ribbon needs to be replaced.	Install a replacement ribbon cartridge. See "Installing the Cartridge Ribbon" on page 46.	
The RPM readings are erratic.	The reflective clip is not registering.	Align the clip with the sensor.	
The RPM sensor is not reading.	The sensor has come loose or is not plugged into the QAII rear panel.	Restore connections.	

Revision K, 11/10	QAII Station	Page 53
P/N 70-9018	Operator's Manual	Maintenance and Troubleshooting

Symptom	Probable Cause	Suggested Solution
The wobble gauge is not reading.	The sensor has come loose or is not plugged into the QAII rear panel.	Restore connections.
	The pointer is not touching the shaft or rim of the basket.	Realign the mounting bracket.
	The wobble gauge is not turned on.	Turn on the wobble gauge.
The level sensor is not reading.	The sensor has come loose or is not plugged into the QAII rear panel.	Restore connections.
The level sensor displays erratic readings.	The sensor is not in the upright position.	Correct positioning.
The vibration sensor is not reading.	The sensor has come loose or is not plugged into the QAII rear panel.	Restore connections.
There is excessive wobble.	The paddle or basket shaft is damaged.	Verify spindle wobble is within specifications using the certified calibration shaft.
The shaft verticality or perpendicularity is excessive.	The paddle or basket shaft is damaged.	Verify the spindle perpendicularity is within specifications using the certified calibration shaft.
	The drive unit is not level.	Contact the Dissolution Systems Service Department.
	The spindle is not centered.	Contact the Dissolution Systems Service Department.

QAII Default Values

The default value prints for any sensor which is not plugged into the unit during testing. The values listed below apply for both standard and metric units (see "Measuring and Recording Wobble" on page 35 and "Main Menu Option 2, Setup" on page 30 for instructions on selecting inches or millimeters for wobble and vibration respectively).



Note If any of the following default values display, ensure all connections are tight.

Variable	Default Values
RPM	000.00
Wobble	00.00
Temp	-1.5
Shaft-X	-6.8
Shaft-Y	-6.8
Table-X	-6.8
Table-Y	-6.8
Vibration-X	0.000
Vibration-Y	0.000
Vibration-Z	0.000
Frequency	000.00
Velocity	00.00
Displacement	.0000

Chapter 6 Service and Warranty

The warranty is provided by Varian, Inc. or one of its authorized representatives.

Service and Warranty Information

Varian Dissolution products carry a one-year warranty on parts and labor. The Dissolution Systems Service Department (or one of its representatives) will, at its option, either repair or replace any mechanical and electrical components in your instrument which prove to be defective. During the first year of warranty coverage, there is no charge for the labor to repair your unit. The Dissolution Systems Service Department (or one of its representatives) will determine the best site to repair the unit, either onsite or returned to Varian, Inc. Any onsite warranty services are provided only at the initial installation point. Installation and onsite warranty services are available only in Dissolution Systems service travel areas.

Page 56	QAII Station	Revision K, 11/10
Service and Warranty	Operator's Manual	P/N 70-9018

Exclusions and Limitations

Excluded from this warranty are expendable or consumable items such as, but not limited to, paddles, baskets, vessels, and acrylic water baths. Also excluded are defects from improper or inadequate maintenance by the customer, user-induced chemical action or contamination, unauthorized modification or misuse, and improper site preparation and maintenance.

Operation of software is not warranted to be uninterrupted or error-free.

Obtaining Warranty Service

To obtain warranty service in the United States, contact the Dissolution Systems Service Department at 800.229.1108 to obtain authorization to return units for repair. At the option of the customer, onsite warranty service is available, but travel charges may be incurred. The customer should prepay all shipping charges for products returned to the Dissolution Systems Service Department (unless otherwise authorized), and Varian, Inc. will pay all charges for return to the customer.

Warranty Limitations

Varian, Inc. makes no other warranty, either express or implied, with respect to this product. Specifically disclaimed are any implied warranties of merchantability and fitness for a particular use. In no event will Varian, Inc. be liable for any indirect, incidental, or consequential damages arising from the use of this product. This warranty gives you specific legal rights which may vary from state to state or province to province, so you may have other rights and some of these exclusions may not apply to you.

Exclusive Remedies

The remedies provided herein are the customer's sole and exclusive remedies. In no event shall Varian, Inc. or its representatives be liable for any direct, indirect, special, incidental, or consequential damages, whether based on contract, tort, or any other legal theory. Some states or provinces do not allow the exclusion or limitation of incidental or consequential damages, so the above limitation or exclusion may not apply to you.

Page 58	QAII Station	Revision K, 11/10
Service and Warranty	Operator's Manual	P/N 70-9018

This page was intentionally left blank, except for this message.

Index

а

apparatus without easealign centering rings 28

С

calibration 45 cartridge ribbon 46 conventions 17

d

default values 54

е

exclusions 56 exclusive remedies 57

f

fuse replacement 51

h

hazards 9

i

installing the printer cartridge ribbon $\,46$ introduction $\,15$

l level sensor 22 limitations 56

m

magnetic clip 21

ο

obtaining warranty service 56 operating the qaii 29

р

periodic maintenance 45 power switch functions 24 printer configuration 49 printer self test 49 printer, toggling online 48 printing results 41

r

reader comment form 61 rpm, measuring and recording 33 running a test 31 Index

s

safety practices 9 self test, printer 49 serial numbers 32 service 55 setup 19 shaft verticality, measuring and recording 38 storage numbers 32 storage table 42 system set up 30

t

tachometer sensor 21 temperature probe 23 temperature, measuring and recording 35 troubleshooting 52

u

universal adapter 22, 28 unpacking your qaii 19

v

vessel plate level 39 vibration 30 vibration sensor 23 vibration, measuring and recording 40 viewing data 42

w

warranty 55 warranty limitations 56 wobble gauge 20 wobble, measuring and recording 35



Tell Us How We Are Doing

We listen to our customers. We work hard to make our technical documentation user friendly, and to make the information in our manuals easy to retrieve and use. We'd like you to tell us the kinds of additional information you'd find helpful in our documentation. Your feedback will be carefully considered when we prepare future editions of this manual.

This manual should contain the following additional information:

The most useful thing about this book is:

This manual would be more helpful if:

My general impressions of this book are:

May we contact you regarding your comments? ____ YES ____ NO (If yes, please write your name, address, and telephone number here.)

Please return this form via mail to: Technical Writing / Dissolution Systems, Varian, Inc., 13000 Weston Parkway, Cary, North Carolina 27513-2250 USA. Optionally, you can return this form via fax at 1.919.677.1550. Always, feel free to telephone us to discuss your comments at 1.800.229.1108.